

# SoftMax<sup>®</sup> Pro Software

Version 5 for Mac and Windows User Guide



Document Number: 5000938 Rev A  
August 2009

---

This document is provided to customers who have purchased MDS Analytical Technologies (US) Inc. ("MDS AT") equipment, software, reagents, and consumables to use in the operation of such MDS AT equipment, software, reagents, and consumables. This document is copyright protected and any reproduction of this document, in whole or any part, is strictly prohibited, except as MDS AT may authorize in writing.

Equipment, software, reagents, and consumables that may be described in this document are protected under one or more patents filed in the United States, Canada, and other countries. Additional patents are pending.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such usage is intended only to designate those manufacturers' products as supplied by MDS AT for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

MDS AT makes no warranties or representations as to the fitness of this equipment for any particular purpose and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use to which the purchaser may put the equipment described herein, or for any adverse circumstances arising therefrom.

For Research Use Only. Not for use in diagnostic procedures.



SOFTMAX, SPECTRAMAX, VMAX, EMAX, FLEXSTATION, PATHCHECK, ANALYST, and FLIPR are registered trademarks of MDS Analytical Technologies (US) Inc. These trademarks may not be used in any type of promotion or advertising without the prior written permission of MDS Analytical Technologies (US) Inc.

Equipment built by MDS Analytical Technologies (US) Inc.  
1311 Orleans Drive, Sunnyvale, California, United States of America 94089.  
MDS Analytical Technologies (US) Inc. is ISO 9001 registered.  
© 2009 MDS Analytical Technologies (US) Inc.  
All rights reserved.  
Printed in the USA.

---

# Contents

<b>Chapter 1 Introduction</b> . . . . .	<b>9</b>
Major Functions of SoftMax® Pro Software . . . . .	10
Instrument Control . . . . .	10
Data Collection and Display . . . . .	10
Data Reduction and Plotting . . . . .	10
Immediate Results Reporting and Analysis . . . . .	10
Two Available Editions: Standard and GxP . . . . .	11
<b>Chapter 2 Installation and Setup</b> . . . . .	<b>13</b>
Installation . . . . .	13
Installing an Instrument . . . . .	13
Computer System Requirements . . . . .	13
Installing SoftMax Pro Software . . . . .	13
Uninstalling SoftMax Pro Software . . . . .	14
Running the Software for the First Time . . . . .	14
Communicating with an Instrument . . . . .	14
Starting SoftMax Pro Software . . . . .	14
Entering Registration Information . . . . .	14
Connecting to the Instrument Via USB . . . . .	15
Installing Single Port Adapters . . . . .	15
Installing Multiple-port Adapters . . . . .	16
Troubleshooting Instrument Connections . . . . .	17
LMax II . . . . .	17
Instrument Preferences . . . . .	17
Setting Preferences . . . . .	18
Reader Settings . . . . .	18
Manual Export Format . . . . .	19
Save Data After Read (AutoSave) . . . . .	21
Print Document After Read . . . . .	23
The Protocols Menu . . . . .	24
<b>Chapter 3 SoftMax® Pro Software Interface</b> . . . . .	<b>25</b>
The SoftMax® Pro Application Window . . . . .	25
WYSIWYG Display . . . . .	25
Status Bar . . . . .	26
Instrument Icon . . . . .	26

Temperature Display . . . . .	26
Read . . . . .	26
Reference Button . . . . .	27
StakMax® Software Interface Button . . . . .	27
Incubator Button . . . . .	27
Automix Button . . . . .	28
Drawer Button . . . . .	28
Instrument Settings and Control . . . . .	28
Instrument Settings . . . . .	28
Read . . . . .	29
Close Drawer, Open Drawer . . . . .	29
Incubator . . . . .	29
Ref . . . . .	29
Experiments . . . . .	30
Selecting an Experiment . . . . .	30
Manipulating Experiments . . . . .	30
Sections . . . . .	31
Active Section Menus . . . . .	31
Manipulating Sections . . . . .	31
Notes Section . . . . .	33
Plate Section . . . . .	34
Group Section . . . . .	37
CuvetteSet Section . . . . .	38
Graph Section . . . . .	39
Creating a Default Protocol . . . . .	40
<b>Chapter 4 Data Collection . . . . .</b>	<b>41</b>
Introduction to Data Collection . . . . .	41
Instrument Settings . . . . .	41
Instrument Operation . . . . .	41
Instrument Settings Vary for Different Models. . . . .	42
Read Type . . . . .	42
Read Mode . . . . .	44
Wavelengths . . . . .	46
Shake . . . . .	47
Integration . . . . .	47
Injection and Delay . . . . .	48
Injection Wells . . . . .	48
Sensitivity . . . . .	49

---

Automix . . . . .	57
Blanking (Pre-Read Plate). . . . .	57
Autocalibrate . . . . .	58
AutoCalibrate for Fluorescence and Luminescence Reads . . . . .	59
Well Scan Editor. . . . .	59
Assay Plate Type . . . . .	60
Strips/Wells to Read. . . . .	61
Compound Source. . . . .	62
Compound Transfer. . . . .	62
Triturate . . . . .	64
Pipette Tips Layout . . . . .	64
Compound & Tips Columns. . . . .	65
Settling Time . . . . .	67
Speed Read . . . . .	67
Column Wavelength Priority . . . . .	67
AutoRead. . . . .	68
Template Editor . . . . .	68
Selecting Wells or Cuvettes in the Template Editor . . . . .	69
Group. . . . .	69
Sample. . . . .	72
Assign. . . . .	72
Clear. . . . .	72
Series . . . . .	73
Blanking . . . . .	74
Plate Blanks . . . . .	75
Copying and Pasting Templates . . . . .	77
Exporting and Importing Templates . . . . .	78
Reading a Microplate or Cuvette . . . . .	78
Data Collection from a Microplate. . . . .	78
Data Collection from a Cuvette . . . . .	79
Calibration . . . . .	81
Data Display During a Reading . . . . .	82
<b>Chapter 5 Data Analysis . . . . .</b>	<b>83</b>
Data Display . . . . .	83
Display Dialog Box. . . . .	83
Graphing Wells . . . . .	86
Masking Wells or Cuvettes. . . . .	88
Data Reduction . . . . .	88

EndPoint Reads . . . . .	89
Kinetic Reads . . . . .	90
Spectrum Reads . . . . .	94
Well Scan . . . . .	95
Flex Reads . . . . .	95
Data Mode (%Transmittance/Absorbance) . . . . .	97
Custom Reduction Formulas . . . . .	97
Recalculation Options . . . . .	99
Group and Graph Sections . . . . .	100
Group Sections . . . . .	100
Graph Section . . . . .	103
<b>Chapter 6 Tutorials . . . . .</b>	<b>115</b>
Tutorial 1: Quantitative EndPoint Protocol . . . . .	116
Step 1: Create a New Data Base . . . . .	116
Step 2: Define Instrument Settings . . . . .	117
Step 3: Define the Template . . . . .	118
Step 4: Set Reduction Parameters . . . . .	121
Step 5: Set Display Parameters . . . . .	121
Step 6: Save the Protocol . . . . .	121
Step 7: Read the Plate . . . . .	122
Step 8: Data Analysis—Group Sections. . . . .	122
Step 9: Data Analysis—Standard Curve. . . . .	125
Step 10: Print a Report. . . . .	125
Tutorial 2: EC 50 Protocol Tutorial. . . . .	126
Step 1: Create a New Data File . . . . .	126
Step 2: Define Instrument Settings . . . . .	127
Step 3: Define a Template. . . . .	128
Step 4: Set Reduction Parameters . . . . .	129
Step 5: Set Display Parameters . . . . .	129
Step 6: Save the Protocol . . . . .	130
Step 7: Read the Plate . . . . .	130
Step 8: Data Analysis—Group Sections. . . . .	131
Step 9: Data Analysis—Graph . . . . .	133
Step 10: Data Analysis—Summary Formulas. . . . .	133
Step 11: Print a Report. . . . .	134
<b>Chapter 7 File Management and Printing . . . . .</b>	<b>135</b>
File Creation and Management . . . . .	135

---

Data Files . . . . .	135
Protocol Files . . . . .	135
Organizing Protocol and Data Files . . . . .	136
Creating New Files. . . . .	136
Opening Files. . . . .	136
Saving Files Manually. . . . .	137
AutoSave . . . . .	137
Printing . . . . .	137
Printing Reports . . . . .	137
Including or Excluding Sections. . . . .	137
Changing Section Order . . . . .	138
Changing Text Format. . . . .	138
Import/Export. . . . .	138
Export. . . . .	138
Import . . . . .	141
Importing and Exporting Templates . . . . .	142
Exporting Graphs. . . . .	144
Copying and Pasting . . . . .	144
<b>Chapter 8 SoftMax® Pro GxP Software . . . . .</b>	<b>145</b>
How the Programs Work Together . . . . .	146
GxP Admin Features . . . . .	146
Using SoftMax Pro GxP Software . . . . .	147
User Log On. . . . .	147
Files . . . . .	148
GxP Menu . . . . .	148
Statements. . . . .	149
Locking and Unlocking Sections . . . . .	150
<b>Chapter 9 Robotics and LIMS Integration . . . . .</b>	<b>153</b>
Introduction . . . . .	153
Sample Source Code and Applications . . . . .	153
Remote Control . . . . .	153
Send/Receive Remote Commands. . . . .	153
The WM_SETTEXT Approach . . . . .	154
The WM_COPYDATA Approach . . . . .	155
Summarizing Remote Commanding . . . . .	157
Programmer Notes . . . . .	158
Remote Control in SoftMax Pro GxP Software. . . . .	158

SoftMax Pro Commands .....	159
Visual Basic (Excel Macro) Example .....	166
MFC C++ Interface to SoftMax Pro .....	169
MFC Example Code. ....	169
<b>Appendix A Glossary of Terms .....</b>	<b>173</b>
<b>Index.....</b>	<b>185</b>



SoftMax® Pro software controls MDS Analytical Technologies spectrophotometers, absorbance, luminescence, and fluorescence microplate readers—providing extensive data calculation and analysis capabilities under a Good Manufacturing Practices (GMP), Good Laboratory Practices (GLP) work environment for pharmaceutical, biotechnology, academic, hospital, and government customers.

Over 120 assay protocols are included in the software to speed life science research and drug discovery assay development and screening. Researchers can customize experiment protocols, analyze and display data, and create meaningful reports. The straightforward yet powerful programming capabilities of SoftMax Pro software can further enhance any specialized data collection and analysis needs through custom assay development.

This industry-leading software package is also widely integrated with industry-leading robotics systems. Optional Validation Tools are also available to reduce the time, effort and cost of certifying and revalidating MDS Analytical Technologies laboratory tools.

Two editions of the software are available: SoftMax Pro Standard software and SoftMax Pro GxP software. Both support all MDS Analytical Technologies instruments including: SpectraMax®, VersaMax, VMax®, EMax®, Gemini, ThermoMax, and FlexStation® readers. Data from Analyst®, and FLIPR® systems can also be easily imported and analyzed.

SoftMax Pro GxP software is a special release that includes full user and permission control and audit trail generation in addition to all the data collection and analysis features found in the Standard edition. With the audit trail features in SoftMax Pro GxP, all experiments are fully traceable and can be completely reconstructed.

This user guide describes features available in SoftMax Pro software release 5.4 and above. SoftMax Pro software release 5.4 is available for PC-compatible computers using Windows XP and Vista (32-bit only), and Mac computers using OSX 10.4 or greater.

## Major Functions of SoftMax® Pro Software

### Instrument Control

SoftMax Pro software allows you to set up and run a complete protocol for all instruments. Instrument settings can be saved as a protocol file and used repeatedly for reading different microplates or cuvettes. All standalone instrument functions can be controlled using the software. In addition, SoftMax Pro software provides capabilities that are not available when using an instrument in standalone mode such as user-defined Kinetic run times, read intervals, and Automix parameters

### Data Collection and Display

SoftMax Pro software collects and stores all raw data received from the instrument. Data is displayed in a grid format that corresponds to the wells in a microplate (all instruments) or individual cuvettes (using SpectraMax Plus, Plus384, M2, M2e, M5<sup>e</sup> or M5 readers).

SoftMax Pro software can collect data from one or more microplates or cuvettes and store it in a single data file, using the same or different instrument settings for different microplates or cuvettes. For example, microplates containing different samples can be read using the same or different modes, all within the same experiment.

In addition to displaying data collected from an MDS Analytical Technologies microplate system instrument, you can import, display, and analyze data collected from either a FLIPR® or Analyst instrument.

### Data Reduction and Plotting

You can manipulate or reduce the raw data using dozens of built-in formulas or define your own analysis structure to quickly and easily summarize the raw data. More than one reduction can be shown, and results from different microplates and cuvettes can be compared within the same experiment.

### Immediate Results Reporting and Analysis

Once you have defined instrument settings, and have customized a SoftMax Pro software data file with assay information, reduction settings, custom columns in Group sections, and summary objects, you can “Save” a Protocol file type to create an assay template that can then be used and

distributed throughout a department or company for highly repeatable data collection and analysis that is completed the second the plate read has completed.

## **Two Available Editions: Standard and GxP**

Two editions of SoftMax Pro software are available: Standard and GxP.

SoftMax Pro software provides instrument control and analysis functions for all SpectraMax, FlexStation, Gemini, and earlier instruments as well as import and analysis of data files from FLIPR and Analyst instruments.

SoftMax Pro GxP software, together with an MDS Analytical Technologies plate reader and MDS Analytical Technologies GxP Admin software, provides a fully integrated hardware and software solution for adherence to Good Manufacturing Practices (GMP), Good Laboratory Practices (GLP), and many other regulatory compliance requirements including 21 CFR Part 11.



## Installation

### Installing an Instrument

Please consult your instrument manual for instructions on connecting your instrument to a computer.

### Computer System Requirements

Recommended computer hardware and operating system software specifications that will ensure proper operation of SoftMax® Pro software can be found online at:

<http://www.moleculardevices.com/softmax>.

MDS Analytical Technologies recommends that you increase the amount of RAM in your computer to a minimum of 2 GB in the following cases:

- If you plan to use SoftMax Pro software with a FlexStation® instrument.
- If you plan on using your microplate reader with the StakMax® integrated plate handling system.
- If you use large (more than 40 Plate sections per file) or complicated (for example, 384 well ADME protocols) files.
- If you plan to use robotic applications.

### Installing SoftMax Pro Software

To install SoftMax Pro software

1. Insert the SoftMax Pro CD into the CD-ROM drive.
2. Windows: Open My Computer and then double-click the CD-ROM drive icon to display its contents.  
Mac: Double-click the CD icon to display its contents.
3. Double-click the SoftMax Pro software version 5.4 Installer icon and follow the screen prompts.

## Uninstalling SoftMax Pro Software

Before uninstalling the program, make sure to backup your data and any saved files to a folder outside of the SoftMax Pro software folder, because uninstalling deletes the entire SoftMax Pro software directory.

Windows: To uninstall SoftMax Pro software

1. Click **Start > Settings > Control Panel**.
2. Double-click **Add or Remove Programs**.
3. In the list that appears, click **SoftMax Pro version 5.4** and then click **Remove**.



---

**Note:** This is the recommended method of removing SoftMax Pro software from a Windows-based computer since it also removes related information from the Windows Registry.

---

Mac: To remove SoftMax Pro software

1. Drag the SoftMax Pro software folder to the **Trash**, then empty the Trash.

## Running the Software for the First Time

### Communicating with an Instrument

To ensure that SoftMax Pro software can communicate with your instrument, turn on the instrument before starting SoftMax Pro software.

If using a USB-to-serial adapter, be sure to obtain the latest software driver from the manufacturer before connecting to your instrument.

### Starting SoftMax Pro Software

After the instrument is powered on and has completed its start-up sequence, double-click the SoftMax Pro icon on your desktop to start the program.

### Entering Registration Information

The first time you run SoftMax Pro software after installation, a first-launch wizard will guide you through the setup process. In particular, it requires

you to enter your name and company, as well as the serial numbers of SoftMax Pro software and the connected instrument.

The software serial number can be found in the SoftMax Pro software box, and the instrument serial number is located on a label affixed to the back of the instrument.

You will also be prompted to register your software with MDS Analytical Technologies over an internet connection. Registration of SoftMax Pro software activates the product warranty and entitles you to receive software updates when available.

## Connecting to the Instrument Via USB

Many Windows and Mac-based computers no longer have a standard RS232 serial port. If you have a computer without a standard serial port, you will need to use a USB-to-serial adapter.

These adapters are available at many computer supply stores as well as through online merchants. MDS Analytical Technologies recommends serial adapters made by Keyspan ([www.keyspan.com](http://www.keyspan.com)), as they have been proven repeatedly in field operations.

Please refer to the instrument manuals for additional instructions related to the USB-only instruments from MDS Analytical Technologies: the SpectraMax® L Luminometer and StakMax® plate handling system.

## Installing Single Port Adapters

To install single port adapters

1. Connect the MDS Analytical Technologies serial cable to the USB-to-serial adapter.
2. Connect the USB-to-serial adapter to an open USB port on your computer.
3. Press the black toggle switch on the back of the microplate instrument to the x position to turn the instrument on.
4. Start the SoftMax Pro software and click **Edit > Preferences**.
5. Select the appropriate serial port (most often COM1) from the Serial Port pull-down menu.

You should now see a three-dimensional image of your instrument in the top-left corner of SoftMax Pro software immediately beneath the File menu.

6. Click on the instrument image to confirm that your instrument is communicating properly. Your specific reader should be automatically selected in the Reader pull-down menu.

## Installing Multiple-port Adapters

Before attempting to use a multi-port adapter, be sure you have downloaded and installed the very latest software driver from the manufacturer's website. If you are using a two-port adapter from Keyspan such as their USA-28x Twin Serial Adapter, you **MUST** be sure to connect your MDS Analytical Technologies microplate reader to PORT 1. Ports are usually identified on the side or top of the adapter itself. With the 28-x, a small "1" is embossed in the plastic of the adapter.

To install multiple-port adapters

1. Connect the MDS Analytical Technologies serial cable to PORT 1 on your USB-to-serial adapter.
2. Connect the USB-to-serial adapter to an open USB port on your computer.
3. Press the black toggle switch on the back of the microplate instrument to the x position to turn the instrument on.
4. Launch SoftMax Pro software.
5. Click **Edit > Preferences**, and make sure the correct serial port is selected.



---

**Note:** If you are using a multiple-port adapter with more than one MDS Analytical Technologies microplate reader, be sure to use the serial cables that shipped with your MDS Analytical Technologies microplate instrument. Non-MDS Analytical Technologies serial cables will **NOT** work with MDS Analytical Technologies microplate instruments.

---



---

**Note:** These USB-to-serial adapters do **NOT** work without installing the correct software driver. If you are not sure you have or are using the latest driver available, check with the manufacturer directly. Drivers can normally be downloaded from the manufacturer's website. Be sure you have installed the latest software drivers before using a USB-to-Serial adapter with your MDS Analytical Technologies microplate instrument.

---



## Troubleshooting Instrument Connections

If the instrument is properly connected and turned on, you will see the instrument's icon in the upper-left corner of the SoftMax Pro software window below the menus, and the correct instrument is selected in the Edit > Preferences dialog box.

If the instrument is not properly connected or is turned off, the instrument icon in the upper left-hand corner of the window will appear with a question mark. Check that the connection is secure and that the instrument is turned on.

### LMax II

The LMax II and LMax II384 instruments require manual selection of the instrument in the Preference dialog box in order to establish communication. This is due to the different baud rate requirement for LMax II instruments.

### Instrument Preferences

If you have problems communicating with a connected instrument, open the Edit > Preferences dialog box and select the appropriate communications port (ports that are not available are disabled):

- Windows: use COM1 through COM9.
- Mac: use a USB-to-serial adapter with an available USB port.
- If you do not have an instrument connected to the computer but want to use SoftMax Pro software, select None.
- Click **OK** to close the Preferences dialog box and the instrument icon should appear without a question mark.

If the serial port setting is correct and the instrument is on, yet communication is still failing, close SoftMax Pro software, turn off power to both the computer and the instrument, and check that the cable connections between the instrument and computer are secure. Turn both machines on again and restart SoftMax Pro software.

For more on instrument preferences, see [Setting Preferences on page 18](#).

## Setting Preferences

Before using SoftMax Pro software we recommend that you check the global application settings in the Edit > Preferences dialog box. Here you will find:

- Reader settings, such as the port used for the connection to the instrument, supported readers, and filter settings for use with Series I microplate readers.
- Automatic file saving and printing options.
- File export settings when data is exported using the File > Import > Export menu item.

## Reader Settings

### Reader

When a reader is connected and is powered on, the Reader list automatically reports the instrument connected to the computer.

If a reader is not connected to the computer, the Reader list displays the model of the last reader connected.

A protocol file can be written if an instrument is neither connected to the computer nor selected in the preferences. Choosing an instrument can be useful, however, since the options available in the Plate > Settings dialog box depend on the choice of instrument here. For example, with a SpectraMax® reader you can configure up to six wavelengths, whereas you cannot do this with a VMax® reader.

### Filters

Filter settings apply only to EMax®, VMax®, UVMax, and ThermoMax instruments.

Check that the filter wavelength displayed for each position matches the wavelength of the filter physically installed in each position of the filter wheel. If the settings in this dialog box are not correct, subsequent readings will generate incorrect data, because SoftMax Pro software has no way to verify the identity of the filters in the filter wheel.

The SpectraMax 250 reader accepts two optional filters, but the wavelengths for these filters are specified using the instrument's front panel. For specific information regarding how to install optional filters in the SpectraMax 250 reader, refer to the instrument manual.

## Serial Port

Check that the setting for the serial port matches the physical port used to connect the computer and the instrument.

If you are using a Keyspan serial adapter with a Mac computer, set the port to Printer.

### None

Select None to use the serial ports on your computer for other purposes, such as a modem connection, a network connection, or printing.

When None is selected, the Status Bar in SoftMax Pro software shows an instrument icon with a ? through it, along with the words No port selected. You cannot communicate with your instrument.

## Port Speed

When a FlexStation instrument is selected as the reader, a Port Speed option is available.

SoftMax Pro software automatically detects the baud rate of the instrument to which the computer is connected and establishes a communication rate accordingly. This option allows you to view the baud rate that is being used and to manually set a different rate. This should not be done unless a serial communications problem occurs.

## Manual Export Format

Data collected by SoftMax Pro software can be exported for use with other software or for archiving.

### Time / Plate / XML

For Plate sections, you can choose one of the following three formats:

- Time exports data in single column text for each well.
- Plate exports data in a text matrix corresponding to a microplate grid.
- XML exports data in an XML file format.

CuvetteSet sections are automatically exported in the time format.

### **Include Labels**

Includes microplate column labels for all exported data.

This is most useful for Group section data, since the column headings in the Group table are shown, including any other calculated data in Summaries.

Labels included with Plate and CuvetteSet sections show "1-12" headers for 96-well Plate section data exported in Plate format; 384-well plate data have "1-24" headers; CuvetteSet section labels show "A1," "A2," representing the cuvettes shown in the CuvetteSet section.

### **Include Filename and Date**

Includes the file name and date in the file with the exported data.

### **Include Account Info**

Includes the name of the user who generated the file. This option is available only if you are using SoftMax Pro GxP software.

### **Interpolate Wells (FlexStation Only)**

Interpolation is used to normalize the Kinetic read times of wells against the read time of the first well in each column. Because injection occurs at the same time in all wells of a column but the wells are read sequentially, the reaction advances in the other wells while the first is being read. Consequently, the data in the other wells must be interpolated against the amount of time that has elapsed between injection and reading in order to compare data from one well to another, and to compare data from one run to another.

Once exported, well-interpolated data cannot be re-imported into SoftMax Pro software.

This interpolation is different from the interpolation option in the **Plate > Reduction** dialog box.

### **384 to 4 × 96 time format**

This is used when you are using 384 well plates and conducting extended Kinetic reads with more than 255 time points. This option enables you to work around architectural issues in some versions of Microsoft Excel that prevent the import of data files with more than 255 columns.

## Save Data After Read (AutoSave)

When this option is enabled, the collected data is saved automatically to a user-defined location immediately after each plate read is completed. This greatly reduces the likelihood of lost data, particularly when AutoSave is set to save files to corporate network volumes that are backed up on a regular basis.

You can add as many AutoSave instances as you wish, each with its own settings. After clicking Add, you are presented with the Add AutoSave Entry dialog box that allows you to specify the location, file naming conventions, and file format. Once specified, simply click OK to save these settings.

To modify the AutoSave settings, in the Preferences dialog box, select the instance you want to modify and then click Edit.

### Location

You can control where files are AutoSaved by modifying the options listed in the Location section. The default location set in SoftMax Pro software is an invalid dummy path that should be deleted before saving another location. To save data files to an assigned folder, click Assigned Folder and choose a different folder.

If the AutoSave function reports an error or if files are saved in the wrong location, it may be that a previous version of SoftMax Pro software was not uninstalled before a newer version was installed.

### Name

You can control the automatic naming of all AutoSaved files by modifying the options listed in the Name section.

#### Protocol's Name

If selected, each AutoSaved file begins with the name of the protocol used to generate the data that has been AutoSaved.

#### Assigned Name

If selected, any text in this field is automatically inserted at the beginning of each AutoSaved file name. Files are also numbered sequentially, for example "Data 1", "Data 2".

If the Assigned Name field is left blank, the date and time of each plate read is used as the file name for each AutoSaved file. For example, the first file AutoSaved on November 18th, 2005 at 9:01am and 58 seconds would be AutoSaved as "11-18-05-090158".

### **Append UserID (GxP ONLY)**

If selected, Append UserID adds the UserID for the user logged on to SoftMax Pro GxP to the Assigned Name or Protocol Name specified.

### **Append Date**

If selected, Append Date adds the date to the Assigned Name or Protocol Name specified.

If Append Time is not also selected, SoftMax Pro software indexes the runs for that day starting with "1" and increments with each successive run. For example, the first file AutoSaved on November 18th, 2005 at 9:01am and 58 seconds would be named as "Data-11-18-05 1", while the second file AutoSaved would be named "Data-11-18-05 2", and so on.

### **Append Time**

If selected, Append Time adds the time in the format "hhmmss" to the Assigned Name or Protocol Name specified. For example, the first file AutoSaved on November 18th, 2005 at 9:01am and 58 seconds would be named as "Data-090158 1", while the second file AutoSaved would be named at the exact time the second read completed, "Data-090744 1", and so on.

### **Format**

Format options allow you to specify what type of file is AutoSaved. Version 5 supports three file formats: SoftMax Pro file format, XML, and Tab-Delimited.

After each Plate section or CuvetteSet section is read, the data is saved to a new file. For protocols with more than one Plate section, each subsequent AutoSave file contains all the data from the just-completed plate read, as well as all previous readings. For example, if you open a protocol containing two Plate sections and a CuvetteSet section and read all three with AutoSave enabled:

- File1 contains data for PlateSection1.
- File2 contains data for PlateSection1 AND PlateSection2.
- File3 contains data for both Plate sections AND the CuvetteSet section.

This functionality applies to both SoftMax Pro and text file formats.

### **SoftMax Pro Program File**

When selected, your data will be AutoSaved as a proprietary SoftMax Pro program file format: .pda (standard) or .eda (GxP) file format. If your protocol contains more than one experiment section, only the Experiment containing the plate/cuvette being read is automatically saved.

### **XML File**

XML is supported for data export and AutoSave. XML is the optimal file format if you plan on importing all read data into other data collection and storage applications—specifically, LIMS (Laboratory Information Management System) or SDMS (Scientific Data Management System) packages.

### **Tab-Delimited File**

When selected, your data will be AutoSaved as a .txt file format, which can be opened by any word processor, spreadsheet, or database program. If you are saving data to a text file, use Append to File to save all data for all Plate and CuvetteSet sections to a single .txt file, or use Create New File to save the data from each plate to a new file.

## **Print Document After Read**

If selected, this box enables automatic printing of a report after each plate is read. The report contains information from each section that is designated to be included in the report. For more information on reports, and including or excluding sections, see [Printing on page 137](#).

## The Protocols Menu

The Protocols menu is a dynamically generated menu, listing all top-level folders and the protocol files they contain. When SoftMax Pro software is installed, several folders—each containing predefined protocol files for use with SoftMax Pro software—are automatically created within the Program Files / Molecular Devices / SoftMax Pro folder.

To change the folder used to generate the Protocols menu, click Protocols > Set Folder and navigate to the appropriate folder. This allows you to display a custom set of protocols that your company may be using instead of the set that ships with SoftMax Pro software.

Additionally, custom folders can be created within Program Files / Molecular Devices / SoftMax Pro. Once a protocol file (.ppr or .epr file) is placed in this custom folder, both the folder name and the protocol name are listed in the Protocols menu.

Some protocols will not be appropriate for your particular instrument (they might have reductions that are appropriate only for absorbance or fluorescence instruments, for example), so check the Revision Notes section located at the top of each MDS Analytical Technologies-authored protocol file for reader suitability before using the protocol.

All protocol files can be opened, run, and customized for any instrument, whether or not SoftMax Pro software is connected to an actual instrument see [Setting Preferences on page 18](#).

This can be useful in the following circumstances:

- If you want to create protocol files on a computer that is running SoftMax Pro software but is not connected to an instrument.
- If you want to create protocols for an instrument other than the one connected to the computer.



This chapter describes some of the main elements of the SoftMax® Pro software interface, and their logical relationships to each other.

For detailed instructions on setting up an experiment, and acquiring and analyzing data, see Chapters 4 through 6.

## The SoftMax® Pro Application Window

The SoftMax Pro software application window consists of two elements:

- A Menu Bar and Status Bar at the top of the application window.
- An untitled document showing an experiment data file divided into a number of sections. The displayed sections depend upon the configuration of the default protocol, and the type of instrument chosen in the Edit > Preferences or connected to the computer.

## WYSIWYG Display

SoftMax Pro software has been designed to provide you with a WYSIWYG (What You See Is What You Get) preview of printed reports.

Consequently, the width of a document is constrained by the width set for the printed page size.

You can change the dimensions and orientation of the printed page and consequently the display size in the File > Print Setup (Windows) or File > Page Setup (Mac) dialog boxes.

## Status Bar

Below the menus at the top of the SoftMax Pro program window is a Status Bar that displays instrument controls. The Status Bar can be shown and hidden using View > Show Status and View > Hide Status respectively.

Depending on the type of instrument that is connected to the computer or selected in the Edit > Preferences dialog box, some or all of the following items are in the Status Bar:

### Instrument Icon

Indicates that the instrument is communicating properly with the computer. If an "?" appears on top of the icon, the computer is not making proper contact with the instrument.

Click the instrument icon to open the Edit > Preferences dialog box.

### Temperature Display

Reports the current temperature within the microplate chamber in degrees Celsius.

With the SpectraMax® M2, M2e, M5, M5e, Plus, and Plus<sup>384</sup> readers, the SoftMax Pro Status Bar shows the temperature within the microplate chamber. This may be different from the front panel of the instrument, which displays the temperature within the cuvette chamber. The readings should be very similar to one another after both chambers have reached equilibrium. During warm-up, however, you may notice a discrepancy in temperatures, which is normal.

### Read

If a Plate or CuvetteSet section is active or if only a single Plate or CuvetteSet section has been created in the Experiment section, clicking the Read button starts the reading. This reading is based on the settings chosen in the Instrument Settings dialog box.

If more than one Plate or CuvetteSet section has been created in the Experiment section and no Plate or CuvetteSet section is currently active, click the Read button to open a dialog box that asks you to select the plate or cuvette to read.

If AutoRead is selected in the Instrument Settings, all of the plates/cuvettes will be read sequentially.

## **Reference Button**

This feature is active only when a CuvetteSet is selected. Use the Reference button to measure a reagent blank. The Reference value is stored and subtracted automatically from the Read values. The reference reading may be taken before or after the sample reading.

This setting applies only to the SpectraMax® M2, M2e, M5, M5e, Plus, and Plus384 readers.

## **StakMax® Software Interface Button**

Displays the StakMax® software interface, allowing you to control the StakMax microplate handling system. The StakMax software interface accessed through SoftMax Pro software is the only method of interacting with the StakMax instrument.

When the StakMax software is started it automatically connects to the StakMax instrument, and the current status of the instrument is displayed in the Status line in the StakMax software.

The StakMax system is compatible with SpectraMax L, M2, M2e, M5, M5e, Gemini EM and XPS readers, and the new models of the following instruments (compatible models are identifiable by their serial numbers): SpectraMax Plus384 (Serial#-MNR-xxxxx), 340PC384 (Serial#- LNR-xxxxx), 190 (Serial#-NNR-xxxxx), VersaMax (Serial#-BNR-xxxxx).

Click the Help button in the StakMax software user interface to open the StakMax User Guide, which provides complete documentation on the use of the StakMax system.

## **Incubator Button**

Displays the Control > Incubator dialog box, allowing you to set and regulate the temperature of the microplate/cuvette chamber. The temperature setting can be left at the default or can be specified by typing a different value.

The incubator setting is independent of the protocol being run. Running an assay does not automatically set the temperature set point. After a reading, however, the temperature set point, range, and average temperature are recorded in the saved file.

## **Automix Button**

Shakes the microplate. This is a manual shaking of the plate as opposed to an automatic shaking that can be chosen in the instrument settings and is not available for the EMax® reader.

## **Drawer Button**

Opens or closes the microplate drawer.

# **Instrument Settings and Control**

In addition to the buttons on the Status Bar, most commands for instrument control are in the Control menu.

## **Instrument Settings**

Use the Settings dialog box to configure all parameters for a plate read, such as the read type (End-point, Kinetic, Fast Kinetic, Spectrum, Well Scan, or Flex), read mode (Absorbance, Fluorescence, Luminescence, Fluorescence Polarization, Time Resolved Fluorescence), wavelength settings, Automix parameters, pre-read blanking, strips, PathCheck, Auto-Read, timing for Kinetic runs, Speed Read for Spectrum scans, and plate type.

Each Plate or CuvetteSet section can have a different instrument setup.

The available options in the Settings dialog box depend on the type of instrument that is connected to the computer, or selected in the Preferences dialog box see [Setting Preferences on page 18](#).

The Settings dialog box is only available when a Plate section is selected in the data file.

## Read

In a single plate experiment, the Read command reads the plate or cuvette immediately.

If no Plate section or CuvetteSet section is active, the Read command displays a dialog box in which you select the Plate section or CuvetteSet section to read.

The default Plate section or CuvetteSet section, selected by pressing the Return key, is chosen on these criteria:

- The selected Plate or CuvetteSet section, if one is selected.
- The next unread Plate or CuvetteSet section.
- The first plate or cuvette, if all have been read.

Menus are disabled while reading is in process.

## Close Drawer, Open Drawer

Use these commands to close and open the instrument drawer so that a plate can be inserted or removed.

## Incubator

The incubator command allows you to turn the incubator on or off. When on, you can enter a setting to regulate the temperature in the microplate chamber.

This command is available for all instruments except Emax, VMax, UVMax and LMax II readers.

## Ref

The Ref command causes the SpectraMax M2, M2<sup>e</sup>, M5, M5<sup>e</sup>, Plus, or Plus<sup>384</sup> readers to read the cuvette in the cuvette port and applies the reading as a reference to all cuvettes in the active CuvetteSet section.

This command is available only when a CuvetteSet section is active and a SpectraMax M2, M2<sup>e</sup>, M5, M5<sup>e</sup>, Plus, or Plus<sup>384</sup> readers is connected or when you open the Preferences dialog box.

## Experiments

SoftMax Pro software collects data from one or more microplates or cuvettes and stores it in a single data file, using the same or different instrument settings for different microplates or cuvettes. For example, microplates containing different samples can be read using the same or different modes, all within the same experiment.

Each SoftMax Pro program file contains at least one experiment. Within an experiment there are one or more sections.

Experiments provide quick access to different types of information without the need to open or refer to multiple files.

### Selecting an Experiment

An experiment can be selected by clicking the Experiment title bar (the area that shows the experiment number) with the mouse cursor. Selecting an Experiment changes its tool bar from light gray to dark gray, and allows you to manipulate the whole Experiment, such as deleting or duplicating it, or adding new sections.

Multiple Experiments can be selected by holding down the Shift key while clicking the individual title bars of the Experiments.

### Manipulating Experiments

#### New Experiment

To reate a new Experiment click the **Experiment > New Experiment**.

If an existing Experiment is active when this command is chosen, all of the information from the previous Experiment except for the data is copied to the new Experiment.

#### Rename Experiment

When an Experiment is created it is given a placeholder name, like "Experiment#1".

To rename an Experiment, double-click on its title bar to open an edit dialog box.

#### Duplicate Experiment

Trp copy of an experiment (excluding data) in the same file, clkick **Edit > Duplicate** [Experiment Name].

To duplicate multiple experiments , shift-click on them and then click **Edit > Duplicate Selection**.

### **Delete Experiment**

To delete a selected experiment, click the **Edit > Delete** [Experiment Name].

To delete multiple experiments, by shift-click on them and then click **Edit > Delete Selection**.

### **Reorder Experiments**

To change the order that experiments appear in a file, drag and drop. Click the title bar of the experiment and drag it to its new position in the file, marked by a thick black line.

## **Sections**

A section is a part of the SoftMax Pro program Experiment window intended to perform a specific set of functions. There are five types of sections—Notes, Plate, CuvetteSet, Group, and Graph—and each type of section has a specific icon that identifies it.

### **Active Section Menus**

When a section is active, a section-specific menu is added to the SoftMax Pro program main application menus between the View menu and the Protocols menu: Notes, Plate, Cuvette, Group, or Graph.

### **Manipulating Sections**

Like Experiments, sections are either active or inactive. Make a section active by clicking it with the mouse, or by selecting it from the **View > Experiment** sub-menu.

When active, the section title bar is colored dark gray. Multiple sections can be made active at the same time by Shift-selecting them.

### **New Section**

You can create a new section (except Group sections) by selecting the appropriate New command from the Experiment menu: New Notes, New Plate, New CuvetteSet, New Graph.

If an existing section is active when this command is chosen, all of the information in the section except data is copied to the new section.

Group sections are created automatically when you define new groups in the Template Editor, and can also be created with the Duplicate command.

### **Section Name**

When a section is created it is given a name based on its section type, like "Notes#1".

To rename a section by double-click on its title bar to open the Section Name dialog box, or select Section Name from the Section (Notes, Plate, Cuvette, Graph) menu.

### **Duplicate Section**

To make a copy of a selected section (excluding data) in the same file select **Edit > Duplicate** [Section Name].

To duplicate multiple sections, by shift-click them and then select **Edit > Duplicate Selection**.

Duplicating a CuvetteSet section creates a copy of the previous CuvetteSet section, including the last reference that was read.

### **Delete Section**

To delete a selected section, select the **Edit > Delete** [Section Name].

To delete multiple sections, shift-click on them and then select **Edit > Delete Selection**.

Deleting a Group section automatically deletes references to that group from the Template.

### **View Section**

Each experiment in a file has its name added to the View menu as a command, and all its sections are available in cascading sub-menu commands. Selecting a section from one of these sub-menus opens the section in the current window.

### **Expand and Collapse Sections**

Newly created experiments and sections are initially shown expanded. To simplify navigation within an Experiment data file, you can collapse sections to show just their tool bars.



To expand or collapse a single section, click its triangular indicator located in the top left corner of the tool bar.

To expand or collapse all sections in an experiment, or multiple experiments in a file, use the **View > Expand** and **View > Minimize** commands, or their associated hot keys

Ctrl+[ and Ctrl+].

### **Reorder Sections**

To change the order that sections appear in an Experiment, click on them and then drag and drop them to where you want them to be. Select the title bar of the section and drag it to its new position in the experiment, marked by a thick black line.

### **Open Section in New Window**

To open an active section and edit it in a separate window, select New Window from the active section menu (Plate, Cuvette, Graph, Notes, or Group).

Closing the new section window or expanding the section in the original window saves changes made to the section.

### **Print Section**

The **File > Print** [Section Name] commands enables you to print just the active sections.

### **Notes Section**

Notes sections are used to record text or to report Summary data pertaining to the Experiment. You can enter text in this section by typing directly, or by creating Summaries containing formulas for displaying reduced data.

A Summary is part of a Notes section that is generated from formulas applied to data in the Experiment.

Text in Notes sections can be formatted (font face, font size, and font style) using the Edit > Text Style dialog box.

You can create multiple Notes sections within the same Experiment.

### **New Summary**

Notes > New Summary opens a dialog box allowing you to:

- Name a new Summary.
- Enter a formula to be used for the Summary.

### **Edit Summary**

Use Notes > Edit Summary to edit the name and formula of an existing Summary.

See the *SoftMax Pro Software Formula Reference Guide* for a complete guide to formulas.

### **Plate Section**

Plate sections are used to collect data from the instrument, and to define data display and data reduction. If you read the same physical plate twice with different instrument settings, you would have two Plate sections.

The Template Editor in the Plate section is used to create a map of the contents of the microplate.

Plate sections are divided into:

- Tool bar.
- Data display (shown in a microplate grid format).
- Instrument settings, shown to the right of or below the data display, depending on the plate type and printer settings.
- Reduction settings, shown to the right of or below the data display, depending on the plate type and printer settings.
- Fluid transfer and checksum status for the FlexStation instrument.

If the plate grid in the Plate section is colored, then a template has been defined for the Plate section. Each group defined in the template has a different color (the icon of the corresponding group table has the same color).

If your experiment requires multiple plates, you can create as many Plate sections as needed.

### **Template**

The Template button (or **Plate > Template**) opens the Template Editor, which allows you to describe the samples in the wells, providing the link between the raw data and analysis groups. You can create new groups in the template or assign wells to already established groups.

A group is a set of associated samples that has been created in the Template. Some group types (such as Standard or Unknown) can be created as part of the default protocol, and you can create others as required. When you create a group and assign samples to it, a corresponding group table is automatically created.

Each Plate section in your Experiment requires a template to be created to identify groups within that plate. However, two or more Plate sections may use the same template, or the same groups may be used in different templates.

### **Reduction**

The Reduction button (or **Plate > Reduction**) opens the Reduction dialog box, which is used to specify the reduction of the data collected from the microplate reader, typically to a single number per well.

Options available in the Reduction dialog box depend on the type of reading performed (Endpoint, Kinetic, Spectrum, Well Scan, or Flex) and the instrument used to collect the data.

See [Chapter 5, Data Analysis](#) for more details on Data Reduction.

### **Graph**

The Graph button (or **Plate > Graph**) creates a graph from the data in the selected wells in the active Plate section. It is enabled both during data collection and after data has been acquired for a Kinetic, Spectrum, Well Scan, or Flex reading.

Double-clicking a well also creates a graph of the data in that well. Graph is disabled if the instrument settings are for an Endpoint reading.

### **Mask**

The Mask button (or **Plate > Mask**) allows you to remove selected wells from the data analysis.

Mask is a toggle, so that selecting masked wells and choosing the Mask command unmask them.

### **Copy and Paste Data**

Data can be copied from one Plate section to another or from one CuvetteSet section to another within SoftMax Pro software:

1. Make the source section active by clicking on it.
2. Select **Edit > Copy** (**⌘C**; Ctrl+C).

3. Create a new section, if necessary, or make the destination section active.
4. Select **Edit > Paste** (⌘V; Ctrl+V).

When copying Plate sections, all wells are copied, even if only a subset of wells is selected. It is not possible to copy and paste the data from selected wells only.

It is possible, however, to copy individual cuvettes or groups of cuvettes and paste these into other CuvetteSet sections. If you are copying cuvette data and the target CuvetteSet section does not contain the appropriate cuvettes for the data that is to be pasted, new cuvettes are created automatically. Note that pasting cuvette data for only some cuvettes clears the existing data from all other cuvettes in that section.

If the Instrument Settings of the target section do not match the settings from the section being copied, the Instrument Settings for the target section are updated to match.

### **Copy and Paste Template**

There are two ways to copy and paste a template, depending on whether or not you want to keep the same group names.

#### **Keep Group Names**

**Plate > Copy Template** copies the template from the active Plate section to the Clipboard, so that it can be pasted into a different Plate section using **Plate > Paste Template**, either in the same or in a different Experiment, or into a Plate section in another SoftMax Pro file.

Pasting a template into an Experiment where the groups from the source template do not exist creates the groups. Existing groups of the same name are not changed by the pasting process.

For example, if the source Experiment contains the groups "Standards" and "Unknown," and the destination Experiment already contains the group "Standards", the group "Unknown" is created but the group "Standards" in the destination Experiment is unchanged.

If you copy and paste the template within the same Experiment, no new groups are created. Group names can be edited in the pasted template, but these changes are made in the original template also.

#### **New Group Names**

The second way to copy and paste a template allows you to keep the same sample lay-out and sample names and concentrations, but also

allows you to change the Group names independent of the source template:

1. Open the Template Editor and select all wells.
2. Select ⌘C or Ctrl+C to copy the layout.
3. Open the Template Editor for a new Plate section and select all wells.
4. Select ⌘V or Ctrl+V to paste the layout.

### **Export Template**

Use Plate > Export Template to export a template to a tab-delimited text or XML file. The file format must be specified first in the global Preferences dialog.

### **Import Template**

Use Plate > Import Template to import a template from a tab-delimited text or XML file.

### **Group Section**

A Group section is a table that shows all of the information for a particular sample group that has been defined in a Plate template.

Unlike other section types, Group sections are created automatically when you create or select groups for an Experiment in the Template Editor.

Group sections can also be created by duplicating existing Group sections.

Group sections are divided into:

- Tool bar.
- The body of the Group table.
- Summaries.

### **Group Settings**

Use the Group > Group Settings dialog box to edit the Group name, descriptors, units and format.

The Column Format field in the Group Settings dialog box specifies the default columns for the data calculated and reported in the active group.

## Delete Group

You can delete a selected group with the **Edit > Delete [Group Name]** command.

Deleting a Group section automatically deletes references to that group from the Tem-plate.

## Duplicate Group

You can duplicate a selected group with the **Edit > Duplicate [Group Name]** command.

Group section tables can also be duplicated. When a Group table is duplicated, its name is added to the Template Editor's list of group names (but the group is not assigned to any wells).

## CuvetteSet Section

CuvetteSet sections are used to collect data from the cuvette port of the SpectraMax M2, M2e, M5, M5e, Plus or Plus384, readers to define an analysis template, and to define data display and data reduction. The Template Editor in the CuvetteSet section is used to describe the contents of the Cuvette.

CuvetteSet sections are divided into:

- Tool bar.
- Data display (can be shown in three different ways: three samples per row, one sample per row, and in a grid of cells similar to a microplate format).
- Instrument settings, shown to the right of the data display.
- Reduction settings, shown at the bottom of the data display.

When colors appear in the name labels above individual cuvettes (when displayed as one or three cuvettes per row) or when colored areas appear when the CuvetteSet section is displayed as a grid, this means a template has been defined for the CuvetteSet section. Each group defined in the template has a different color (the icon of the corresponding group table has the same color).

If your experiment requires multiple cuvettes, you can add up to 96 cuvettes in a single CuvetteSet section or create more than one CuvetteSet section.

The commands described in the Plate section above also apply to CuvetteSet sections.

## Graph Section

Graph sections are used to plot information from groups as scatter plots or bar graphs.

Once a graph has been created, new plots can be added and deleted, the axes can be customized, and the size of the graph can be changed. Grid lines can be enabled or disabled (default is enabled).

Graph sections are divided into:

- Tool bar.
- The body of the Graph section.
- Plot information below the graph.

You can create more than one Graph section within an Experiment data file, and plots in the Graph section can be created from any Experiment in the file.

## Graph Options

Use the Graph Options button (or **Graph > Graph Options**) to configure the style and content of a graph.

You can select from three types of graphs:

- Scatter (with symbols plotted and the points connected). - Cluster Bar.
- Stack Bar.

Select New to add a new plot to the graph from any of the Group sections in the Experiment.

## X-Axis

Use the X-Axis button (or **Graph > X-Axis**) to edit X-axis parameters (tick marks, scale, label, gridlines, and auto-range).

## Y-Axis

Use the Y-Axis button (or **Graph > Y-Axis**) to edit Y-axis parameters (tick marks, scale, label, gridlines, and auto-range).

## Export Graph

Exports a graph to file.

On the Mac, the graph is saved in PICT format; on Windows, it is saved in EMF (Enhanced Metafile) format.

## Creating a Default Protocol

While earlier versions of SoftMax Pro software opened to a specific Tutorial template, SoftMax Pro software versions 4.8 and later open to a blank protocol that matches the instrument connected to the computer.

If no instrument is connected to the computer and if no default protocol file exists, SoftMax Pro software uses the instrument chosen in **Edit > Preferences** to create a new protocol.

To use a customized protocol file as your default template every time SoftMax Pro software is

launched, you must specifically save a file as your default protocol:

1. Create a new file using the **File > New** menu command.
2. Modify all preferences and settings as necessary.
3. Add any desired text information in Notes sections.
4. Customize the columns, formulas, and reductions to suit your research needs.
5. When the file is to your liking, select the **Protocols > Save As Default Protocol** menu item to save this as your new default protocol that will be displayed every time SoftMax Pro software is launched.

It is not mandatory to create a default protocol because the software detects the instrument connected to the computer and automatically creates a new file based on instrument. For example, if connected to a Fluorescence instrument with a cuvette port, the software displays fluorescence units on both a Plate section and a CuvetteSet section; if connected to an Absorbance instrument with no cuvette port, the software displays absorbance units on a Plate section only. In addition to these sections, a blank Notes section is created.

Finally, you can change your default protocol at any time. To remove the default protocol used, simply delete the file "Default Protocol" from Program Files / Molecular Devices / SoftMax Pro v5.4.



## Introduction to Data Collection

A typical process for preparing, collecting, and analyzing data is as follows:

1. Launch SoftMax® Pro software.
2. Create Plate and CuvetteSet sections as needed.
3. Specify instrument settings using **Plate > Settings**.
4. Define templates, reduction parameters, and display parameters.
5. Prepare the microplates or cuvettes to be read and place them in the instrument.
6. Start the reading.
7. Save the data file.
8. If required, modify templates, reduction parameters, or display parameters.
9. Analyze the data using Group section tables and Graph sections.
10. Save the final data file.

## Instrument Settings

Instrument settings are associated with a Plate or CuvetteSet section in a SoftMax Pro Experiment. You can create more than one Plate or CuvetteSet section in a single Experiment, and the instrument settings associated with each Plate section can be different.

You must specify instrument settings (wavelengths and read mode) prior to reading a microplate or cuvette.

When you have defined instrument settings, a summary is displayed to the right of the data display in the Plate or CuvetteSet section.

## Instrument Operation

Instrument settings are made in the Plate > Settings dialog box. You must select a Plate section to make this menu available.

The operation of individual instruments is only described here to illustrate the functionality of the software. For complete information about the function of the instrument or how to set it up for a reading, consult the instrument manual.

## **Instrument Settings Vary for Different Models**

Only the settings relevant to the instrument connected to the computer or chosen in Edit > Preferences are available in the Plate > Settings dialog box. However, this guide lists many of the common settings across instrument models.

### **Read Type**

The Read type specifies the spatial, temporal and chromatic properties of the read.

#### **Endpoint**

In an Endpoint read, a reading of each microplate well is taken at a single or at multiple wavelengths.

Depending on the read type selected, values can be reported as optical density, %Transmittance (if this is selected in the Reduction dialog box), relative fluorescence units (RFU), or relative luminescence units (RLU).

For multiple reads per well, see [Kinetic Reads on page 90](#) and [Well Scan Editor on page 59](#).

#### **LMax II and LMax II384 and SpectraMax® L Readers Only**

Endpoint on an LMax® II or SpectraMax® L instrument line means:

- ◆ A single integration of samples in 96- or 384-well plates.
- ◆ Two injections, if desired, in selected wells of a 96-or 384-well plate (only one injection is possible in 384-well plates with the LMax II instrument line).

A pre-read of the microplate can be made before injection of buffer or reagent, or reading of the samples. Normal pre-read uses the same integration time as sample read, while Special pre-read allows you to enter a different integration time.

Values are reported as relative luminescence units (RLU); the default reduced value is !Lm1.

## **Kinetic**

In a Kinetic read the data are collected over time with multiple readings taken at regular intervals. To achieve the shortest possible interval for Kinetic readings, choose wavelengths in ascending order.

The default reduced values calculated for Kinetic data are VMax (units/sec or units/ min), Time to VMax (seconds), Onset Time (seconds), Time at Minimum, Time at Maximum, Time at 1/2 Maximum, Slope, and Area Under Curve.

## **Spectrum**

Depending on the read mode selected, a Spectrum read measures optical density, %Transmittance (if this has been selected in the Reduction dialog box), relative fluorescence units (RFU), or relative luminescence units (RLU) across a spectrum of wavelengths.

All Spectrum readings are made using the scanning monochromator of the instrument.

## **Well Scan**

A Well Scan read takes one or more readings of a single well of a microplate at single or multiple wavelengths. Every option available for Endpoint reads is available for Well Scans.

Values reported are optical density, %Transmittance, relative fluorescence units (RFU), or relative luminescence units (RLU).

## **Flex**

Flex is used to set up reads on FlexStation® instruments, where one has to configure compound additions as well as the usual plate reading parameters.

## **Dual Read**

### **(SpectraMax L and LMax II384 Readers Only)**

Also an Endpoint reading, Dual Read is designed to perform two separate integrations, one after each of two possible injections in any or all wells of a 96-well plate or 384-well plate (LMax II384 can inject twice in 96-well plates only). Integration times for both readings are the same, unless otherwise specified.

You can program a delay after each injection, and you can read all or only some of the wells in the plate. Pre-reading a microplate is not possible with this read type.

Default values reported are relative luminescence units (RLU). The default reduced value is !Lm1.

### **Fast Kinetic**

#### **(SpectraMax L, LMax II and LMax II384 Readers Only)**

Fast Kinetic performs repeated readings of one or more wells of a 96- or 384-well microplate up to a 100-point maximum integration. All readings of a single well are made before the next well is read. Pre-reading a microplate is not possible in Fast Kinetic reads.

One or two injections can be made in each well of a 96-well microplate (and 384-well plate for the SpectraMax L reader) at the start of the reading. Integration time can be set from 0.1 to 100 seconds.

Default data reductions are VMax per Sec, VMax, Time to VMax, or Onset Time.

## **Read Mode**

The Read mode is the form of detection used by the instrument.

All read types except Well Scan are available for all read modes. Not all read modes are supported by all instruments.

### **Fluorescence (RFUs)**

Fluorescence is light emitted by a substance resulting from the absorption of incident radiation. The governing equation for fluorescence is:

Fluorescence = (extinction coefficient) × (concentration) × (quantum yield) × (excitation intensity) × (pathlength) × (emission collection efficiency)

### **Absorbance (ODs)**

Absorbance is the amount of light absorbed by a solution.

In the absence of turbidity:

Absorbance = optical density – reflection

### **Luminescence (RLUs)**

Luminescence is the emission of light by processes that derive energy from essentially non-thermal changes, the motion of subatomic particles, or the excitation of an atomic system by radiation.

With the SpectraMax L, M5 and M5e readers, selecting the Luminescence read mode displays the integration time box next to the Read Type selection.

### **Time Resolved Fluorescence (RFUs)**

Most fluorescent substances are not suitable for this type of reading. However, the fluorescence emitted by lanthanide dyes is delayed long enough to measure fluorescence after the lamp is turned off. Time resolved fluorescence is used to reduce the amount of background noise that interferes with fluorescence. The excitation lamp flashes, and after it is off, the delayed emission is collected for a set period of time before the lamp is flashed again.

With the SpectraMax M5 and M5e readers, Time Resolved Fluorescence read mode displays Integration Delay and Integration next to the read type selection instead of Start and End.

### **Fluorescence Polarization (RFUs)**

By using a fluorescent dye to label a small molecule, its binding to a large molecule can be monitored through its speed of rotation.

Fluorescence Polarization mode returns two sets of data: one for fluorescence intensity parallel (P) to the excitation plane, and the other for fluorescence intensity perpendicular (S) to the excitation plane. These S and P values are used to calculate the Polarization (mP) and Anisotropy (r) values in SoftMax Pro software. Although the Raw S&P value is the true raw data returned from the instrument, the calculated Polarization (mP) and Anisotropy (r) values are treated as the raw data, and these values become the basis for further reduction calculations. You can choose to display any of these data types on the Plate section.

When Raw (S&P) is displayed, the mP value is by default used for all reduction calculations. When exporting, the displayed raw values are exported, whether mP, r, or Raw (S&P).

Polarization (mP) is calculated as follows:

$$mP = 1000 \times \frac{(\text{parallel} - (G \times \text{perpendicular}))}{(\text{parallel} + (G \times \text{perpendicular}))}$$

Anisotropy (r) is calculated as follows:

$$r = \frac{(\text{parallel} - (G \times \text{perpendicular}))}{(\text{parallel} + (2G \times \text{perpendicular}))}$$

## Wavelengths

With certain instruments, when reading using Fluorescence and Time Resolved Fluorescence, you must also choose whether to read excitation at a fixed wavelength and scan emission wavelengths, or vice versa, within the range of the instrument (250 nm – 850 nm). A procedure to optimize excitation and emission wavelengths for a given protocol is outlined in the instrument manuals for the readers that support this function. Please see those manuals for more information.

### For Endpoint or Kinetic Readings

- Choose the number of wavelengths to read.
- Specify the wavelengths from the entries in the lists, or type any wavelength within the available range of the instrument over an existing setting. The lists show only common monochromator settings.

### For Spectrum Readings

Choose a Start and a Stop wavelength, and a wavelength step for each increment between reads.

The minimum selectable increment is 1 nm.

### Cutoffs

The choices in this portion of the Wavelengths setting depend upon the selections for read type (Endpoint, Kinetic, Spectrum, or Well Scan) and read mode (Fluorescence or Time Resolved Fluorescence). The term cutoff refers to filters used to block unwanted residual excitation light and minimize background interference.

No emission cutoff filter is used when Luminescence is selected. With other read types, choices are to enable or disable Auto Cutoff with specific filter settings available if it is disabled (manual cutoff selection). When Auto Cutoff is enabled, the instrument sets the cutoffs based upon the wavelengths chosen for reading; when Spectrum is selected as the read type, however, a manual setting for the emission monochromator is required, with the default being no cutoff filter.

Determining a manual setting for a cutoff filter is based upon the known value of the Stokes shift, which is the difference between the wavelengths of the excitation and emission maxima. If the Stokes shift is small, it may be advisable to choose an excitation wavelength that is as far as possible from the emission maximum while still being capable of exciting the fluorophore. Doing this causes less of the excited light to overlap the emission spectrum, allowing better selection and quantitation of the emitted light. See the individual instrument manuals for more information regarding cutoff filter settings.

## **Shake**

### **(LMax II AND LMax II384 Readers Only)**

This function shakes the plate to mix the injected reagent with the solution in the microplate. You can specify a duration in seconds.

Shaking begins upon clicking the Read button.

## **Integration**

### **(SpectraMax L, LMax II AND LMax II384 Readers Only)**

Enter the amount of time to perform a single integration.

Dual Read type allows the choice of Special, where you can specify a separate integration time for P and M injections.

For Kinetic, enter the total Run Time and the desired Interval between readings. The resulting number of Kinetic reads is reported below the run time.

For Fast Kinetic, enter the integration time followed by the number of times this integration should be repeated for each well before moving to the next well.

## **Injection and Delay**

### **(SpectraMax L AND LMax II384 Readers Only)**

The SpectraMax L and LMax II384 readers are equipped with two injectors having variable volumes. One (labeled M for measurement) injects directly in the reading position, and the other (labeled P for preinjection) injects into the preceding position.

High-precision pumps are used for the injections, ensuring both instantaneous sample mixing after injection and precise injection volumes.

The injection system can only be used with 96- and 384- well plates. SoftMax Pro software allows selection of one or both injectors for any well of a 96-well or 384-well plate (the LMax II384 can use both injectors in a 96-well plate only).

For the LMax II384 reader you can define the injection volume and the post-injection delay, which is the time between the last injection and the start of the reading.

For the SpectraMax L reader you can define a large range of injection parameters.

## **Injection Wells**

### **(SpectraMax L AND LMax II384 Readers Only)**

Use the Injection Wells options to define the wells to which P and M injections are applied. To apply injections:

1. Enable P-injection or M-injection on the Injection and Delay section of the Settings dialog box.
2. Select the Injection type on the Injection Wells section of the Settings dialog box by clicking the P injector or M injector icon. Only the wells specified in Wells to Read are displayed.
3. Click the wells to which you want the injections to apply.



## **Sensitivity**

(Gemini; SpectraMax L, M2, M2e, M5, M5e; FlexStation® Readers; Fluorescence/Luminescence Only)

### **Readings**

This setting determines the number of readings made in each well of the microplate, or each spot in a well for Well Scan reads. Readings are averaged, and the average reading is displayed.

The default number of readings per well varies with the read mode: for Fluorescence, the default is 6; for Luminescence, the default is 30; for Time Resolved Fluorescence, the default is 20. The number of readings can vary from 1 to 30.

### **PMT Sensitivity**

The PMT (photomultiplier tube) is a vacuum tube that detects light, especially from dim sources, through the use of photo emission and successive instances of secondary emission to produce enough electrons to generate a useful current.

PMT sensitivity can be set to low (for samples having a higher concentration), medium, or high (samples having lower concentration).

In Endpoint reads, an Automatic setting is available, allowing the instrument to adjust the PMT voltage automatically for varying concentrations of sample in the plate. Automatic is not available for Fluorescence Polarization reads; if Automatic was previously selected, Fluorescence Polarization defaults to Medium.

For the SpectraMax M5 and M5e readers, the PMT Sensitivity is always set to Automatic for Luminescence and Time Resolved Fluorescence read modes.

For the SpectraMax L reader, sensitivity options are Auto Range, Analog Only Photon Counting, and MaxRange.

### **Target Calibration Wavelength**

The SpectraMax L reader supplies the option to choose a target wavelength for your experiment, based on the assay you are using.

### **Timing**

For Kinetic reads you can configure the total Kinetic run time and the time interval between readings.

After entering a total run time and an interval between readings, the total number of readings is calculated and reported. The minimum read interval is determined by the instrument, and depends on many factors including the number of wavelengths, Automix status, the number of strips, and the distance the instrument's filter wheel must move.

If you are writing a protocol to be used with a different model instrument:

1. Select **Edit > Preferences**.
2. Set the Serial Port to None.
3. Select the other instrument from the Reader list.

Failure to do so could cause instrument errors when you attempt to read.

To achieve the shortest possible interval for Kinetic readings, choose wavelengths in ascending order.

The minimum interval for Flex reads is dependent on the Compound Transfer settings and the number of wells to be read per column.

### **PathCheck® Pathlength Correction**

**(SpectraMaxPLUS, Plus384, 190, 340PC, M2, M2e, M5, M5e Readers)**

The Beer–Lambert law states that absorbance is proportional to the distance that light travels through the sample:


$$A = \epsilon \cdot b \cdot c$$

where  $A$  is the absorbance,  $\epsilon$  is the molar absorptivity of the sample,  $b$  is the pathlength and  $c$  is the concentration of the sample. In short, the longer the pathlength, the higher the absorbance.

Microplate readers use a vertical light path so the distance of the light through the sample depends on the volume. This variable pathlength makes it difficult to perform extinction- based assays and also makes it confusing to compare results between microplate readers and spectrophotometers.

The standard pathlength of a cuvette is the conventional basis for quantifying the unique absorptivity properties of compounds in solution. Quantitative analyses can be performed on the basis of extinction coefficients, without standard curves (for example, NADH-based enzyme assays). When using a cuvette, the pathlength is known and is independent of sample volume, so absorbance is proportional to concentration.

In a microplate, pathlength is dependent on the liquid volume, so absorbance is proportional to both the concentration and the pathlength of the sample. Standard curves are often used to determine analyte concentrations in vertical-beam photometry of unknowns, yet errors can still arise from pipetting the samples and standards. The PathCheck feature automatically determines the pathlength of aqueous samples in the micro-plate and normalizes the absorbance in each well to a pathlength of 1 cm. This novel, patented approach to correcting the microwell absorbance values is accurate to within 2.5% of the values obtained directly in a 1-cm cuvette.

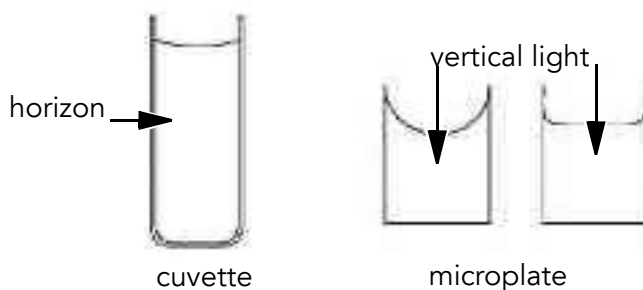

#### Figure 4-1 Cuvette and Microwell light paths

Reference measurements made by reading the cuvette or using factory-stored values

derived from deionized water can be used to normalize the OD data for microplate wells. PathCheck software is used to normalize the data acquired from absorbance endpoint microplate readings to a 1 cm pathlength, correcting the OD for each well to the value expected if the sample were read in a 1 cm cuvette.

The SpectraMax Plus, Plus384, M2, M2e, M5, and M5e readers offer two methods of pathlength correction: the first uses a cuvette reference, and the second uses a water constant that is stored in the instrument.

The SpectraMax 190 and 340PC use the water constant only.

The actual pathlength,  $d$ , of a solvent is found from the following equation:

$$d(\text{cm}) = \frac{\text{Sample}(OD_{1000} - OD_{900})}{k}$$

When a cuvette reference is used for pathlength correction, the value of  $k$  is obtained by taking OD measurements on the fluid in the cuvette at two wavelengths, 1000 and 900 nm:

$$k = \text{Cuvette}(OD_{1000} - OD_{900})$$

When the water constant is used for pathlength correction, the value of  $k$  is obtained from the instrument. This constant is saved in the instrument in the factory and may differ slightly from instrument to instrument.

Once the pathlength  $d$  is found, the following equation is used for the pathlength correction:

$$\frac{OD}{\text{cm}} = \frac{OD_{\text{sample}}}{d(\text{cm})}$$

### **PathCheck Software: Cuvette Reference or Water Constant (For SpectraMax Plus, Plus384, M2, M2e, M5 and M5e Readers)**

To enable PathCheck for Absorbance readings:

1. Open **Plate > Settings** and select PathCheck.
2. Check the PathCheck option.
3. Select either Water Constant or Cuvette Reference (see below).
4. If you have determined a plate background constant for your microplate, enter it in the designated field (see ["Use Plate Background Constant on page 55"](#)).

The PathCheck program measurement is based on the absorbance of water in the near infrared region (between 900 nm and 1000 nm). If the sample is completely aqueous, has no turbidity and has a low salt concentration (less than 0.5 M), the Water Constant is adequate. The Water Constant is determined during manufacture and is stored in the instrument.

If the sample contains an organic solvent such as ethanol or methanol, we recommend using the cuvette reference. It is important that the solvent does not absorb in the 900 nm – 1000 nm range (to determine whether or not a given solvent would interfere, see the discussion of interfering substances below). When a non-interfering solvent is added to the aqueous sample, the water absorbance decreases proportionally to the percentage of organic solvent present. For example, 5% ethanol decreases the water absorbance by 5% and results in a 5% underestimation of the pathlength. You can avoid the error by putting the same water/solvent mixture in a cuvette and using the Cuvette Reference.

To use the Cuvette Reference, place into the cuvette port a standard 1 cm cuvette containing the aqueous/solvent mixture that is used for the samples in the microplate. The cuvette must be in place when you read the microplate. When you click the Read button, the instrument first makes the 900 nm and 1000 nm measurements in the cuvette, and then makes the designated measurements in the microplate. The cuvette values are stored temporarily and used in the PathCheck program calculations for the microplate samples.

Use of Cuvette Reference with PathCheck software is different from a reference reading of a cuvette in a CuvetteSet section (by clicking the Ref button in the CuvetteSet section tool bar). The cuvette reference used for PathCheck calculations (measurements at 900 nm and 1000 nm) does not produce data that can be viewed in a CuvetteSet section and is used only with data in microplates, not cuvettes.

Once you have read a plate with the PathCheck program turned on, PathCheck application information is stored permanently in the data file. You have the option of applying, or not applying, the PathCheck application to the absorbance values as you choose. If you did not have the PathCheck program turned on during the plate read, you cannot apply the PathCheck program after the read.

## PathCheck on SpectraMax 190 and 340PC Readers

The PathCheck application on the SpectraMax190 and 340PC readers uses the water constant only. Be aware that if your sample matrix contains an organic solvent such as ethanol or methanol, the estimated pathlengths will be lower than the true values, and PathCheck program normalized values will be higher, than the corresponding 1-cm values (see discussion above):

- Open **Plate > Settings** and select PathCheck.
- Check the PathCheck option.
- If you have determined a plate background constant for your microplate, enter it in the designated field (see [Use Plate Background Constant on page 55](#)).

Once you have read a plate with the PathCheck application turned on, PathCheck program information is stored permanently in the data file. You have the option of applying, or not applying, the PathCheck application to the absorbance values as you choose. If you did not have the PathCheck application turned on during the plate read, you cannot apply the PathCheck program after the read.

## Background Constant Subtraction and Blanking Considerations

Raw OD measurements of microplate samples include both pathlength-dependent components (sample and solvent) and a pathlength-independent component (OD of microplate material). The latter must be eliminated from the PathCheck program calculation in order to get valid results that have been normalized by the PathCheck program. There are 3 ways to accomplish this, described below, beginning with the easiest:

### Use a Plate Blank

This method can only be used if all samples in the microplate are the same volume and you are not depending on the PathCheck program to correct for variability in volumes.

To use this method:

1. Designate at least one well (preferably several) as Plate Blank (see instructions under Template).
2. Pipette buffer (for example, your sample matrix) into those wells and read along with your samples. SoftMax Pro software automatically subtracts the average of the blank wells from each of the samples. The OD of the microplate material is subtracted as part of the blank.
3. Make sure that Apply Plate Blank is checked under Reduction.

### **Use Plate Background Constant**

If your sample volumes are not identical or if you choose not to use a Plate Blank, then you can use the Plate Background Constants.

To determine Plate Background Constants:

1. Fill a clean microplate with water.
2. Read at the wavelengths you will be reading your samples.
3. The average OD value is the Plate Background Constant. Enter it in the designated field under **Plate > Settings > PathCheck**.
4. If you intend to read your samples at more than one wavelength, there should be a corresponding number of Background Constants.

It is important that you put water in the wells and not read a dry plate for the Back-ground Constant. Dry plates have a slightly higher OD value than a water-filled plate because of differences in refractive indices. Using a dry plate results in PathCheck normalized values that are lower than 1-cm cuvette values. Omitting the Background Constant results in values that have been normalized by the PathCheck program and are higher than 1-cm cuvette values.

### **Use Pre-Read**

This method is less desirable and less convenient than the first two. When you fill the microplate with water and pre-read the plate, SoftMax Pro stores the pre-read values. Then you empty the microplate and add your samples and read again. SoftMax Pro software subtracts the pre-read values on a well-by-well basis (A1 from A1, B1 from B1).

MDS Analytical Technologies enabled the Pre-read feature in earlier versions of SoftMax Pro software because the early UV-transparent

microplates had high and variable background OD values. Use of Pre-read dramatically improved the quality of the data at that time. Now, however, commercially available microplates are of such high quality (low and uniform background) that pre-reading is not necessary. In fact, it is less desirable because if there is a speck of dirt or a dust particle in any well, the erroneous value is subtracted from the sample OD in the subsequent read.

### **PathCheck Program and Interfering Substances**

Any material that absorbs in the 900 nm – 1000 nm spectral region could interfere with PathCheck program measurements. Fortunately, there are few materials that do interfere at the concentrations typically used.

Turbidity is the most common interference: if you can detect any turbidity in your sample, you should not use the PathCheck program. Turbidity elevates the 900 nm measurement more than the 1000 nm measurement and causes an erroneously low estimate of pathlength. Using Cuvette Reference does not reliably correct for turbidity.

Samples that are highly colored in the upper visible spectrum may have absorbance extending into the near infrared (NIR) and can interfere with the PathCheck program. Examples include Lowry assays, molybdate-based assays and samples containing hemoglobins or porphyrins. In general, if the sample is distinctly red or purple, you should check for interference before using the PathCheck program.

To determine possible color interference, do the following:

1. Measure the OD at 900 nm and 1000 nm (both measured with air reference).
2. Subtract the 900 nm value from the 1000 nm value.
3. Do the same for pure water.

If the delta OD for the sample differs significantly from the delta OD for water, then it is advisable not to use the PathCheck feature. Use of Cuvette Reference (on SpectraMax Plus, Plus384, M2, M2e, M5, M5e readers) does not correct for the interference with the current calculation scheme in SoftMax Pro software. Currently, Cuvette Reference involves a single (automated) read at 900 nm and 1000 nm and the automated calculations in SoftMax Pro software do not compensate for color or solvent interference. However, you could correct for such interference by taking two cuvette measurements and using a different set of calculations. For further information, contact MDS Analytical Technologies Technical Support.



Organic solvents could interfere with the PathCheck program if they have absorbance in the region of the NIR water peak. Solvents such as ethanol and methanol do not absorb in the NIR region, so they do not interfere, except for causing a decrease in the water absorbance to the extent of their presence in the solution. Their passive interference can be avoided by using the Cuvette Reference. If, however, the solvent absorbs between 900 and 1000 nm, the interference would be similar to the interference of highly colored samples described above. If you are considering adding an organic solvent other than ethanol or methanol, you are advised to run a Spectrum scan between 900 nm and 1000 nm to determine if the solvent would interfere with the PathCheck program.

## **Automix**

Automix is a feature that allows you to set automated shaking of the microplate. The options available for Automix depend on the read type chosen and the type of instrument.

### **Before First Read**

This setting can be used with any read type. Whether you read at a single wavelength or at multiple wavelengths, checking Before First Read shakes the plate for the set amount of time before the first wavelength reading only.

### **Between Reads**

This option is available only when Kinetic read is selected.

For readings at a single wavelength, Between Reads shakes the plate for the specified amount of time prior to each reading at that wavelength.

For readings at multiple wavelengths, enabling Between Reads shakes the plate for the specified amount of time before each reading at the first wavelength only.

### **Blanking (Pre-Read Plate)**

The blanking specified in the Plate > Settings dialog box is known as pre-read plate blanking, but pre-reading the plate is only one of a number of types of blanking available in SoftMax Pro software. Pre-read plate blanking adds considerable time to a read, and is generally not required for modern plates that have high uniformity.

If you choose pre-read plate blanking, SoftMax Pro software prompts you to read the blank plate after you click the Read button.

The pre-read values are stored in the data file and they are subtracted from the individual well values obtained during sample reading. Pre-read blanking can be disabled in the Reduction dialog box so the data can be reviewed with and without pre-read blanking subtracted.

The purpose of pre-read plate blanks is to correct for well-to-well variability over a whole plate. Pre-read blanking is available for all read types but is most useful for Endpoint readings. When used in Spectrum scan reads (SpectraMax instruments only), the pre-read scan is subtracted on a point-by-point basis from the normal scan.

If pre-read is selected using Kinetic read (not generally recommended), a pre-read with the same Kinetic read interval and duration as the normal reading is done and the OD/RFU/ RLU at each time point of the pre-read is subtracted from the OD/RFU/RLU of the corresponding time point of the normal reading.

Background constant subtraction is available for Endpoint reads as part of the PathCheck program correction. If you pre-read a plate, make sure that background constant subtraction is disabled.

## **Autocalibrate**

This checkbox allows you to disable or enable automatic pre-read calibration for all instruments.

### **AutoCalibrate for Absorbance Reads**

For Absorbance reads, an air calibration value is used to provide the I0 value for the  $\text{Log}(I0/I)$  calculation.

Air calibration values for all wavelengths offered by the instrument are stored in nonvolatile RAM and are loaded into volatile RAM when the instrument is turned on. Every time you perform a plate or cuvette read (unless you have turned off AutoCalibration), new air calibration measurements are made (at the wavelengths used for the read) and the volatile RAM values are updated for those wavelengths. If you are doing a Kinetic read, an Auto-Calibrate value is taken before each read.

If you read a plate or cuvette with AutoCalibration turned off, the current air calibration values in volatile RAM are used without updating. Turning off AutoCalibrate allows the instrument to complete readings more quickly and to begin the first reading of a Kinetic plate faster. If Speed Read is turned on, AutoCalibrate is automatically turned off.

MDS Analytical Technologies recommends that you periodically update the air calibration information stored in the instrument's NVRAM. This is particularly important if you are using Speed Read, reading cuvettes without a reference, or running Kinetics in microplates with read intervals that are too short to allow AutoCalibration.



**Note:** NVRAM air calibration update is not supported by VMax® and EMax® readers.

To update the instrument air calibration values, use the **Control > Calibration** command:

- When you select the Calibrate Plate checkbox to update the plate air calibration values, the instrument steps through each wavelength available to the instrument one by one, taking a reading through the air calibration hole on the plate holder drawer. These values are stored in the non-volatile RAM until the next calibration is done.
- When you select Calibrate Cuvette, the instrument steps through each wavelength one by one taking air calibration values for the cuvette. It is therefore important that the cuvette port is empty.

## AutoCalibrate for Fluorescence and Luminescence Reads

In Fluorescence and Luminescence modes, there are no air calibration values. Instead, reference reads are made on the reference chip on the carriage before each plate read. These reference reads are used to set the RFU/RLU scale (at each wavelength to be used in the read) by adjusting internal parameters including PMT gain. When you turn off AutoCalibrate, the instrument uses the setting from the previous read (or from the default settings if there has been no read since the instrument was turned on).

Turning off AutoCalibrate allows the instrument to complete readings more quickly and to begin the first reading of a Kinetic plate faster. However, you should allow the instrument to perform AutoCalibration of at least one plate prior to disabling this function.

## Well Scan Editor

Well Scan reads allow readings to be taken at more than one location within a well. The Well Scan Editor allows the user to specify how a Well Scan is performed.

You can specify the scanning pattern, and the scanning density. Only odd numbers of points are included in density settings.

Four scanning patterns are available: a horizontal line, a vertical line, a cross (which is a combination of a horizontal and vertical line), and a grid pattern. The density setting determines:

- The number of points read in the line patterns.
- The maximum number of horizontal and vertical points included in the grid pattern.

With a grid pattern, all points that land within 2.8 mm of an edge are excluded.

The minimum number of points is 3 for most plate sizes, while the maximum number of points depends on the well diameter; for 384-well plates, only one read per well is allowed.

The SoftMax Pro software export format supports Well Scans. The data from each position is out-put in subsequent rows or plates starting with the top, left point and proceeding with horizontal priority.

## **Assay Plate Type**

This setting determines the display of wells in the microplate—it should be set to match the type and number of wells in the microplate to be read.

Changing the Plate Type within a Plate section causes well assignments stored in the previous template to be discarded. The groups created previously remain, however, so you can select new wells and apply existing groups to these wells.

## **Plate Types**

The Plate Type list includes entries for microplates with standard dimensions and additional entries for plates that work best with custom settings. For both 96- and 384-well plates, the Standard plate should be used unless the microplate you are using is listed.

For Luminescence reads, white, opaque microplates yield optimum results for most applications. The light reflection with white microplates provides a high signal-to-noise ratio with very low crosstalk from adjacent plate positions. Clear-bottom microplates also give low crosstalk values, although the signal efficiency is reduced in comparison to white microplates. Transparent sample plates cannot be used with the LMax II

series instruments due to extremely high crosstalk. The reading efficiency and the signal-to-noise ratio are lower in black plates because of the decreased light reflection, so black plates are not usually recommended for luminescence.

## **Edit Plate Type**

**(Gemini, SpectraMax® M2, M2e, M5, M5e, FlexStation® Readers Only)**

You can customize the plate type with **Plate > Edit Plate Type**.

The Plate Type Editor includes five fields for entering, in millimeters:

- The distance from the left edge of the plate to the center of well A1
- The distance from the top edge of the plate to the center of well A1
- The horizontal distance between wells
- The vertical distance between wells
- The well diameter

The distance from the left edge of the plate to the center of well A1 must also include an offset to accommodate the angle of the optics. Appropriate measurements should be taken in the lab to determine the correct settings for these fields.

Plate information is saved to the file Custom Plates in the SoftMax Pro folder. Because this is specific to the computer on which it is saved, when a file with a custom plate is opened on another computer, the plate type is changed to the standard plate type.

## **Strips/Wells to Read**

You can choose to read a subset of wells or columns of wells in the microplate. Partial-plate reading is available for all read types, although it is most useful for fast Kinetic analyses. The time required for Kinetic readings can be significantly reduced using this setting since the instrument does not have to read the entire plate.

To enable a partial-plate read, highlight the columns to be read. Columns do not need to start with "1" but must be contiguous.

If you enable a partial-plate reading, only the wells to be read are visible in the data display, indicating that no data will be collected for the other wells (all wells are present in the Template Editor, however).

## Compound Source

(Flex Only)

Select the plate used to store fluid compounds that are withdrawn and then injected into plate wells during the run.

The type of plate used should match the actual plate used in the Compound Source drawer, and it should have the same number of wells as the plate selected in the Assay Plate Type.

## Compound Transfer

(Flex Only)

Use the settings in this section to specify the volumes and rates of compound transfers.

### Assay Plate Fluid

The entry in the Initial Volume field should be equal to the largest initial volume for any well in the plate. SoftMax Pro software assumes all wells hold the same initial volume. This setting is used to compute the total volume in each well after all fluids have been dispensed, to check for potential fluid overflow.

### Transfers

You may enable up to three compound transfers during a reading (or choose none).

When transfers are enabled, numbered transfer buttons are displayed to enable you to set the parameters for that transfer.

### Pipette Height

Determines the amount of fluid (in microliters, measured from the bottom of the microplate well) above which the tip of the pipette remains. You can set up to 130  $\mu\text{L}$  for a 384-well plate, and up to 300  $\mu\text{L}$  for a 96-well plate.



---

**Note:** Subsequent transfers require that you calculate the amount of fluid added.

---

**Rate—from 1 to 8**

Determines the rate at which the fluid is pipetted into the well. A setting of 1 is equal to 26 microliters per second and each subsequent number increases in increments of 26 microliters. A setting of 3 to 4 may help to minimize cell damage and splashing.

**Volume**

Determines the volume of material that is pipetted from the source to each well chosen to receive that transfer.

You can set up to 30  $\mu\text{L}$  for a 384-well plate, and up to 200  $\mu\text{L}$  for a 96-well plate.

**Time Point**

Determines the time after the start of the run when the fluid is scheduled to be dispensed.

**Minimum Time**

This is the minimum time required before a pipetting event can occur. The Time Point setting above cannot be set smaller than this value.

The minimum time value is cumulative: it is not an interval between pipetting events. The value is the minimum number of seconds of elapsed time from the beginning of the read, taking into account:

- The mechanical speed of the pipetter head
- The time needed to aspirate and dispense fluids
- Trituration
- Automixing.

The minimum time for the second pipetting event depends on when the first pipetting event occurred. The calculation for the second event starts at the end of the first event and adds to that the total time necessary to perform actions prior to aspirating new fluids from the compound plate and dispense them into the assay plate.

Compound Transfer settings work in conjunction with the Compound & Tip Columns setting described below.

## **Triturate**

(Flex Only)

The available Triturate choices are derived from the number of compound transfers specified. If no transfers are enabled, no triturate settings are enabled.

Trituration mixes the contents of the wells in either or both the compound source and assay plates by withdrawing and re-pipetting from the wells. You can set the volume of fluid to be withdrawn from the well and the number of times (cycles) it is pipetted back into that well.

For assay plates you must enter a value for the height at which the transfer occurs. The height must take into account the amount of fluid currently in the well in order for the pipette to engage the fluid and allow the desired mixing. The tip must be in the fluid in order to triturate.

## **Pipette Tips Layout**

(Flex Only)

Use the Pipette Tips Layout to specify where the pipette tips are physically loaded in the FlexStation instrument.

MDS Analytical Technologies recommends that you use a full rack of tips each time you perform a transfer of fluid for Flex reads, since this prevents potentially serious errors from occurring when partial tip racks are used. For example, if you should mistakenly enable a pipetting function from a tip that is not present, your samples will not receive the intended compound. If you use a partial rack of tips, ensure that the layout described in Pipette Tips Layout matches the tips inserted in the tip drawer.

If you change the Assay Plate Type setting after making selections in the Pipette Tips Layout tab, for example from a 96-well plate to a 384-well plate, the selection of pipette tips is reset to all tips. The Assay Plate setting takes precedence over all other fluidics module settings.



## Compound & Tips Columns

(Flex Only)

When one or more compound transfers are enabled, use these settings to choose the tips and compounds that are used for transfers.

This tab displays:

- One, two, or three transfers matching the setting made in the Compound Transfer tab.
- The Tips Column list contains tip numbers matching the selected columns in the Pipette Tips tab.
- The Compound Column list contains numbers from 1 through the number of columns of the compound plate or the number of troughs selected in the Compound Source tab.

A Tips (T) or Compound (C) row has up to 12 columns to represent the number of columns in a 96-well assay plate. The actual number of columns and their placement further depends on the number of columns and their locations as selected in the Wells to Read tab.

To assign a tip to an assay plate well:

1. Highlight one or more tip cells.
2. Select one of the items available in the Tips list. Choosing Fill fills the selected cells with the tip numbers that correspond to the same column numbers (tip 3 with column 3).



---

**Note:** Fill does not take into account the actual tips that are available if less than a full rack is used; the choices in the Fill list represent the actual tip numbers that are specified as being available in the Pipette Tips Layout tab.

---

Similarly, to assign a compound:

1. Highlight one or more compound cells.
2. Select one of the items from the Compound list.

The wells shown in this tab setting convey graphically the volumes of liquid in the assay plate for each pipetting event. Using the color associated with each event, the compound settings display the dispensed compound volume as a percentage of the total assay well volume. An initial volume entered on the Compound

Transfer tab is shown as a gray fill. As the liquids are cumulative, the first event's volume is shown above the initial liquid volume (if any), the second event's volume is shown above the first, and the third event's volume above the second.

To deselect a tip or compound assignment:

1. Select the appropriate cells.
2. Select Backspace on a Windows PC or Delete on a Mac.

To change an assignment:

1. Select the wells.
2. Choose new values or type a value on the keyboard.



---

**Note:** If you have multiple wells selected and you type a value, that value is shown in the first selected well and subsequent wells increment to the next higher value. Again, this does NOT take into account the actual pipette tips shown to be available in the Pipette Tips Layout selection.

---

SoftMax Pro software automatically enters information in the Compound and Tip Columns tab using a best-fit algorithm that uses the number of transfers, the Wells to Read section, and the Pipette Tips Layout selected area.

The algorithm assumes:

- All columns in the Wells to Read selection should receive fluid.
- The fluids are dispensed from left to right.
- The fluid transfer targets are cumulative from transfer to transfer (that is, the second transfer's targets start with the next available clean tip and untargeted compound column rather than reusing tips and compound columns targeted by the preceding fluid transfer).
- Each fluid transfer uses a new tip.

When the number of columns in the Wells to Read area is the same or greater than the number of columns of both Tips and Columns, the algorithm fills the Tips and Columns targets as before (that is, each assay plate column is targeted with a new tip column and new compound column until all tips and columns have been targeted).

SoftMax Pro software does not fill beyond the limitations of the available Tips and Compound columns. Once all available tips and compound columns have been assigned, SoftMax Pro software stops assigning targets.

## **Settling Time**

### **(SpectraMax M5 and M5e Readers Only)**

In Fluorescence Polarization mode, Settling Time specifies a well-to-well read delay. After the read head moves to the next well, the read is delayed by the settling time to allow the fluid in the microplate wells to settle, thereby minimizing sloshing.

For all other read modes, the selected delay is added between the reading of each column (there is no delay between each well in a column).

## **Speed Read**

### **(Absorbance Only; Endpoint, Kinetic and Spectrum Scan Only)**

Speed Read uses fewer flashes per read in order to decrease read time.

Speed Read turns off AutoCalibration and uses the air calibration values stored in the instrument. As a result, Speed Read may not provide as accurate an absorbance measurement of each wavelength in the scan as with the normal read.

## **Column Wavelength Priority**

Choosing column or wavelength priority determines the order in which a plate is read in multi-wavelength readings:

- In column priority, an instrument reads the entire microplate (or chosen number of strips) at the first wavelength, and then goes back and reads the microplate at the second and subsequent wavelengths.
- In wavelength priority, the instrument reads all wavelengths for the first column of wells in the microplate, and then reads all wavelengths for the second column.

For most instruments there is a choice between column or wavelength priority when doing Endpoint or Kinetic readings, for all read modes.

If Column Wavelength priority choices are not offered, the following priorities are used by default:

- Endpoint, Kinetic: column priority.
- Spectrum, Well Scan, Fast Kinetic, Flex: wavelength priority.

## AutoRead

AutoRead enables the automatic reading of all Plate sections in the order in which they appear in an Experiment. The time in the Delay field specifies the delay time between plate reads.

## Template Editor

The Template Editor is used to describe the location of samples in a microplate or cuvette, and thus provides the link between raw data and analysis groups.

Each template is composed of samples and groups:

- A set of one or more replicate wells makes up a sample.
- A set of related samples forms a group.

For example, you may have a group named "Standards" that consists of seven samples named STD0 1, STD02, ..., STD07, and a group named "Unknowns" that consists of five samples names UNK01, UNK02, ..., UNK05.

Each sample would be typically applied to a column or some other subset of wells on a microplate, or to one or more cuvettes.

Each well designated as being part of a group has associated with it a group name, a sample name (or replicate ID), a sample descriptor (optional), and a column format for the calculations and data reported in the associated Group section.

Samples and groups may exist across multiple Plate sections as well as within a single Plate section.

Deleting the wells assigned to a group from the template does not delete the group. To delete a group name in the Template Editor you must delete the corresponding Group section from the Experiment.

## Selecting Wells or Cuvettes in the Template Editor

Open the Template Editor for a Plate section with the **Plate > Template** command, or the Template button (the same applies to CuvetteSet sections):

- Select one or more wells by clicking and dragging on the grid.
- Non-contiguous selections can be made by Shift-clicking individual wells.
- Clicking and dragging row or column labels selects entire rows or columns.
- Non-contiguous row or column selections can be made by ⌘-clicking (Mac) or Ctrl- clicking (Windows) row or column labels.
- Double-clicking anywhere within a group selects the entire group.
- Clicking in the empty border around the Template deselects all wells.

## Group

Use the groups in the Group list to assign groups to well selections on the plate grid. The list contains two default entries—Blank and New—and any user-defined groups:

- Blank—Creates a plate blank in the active section. The name blank is reserved for use by SoftMax Pro software and should not be used when creating custom groups.
- New—Opens the Group Settings dialog box in which you can create a new group.

## Group Settings

The Group Settings dialog box allows you to define the name for a group of related samples, a descriptor associated with the samples, and the initial column format for the data calculated and reported in the associated Group section.

Whenever a new group is created (whether or not wells are selected in the template), a Group section is created in your Experiment. To delete a group, you must delete the Group section. Clearing a group from the Template Editor removes only the assignment of wells to that group name; it does not delete the group.

### **Name**

Use either the default name, "Group#X," where "X" is an integer that is initially set to "1" and is incremented (for example, "1," "2," and "3,") as other groups are created, or enter a custom name (up to 32 characters in length).

### **Sample Descriptor**

Check Sample Descriptor to apply a descriptor and its units to the group. Select a set of units from the list or type them directly in the Units field.

Sample Descriptor is automatically selected when either the Standards or Unknowns (Dilution) column format (see below) is selected. The Sample Descriptor field can also be used for time or fraction number information.

No sample descriptor is assigned automatically for the Basic or Unknowns column format. You can assign a sample descriptor to these column formats manually but the information does not appear automatically in the Group section. To see it, create a new column containing the formula !Sampledescriptor, !factor or !concentration. For more information about creating columns, [Adding Columns on page 100](#).

## Column Format

The Column Format setting defines the default columns for the data calculated and reported in the associated Group section of the current sample group.

Each of the four column formats—Basic, Standard, Unknowns, and Unknowns (Dilution)—creates a different set of columns in a new or existing group. The table below shows the default columns created with each selection type.

**Table 4-1** Columns Created and Their Formulas

Name	Formula	Basic	Standard	Un-known	Un-known (Dilution)
Sample	!SampleNames	✓	✓	✓	✓
Wells	!WellIDs	✓	✓	✓	✓
Sample#	Index	✓			
Concentration	!Concentration		✓		
Mean Value	Average(Values)	✓	✓		
Values	!WellValues	✓	✓	✓	✓
R (Outside Standard Range)	If(Values)>='MinStd@Standards' and Values <='MaxStd@Standards', "", "R"			✓	✓
Result	InterpX(STD@StandardCurve, Values)			✓	✓
MeanResult	Average(Result)			✓	✓
Std.Dev.	Stdev(Result)		✓	✓	✓
CV%	Cv(Result)		✓	✓	✓
Dilution	!Factor				✓
Adj.Result	MeanResult*!Factor				✓

### **Edit**

The Edit button opens the Group Settings dialog box for a selected group in the Group list.

The name, sample descriptors, and column format can be changed for the group that you are editing.

Changing the column format causes the columns in the group to revert to their defaults. If you have added columns or changed column information in the section for the group you are editing, these changes are removed.

### **Sample**

The Sample area contains the Sample list and sample descriptor values.

#### **Sample List**

The Sample list reports the sample name assigned to the currently selected wells.

When samples are assigned to a group, the default name is the first 3 letters of the group name followed by a number. For a group called "Standard," for example, the default sample names would be Sta01, Sta02, Sta03, and so on.

SoftMax Pro software automatically increments sample names for a group as you successively assign samples to columns, for example.

However, you can type a sample name directly into the Sample field.

#### **Sample Descriptor Field**

Use the Sample Descriptor field to enter sample information, such as the sample dilution or concentration.

The Sample Descriptor field is displayed only if the selected group has associated descriptor information.

### **Assign**

Click this button to apply a sample name to selected wells.

### **Clear**

Clears the group assignment of the selected wells.



## Series

This button is used to define several samples as a series, allowing you to easily enter incremental sample descriptors (for example, dilutions or concentrations) and sample names to the template as long as the increment can be expressed as a mathematical operation.

### Replicate Samples and Series

You can assign the same group or sample definition (same sample name and description within a group) to multiple wells to create replicates. For example, you might want to read standards in replicate to ensure that anomalies can be excluded prior to generating a standard curve. Replicates can be created using the Assign button or with the Series dialog box.

A series is defined in one direction (starting from left, right, top, or bottom) and therefore samples must be arranged sequentially in either ascending or descending order.

To create a series, do the following:

1. In the Template Editor, highlight a rectangular selection of wells to be part of the series.
2. Click the Series button.
3. Specify the name of the first sample in the series in the box. Subsequent replicates use this name as a base and either increment the number within the name or append a number to it.
4. The Sample Descriptor information is derived from the information entered in the Group Settings dialog box. Specify the starting value for the first sample in the series.
5. Choose the operator for the series (+, -, \*, or /) from the Step by list and enter a value for the increment. For example, with a starting value of 0.5, the division operator (/) and the increment 2, the series will start with 0.5 and divide each subsequent sample by 2 to produce the series 0.5, 0.25, and 0.125.
6. Choose a fill direction to describe how the series fills the wells, and a number of replicates. The maximum number of replicates depends on the selection of wells: if you are filling down from the top, for example, and you have chosen a block of wells that is 4 wells wide by 8 wells high, the maximum number of replicates will be four; with the same block of wells selected, filling from the left

would allow eight replicates. If the number of replicates you choose does not divide evenly into the number of rows or columns you select, depending on the filling direction, the remaining wells that cannot contain replicates will be labeled as additional individual wells in the series. To create a series with all 96 samples, highlight all the wells in the plate, select the fill direction, and set the replicates to 1.

7. Click OK to save the series and close the dialog box.

To increment the names of a series of samples without applying a concentration or dilution:

1. Create a series in the Series dialog box.
2. Uncheck the Sample Descriptor field.

To increment the names of a series of samples with a constant concentration or dilution:

1. Create a series in the Series dialog box.
2. Check the Sample Descriptor field.
3. Enter the concentration or dilution, and select the multiply operator with a factor of 1.

If sample names are set to increment automatically using the Series function, be aware that SoftMax Pro software will automatically truncate the sample name to three or four characters, including the incrementing number if the sample name starts with letters. If the sample name is entirely numeric, it is not truncated. If sample names are very long, and they must be kept long for reference, you will have to set up your replicates manually.

Hold down Ctrl+ Shift keys to display the values for all wells.

## Blanking

In addition to the pre-read plate blanking configured in **Plate > Settings**, and a reference reading of a cuvette, in the Template Editor you can configure Plate blanks and group- associated blanks

The different blank types have these features:

- Pre-read plate blank (not in the Template Editor): The pre-read value of each well of a microplate is subtracted from any reading made in the well. It is used to correct for well-to-well variability

across a whole plate; generally not required for current plates that are highly uniform.

- Plate blanks: used to correct readings when all samples on a plate have been prepared in the same buffer or matrix, and therefore can use the same correction value. Plate blanks are subtracted from the raw data (or after pathlength correction if PathCheck is selected).
- Group-associated blanks: used to correct readings when samples on a plate have been prepared in a different buffer or matrix, and therefore have to be individually corrected. Group blanks are subtracted after wavelength and Kinetic reduction.

Blanks can be used in combination and have cumulative effects.

## Plate Blanks

### Plate Sections

Wells can be assigned to a “Blank” group in the Template Editor to create a Plate Blank:

1. Open the Template Editor for the Plate section.
2. Select the wells to be used for the Plate Blank.
3. Select Blank from the Group list.

Unlike pre-read plate blanks, in a plate blank group the average value of all the wells in the Blank group is subtracted from each individual well value on the microplate. If you have an instrument that supports the PathCheck software, the Plate Blank is subtracted after pathlength correction.

Use Plate Blanks when all samples on the plate have been prepared in the same way, and therefore can be corrected from a single blank reading.

Plate Blank subtraction can be turned off in the Reduction dialog box so that the data can be reviewed with or without plate blank subtraction.

For Kinetic and Spectrum data, the blank value is averaged and subtracted from each point in the read.

## CuvetteSet Sections

To perform a similar function to Plate Blanking for CuvetteSet sections, you can assign cuvettes to a Blank group in the Template Editor of the CuvetteSet section and create a Template Blank:

1. Open the Template Editor for the CuvetteSet section.
2. Select the cuvettes to be used for the Template Blank.
3. Select Blank from the Group list.

The average value of all the cuvettes in a Template Blank group is subtracted from individual cuvette readings in the same CuvetteSet section.

For Kinetic and Spectrum data, the blank value is averaged and subtracted from each point in the read.

Template blanks are subtracted after the cuvette reference reading is subtracted.

## Group-Associated Blanks (Plates and Cuvettes)

Blank wells can be assigned within any group (other than the "Blank" group) in the Template Editor:

1. Open the Template Editor for the plate.
2. Select the wells or cuvettes to be used for the group-associated blank. These should be wells or cuvettes that are already assigned to a Group (not the "Blank" group).
3. Select BL from the Sample list.

The average reduced value of the blank wells or cuvettes within that group is subtracted from individual reduced values within that group only.

Use Group-Associated Blanks when samples on the plate have been prepared differently, and therefore require individual blank correction.

This blanking function uses the reduced number (example: !Lm1-!Lm2, VMax reader) only—it does not subtract on a point-by-point basis. When a group blank has been assigned, it is automatically subtracted and you do not have the option of reviewing the uncorrected data. To see the reduced number without the group blank subtracted, you must mask the group blank wells or cuvettes (see [Masking Wells or Cuvettes on page 88](#)).

Group blanks apply to all wells or cuvettes in a group and thus can be subtracted from wells or cuvettes in more than one Plate or CuvetteSet section.

## Copying and Pasting Templates

To copy a template:

1. Activate the source template by clicking in the Plate section containing it.
2. Choose the **Plate > Copy Template** command.
3. Activate the destination Plate section by clicking it.
4. Choose the **Plate > Paste Template** command.

### Copying and Pasting within the Same Experiment Section

When copying and pasting within the same Experiment section, the group names and sample names of the pasted template are identical to those of the source template. In effect, the destination template is an extension of the original template and all wells on the destination plate are considered replicates of the wells on the source plate (same group and sample names).

Changes made to the template on the destination plate are also made on the source plate. Data from the two Plate sections are analyzed together.

### Copying and Pasting between Different Experiment Sections

When copying and pasting between different Experiment sections, the group names and sample names of the pasted template are identical to those of the source template but, since they are in different Experiments, the full name is different (for example: group@experiment#2 instead of group@experiment#1).

Changes made to the template of the destination Plate section do not affect the source section.

Data from the two Plate sections are not analyzed together.

### Copying and Pasting Template Regions

Contiguous regions of templates can be copied and pasted to other regions of templates within the same or to a different Experiment. All template settings are pasted into the new template except the group

name: the new group is given the name "Group#X", where X is the next greater number than that used already for group names.

The real value in this method is the ability to copy a complicated template and then to be able to analyze the data separately from the original Plate.

To copy and paste a region of a template:

1. Open the Template Editor of the source Plate and select the wells to copy.
2. Select ⌘C or Ctrl+C to copy the layout.
3. Open the Template Editor for a new Plate section and select all wells.
4. Select ⌘V or Ctrl+V to paste the layout.

Only the highlighted wells have template information pasted into them. Data is pasted using row priority: all wells in a row of highlighted wells have information pasted into them before proceeding to the next row, from left to right.

If fewer wells are highlighted for pasting than were copied, the data is pasted until all paste-highlighted wells are filled.

If more wells are highlighted for pasting than were highlighted for copying, the extra wells are left empty.

## Exporting and Importing Templates

Templates can be exported or imported by opening a Plate section and **Plate > Import Template** or **Plate > Export Template**.

Template files can be prepared outside SoftMax Pro software and then imported as needed when ready to run samples.

The easiest way to determine the correct format for importing a template is to export a template from SoftMax Pro software.

## Reading a Microplate or Cuvette

### Data Collection from a Microplate

You can start a reading at any time after defining instrument settings. It is not necessary to define groups and assign wells within the Template Editor first. The values received from the instrument are raw absorbance,

fluorescence, luminescence, or fluorescence polarization values and are not affected by settings in the Template Editor.

To read a microplate:

1. Open the drawer of the instrument by clicking the Drawer button in the Status Bar, or selecting the **Control > Open Drawer** command.
2. Insert the prepared microplate.
3. Click the Read button in the Status Bar or choose the **Control > Read** command.
4. The active Plate section is read. If no Plate section is active, a dialog box is displayed in which you can choose a Plate section to read.
5. If you have chosen to pre-read the plate for blanking, you are offered the choice of a Pre-read or a Normal read (for example, no pre-read). After reading the blank plate, you must start the read process again for the normal plate.
6. When a reading begins, the Read button changes to Stop, allowing you to terminate a reading if desired.

## Data Collection from a Cuvette

Two types of data collection are possible using a cuvette: a reference and a sample reading.

You can start a reading at any time after defining instrument settings. It is not necessary to define groups and assign cuvettes within the Template Editor first, but you can use the Template Editor to create the appropriate number of cuvettes in the section. Alternatively, it is easy to create a small number of cuvettes one at a time with the **Cuvette > New Cuvette** command.

Instrument settings must be the same for all cuvettes in a CuvetteSet section.

### Sample Reading

You can highlight any cuvette in the CuvetteSet section and choose to read it. The Read button changes to Stop, allowing you to terminate a reading if desired.

If you select more than one cuvette, the software starts with the left most-selected cuvette that does not contain sample read data.

CuvetteSet sections have a limit of 96 cuvettes. If this limit is reached, create a new CuvetteSet section.

### Reference Reading

As in a spectrophotometer, taking a reference reading can be done either on air or using a cuvette containing the buffer of your sample.

The reference may be read before or after reading samples.

If no reference has been taken in a CuvetteSet section before samples were read and you then read a reference, the following occurs:

- The reference reading is applied to all cuvettes in the CuvetteSet section, unless you have highlighted specific cuvettes (see below).
- The time and date of the reading are displayed under each cuvette in the section (three cuvettes per row and one cuvette per row displays only).
- The reference is applied to any new cuvettes subsequently created in that section.

You can have more than one reference per CuvetteSet section. If you have cuvettes to which you want to apply a different reference:

1. Highlight the cuvettes in the CuvetteSet section.
2. Place the appropriate reference in the cuvette port.
3. Click the Ref button or select the **Control > Ref** command.
4. The time and date stamp of the reference for the selected cuvettes changes.

Any new cuvettes created in the CuvetteSet section after reading a second reference use the new reference reading.

The reference read is automatically applied to the normal read. The reference value is not displayed in SoftMax Pro software. If you wish to see the reference value, you can choose one of the following procedures:

#### Procedure 1:

1. Create a cuvette in the CuvetteSet section and Ref on air (empty cuvette port).
2. Read a cuvette containing the blanking solution.



If you want to have both reference and read values available, do the following:

1. Create a CuvetteSet with up to the maximum of 96 cuvettes.
2. Define a template blank in the CuvetteSet section.
3. Place buffer in the cuvette and select Read.
4. Click the Read button to read the remaining samples.

The optical density of the cuvette designated as the template blank is subtracted from the optical density of all the other cuvettes in the CuvetteSet section.

### Procedure 2:

1. Create a CuvetteSet with up to the maximum of 96 cuvettes.
2. Reference and read the cuvette.
3. Create a template with the appropriate groups.
4. Create a column in each Group table with the formula !wellprereadLm1.

If you have a multiple-wavelength reading and want to see the reference at each wave-length, create additional columns with the formula !wellprereadLmx where "x" is 2, 3, or 4, corresponding to the wavelength for which you want to see a reference.

The OD/RFU/RLU with the reference subtracted is displayed in the Values column. The reference value for each sample is displayed in the custom column you have created.

## Calibration

For some instrument models, **Control > Calibration** performs an air reading at all available wavelengths and stores those values in the instrument.

When you select the Calibrate Plate checkbox to update the plate air calibration values, the instrument steps through each wavelength one by one, taking a reading through the air calibration hole at the edge of the plate holder drawer. Those values are stored in the non-volatile RAM until the next calibration is done.

When you select Calibrate Cuvette, the instrument steps through each wavelength one by one, taking air calibration values for the cuvette. It is therefore important that the cuvette port be empty.

These air calibration values are used for Absorbance reads only, in the following circumstances:

- When the **Instrument Settings > Auto Calibrate** option is disabled.
- When a Spectrum reading is performed using the Speed Read option.

MDS Analytical Technologies recommends that you periodically update the air reference information that is stored in the instrument. This is particularly important if you are using Speed Read, reading cuvettes without a reference, or running Kinetics in microplates with read intervals that are too short to allow AutoCalibration.

## Data Display During a Reading

The values read by the instrument appear in real time in the data display as they are received from the instrument. Depending upon the read mode and type, the number of wavelengths, the plate type, and the display options selected, the data display shows one or more numbers or plots.

During a Kinetic or Spectrum reading you can enlarge the display of one or more wells. To enlarge a single well to a Well Graph, double-click the well.

To display the data from multiple wells in one enlarged Well Graph:

1. Hold the Shift key while selecting more than one well.
2. Choose the **Plate > Graph** command or click the Graph button on the tool bar.

The data in the enlarged Well Graph update to show new data points as they are received from the instrument.

Other choices and settings within SoftMax Pro software affect the way in which data is displayed—the type of reduction, specific time settings, and so on. See [Chapter 5](#) for more information about data analysis.

Any SoftMax® Pro software data file has a minimum of one Plate section or one CuvetteSet section. Most files also have an assigned plate template and associated Group, Graph and Notes sections. However, depending upon the protocol and how you want to report the data, only some of these sections may be present in every file.

Some simple protocols, such as reading the absorbance of proteins at 280 nm in a cuvette (or in a microplate with PathCheck® software), allow you to calculate concentration in the Plate or CuvetteSet sections. In such an example, Group and Graph sections are necessary only if you wish to display the data in tables or graphically.

Many common protocols consist of data acquired in a Plate section or CuvetteSet section. Standards and unknowns are defined in a template. Unknowns are then interpolated from a standard curve.

## Data Display

### Display Dialog Box

At any time in an active Plate section you can click the Display button or select Plate > Display to change how the data is presented.

Choices available in this dialog box depend on whether a section is active, which instrument is chosen in the Preferences or is connected to the computer, and which read type you have chosen (for example, the choice for Plots is not shown if Endpoint is selected).

### Raw

Selecting Raw displays the default data type for the selected read type:

- Endpoint: Raw absorbance, fluorescence or luminescence values.
- Kinetic: The change in raw OD/RFU/RLU values over time, displayed as a plot.
- Spectrum: Raw OD/RFU/RLU values for the range of wavelengths, displayed as a plot.
- Well Scan: Raw OD/RFU/RLU values as shades of gray.
- Flex: Raw RFU/RLU values over time, displayed as a plot.

## **Reduced**

The reduced number is a combination of plate blank subtraction in the Plate section, wavelength reduction, and if applicable, a Spectrum, Well Scan, Kinetic or Flex reduction.

The reduced number is reported in the Group section when a template has been defined.

### **Number**

To view by reduced number alone, select Number.

To see a reduced number as well as another type of display, check With reduced number.

### **Threshold**

Threshold shows the raw data as a plus (+) for values above the high limit, an asterisk (\*) for values within the limits, and a minus (-) for values below the low limit.

### **Ranged**

When Ranged is chosen, raw data that falls between the high and low limits is scaled on an integer scale from 0 to 9. Values above the high limit are displayed as a plus (+), and values below the low limit are displayed as a minus (-).

### **Grayscale**

Grayscale presents the raw data in eight shades of gray, changing from light (for values less than or equal to the low limit) to dark (for values greater than or equal to the high limit).

### **Plots (All Read Types Except Endpoint)**

The Plots option presents Kinetic or Spectrum data as plots of OD/RFU/RLU versus time or wavelength.

## **Show**

### **With Reduced Number**

When you select this option you can see both the raw and reduced numbers displayed.

### Raw Data in 1000's

Use this option to display Luminescence readings, for example, as multiples of 1000 to make the numbers in the wells easier to read (e.g. "20,000" would be shown as "20").

### 384 Well Display

Data in the Plate section is displayed in a grid array (microplate format). If you are using a 384 well plate, you have the following additional options for displaying the data:

- Normal: Shows the data in standard left-to-right format, with well A1 in the upper left corner. Columns 1 through 24 are along the top and rows A through P are shown from top to bottom. The wells are smaller and data is shown with fewer digits (less precision).
- Vertical: The plate is shown in the same orientation as Normal except that columns 1 through 12 (all rows) are shown above columns 13 through 24 (all rows).
- Rotated: Same as the Normal view except that it is rotated 90 degrees clockwise (well A1 is in the upper right corner).
- Large: Same as the Normal view except larger. This display can be seen in its entirety only if you have the page size set for legal paper (8.5 x 14 inches) in Landscape orientation (set this in **File > Print Setup**).
- Interleaved: The wells are shown in a format that skips every other well as follows: all odd columns and rows begin in the upper left corner of the plate display and are followed by all even columns and rows. So well A1 is in the upper left corner, but it is followed horizontally by well A3, A5, ... A23, A2, A4, and vertically by well C1, E1, ... O1, B1, D1. This display is most useful when the 384-well plate is composed of 4 daughter plates of 96 wells each.

### Polarization Display

Polarization Display options are available in the Display dialog box when the read mode is Fluorescence Polarization.

You can choose to display Polarization (mP), Anisotropy (r) or Raw (S&P) values. The G-factor in the Polarization and Anisotropy calculations (default is 1) can be manually adjusted within the range of 0.01 to 99.

## Graphing Wells

You can enlarge the well plots in Kinetic, Spectrum, Well Scan or Flex reads to get a close-up view of the data:

- Double-click a well or cuvette in the data display, or
- Select the wells or cuvettes, and select **Plate > Graph**.

### Kinetic and Spectrum Well Graphs

Well Graphs plot the data for the individual wells in a microplate or for individual cuvettes and show the well ID and reduction information, including the goodness of fit (R2 value), if appropriate. The default display for the X-axis shows time for Kinetic plots and wavelength for Spectrum plots; OD, %T, RFU/RLU, or mP is displayed on the Y-axis for both plots.

### Well Scan Well Graphs

Well Graphs for Well Scan data default to displaying raw data in a grayscale plot.

### Well Graph Options

#### Print

Opens the Print dialog box.

#### Reduction

Opens the Reduction dialog box.

By default, the ranges on the plot are set to the limits defined in the Reduction dialog box, and the information included near the bottom of the window is based on the reduction settings currently in use. You can change these settings, but changes made within well graphs are applied to all wells or cuvettes in the Plate or CuvetteSet section.

When viewing Raw data, you can scale the graph to see data outside the reduction limits (lines are included on the graph indicating the limits and the area outside of them will be shaded slightly). This view should assist you in determining the best limits for reduced data.

#### Show Raw/Show Reduced

Toggles the Well Graph between views of raw and reduced data.

When the Well Graph is opened, the initial view depends on the Display settings for the Plate section.

- ◆ If raw data is being displayed in the Plate, the initial view in the Well Graph is raw data, and the button is set to Show Reduced.
- ◆ If reduced data is being displayed in the Plate section, the initial view in the Well Graph is reduced data, and the button is set to Show Raw.

### **Scale to Data/Scale to Limits**

Toggles between a view of the data within the reduction limits and a view of the entire data set.

Switching between these views helps to visualize how the limit settings affect data reduction:

- ◆ If raw data is being viewed in the Plate section, the display in the Well Graph is set initially to show the plot scaled to the reduction limits.
- ◆ If reduced data is being viewed in the Plate section, the Well Graph display is set initially to show the plot scaled to the data.

If you are viewing raw data and click the Scale to Limits or Scale to Data button, only the display of raw data in the Well Graph is affected (the button will not affect the view of reduced data). The same is true if you click this button while viewing reduced data. To produce comparable scaling of both raw and reduced displays, this button must be set to either Scale to Limits or Scale to Data for both raw and reduced views of the Well Graph.

### **Graph Options**

Opens the Well Graph Options dialog box, which allows you to enable or disable connected points or plotted symbols on the graph.

### **X-Axis Settings/Y-Axis Settings**

Opens the Well Graph Axis Settings dialog box, allowing you to set Auto-range parameters, choose the minimum and maximum values for the selected axis, cause the axis to be auto-ranged, add or remove gridlines, and add or remove axis tick marks.

Only the Well Graph being viewed is affected by changes in these dialog boxes.

## Masking Wells or Cuvettes

Masking is the process of excluding outliers in certain wells or cuvettes from data reduction calculations. To mask wells or cuvettes:

1. Select the wells or cuvettes to be masked.
2. Click the Mask button, or select **Plate > Mask** or **Cuvette > Mask**.

Masking can be used as a “what if?” tool. Suppose you have included a group blank in the template and want to see the data with and without the blank. Masking the group blank wells suppresses the blanking function, while unmasking them enables it again.

## Data Reduction

Because this chapter deals with more than one instrument model, not all of the information found here may apply to your instrument.

The reduction process in SoftMax Pro software is based on formulas that reduce the raw data to show a single number for each well or cuvette. Further analysis of this reduced number is done in Group and Graph sections.

SoftMax Pro software performs calculations hierarchically when reducing raw information collected from the instrument. Calculations in the Plate section or CuvetteSet section are performed in the order shown below. If an option has either not been selected in the Instrument Settings dialog box, is not available for the instrument you are using, or has not been defined in the template or Reduction dialog box, SoftMax Pro software continues with the next listed calculation.

### Plate Section

1. Convert OD to %T.
2. Pre-read plate or plate background constant subtraction (or division for %T).
3. PathCheck pathlength correction if available (different equations for OD and %T).
4. Plate blank subtraction.
5. Wavelength reduction.



6. Kinetic or Spectrum reduction.
7. Group blank subtraction.

### **CuvetteSet Section**

1. Convert OD to %T.
2. Reference is subtracted from the reading (optical density mode) or the read data is divided by the reference (%T mode—but see below).
3. Template blank subtraction.
4. Wavelength reduction.
5. Kinetic or Spectrum reduction.
6. Group blank subtraction.

### **EndPoint Reads**

#### **Wavelength Combination**

The default reductions for Endpoint readings are !Lm1, !Pathlength, and Custom for all instruments when you read at a single wavelength or if you choose 1-Ref. The !Pathlength reduction reports the pathlength in each well. With all SpectraMax®, Gemini and FlexStation® instruments, additional wavelength combinations become available depending on the number of wavelengths chosen.

For any instrument, you can choose Custom and create different reduction formulas; see the *SoftMax Pro Software Formula Reference Guide* for more on custom formulas.

The !Pathlength option is only applicable to Plate sections when using the SpectraMax 190, 340PC, 340PC384, Plus, Plus384, M2, M2e, M5 or M5e. However, the wavelength reduction display is not instrument-dependent in SoftMax Pro software, so this choice is always shown, regardless of the instrument.

#### **PathCheck® Software**

Pathlength correction allows microplate data in Endpoint reads (for those instruments that support PathCheck) to be normalized to a 1-cm pathlength. When PathCheck software is turned on in Plate > Settings for the plate, Apply PathCheck is displayed as an option in the Reduction dialog box.

## Kinetic Reads

### Wavelength Combination

The default wavelength combination for Kinetic readings is !Lm1 and is the only choice (besides Custom) for all instruments except the SpectraMax reader.

For the SpectraMax reader, the choices depend on the number of wavelengths chosen. For example, if you read at two wavelengths, choices default to !Lm1 and !Lm2, along with Custom.

### Kinetic Reduction

Kinetic reductions are applied to the value at each time point after the wavelength combination formula is applied.

VMax per min and VMax per sec reductions are available for all instruments that are capable of Kinetics; the default Kinetic reduction is VMax.

#### VMax

VMax is the maximum slope of the Kinetic display of mOD/min or RFU/RLU per second. VMax is calculated by measuring the slopes of a number of straight lines, where VMax Points (see below) determines the number of contiguous points over which each straight line is defined.

#### VMax Points

This setting defines the maximum size of the line segment used to determine the slope of the line used in calculating the rate of the reaction. The default is the total number of points taken in the reading (reading except on the FlexStation instrument, where the default is 5).

The first slope is calculated for a line drawn beginning at the first reading as defined by Lag Time (described below) and ending at a total number of readings equal to the VMax Points setting. The second and any subsequent slopes are calculated beginning at the second time point and ending at a total number of readings equal to VMax Points. The steepest positive or negative slope is reported as VMax.

If the data plot displays fewer time points (data points) than VMax Points, all of the time points are used to determine the slope of the data.

**Time to VMax**

This is an alternative method for analyzing non-linear Kinetic reactions that reports the elapsed time until the maximum reaction rate is reached, rather than reporting the maximum rate itself. Used in conjunction with VMax Points, Time to VMax is the time to the midpoint of the line defined by VMax Points and used to calculate VMax.

This elapsed time data is useful for applications including coagulation chemistry where the changing concentration of the reagents does not change VMax, but rather changes the time at which the reaction reaches the maximum rate.

**Onset Time**

This is another method for analyzing non-linear Kinetic reactions. Onset Time reports the time required for a Kinetic reaction to reach a specified OD or RFU/RLU (onset OD/RFU/FLU).

This elapsed time data is useful for cascade reactions including clot formation (endotoxin testing, for example) and clot lysis applications where the change in reagent concentration does not affect the maximum optical density change but changes the time required for the reaction to reach completion.

**Time at Minimum**

This setting reports the time at the minimum OD, RFU/RLU, or %T that falls within the reduction limits.

**Time at Maximum**

This setting reports the time at the maximum OD, RFU/RLU, or %T that falls within the reduction limits.

**Time at 1/2 Maximum**

This setting reports the time at the half of the maximum OD, RFU/RLU, or %T that falls within the reduction limits.

To calculate this reduction, SoftMax Pro software determines the Kinetic point (within the reduction limits) that has the maximum signal level (OD or %T) and divides it by 2 to get the 1/2 Maximum value. Then it finds the time value at the 1/2 Maximum.

**Slope**

This setting determines the slope of the combined plot (for example, the slope of the line using linear regression after the

wavelength combination reduction). It is calculated using all visible time points in the reduction window.

Slope is the same as VMax Rate when VMax Rate is set to the same number of points as the run, but is different if you have modified VMax Points.

### **Area Under Curve**

This reduction estimates the area under the curve as defined by the data plots within the reduction limits. The data plots are treated as a series of trapezoids with vertices at successive data points and at the X-axis coordinates of the data points. The areas defined by each of the trapezoids are then computed and summed.

### **Kinetic Limits**

Limits define the data that are viewed and included in data reduction. If you alter a limit to show less data, you can always display the excluded data again by changing the limit.

Unless you have selected the Absolute Values option, the display of OD/RFU/RLU values is relative to the first point measured for each well for example, the first point is always set to (0,0), and all other data are shifted accordingly. The Absolute Values option allows you to see the unshifted data (see below).

Negative Kinetic values decrease with time, and limits should be set accordingly (below 0) to view negative Kinetic data.

### **MaxOD/RFU/RLU**

The limit for the maximum value to report. Any values from the reading that are above this limit are excluded from data reduction. The default for MaxOD is 1.0, while it is 20,000 for RFU/RLU.

The MaxOD/RFU/RLU reduction parameter can be used to exclude the non-linear portion of the reaction from data analysis. This type of data reduction is most useful in reactions where the initial portion of the data is linear. You might also adjust the End Time setting to remove a non-linear portion after a certain time point in the reaction. Using MaxOD/RFU/RLU allows the use of the maximum number of linear points to calculate the slope of the line and, thus, to determine the rate for each well.

**MinOD/RFU/RLU**

The limit for the minimum value to report. Any values from the reading that are under this limit are not shown and are excluded from data reduction. The default is 0 OD or RFU/RLU. To display negative Kinetics, the value should be set below 0 (zero).

**Lag Time**

Specifies how many initial data points are excluded from the calculation of VMax Rate.

Lag Time truncates the data used in the calculation, but does not prevent data from being collected. Kinetic plots do not display the data collected prior to the Lag Time. The default is 0.

**End Time**

Specifies the time at which to stop showing data in the display. Any values occurring after this limit are not reported in the display. The default setting is the total assay time.

**Absolute Values**

This setting displays OD/RFU/RLU values unshifted, as opposed to the default that offsets the first point to (0,0) and shifts all other data accordingly. Enabling Absolute Values can cause some data points that do not fall within the default limits to disappear.

To see more data, yet still display absolute values, increase the limits for MinOD/ RLU/RFU and MaxOD/RLU/RFU.

## Spectrum Reads

### Spectrum Reduction

The Spectrum reduction formula is applied to the list of numbers in each well (values at each wavelength) after the wavelength combination formula is applied. The default reduction for Spectrum readings is Lambda at Maximum.

#### Maximum

Reports the maximum absorbance (optical density) or percent transmittance (%T), RFU or RLU within the reduction limits.

#### Minimum

Reports the minimum absorbance (optical density) or percent transmittance (%T), RFU or RLU within the reduction limits.

#### Lambda at Maximum

Reports the wavelength at which maximum absorbance (optical density) or percent transmittance (%T), RFU or RLU within the reduction limits.

#### Lambda at Minimum

Reports the wavelength of minimum absorbance (optical density) or percent transmittance (%T), RFU or RLU within the reduction limits.

#### Area Under Curve

Estimates the area under the curve as defined by the data plots (within the reduction limits). The data plots are treated as a series of trapezoids with vertices at successive data points and at the X-axis coordinates of the data points. The areas defined by each of the trapezoids are then computed and summed.

### Spectrum Limits

Limits define the data that are viewed and included in data reduction. If you alter a limit to show less data, you can always display the excluded data again by changing the limit.

#### Start (Wavelength) nm

Specifies the limit for the minimum wavelength setting to report. Any values from the reading that are under this limit are not shown and are excluded from data reduction.

### **End (Wavelength) nm**

Specifies the limit for the maximum wavelength setting to report. Any values from the reading that are above this limit are not shown and are excluded from data reduction.

### **Well Scan**

#### **Wavelength Combination**

!Lm1 is the only choice besides Custom.

#### **Well Scan Reduction**

Choices are Maximum, Minimum, Average (default), and Custom. Average provides the average value for all points in the Well Scan.

### **Flex Reads**

#### **Smoothing**

Smoothing is used to reduce noise in Flex readings. This setting takes the moving average of the raw data and displays the Flex plot using the averaged values.

The default is no smoothing (setting of 1 "averaged" point). A setting of "3" provides the least amount of smoothing: the value of the center point is added to the values of the points on either side; the sum of these points is then divided by three and the resulting value is substituted for the center point. A setting of "9" averages the center point with the four points on each side.

When multiple wavelengths are read, smoothing is performed before performing wavelength combination calculations such as ratios or multiplication. For example, if you read at two wavelengths and choose to take a ratio of those wavelengths, smoothing is performed on each individual wavelength and the wavelength ratio is calculated using smoothed data.

#### **Baseline Options**

Baseline options affect the way that the data is graphed.

##### **Absolute**

Shows the raw data without any change in the way it is graphed. Raw data is used in subsequent reduction calculations.

### **Zero Baseline**

Offsets the graph slightly.

If you enter "5" for the number of points, for example, the average of the first 5 points is used as the baseline value and all raw data are subtracted from this value.

Baseline subtraction is applied only to the reduced data—raw data plots do not change. To view the plot with baseline subtraction, choose Reduced Plot in the Display dialog box or select Show Reduced in a zoomed well Graph window.

If two or more wavelengths are present in a reading and you choose to combine these wavelengths while Zero Baseline is selected, the zero baseline applies only after the wavelength combination.

### **% Baseline with Multiplier**

This option makes Flex data look more like FLIPR® data. FLIPR data can be imported into SoftMax Pro software and this option modifies the plot and reduction calculations to make Flex data comparable to the FLIPR data.

This method averages the first number of points you specify (the upper box), divides the raw data by this average value, and then multiplies by 100 (%). The Multiplier field allows you to enter a number from 0.000 1 to 9999 that is used to multiply the % Baseline value.

This calculation is performed only on the reduced data and only after any wavelength combinations are calculated.

To see the modified plot, choose Reduced Plot in the Display dialog box or click Show Reduced in a zoomed well Graph window.

### **Interpolate Raw Data**

For Flex reads each data value is logged with its own read time, which can make it difficult to perform wavelength combination calculations correctly. For example, to take a ratio of two different wavelengths in a single well, it is necessary to take a ratio ( $Lm1/Lm2$ ) of values that have been read at two different points in time.



To solve this problem, select the Interpolate Raw Data checkbox in the Reduction dialog box. This checkbox is shown only when two or more wavelengths are present in the reading.

Enabling this checkbox causes SoftMax Pro software to use quadratic interpolation to calculate the interpolated values. This results in the loss of one point at each end of the plot. Prior to beginning the calculation, SoftMax Pro software determines the starting point by averaging the initial data points of the separate wavelengths. For example, if the first data points of the two wavelengths have time values of 0.3 and 0.5 seconds, SoftMax Pro software averages these and uses 0.4 seconds as its initial point for interpolation.

### **Data Mode (%Transmittance/Absorbance)**

For Absorbance reads only, you can choose whether to display absorbance data as %Transmittance or Absorbance (OD).

Separate mathematical calculations are used for handling OD and %T calculations for Pre-read plate blanking, PathCheck, and Reference, because OD calculations are performed on a linear scale, whereas %T calculations are performed on a logarithmic scale. However, SoftMax Pro does not perform other calculations differently for absorbance and %T modes. Because of this, MDS Analytical Technologies recommends that %T be used only for raw OD, raw OD with Pre-read plate blank subtraction, or raw OD readings which are normalized by the PathCheck program. Use caution when using %T on reduced numbers or any readings that apply other calculations since the data may not be calculated correctly.

### **Custom Reduction Formulas**

If the predefined reduction formulas do not meet your needs, you can create one or more different reduction formulas for any read type (Endpoint, Kinetic, Spectrum, or Well Scan). Choosing Custom from any of the menus or dialog boxes displays a Formula button with which you can open a Formula dialog box.

Examples of some formulas that can be used when combining multiple wavelengths in a Formula dialog box are given in the table below.

**Table 5-1 Wavelength Reduction Formula Examples.**

<b>2</b>	<b>3</b>	<b>4 to 6</b>
$!Lm1 + !Lm2!$	$Lm1 + !Lm2 + !Lm3!$	$Lm1 + !Lm2 + \dots + !Lmn$
$!Lm1 - !Lm2$	$(!Lm1 - !Lm3) / (!Lm2 - !Lm3)$	$(!Lm1 - !Lm6) / (!Lm2 - !Lm5) / (!Lm3 - !Lm4)$
$!Lm1 / !Lm2$	$!Lm1 / !Lm3$	$!Lm1 / !Lmn$
$!Lm1 * !Lm2$	$!Lm1 * !Lm3$	$!Lm1 * !Lmn$
$\text{Log}10(!Lm1/!Lm2)$	$\text{Log}10(!Lm1/!Lm3)$	$\text{Log}10(!Lm1/!Lmn)$
$!Pathlength$	$!Pathlength$	$!Pathlength$
$!Lmx/\text{constant}$	For example, $Lm1/1.44$ for quantitating a polyclonal antibody by measuring the absorbance at A280 with PathCheck on.	
$\text{Average}(!allvalues)$	Averages together the optical densities for multiple readings at the same wavelength (for example, if you read the well six times at 280 nm).	
$\text{Min}(!allvalues)$	Reports the minimum OD/RFU/RLU recorded for multiple wavelength readings in each well.	
$\text{Max}(!allvalues)$	Reports the maximum OD/RFU/RLU recorded for multiple wavelength readings in each well.	
$\text{If}(!Lmz < A, \text{makeerr}(118), (\text{if}(!Lmx > B, \text{makeerr}(117), !Lm1)))$	Reports "low" for any well with an OD/RFU/RLU less than A, "high" for any OD/ RFU/RLU greater than B, and the OD/ RFU/RLU of any well that lies between A and B.	

For EMax®, VMax®, UVmax, or ThermoMax instruments, the choice available for reading at two wavelengths is designated "1-Ref." While the instrument performs a reading at the two wavelengths you set (one for the measurement and one for the reference), data from these readings is automatically reduced to a single value in the display, and you cannot view the data from the individual wavelengths.

Custom reduction formulas utilizing mathematical operators or terms can be used to obtain specific types of data. The table below provides some examples of such formulas for Kinetic and Spectrum readings.

**Table 5-2** Kinetic and Spectrum Reduction Formula Examples.

<b>Kinetic</b>	
Vmaxcorr(!combinedplot, !Vmaxpoints,!readinterval)	Reports VMax correlation coefficient for plots in all 96 wells.
Vmax(Delta(!combinedplot), !Vmaxpoints,!readinterval)	Reports the VMax Rate of the delta between each time point.
<b>Spectrum or Kinetic</b>	
Nthitem(!Lm1,X)	Reports the optical density at item X in the list of readings. For example, if you have a Kinetic run with 20 time points and X is 10, it reports the OD/RFU/RLU of the 10th time point. Similarly, if you have a Spectrum scan with 20 measurements and X is 10, it reports the OD/RFU/RLU of the 10th wavelength measured.

For a full discussion of custom formulas, see the *SoftMax Pro Software Formula Reference Guide*.

When you finish your entry in the Calculation dialog box and click OK, the formula is displayed and becomes the default selection for the Custom option.

## Recalculation Options

By default, SoftMax Pro software performs continuous recalculation of the data when you read plates, create or change formulas, and so on. At certain times, it is useful to disable recalculations so that you can edit the Experiment without waiting for recalculations to be completed.

**Edit > Suspend Recalculation** disables automatic recalculation. While suspended, no recalculation occurs regardless of what you may change, add, or delete from the Experiment. This is useful when creating or changing column formulas within Group sections. Select **Edit > Suspend Recalculation** again to re-enable automatic recalculation.

When recalculation has been suspended, you may want to see the results of the changes you have made but may still not wish to enable continuous recalculation. Choose **Edit > Recalculate Now** to recalculate all data once without enabling continuous recalculation.

## Group and Graph Sections

### Group Sections

Unlike other section types, Group sections are created automatically when you create and assign a Group to a Plate section within the Template dialog box (after clicking the Template button in a Plate section).

Group sections can also be created by duplicating existing Group sections within the Experiment file itself, which is often an easier way to replicate Groups if many custom columns or Summary objects have been added.

Reduced numbers from the Plate section are displayed in the Values column of the Group section tables by default.

Custom reduction formulas require the use of accessors and operators that are understood by SoftMax Pro software. For a list of these accessors and operators see the *SoftMax Pro Software Formula Reference Guide*.

### Adding Columns

Click the New Column button or use **Group > New Column** to add a column to the active Group table:

- If no columns are highlighted, the column is added at the end of the group.
- If you select a column first, the new column is added to the right of the highlighted column.

New columns can contain references to other columns, either in the current section or in a different one. For example, if you wanted to subtract the mean values in one group from those in another, you could create a new column (in either group) to do this. If the two groups were named "Group 1" and "Group 2," for example, and both groups contained a column entitled "Mean," you could create a column within "Group 1" that would subliterate the mean values in "Group 2" from those in "Group 1." The column formula needed to do this would be:

Mean-Mean@Group2

Many possibilities exist for creating and editing column formulas: see the *SoftMax Pro Software Formula Reference Guide* for a complete discussion.

## Copy and Paste Column

You can copy a Group column and paste it within another Group in the same Experiment or into a Group of another Experiment:

1. Highlight the column to be pasted.
2. Select **Edit > Copy** (⌘C; Ctrl+C).
3. Activate the destination section. If you highlight a column in the destination section, the column you are pasting is placed to the right of the highlighted column; otherwise, it is appended to the right of all existing columns.
4. Select **Edit > Paste** (⌘V; Ctrl+V).

Columns and Summaries in Group sections can also be copied and pasted to other programs.

## Showing Existing Column Formulas

To show Group formulas in a Group table, select the **Group > Show Formulas** command.

To see formulas briefly, hold down the Shift+Ctrl keys.

## Modifying Column Formulas

When a column is selected, you can use **Group > Edit Column** (or double-click the column) to edit the formula associated with the column, and the column name.

## Adding or Editing a Summary in a Group Section

Click the Add Summary button or use **Group > New Summary** to create a Summary (or double-click an existing Summary) and enter a formula for this Summary for the Group.

Information in the New Summary dialog box is placed at the bottom of the Group section as follows:

Name = Description Formula Result

You can check the Hide Name option to display the Summary as:

Description Formula Result

The name assigned to a Summary is important since you can reference the Summary by name in other formulas (building one upon another).

Move Summaries horizontally by dragging them left or right. If you do this while holding down the Ctrl key, the Summary snaps in line with a column.

### **Copy and Paste Formulas and Summaries**

Formulas are used in Group columns, custom reductions, and Summaries. Sometimes it is useful to copy an existing formula, column, or Summary from one area of the program to another, or from one Experiment to another, rather than recreating it. Summaries can also be copied and pasted within SoftMax Pro software. After pasting, both formulas and Summaries can be edited.

Contiguous column formulas may be copied and pasted together. Multiple Summaries may also be copied and pasted together if the Shift key is held down while selecting the Summaries.

To copy a formula or Summary:

1. Highlight the formula or Summary.
2. Select **Edit > Copy** (⌘C; CTRL+C).
3. Activate the destination; if necessary with a formula, access the dialog box into which it will be pasted.
4. Select **Edit > Paste** (⌘V; CTRL+V).

### **Autosize**

Use Group > Autosize (or double-click the column dividers) to set the width of selected columns to accommodate the longest text in the column.

### **Hide**

Group > Hide hides selected columns from display.

You can show hidden columns with Group > Show All.

### **Hide Replicates**

Group > Hide Replicates hides data for all replicates.

You can show hidden data with Group > Show All.

### **Show All**

**Group > Show All** shows all hidden columns or replicates.

## Graph Section

### The Plot Dialog Box

The Plot# dialog box opens when you create a new Graph section with Experiment > New Graph, or when you add a plot to an existing graph from the Graph > Graph Options dialog box. In this dialog box, you can name the plot, assign specific information to be plotted on the X- and Y-axes, and choose which icon and color will be used.

All groups that have been created or assigned within the file are listed on the left side in the Source field. To create a new plot:

1. Select a group in the Source field.
2. Select group columns to assign to the plot in the X-Axis and Y-Axis lists.

Summaries and user-defined formulas can also be graphed:

- ◆ Click the Formula button for either the X-axis or Y-axis to open a Formula dialog box.
- ◆ To plot a Summary, enter the Summary name in this dialog box.
- ◆ Formulas can be plotted for comparison with other data. For example, if the data is assumed to follow a specific model, the model formula can be entered as a plot and then the data can be chosen for another plot within the same graph.
- ◆ You can use this feature to plot Kinetic or Spectrum data in Graph sections. For example, to plot Kinetic data from well A1 in a Graph section, you can enter the formula !A1Lm1 for the Y-Axis and the !TimeRun accessor for the X-axis.
- ◆ The entered formulas are shown in the Source field of this dialog box.

The size, color, and style of the icon used for a plot can be selected on the right side of the dialog box. Four sizes, nine colors, and nine styles are available from drop-down menus.

### The Graph Options Dialog Box

The Graph > Graph Options dialog box is displayed when you are editing or creating a graph. It allows you to:

- Change the name and size of the Graph section.

- Select the type of graph to be displayed (scatter, cluster bar, stack bar).
- Change the fonts used on the title, legend, and axes.
- Access the Plot# dialog box for creating, editing and deleting plots.

Each of the plots that is created for the Graph section is listed in the Plots field. Choosing any of these plots provides an indication of the data to be used for the X- and Y- axis data points and error bars on the right side of this field. Plots can be deleted by highlighting the plot name and clicking the Delete button.

Clicking the New button opens the Plot# dialog box, allowing you to create a new plot for the graph. Existing plots can be edited by highlighting the plot name and clicking the Edit button which opens the Plot# dialog box containing all the information for the selected plot.

### **Creating Multiple Plots**

You can plot data from any group or from several groups on a graph, even if the groups are in separate Experiments. It is possible, for example, to plot multiple dilution series of an unknown, to graph together a series of patient samples, or to plot several controls that have been run over time.

This example shows how three groups, all of which are Standards, can be plotted on the same graph.

1. Groups are created in the Template Editor. For this example, four groups are created: three for standards (Standard 1, Standard2, and Standard3) and one for Unknowns (Unknown 1).
2. The associated Group sections are created using the Standards column format that, by default, generates columns for concentration and OD/RFU/RLU values.
3. Create a new Graph section with Experiment > New Graph. Create a new plot named "Standard#1 ". Click OK to open the Graph Options dialog box.
4. Select Standard1 for the group to be plotted, the concentration of Standard1 for the X- axis, and values of Standard 1 for the Y-axis.
5. Add a second plot to this graph by clicking New in the Graph Options dialog box to open the Plot#2 dialog box. Change the second plot name to "Standard#2," choose Standard2 for the



group, and change the X- and Y-axes to show concentration and values, respectively, for Standard2.

6. Repeat the process for a third plot, labeled "Standard#3."
7. Now that all plot information is complete, click OK in the Graph Options dialog box to close it, and the resulting graph shows the three plots created.

### **Error Bars**

When data is displayed as a scatter or bar graph, you can choose to display error bars for the plot of the data for the X- and Y-axis. Error bars are lines that extend beyond a plotted value in either or both directions and graphically represent some amount of error in plotted data.

### **Curve Fitting**

When you first create a graph of the data, it does not have a fit associated with it. You can fit any plot to one of ten curve-fitting functions: Linear, Semi-Log, Log-Log, Quadratic, 4-Parameter (logistic), 5-Parameter (logistic), Log-Logit, Point-to-Point, Exponential, and Cubic Spline. These selections are shown in the Fit list in the Graph section tool bar. All plots on a graph must have the same type of fit.

Typically, a standard curve refers to the curve fitted to the plot of concentration versus mean value for the Standard group.

Once you have selected a fit type, SoftMax Pro software determines the parameter values that best fit the data. The function with these parameters is drawn on the graph. Ideally, the type of fit used should be determined by the underlying chemistry of the assay and could be set before data is read.

When a fit is performed, the parameter values are displayed in the legend at the bottom of the graph. The correlation coefficient describes how well a change in x values correlates with a change in the y values. The R<sup>2</sup> value should only be used for linear curve fits.

A good discussion of curve fitting appears in "Data Analysis and Quality Control of Assays: A Practical Primer" by R. P. Channing Rogers in Practical Immuno Assay, edited by Wilfrid R. Butt (published by Marcel Dekker, Inc., New York, 1984).

#### **Linear**

The linear function fits the best straight line to the data. The equation for this fit has the form of:

$$y = A + Bx$$

where A is the y-intercept of the line and B is the slope. A linear fit should be used whenever the values appear to lie on or scattered around a straight line.

### **Semi-Log**

The semi-log function fits the best straight line to a set of data for  $\log(x)$  plotted against y. The resulting curve displayed will be a straight line with the X-axis drawn in logarithmic scale. The equation for a semi-log fit is:

$$y = A + B \log(x)$$

where A is the y-intercept of the line and B is the slope (log in this equation is the common or base-10 logarithm).

### **Log-Log**

The log-log function fits the best straight line to the set of data which consists of the logarithm of the readings on the Y-axis (the response) and the logarithm of the dose on the X-axis. The resulting display is a straight line with both axes drawn in logarithmic scale. The equation for the log-log fit is:

$$\log_{10}(y) = A + B \log_{10}(x)$$

where A is the  $\log_{10}$  Y-intercept of the line for  $\log_{10}(x) = 0$ , and B is the slope (log in this equation is the common or base-10 logarithm).

### **Quadratic**

The quadratic function fits the best parabola to the data. The parabola is a curved line based on the equation:

$$y = A + Bx + Cx^2$$

which is shown on the screen as:

$$y = A + B*x + C*x^2$$

where A is the intercept, B is the slope of the curve at the intercept, and C is the measure of the curvature of the parabola.

The quadratic fit is most appropriate when the standard curve has a tendency to curve up or down.

### 4-Parameter Logistic, 5-Parameter Logistic and Log-Logit

If the standard curve has a sigmoidal shape when plotted on the semi-log axes, it may be appropriate to use either the log-logit, 4-parameter, or 5-parameter fit. Both logistic and log-logit fits are based on the equation:

$$y = \frac{A - D}{1 + (x/C)^B} + D$$

where D is the Y-value corresponding to the asymptote (for example, the flat part of the curve) at high values of the X-axis, and A is the Y-value corresponding to the asymptote at low values of the X-axis. The coefficient C is the X-value corresponding to the midpoint between A and D, commonly called the IC50 or EC50. The coefficient B describes how rapidly the curve makes its transition from the asymptotes in the center of the curve, and is commonly known as the slope factor. A large value of B describes a sharper transition. Typically, B has a magnitude of about 1 for inhibitory experiments and -1 for excitatory experiments.

Both the log-logit and 4-parameter equations are displayed as:

$$y = (A - D) / (1 + (x / C) ^ B) + D$$

The difference between the log-logit fit and the 4-parameter fit are in the way the coefficients A and D are calculated. For the log-logit method, the standard values for the lowest and highest values of x are used. The corresponding y values are assigned to A and D, respectively. Based on these fixed values for A and D, the algorithm then computes values for B and C. This technique works well if there are standard points along the upper and lower asymptotes. If this is not true, the log-logit fit should be avoided in favor of the 4-parameter logistic.

The 5-parameter logistic equation has an extra exponential term G that is sometimes described as an asymmetry factor:

$$y = (A - D) / ((1 + (x / C) ^ B) ^ G) + D$$

As with the 4-parameter logistic equation, D is the high asymptote, A is the low asymptote and B is the slope factor. Unlike the 4-parameter logistic equation, however, C is not the IC50 value. The IC50 value needs to be calculated separately. This is described in Chapter 3 of the Formula Reference Guide.

Use the 5-parameter logistic equation instead of the 4-parameter logistic equation when you need to fit to a non-symmetrical sigmoidal shape. The non-symmetry is typically seen in the rate of change of the slope of the curve as it approaches the asymptote. With 4-parameter fits, the rate of change of the slope is the same at the lower asymptote as at the upper asymptote. With 5-parameter fits, the upper and lower portions of the curve can have very different shapes.

The minimization method for the 4-parameter and 5-parameter equations is based on the Levenberg-Marquardt Method.

Discussion of these methods can be found in Numerical Recipes in C: The Art of Scientific Computing by William H. Press, Brian P. Flannery, Saul A Teukolski, and William T. Vetterling, published by Cambridge University Press, New York, 1988.

### **Point-to-Point**

A linear equation is fit to each pair of data points. Fit parameters are not given. The point-to-point curve fit is a linear fit composed of pieces that assume a linear relationship between each pair of data points. The line segment defined by each pair is used to interpolate data between those points. Since there are multiple line segments, fit parameters are not shown on the graph.

### **Cubic Spline**

This curve fit generates a fit to a cubic equation between each pair of data points. The general form of a cubic equation is:

$$y = A + Bx + Cx^2 + Dx^3$$

The equations are computed with the requirement that the first and second derivatives of the equations are continuous throughout the range of the data. As multiple curve fits are performed with this routine, fit parameters are not shown with the graph.

## Exponential

The exponential function used to generate this curve fit is:

$$y = A + B(1 - \exp^{-x / C})$$

Previous versions of SoftMax Pro software performed a linearization of the exponential curve fit. Starting with the version 5.3 release, SoftMax Pro software uses a non-linear fitting method whenever the exponential curve fit is selected. For previously saved files that use the exponential function, when opened in SoftMax Pro software version 5.3 and above, the linear fitting will still be retained. You may freely switch between the two fitting methods by toggling between them in the *Fit Setting* dialog box window, under the *Fit Methods* heading.

## Minimum Number of Standards

The number of standards used to generate a standard curve depends on the particular curve fit selected and on the judgment of the operator. The table below gives the minimum number of standards required mathematically by each of the curve fitting algorithms.

**Table 5-3** Minimum Number of Standards Required for Curve Fits.

Curve	Minimum Number of Standards
Linear	2
Quadratic	3
Semi-Log	2
Log-Log	2
Log-Logit	4
4-Parameter	4
5-Parameter	5
Point-to-Point	2
Exponential	3
Cubic Spline	4

For optimal results, you should always exceed the minimum number of standards required for any given fit.

There are other requirements on the standards for the log-logit and 4-parameter and 5-parameter logistic curves. The fits for the logistic curves are based on the assumption that the curve has a high and a low

asymptote, and a certain steepness for the linear portion around the inflection point between the asymptotes. If a sufficient number of points do not define the inflection point and the asymptotes, you will receive a "no fit" error message or an inappropriate fit. For best results, make sure that the logistic fit is appropriate for your assay and that the entire range of the assay is represented in your standards.

If your standards do not clearly define both a high and low asymptote, the spline fit may be more appropriate for your assay.

### Fit Settings

#### 4- or 5-Parameter Curve Fit Weighting

Curve fit weighting is useful when uncertainties in the data are not random, and may be related to the time, the concentration or the number of data points, for example. In such cases, you can up-weight or down-weight points on the curve using a weighted sum of squared errors:

$$wSSE = \sum w_i |y_i - f(x_i)|^2$$

for weights  $w_i$ , observed values  $y_i$  and function values  $f(x_i)$ .

By default SoftMax Pro software does not apply weighting (referred to as "Fixed Weight" in the software), but you may choose a variable weight for each point and enter it as a function in the Parameter Settings dialog box. The formula must be an array of numbers that has the same number of points as the curve. For example, you could use a formula `Weight@Group#1` to use a column of weights from the Group section `Group#1`.

In SoftMax Pro software version 5.4, the weighting feature has been expanded to allow independent weighting of individual plots within a Graph section. Prior to this software release, the same weighting scheme was applied to all plots. You may select the specific plot for the weights to be applied to by using the drop-down menu in the *Fit Settings* dialog box.

#### Manually Set Parameters

SoftMax Pro software offers the option of manually assigning individual fit parameters for 4 and 5-parameter curves.

Parameters can be fixed to a constant value (for example, a number), or they can reference by formula other information in the SoftMax Pro file (for example, a curve fit parameter from another Graph section).

This option can be used independently of weighting and parallel line analysis selection.

### **Parallel Line Analysis**

This feature, accessible through the Parameter Settings dialog box, is only available for the 4 and 5-parameter curve fits.

Determining parallelism between a standard and test compound, and subsequent estimation of relative potency, are fundamentals of bioassay use in biopharmaceutical development and testing. For linear data, parallelism is evaluated by examining the similarity between the slopes of straight lines.

Bioassay data, and immunoassays in particular, has been demonstrated to fit well with a sigmoidal shape defined by a 4- or 5-parameter curve fit. These S-shaped curves are not parallel within the classic definition of never intersecting. However, the basic question of whether a curve with a specified shape can fit both standard and test data sets when shifted along the X-axis can still be addressed, following a method first detailed by Gottschalk and Dunn (Journal of Biopharmaceutical Statistics, 15:437- 463, 2005).

SoftMax Pro software begins by fitting the standard and test samples to non-linear curves, minimizing the weighted sum-of-squares error. The possible fits evaluated by the software algorithms are constrained by assuming they are parallel functions related by a relative potency term.

The Graph section where the parallel line analysis feature has been enabled then displays the curve fit parameters, relative potency (relative to the selected standard curve), the degrees of freedom from both the independent and constrained curve fits, and the probabilities obtained from the F-test and Chi-squared test for parallelism (if applicable).

New function accessors have been added to SoftMax Pro software to quickly summarize this information in a Notes section. Please see the *SoftMax Pro Software Formula Reference Guide* for additional details.

To use parallel line analysis with a standard curve:

1. You need to have a Graph section containing at least two plots, one a Standard and one a Test sample.
2. Select 4-Parameter or 5-Parameter from the Fit list on the Graph section tool bar.
3. Open the Parameter Settings dialog box.
4. Check the Use Parallel Line Analysis check box and select your standard curve from the Standard Curve menu.
5. Click OK.

### **Confidence Intervals**

For the 4 and 5-parameter curve fits with and without the parallel line analysis feature enabled, confidence intervals for the curve fit parameters can be calculated. Due to the computing power that may be required, these options are not enabled by default.

In the *Fit Settings* dialog box, you can select for which parameters (A-G and relative potency) confidence intervals are desired. The confidence level should be specified as a percentage. There are three methods by which the confidence intervals can be calculated: asymptotic approximation, profile with a F distribution, and profile with a Chi-squared distribution.

To report these values, specific formula accessors must be used as described in the Formula Reference Guide. The general format to report the lower and upper confidence intervals for a parameter are as follows:

ParmXCILower(Plot#1@Graph#1)

ParmXCIOpper(Plot#1@Graph#1)

Where X stands for any of the curve fit parameters, and Plot#1@Graph#1 refers to the specific plot and graph.

### **Fit Method**

As previously described for the exponential curve fit, you may toggle between the "Linearized" and "Nonlinear" fit methods in the *Fit Settings* dialog box.



### **Axis Settings**

Selecting Auto-range draws the graph so that all data points are displayed by automatically adjusting the minimum and maximum range settings to encompass the data.

### **Judging a Good Curve Fit**

The correlation coefficient is also known as the goodness of fit value because it shows how good a fit exists between the calculated curve and the data points. It measures the degree to which the points fall on the calculated curve and may give an indication as to what residuals are to be expected. The correlation coefficient is popularly known as R and is between  $-1$  and  $1$ , inclusive. When the data points lie on a perfectly straight line with negative slope, then  $R = -1$ . If the correlation coefficient is not suitably close to  $1$  or  $-1$  (a perfect fit), you can simply apply one of the other types of curves.

For non-linear curve fits, SoftMax Pro software performs transformations that cause the curve to become more linear. The correlation coefficients are derived from that transformed data. Because of this, you should be cautious about using correlation coefficients as an indicator of goodness of fit. Do not let this single value for goodness of fit sway your intuition about which fit is best. Some curve fits may seemingly give a good fit (e.g., fit values close to  $1$ ) but, when inspected by eye, show a poor fit along a major portion of the plot.



The hands-on tutorials included here are designed to take you through the basic features of SoftMax® Pro software:

- Tutorial 1 leads you through the steps required to create and run a quantitative protocol using a standard curve. It is designed to be used by both Absorbance and Fluorescence instruments.
- Tutorial 2 leads you through the steps required to create and run an EC50 protocol using a graph of concentration vs. RFU on a FlexStation® instrument using a Flex read.

The source files for these tutorials are installed with SoftMax® Pro software, and can be found in Program Files / Molecular Devices / SoftMax Pro v5.4 / Tutorials.

If you are new to SoftMax Pro software, we suggest that you read and follow the steps in one of the tutorials before using SoftMax Pro software for the first time, since it will help you learn the basic features quickly.

More information about any of the operations performed in this tutorial can be found in Chapters 4 and 5.

## Hard and Soft Parameters

The only settings that must be defined before reading a microplate are the instrument settings in **Plate > Settings**. These settings tell the instrument the conditions that should be in place in order to collect data. They include the read type (Endpoint, Kinetic, Spectrum, Well Scan or Flex), the number of wavelengths, specific wavelength settings, and so on. These settings, called hard parameters, define the conditions under which the data is obtained. Once a microplate has been read, the instrument settings cannot be changed without deleting the data.

Soft parameters include the template definition and data reduction settings. These parameters can be defined prior to reading a microplate, but they are applied only after the plate is read and can be changed at any time.

In these tutorials, you will build a complete protocol (based on the settings found in the Tutorial file), which will include both hard and soft parameters. You will save the completed protocol file prior to collecting data (this file could then be used to generate future data files).

Before starting, ensure that the instrument that is connected to the computer is chosen in the Edit > Preferences dialog box:

1. Choose your instrument from the Reader list (if it is not already chosen).
2. Click OK.



---

**Note:** Some settings mentioned in the tutorials are not available for some instruments.

---

## Tutorial 1: Quantitative EndPoint Protocol

Tutorial 1 is designed for both Absorbance and Fluorescence instruments. It assumes that you have read the previous chapters in the manual and have a basic understanding of the logic and operation of SoftMax Pro software.

The following steps are required to create and run a quantitative protocol using a standard curve:

1. Create a new data file.
2. Define instrument settings.
3. Define template.
4. Set reduction parameters.
5. Set display parameters.
6. Save the protocol.
7. Read the plate (start the simulator).
8. Data analysis—Group sections.
9. Data analysis—Standard curve.
10. Print a report.

### Step 1: Create a New Data Base

1. Double-click the SoftMax Pro icon to start the program.
2. Choose the **File > Open** command, and open the file "Tutorial1.pda".

If you are using an Absorbance instrument, the CuvetteSet#1 section can be deleted, as the use of a cuvette is not covered in this tutorial. To delete this section:

1. Click the title bar of the CuvetteSet#1 section.
2. Select **Edit > Delete** CuvetteSet#1.

The Experiment should now have a Notes section (Introduction), Plate section (Plate#1), a Graph section (Graph#1), and two other group sections (Controls and Unknowns) that will be used in this tutorial.

## Step 2: Define Instrument Settings

1. Click the Plate#1 section to make it active and then click the Settings button in the tool bar, or select **Plate > Settings**.
2. At the top of the dialog box, Endpoint is selected by default as the read type; no change is required.
3. Clicking the setting name (bold heading) on the left side of the Instrument Settings dialog box shows the choices available for that setting in the right pane, allowing you to make changes.
4. Following are the details about the quantitative protocol settings we will use to create a protocol file using SoftMax Pro software. Select Fluorescence as the read type, and make sure all other settings match those in the table:

**Table 6-1** Example Settings

Read Type	Endpoint
Automix	Off
AutoCalibrate	On
Instrument Settings Specific to Absorbance Instruments:	
Wavelengths	One: 405 nm
Strips	Read entire plate
Instrument Settings Specific to Fluorescence Instruments:	
Wavelengths	One: Excitation: 485; Emission: 538; Auto Cutoff
Sensitivity	6 readings; Automatic PMT
Assay Plate Type	96 Well Standard

**Table 6-1** Example Settings

Read Type	Endpoint
Wells to Read	Read entire plate
Autoread	Off
Standards in $\mu\text{g/ml}$	0, 0.1, 0.25, 0.5, 0.75, 1.0, and 1.4 (each standard run in duplicate)
Standard	Curve Fit: Linear
Blank	Plate blank run in duplicate

The following steps guide you through creating a protocol file for a quantitative protocol. Reading an actual plate is not required as SoftMax Pro software can generate simulated data.

### Step 3: Define the Template

The Template Editor is used to create a map of the contents of the microplate.

1. Click the Template button in the Plate Section tool bar. This opens the Template Editor, showing all wells empty.
2. Click and drag the mouse to select the first two wells, A1 and A2.
3. Click the Group list on the Template Editor to see group selections. Choose Blank. The two selected wells in the template are now defined as the plate blank.
4. Select the two wells B1 and B2 and select Standards from the Group list.
5. The selected wells B1 and B2 now show the default first sample name for the Standards group, St01. The area in the top center of the Template Editor is now active, showing the default values for concentration and units.
6. In this tutorial, the concentration for the first Standard should be set to 0  $\mu\text{g/ml}$ . If a value other than 0 is shown in the Concentration field, highlight the existing value and enter 0. Click the Assign button or press either the Tab or Return key on the keyboard to assign this concentration to the selected wells. You can check the concentrations you have assigned to the template by pressing and holding the Ctrl + Shift keys.

7. Select wells C1 and C2. Standards is still showing as the default group name, but a new sample descriptor is created automatically (St02), so all you need to do is assign a concentration value of 0.1 µg/ml to these wells. After entering the value, click the Assign button.
8. A total of seven standards will be assigned in duplicate in this template. Select wells D1/D2 through G1/G2 and click the Series button.
9. In the Series dialog box, leave the default values for the sample descriptor (Std03), starting from the top, and the concentration in µg/ml. Enter 0.25 for the starting value, choose the Step by operator + and enter 0.25 for the increment of the series.
10. Click OK.
11. These choices assign concentrations of 0.25, 0.5, 0.75, and 1.0 µg/ml, respectively, to the selected wells, which you can check by holding down Ctrl + Shift.
12. Select wells H 1 and H2. A new sample descriptor is created automatically (St07), so all you need to do is assign a concentration value of 1.4 µg/ml to these wells. After entering the value, click Assign (or press the Tab or Return key).
13. 13 This tutorial example uses seven separate unknown samples, with each unknown being placed in four wells. The unknown samples should be entered on the template as follows:
  - Un01: wells A5 through D5
  - Un2:wells A6 through D6
  - Un3:wells E6 through H6
  - Un4:wells A7 through D7
  - Un5:wells E7 through H7
  - Un6:wells A8 through D8
  - Un7:wells E8 through H8
14. Start by highlighting wells A5 through D5. Select Unknowns from the Group menu.
15. The selected wells will now show the default first sample name for the Samples group, Un01.

16. Continue highlighting the wells as described above, assigning new unknown sample descriptors to each subsequent group.
17. This tutorial example uses two controls, with each control being placed in three wells.

The controls should be entered on the template as follows:  
Co0 1: wells A9 through C9  
Co02: wells D9 through F9
18. Start by highlighting wells A9 through C9. Select Controls from the Group menu. The selected wells now show the default first sample name for the Controls group, Co01.
19. Do the same for wells D9 through F9, assigning the default second sample name, Co02 to these wells.
20. The last items to be added to the template are samples with dilution. To do this, you need to create a new group. Choose the wells of the template that will contain these samples: A11 through G12.
21. Click the Group menu and choose New. In the Name field enter "Sam w/dil" and choose Unknowns (dilution) for the Column Format. Choose  $\mu\text{g/ml}$  in the Sample Descriptor Units field.
22. Click OK.
23. The samples with dilution need to be changed to be part of a series, however, since we really want to include six different samples with replicates. Highlight wells A11 through G12 again and click the Series button.
24. Leave the default direction from the top down and leave the default operator of divided by 1. Enter 10 for the starting value.
25. The template is now complete.
26. Click OK to accept the entries and close the Template Editor. The outlines of the groups you have defined in the Template Editor are now shown in the Plate section.
27. An associated Group section is created each time a sample group is created in the template, so we now have Group sections named Standards, Unknowns, Controls and Sam w/dil.



---

## Step 4: Set Reduction Parameters

1. Click the Reduction button in the Plate section tool bar to open the Reduction dialog box.
2. The default reduction setting for a single measurement wavelength, !Lm1, is fine.
3. For Absorbance instruments, make sure that Absorbance is chosen as the Data Mode and that Use plate blank is not checked.
4. Click OK to close the dialog box.

## Step 5: Set Display Parameters

1. Click the Display button to open the Display dialog box.
2. The default display for Endpoint readings is raw OD for Absorbance instruments and raw RFU for Fluorescence instruments. It is not necessary for the purposes of this tutorial to show a reduced number, so make sure the check boxes next to the options under Show are not checked. If you decide later that you would like to view the reduced data, you can change the display options after the reading since this does not affect the actual data—only how it appears on the screen and printout.
3. Click OK to close the dialog box.

## Step 6: Save the Protocol

1. Choose the **File > Save As** command.
  - ◆ On the Mac, select the popup menu command SoftMax Pro Protocol.
  - ◆ On Windows, select Pro Protocol Files (\*.ppr) from the Save as type list and give the file the name "Endpoint Tutorial" in the File name field.
2. Click Save to save a copy of the protocol file to the hard disk.
3. Close "Tutorial1 .pda". You are asked whether or not to save changes; say Don't Save.
4. Use **File > Open** to open the .ppr file you just created. This creates an untitled data file.

Congratulations! You have just created a SoftMax Pro protocol file and data file.

## Step 7: Read the Plate

If you were reading an actual microplate at this time, you would:

1. Prepare the samples in the microplate as it is defined in the Template Editor. (Before an actual reading, you might find it helpful to print a copy of the template to assist in properly filling the plate.)
2. Place the filled microplate in the drawer of the instrument.

Since we are using a simulator, you don't need to load the instrument.

1. To enable the simulator, select the Plate#1 group, hold down Ctrl+Shift and choose **Control > Enable Simulator**.
2. To read a plate selected **Control > Read**, or click the Read button in the Status Bar at the top of the window.
3. Values appear in the data display in the Plate section as if they were being read by the instrument, although faster than in real time.
4. The display for each well defined by a group boundary shows a raw value. The reduction formula is shown at the bottom of the display: Wavelength Combination: !Lm1.
5. Click the Display button in the Plate section tool bar. Select Reduced. Choose Grayscale from the **Options > Display** list. The dialog box updates to show the low and high limits for grayscale display values (which are automatically set to the minimum and maximum data values).
6. Click OK.

The data display in the Plate section updates to show the data represented by a heat map.

## Step 8: Data Analysis—Group Sections

SoftMax Pro software creates Group sections automatically when you define groups in the Template Editor.

The content of Group sections depends on the wells that are assigned to the groups in the Template Editor. Before a plate is read, these sections show no data for the wells. After a reading, Group sections contain information about the values for the wells contained within the groups.

Scroll down until the Standards group is visible.

The columns in this group are set up as follows, from left to right: - Sample: the sample names assigned for the Standards group.

- Concentration: the concentrations assigned to each sample in the template.
- BackCalcConc: Returns the curve fit associated with the plot of Standards from the
- Values: raw OD values obtained for each sample.
- MeanValue: the mean of data from the Values column for each sample. - StdDev: standard deviation for each sample value.
- CV%: coefficient of variation for each sample value.

In this part of this tutorial, you will view formulas associated with columns and Summaries, resize individual columns, Autosize all the columns, add a column, and hide replicates.

### Viewing Formulas

- To view the formulas associated with all columns in the Standards group section, hold down Ctrl+ Shift while the group is active, or select **Group > Show Formulas**. A row of formulas is displayed beneath the column titles.
- If formulas are long they may not fit in the columns. To see the complete formula for a particular column, click once on a column and the formula is reported in the Group section's tool bar.
- To modify or change a formula for a selected column, select the **Group > Edit Column** command or click the Formula button on the section tool bar.

### Resizing and Autosizing Columns

- You can resize columns by double-clicking or dragging column dividers.
- To Autosize columns, select the columns and use the **Group > Autosize** command.

### Adding a Column

You can add a column to any Group section and define the purpose of this column in your data analysis. By default, new columns are added to the right of any selected (highlighted) column or, if no columns are selected, after the last column in the group.

**To add a new column to the Standards group:**

1. Select the group's tool bar to make it active.
2. Deselect any highlighted columns by clicking outside the column area.
3. Choose **Group > New Column** or click the New Column button.
4. We are creating a new column that displays the maximum value obtained for each sample. Enter "MaxValue" in the Name field.
5. Enter "Max(Values)" in the Formula field and click OK.

A new column is displayed at the right of the previous columns, showing the maximum value obtained for each sample.

**Viewing an If Statement in a Column Formula**

The Unknowns Group section contains a column called Outliers. The formula for this column uses an If statement that uses data from another Group section to report when data points lie outside the range of the minimum and maximum Standard values.

An If statement in a formula defines a condition that, if found to be true, gives a certain result and, if not true, gives a different result.

The basic structure of an If statement is:

If (condition, result-if-true, result-if-false)

The Outliers formula is as follows:

```
If(Values>=MinStd@Standards and  
Values<=MaxStd@Standards,"","Outlier")
```

This statement says that, if the data shown in the Values column is greater than or equal to the calculation of MinStd found in the Standards group section and less than or equal to the calculation of MaxStd found in the Standards Group section, then report nothing (signified by two quotation marks with nothing between them); otherwise report the word "Outlier".

The MinStd or MaxStd calculations are used in the Summaries at the bottom of the Standards section, but the names are not displayed by default. To display them:

1. Double-click the first Summary, "Smallest standard value", to open the **Group > Edit Summary** dialog box.

2. Uncheck the Hide Name option and click OK. MinStd and MaxStd are now displayed next to their Summaries at the bottom of the group.
3. The formulas for MinStd and MaxStd are:  
Min(MeanValue)  
Max(MeanValue)

The MeanValue column reports the average of the well values in the Values column (the Standard replicates), and these are used in the Outliers formula. Such a formula might be useful in finding points that might be better masked and thus not used in the calculations.

For a complete discussion of writing formulas in SoftMax Pro, as well as detailed examples of formulas and customized protocols, see the *SoftMax Pro Software Formula Reference Guide*.

### Showing and Hiding Replicates

- Use the **Group > Hide Replicates** command to hide individual replicate values. Where appropriate, average values are shown instead.
- Show replicates again with the **Group > Show All** command.

### Step 9: Data Analysis—Standard Curve

1. Scroll down the SoftMax Pro window until you see the section tool bar for Graph #1.
2. Click the indicator on the left side of the tool bar to open it.
3. This graph is set to show concentration on the X-axis and mean value on the Y-axis.
4. You can edit graph properties in the **Graph > Graph Options** dialog box.

### Step 10: Print a Report

1. SoftMax Pro software has very flexible report printing. The default setting in SoftMax Pro software is for all Experiment sections to be included in the printed report. You can exclude sections from printing as described in Chapter 7, "File Management & Printing".
2. Choose **File > Print** to print the whole Experiment.

3. Select the appropriate settings for your printer and then click Print or OK to print the Experiment report.

Congratulations! You have completed this tutorial.

## Tutorial 2: EC 50 Protocol Tutorial

### (FlexStation Only)

Tutorial 2 is designed for a FlexStation instrument using a Flex read. It assumes that you have read the previous chapters in the manual and have a basic understanding of the logic and operation of SoftMax Pro.

The following steps are required to create and run an EC50 protocol using a graph of concentration vs. RFU:

1. Create a new data file.
2. Define instrument settings.
3. Define template.
4. Set reduction parameters.
5. Set display parameters.
6. Save the protocol.
7. Read the plate (you won't actually do this—instead, open a saved file).
8. Data analysis—Group sections.
9. Data analysis—Graph.
10. Data analysis—Summary formulas.
11. Print a report.

The following steps guide you through creating a protocol file for an EC50 protocol using a Flex read. Reading an actual plate is not required as SoftMax Pro software can generate simulated data.

### Step 1: Create a New Data File

1. Double-click the SoftMax Pro icon to start the program.
2. Choose the **File > New** command to create an untitled document. This will select your default Protocol format if you had previously selected one.

- If you are not connected to an instrument and want to emulate a FlexStation® instrument, choose **Edit > Preferences**, and choose a FlexStation instrument from the Reader list.

The Experiment has a Notes section (Notes#1) and a Plate section (Plate#1).

## Step 2: Define Instrument Settings

- Select the tool bar of Plate#1 to make it active.
- Click the Settings button in the tool bar or select **Plate > Settings**. Endpoint is selected by default. This tutorial uses a Flex read, so select Flex.
- Clicking the setting name (bold heading) on the left side of the Instrument Settings dialog box shows the choices available for that setting in the right pane, allowing you to make changes. Following are the details about the EC50 protocol settings we will use to create a protocol file using SoftMax Pro software. Make sure all settings match those in the table:

**Table 6-2** Example Settings

<b>Read Type</b>	<b>Flex</b>
Wavelengths	One: Excitation: 485; Emission: 525; Auto Cutoff
Sensitivity	3 readings; PMT set to High
Timing	Time: 200 seconds; Interval: 2 seconds; Total number of readings: 101; Minimum interval: 1.28 seconds
Assay Plate Type	96 Well Standard clrbtm
Wells to Read	A1–H10
Compound Source	Beckman 96 2.3 mL
Compound Transfer	None (initial volume 0 µL)
Triturate	N/a
Pipette Tips Layout	Full Rack

**Table 6-2** Example Settings

Read Type	Flex
Compound & Tip Columns	None
Automix	Off
AutoCalibrate	Once

### Final Check of Instrument Settings

Click OK to accept the settings and close the dialog box.

### Step 3: Define a Template

1. Click the Template button in the Plate section tool bar. This opens the Template Editor, showing all wells empty.
2. Select the first column of wells, A1 through H1.
3. Select New from the Group list to create a new group.
4. Enter "Data" in the Name field. Select Unknowns[Dilution] from the Column Format list. Select  $\mu\text{g}/\text{ml}$  as the units for the dilution factor. Click OK to close the dialog box.
5. The selected wells now show the default first sample name for the Data group, Data01. The area in the top center of the Template Editor is now active, showing the default values for Sample name and Dilution Factor.
6. In this tutorial, the default value of 0  $\mu\text{g}/\text{ml}$  for the dilution factor is correct. You can check the concentrations you have assigned to the template by pressing and holding the Ctrl + Shift keys. Click the Assign button or press the Return key on the keyboard to assign this dilution factor if this is not the case.
7. Select wells A2 through H2. Assign a dilution factor of 0.01  $\mu\text{g}/\text{ml}$  to these wells by entering 0.01 in the Dilution Factor field and clicking the Assign button.
8. A total of eight samples will be assigned to columns in this template. Continue to highlight subsequent columns of wells and enter the following dilution factors:  
Column A3 through H4:0.05  
Column A5 through H5:0.5



Column A6 through H6:1

Column A7 through H7:5

Column A8 through H9:10

Column A10 through H10:100

9. The template is now complete. Click OK to accept the entries and close the Template Editor. The outlines of the groups are now shown in the Plate section.
10. An associated Group section is created each time a group is created in a template, so we now have a Group section called Data.

#### **Step 4: Set Reduction Parameters**

1. Click the Reduction button in the Plate section tool bar to open the Reduction dialog box.
2. The default setting for Wavelength Combination, !Lm1, is fine.
3. For Kinetic Reduction change the setting to Max–Min.
4. Set Limits as follows:  
Min RFU:0  
Max RFU:50000  
Lag Time:0  
End Time:200
5. Baseline Options: Zero Baseline: 3 Pts.
6. Smoothing: Moving Average 1 point (no smoothing).
7. Click OK to close the dialog box.

#### **Step 5: Set Display Parameters**

1. Click the Display button to open the Display dialog box.
2. The default display for Flex readings is raw RFU. This tutorial example shows a reduced number, so check the With reduced number option.
3. If you decide later that you would like to view reduced data, you can change the display options after the reading since this does

not affect the actual data—only how it appears on the screen and when printed.

4. Click OK to close the dialog box.

### **Step 6: Save the Protocol**

1. Choose the **File > Save** command.
  - ♦ On the Mac, select the popup menu command SoftMax Pro Protocol.
  - ♦ On Windows, select Pro Protocol Files (\*.ppr) from the Save as type list and give the file the name “Flex Tutorial” in the File name field.
2. Click Save to save a copy of the protocol file to the hard disk.
3. Close “Untitled 1”. You are asked whether or not to save changes; say Don’t Save.
4. Use **File > Open** to open the .ppr file you just created.

Congratulations! You have just created a SoftMax Pro protocol file.

### **Step 7: Read the Plate**

If you were reading an actual microplate at this time, you would:

1. Prepare the microplate as it is defined in the Template Editor. (Before an actual reading, you might find it helpful to print a copy of the template to assist in properly filling the plate.)
2. Place the filled microplate in the appropriate drawer of the FlexStation. A new set of tips would also be loaded in the tip drawer.

Since we are using a saved file, it is not necessary to load anything in the actual instrument.

1. Choose **File > Open** and navigate to the Tutorial folder.
2. Open the file named Tutorial2.pda. The data in this file was generated using the parameters you have just created in this protocol file.
3. The display for each well that is defined by a group boundary shows a raw value with a reduced number.

4. The wavelength reduction formula is shown at the bottom of the display: Wavelength Combination: !Lm1.
5. The Kinetic reduction formula is shown at the top of the Plate section: Max – Min.
6. Click the Display button in the Plate section tool bar to open the Display dialog box.
7. Click the Reduced button.
8. Choose Grayscale from the drop-down menu. The dialog box updates to show the low and high limits for grayscale display values, which are automatically set to the minimum and maximum data values.
9. Click OK to close the dialog box.

The data display in the Plate section updates to show the data represented by a heat map.

## Step 8: Data Analysis—Group Sections

SoftMax Pro software creates group sections automatically when you define groups in the Template Editor.

The content of Group sections depends on the wells that are assigned to the groups in the Template Editor. Before a plate is read, these sections show no data for the wells. After a reading, Group sections contain information about the values for the wells contained within the groups.

Scroll down and click the small triangle on the tool bar of the Data group to open it and make it active.

The columns in this group are set up as follows, from left to right:

- Sample: the sample names assigned for the Data group.
- Wells: the wells that are assigned to the individual sample names (these correspond to the individual wells in the template).
- Concentration: the concentration assigned to each sample in the template.
- WellValues: raw RFU values obtained for each sample.
- MeanValue: the mean of data from the Values column for each sample.
- % Activation: sample activation relative to selected control group.
- CV%: coefficient of variation percent for each sample value.

## Viewing Formulas

- To view the formulas associated with all columns in the Data group section, hold down Ctrl + Shift while the group is active, or select **Group > Show Formulas**. A row of formulas is displayed beneath the column titles.
- If formulas are long, they may not fit in the columns. To see the complete formula for a particular column, click once on a column and the formula is reported in the tool bar.
- To modify or change a formula for a selected column, select the **Group > Edit Column** command or click the Formula button on the section tool bar.

Let's look more closely at the formula for % Activation:

- Double-click the % Activation column. The Calculation dialog box is displayed showing the formula used to derive the data for this column:

MeanValue /(Summary#1)\*1 00

This formula divides the numbers from the MeanValue column by the value of Summary#1, and then multiplies the result by 100 to construct a percentage.

- Close this dialog box by clicking OK.

## Resizing and Autosizing Columns

- You can resize columns by double-clicking or dragging column dividers.
- To Autosize columns, select the columns and use the **Group > Autosize** command.

## Adding a Column

You can add a column to any Group section and define the purpose of this column in your data analysis. By default, new columns are added to the right of any selected (highlighted) column or, if no columns are selected, after the last column in the group.

To add a new column to the Data group:

1. Select the group's tool bar to make it active.
2. Deselect any highlighted columns by clicking outside the column area.

3. Choose **Group > New Column** or click the New Column button.
4. We are creating a new column that displays the maximum value obtained for each sample. Enter "MaxValue" in the Name field.
5. Enter "Max(WellValues)" in the Formula field and click OK.

A new column is displayed at the right of the previous columns, showing the maximum value obtained for each sample.

### Step 9: Data Analysis—Graph

1. Scroll down the SoftMax Pro window until you see the section tool bar for Graph #1.
2. If this section is closed, click the triangle indicator on the left side of the tool bar to open it. The graph contains a single plot.
3. The graph is configured to show Concentration on the X-axis and % Activation on the Y-axis. A 4-Parameter curve fit has been applied. To apply a different fit, select it from the Fit list in the Graph#1 section tool bar.
4. To display the data in a bar graph or another format, open the **Graph > Graph Options** dialog box, or click the Graph Options button in the Graph#1 section tool bar.
5. This dialog box allows you to rename the graph, choose the type of graph that is displayed, and to format the text used in the graph. You can add or remove plots (click the New or Delete) and you can edit the properties of existing plots.
6. Change the type of graph to Cluster bar and click OK.

### Step 10: Data Analysis—Summary Formulas

- Two Summaries are displayed at the bottom of the Data group section. The second is Summary #2 = EC 50% 0.044. Click this Summary to highlight it.
- When selected, the formula used to create this Summary is displayed in the Data group section tool bar: "Summary #2: InterpX(Plot#1 @Graph#1 ,50)".
- Formulas can be simple or quite complex and are composed of elements that perform mathematical operations, statistical analysis functions, or refer to other elements contained within SoftMax Pro

software. Complete information about the possible choices for creating formulas can be found in the *Formula Reference Guide*.

- The formula from the Summary in the Data section of this Tutorial file uses a function, InterpX, that interpolates data from the curve fit in the Graph section. The InterpX function returns the X-values of the specified plot interpolated using the Y-values and the fitted curve. The complete formula is written InterpX(Plot@Graph,Yvalues). In this summary formula, Plot#1@Graph#1 describes the plot from which the data is derived. Yvalues in this example is 50, which describes the number to be interpolated. So, looking at the graph, the Y-value at 50% activation appears at an approximate concentration of 0.044.
- You can create a Summary formula either by clicking the Summary button in the tool bar or by choosing **Group > New Summary** to open the Calculation dialog box.
- The same Calculation dialog box appears when you edit an existing formula. For example, double-click Summary#2 at the bottom of the Data section to see the Calculation dialog box.
- If you wanted to interpolate a value at a different percentage, you could enter a different value in place of 50. You could also reverse the formula to show the percent activation that is interpolated for a certain concentration (0.1 in this example) by writing the formula this way:  
  
InterpY(Plot#1 @Graph#1 ,0. 1)
- The result is 62.867 (to three decimal places). The number of decimal places is configurable in the Calculation dialog box.

## Step 11: Print a Report

1. SoftMax Pro software has very flexible report printing. The default setting in SoftMax Pro software is for all Experiment sections to be included in the printed report. You can exclude sections from printing as described in Chapter 7, "File Management & Printing".
2. Choose **File > Print** to print the whole Experiment.
3. Select the appropriate settings for your printer and then click Print or OK to print the Experiment report.

Congratulations! You have completed this tutorial.

## File Creation and Management

SoftMax® Pro software uses two basic file types: data files and protocol files.

### Data Files

Data files contain the raw data collected by your reader along with any data analysis you may have completed.

SoftMax Pro software automatically creates data files based on the capabilities of the connected reader, or if a reader is not present, based on the reader type specified in **Edit > Preferences**.

### Protocol Files

Protocol files are Experiment template files that contain microplate well layout assignments and all other Reader configuration data, but no microplate data. A large number of predefined protocols are installed with SoftMax Pro software.

Saving a data file as a protocol file removes any data in the file, leaving only the configuration information.

### Creating a Default Protocol

If you frequently find yourself adding the same sections (Graph, Group) to the majority of your Experiment data files, or if you want to customize the document window that is opened every time you launch SoftMax Pro software (for example, to include Summaries or other information in the Notes section), you can create your own default protocol file.

Create the protocol by opening the document you want to be the default protocol and select **Protocols > Save As Default Protocol**. This file resides in the Protocols folder.

Whenever you launch SoftMax Pro software or use the **File > New** command, the document created will be based on your default protocol file.

## Resetting the Default Protocol Folder

The Protocols menu is a dynamically generated menu listing all the folders (one level deep) and protocol files in a selected folder.

To change the folder used to generate the menu, choose the **Protocols > Set Folder** command and navigate to the desired folder.

## Organizing Protocol and Data Files

In a multi-user environment, it is usually helpful to:

- Create a protocol directory for each SoftMax Pro user, or
- Create separate subdirectories for different types of experiments.

## Creating New Files

The File > New command opens a new Experiment file named "Untitled", based on the default protocol, or the instrument selected in the Edit > Preferences dialog box.

## Opening Files

To open a supported protocol or data file, use the File > Open command. SoftMax Pro software allows you to open files from prior versions of the program. However, once you save a file, it is readable only by the current version.

If you need to continue using a data file with an older version of SoftMax Pro software, do not overwrite the original when re-saving it.

While SoftMax Pro software opens both SoftMax and SoftMax Pro data files, two features are not forwards compatible: sample blanks and Auto-range limits in Graph sections. The data for sample blanks obtained using SoftMax software is read when these files are opened in SoftMax Pro software, but no calculations are performed automatically using this data. Auto-range limits are not available when SoftMax Pro software files are opened using SoftMax Pro software, so limits for the display of information in graphs must be re-entered manually.

## Cross-Platform Files

SoftMax Pro files created by either Windows or Mac users can be opened on either platform. The Mac honors SoftMax Pro software 3-letter file extensions as well.



**Table 7-1** File Extension Types

File Type	File Extension	Mac Document Type
SoftMax Pro Protocol File	.ppr	sPFd
SoftMax Pro Data File	.pda	sPFd
SoftMax Protocol File	.pro	sPFd
SoftMax Data File	.dat	sPFd

## Saving Files Manually

Use the **File > Save** command to save data or protocol files.

## AutoSave

SoftMax Pro software can be configured to save one or more copies of your data automatically—immediately after each plate read has completed. See [Setting Preferences on page 18](#) for more information.

## Printing

As explained in [Chapter 3](#), SoftMax Pro software is designed for WYSIWYG report printing: the software interface is identical to the report.

## Printing Reports

You can create customized reports by including or excluding Experiments and individual sections, changing the order in which sections print, and adding text describing each Experiment in Notes sections. You can also create Summaries (consisting of reduction formulas and associated text) in Notes and Group sections. The size and content of Group sections can also be changed.

## Including or Excluding Sections

When sections are created, they default to being included in the printed report. You can exclude sections from printing by clicking the print setting button on the tool bar of that

section (it is a toggle, so it will change from enabled to disabled and vice versa).

Changing the print setting option for the Experiment changes the state of all sections within that Experiment.

Changing the print setting option for one or more sections within an Experiment causes the print setting button for the Experiment to change to undecided. Toggling the Experiment's print setting button overrides any changes made to individual sections.

The File > Print dialog box contains a Print All option that overrides the print settings made to sections or Experiments.

## Changing Section Order

You can change the order that sections appear in an Experiment by drag and drop. Select the title bar of the section and drag it to its new position in the Experiment, marked by a thick black line.

## Changing Text Format

Use the Edit > Text Style dialog box to format text in Notes and Group sections.

# Import/Export

## Export

The File > Import/Export > Export command creates a data file from the data in one or more Plate sections, CuvetteSet sections, or Group table sections within a SoftMax Pro file. The exported data and format depends on the options in the Setting Preferences dialog box. Please contact MDS Analytical Technologies Technical Support for a copy of the SoftMax Pro XML schema.

Data is exported as it is displayed: if raw data is displayed, raw data is exported; if reduced data is displayed, reduced data is exported.

The File > Import/Export > Export dialog box contains additional export options that allow you to export selected plates and cuvettes or all plates and cuvettes, and selected groups or all groups.

## Export Format

The first line of an exported text file includes one field indicating the number of data sets in the file. This single field includes the text

"##BLOCKS=" followed by a number that equals the number of data sets.

The second line of a file for a Plate or CuvetteSet section includes 19 fields that provide all of the information needed to determine the type of test that was run to measure the first data set. The following table describes each of these fields.

**Table 7-2** Plate or CuvetteSet Field Descriptions

Field	Field Entry	Field Description
1	Section kind	Either Plate or Cuvette.
2	Section name	Defined in Section dialog box.
3	Export version	1.3 in SoftMax® Pro software version 5.x.
4	Export format	PlateFormat or TimeFormat, set in Preferences.
5	Read type	Endpoint, Kinetic, Spectrum, Well Scan, or Flex.
6	Data mode	For absorbance plates: Absorbance or %Transmittance. For others: Fluorescence, Luminescence, or Time Resolved Fluorescence.
7	Data type	Raw or Reduced.
8	Pre-read included	TRUE or FALSE; always FALSE for Cuvettes.
9	Kinetic points	Reduced plates and Endpoint plates = 1Kinetic / Spectrum plates: number of reads.
10	Kinetic or Flex read time; Well Scan read pattern	Kinetic / Flex read time in seconds or Well Scan read pattern (horizontal, vertical, X or fill); otherwise blank field.
11	Kinetic or Flex interval; Well Scan density	Kinetic or Flex read interval in seconds or Well Scan density; otherwise blank field.
12	Start wavelength	Spectrum start wavelength; otherwise blank field.
13	End wavelength	Spectrum end wavelength; otherwise blank field.
14	Wavelength step	Spectrum wavelength step; otherwise blank field.
15	# of wavelengths	# of wavelengths in Endpoint, Kinetic, Well Scan, Flex reads; for Spectrum; blank field.

**Table 7-2** Plate or CuvetteSet Field Descriptions (cont'd)

<b>Field</b>	<b>Field Entry</b>	<b>Field Description</b>
16	Read wavelengths	For Absorbance / Luminescence; wavelengths read, separated by a space; for Spectrum; blank field; for Fluorescence: emission wavelengths; for Ex / Em sweep: blank field.
17	First column read	First column read (1 to max), always 1 for cuvettes.
18	Number of columns	Last column read (1 to max), always 12 for cuvettes.
19	Number of wells or cuvettes	Number of wells in a plate (24, 96, 384) or number of cuvettes read for a CuvetteSet.
After Field 19, the format splits, depending on the read mode. Fields 20–28 are not present for Absorbance or Transmittance but are included for Fluorescence and Luminescence.		
20	Excitation wavelengths	Excitation wavelengths, separated by spaces; For Luminescence or sweeps, this is a blank field (the emSweep excitation wavelength is in field 16).
21	Cutoff option	Manual or Automatic.
22	Cutoff filters	Cutoff filter wavelength values, separated by spaces.
23	Sweep wave option	EmSweep or ExSweep or blank field.
24	Sweep fixed wave	The value for the fixed (em or ex) wavelength.
25	Reads per well	Number of times a well is read for a single reading.
26	PMT setting	Automatic, High, Medium, or Low.
27	Start integration time	Time to start integration for Time Resolved Fluorescence; otherwise, blank field.
28	End integration time	Time to end integration for Time Resolved Fluorescence; otherwise, blank field.
29	First row read	First row read (1 to max), always 1 for cuvettes.
30	Number of rows	Last row read (1 to max), always 8 for cuvettes.
31	Time tags	Indicate whether the exported data has read time tags.

The information for the first Plate or CuvetteSet section is provided in the lines following the descriptors. The number of data lines depends on the export format and the type of read that was run. A minimum of one data line is exported when the Time format is used with a single wavelength Endpoint read.

If your instrument has a temperature associated with the data (all readers except EMax®, VMax®, and UVMMax readers), the temperature at the time of the reading appears before the first data point in the reading.

The second line of a file for a Group table section includes fields that correspond to the number and types of columns found in that section (they are labeled as the table is labeled and look very much like the table itself).

Following the data, a blank line is included and then a line with a single termination field: ~End.

Each subsequent data set in the file includes the lines described above, a blank line, and a line with the termination field.

## **Import**

SoftMax Pro software imports data in tab-delimited ASCII text format, allowing you to import data collected by Criterion Host (Analyst AD/HT/Acquest), Analyst Host (Analyst GT), ScreenPlay (ScreenStation), or FLIPR® operating software (FLIPR) into SoftMax Pro software.

Only data that can be displayed in Plate or CuvetteSet sections can be imported.

Import supports up to 384-well plates; if you import 1536-well plate data only the top left quadrant of data is imported.

## **Import**

When data is imported a new Experiment is created in the open SoftMax Pro file for the imported data. New Plate sections and CuvetteSet sections are created for each data set within the file. Plate pre-read and CuvetteSet reference data can also be imported, but data for pathlength correction cannot.

The time/date stamp for a section containing imported data shows imported data rather than a date or time.

**File > Import/Export > Import Analyst** allows you to import data files generated by Analyst instruments (including Analyst, AD/HT/GT, Acquest,

and ScreenStation). Files for import must be generated as verbose reports using Criterion Host 2.0 or ScreenPlay 2.5 for all modes (ABS, FI, FP, Epi ABS, Lum, and the dual method) saved as text (\*.txt).

When importing 1536-well data from the Analyst instrument only, the first quadrant of 1536-well plate data (the top-left section of 24 by 16 wells) can be imported. This data is displayed in the standard 384-well plate format.

Instrument parameters for each plate in the text file are imported into a Notes section as text. Raw data is imported into a Plate section. The plate ID from the text file is used to name both the Plate and the Notes sections and the file name is used to name the Experiment.

More than one text file can be imported into the same SoftMax Pro file by creating or copying another Experiment into an existing file.

All analysis features of SoftMax Pro software can be used to analyze Analyst files.

### **Import FLIPR**

**File > Import/Export > Import FLIPR** allows you to import \*.FPD files that have been generated using FLIPR instrument control software (version 2.0 and above).

Files are exported from the FLIPR software using the **Plug Ins > Import/Export > FLIPR** command. In order to import FLIPR files, they must be exported by the FLIPR software into the specific \*.FPD format. Standard FLIPR files of the \*.FID and \*.FWD formats cannot be imported into SoftMax Pro software.

All analysis features of SoftMax Pro software can be used to analyze FLIPR files.

### **Importing and Exporting Templates**

Template information, saved as a tab-delimited ASCII text file or in XML format, can be imported to or exported from Plate or CuvetteSet sections. A template file consists of seven columns of data separated by tabs. Each line of the file provides information for one well in a plate or one cuvette.

When template information is exported, the information for well/cuvette A1 is written into the first line of the file, information for well/cuvette A2 is written to the second line of the file, and so on, until a line has been created in the export file for each well in a plate or all cuvettes. If no

template information exists for a well in a plate or for a cuvette, that line is left blank. The ASCII file terminates with the string "~End".

The order for importing template information must match the export order. Similarly, blank lines in the file are assigned to a well or cuvette, providing a well/cuvette with no template setting.

**Table 7-3** ASCII Template File Columns and Descriptions

Column	Column Entry	Description
1	Group Name	Assigned in the Group Settings dialog box. If this field is missing, the rest of the line is ignored. Any text string can be entered in this field.
2	Column format	Assigned in the Group Settings dialog box. Five text strings are supported in SoftMax® Pro software version 5.x: Empty, Basic, Standards, Unknowns, and Unknowns[Dilution]. Blank groups have an Empty column format. If this field is left blank, no sample name will be assigned to the well, although the sample name will be included in the group.
3	Sample name	The sample name is set in the Sample field of the Template Editor dialog box. If this field is left blank, no sample name will be assigned to the well, although the sample name will be included in the group.
4	Number of description	0 for Empty and Unknown column formats.
5	Descriptor units	Assigned in the Group Settings dialog box. Supported strings are unit/ml, mg/ml, µg/ml, ng/ml, mg, ng, and ml. If this column is left blank, unit/ml is assigned.
6	Sample descriptor	Assigned in the Group Settings dialog box. Any text string can be entered in this field or it may be left blank.
7	Descriptor value	Assigned in the Series dialog box. Can be set to any number or left blank.

The tab-delimited ASCII files can be opened with spreadsheet programs, such as Microsoft Excel, but if edited in Microsoft Excel they cannot be imported back into SoftMax Pro software.

SoftMax Pro software does not import files in the Microsoft Excel XLS format.

## Exporting Graphs

The **Graph > Export Graph** command exports a graph to file.

On the Mac, the graph is saved in PICT format; on Windows, it is saved in EMF (Enhanced Metafile) format.

Well Graphs can be exported with the **Plots > Export Graph** command.

## Copying and Pasting

Plate section data, CuvetteSet section data, and Notes section text can be copied from external programs and pasted into a Plate section, CuvetteSet section, or Notes section.

For Plate section data, the format of the copied information must match the Instrument Settings of the target Plate section and the settings chosen in the Export Format of **Edit > Preferences**. Pre-read data cannot be pasted into a Plate section. After data is pasted into a Plate section, the information to the right of the display shows imported data.

If you have created a template for a Plate section prior to pasting data, numbers are shown only in the groups you have created in the template. Data in other wells is not reported (but is available if you alter the template). Clearing the template or assigning groups to all wells in the plate will show the complete data set.



MDS Analytical Technologies provides a fully integrated hardware and software solution for adherence to Good Manufacturing Practices (GMP), Good Laboratory Practices (GLP), and many other regulatory compliance requirements including 21 CFR Part 11. This solution consists of:

- An MDS Analytical Technologies microplate reader.
- SoftMax® Pro GxP software for data collection and analysis.
- GxP Admin software for user and permission management.
- MDC FileServer (optional) for TCP/IP-based file serving of one or more multiple User Accounts file.

The GxP Admin application manages all SoftMax® Pro GxP software users and user permissions and creates and stores this information in a proprietary User Accounts (.edb) file. SoftMax Pro GxP software is a version of SoftMax Pro software that includes features to provide regulatory compliance with 21 CFR Part 11 requirements.

SoftMax Pro software connects to and controls all MDS Analytical Technologies microplate readers, and it analyzes all collected microplate and cuvette data. SoftMax Pro GxP software includes full user and permission control and audit trail generation in addition to all the data collection and analysis features found in the Standard edition. With the audit trail features in SoftMax Pro GxP software, all Experiments are fully traceable and can be completely reconstructed.

Before SoftMax Pro GxP software application can be launched and used, a User Accounts file must be created by the administrator using GxP Admin software.

The User Accounts file records all authorized users and their associated information including:

- User names, user login IDs and passwords.
- Assigned user permissions to SoftMax Pro GxP features.
- Global settings pertaining to all listed users (password aging, non-network use, and guest access).

Audit histories of all administrator actions within GxP Admin (user creation, modification).

- Electronic signatures.

If set by the administrator, SoftMax Pro GxP software users do not require continual access to a User Accounts (.edb) file in order to use SoftMax Pro GxP software for data collection and analysis when no network connection is available (corporate network failure, use of SoftMax Pro GxP software while traveling).

## How the Programs Work Together

SoftMax Pro GxP software, GxP Admin, and the MDC FileServer applications work together as follows:

- The administrator installs GxP Admin and MDC FileServer software on a Windows PC.
- The administrator creates a User Accounts (EDB) file with GxP Admin after entering a special Creation Key number.
- Using GxP Admin software, the administrator creates SoftMax Pro GxP users, and assigns the appropriate user permissions.
- SoftMax Pro GxP software is installed on individual computers that are used for data collection or data analysis. Each Windows log on account must be linked to a User Accounts file before SoftMax Pro GxP software can be used. This can be completed by each end-user, by the administrator, or by IT personnel.
- After each Windows log on account is linked to a User Accounts file, users can launch SoftMax Pro GxP software, log on, and use the software to collect and analyze microplate data.
- If enabled by the administrator, Guest users can open, view, and print Experiment data files.

## GxP Admin Features

The GxP Admin software supports the following major functions:

- Template Management: Create templates or groups of SoftMax Pro GxP permissions. This makes managing large numbers of users easy, as the administrator can use templates to quickly and easily assign sets of permissions to new users.
- User Management: Add new users, modify user permissions, and inactivate users. Once created, users cannot be deleted.
- File Reporting: Export user information and permissions to a text file. View the audit trail of administrator activities.

- License Management: Add new site license information, which is used by GxP Admin to control the number of users registered to use MDS Analytical Technologies software programs.
- Password Aging: Users can be required to change their password on a regular basis; users are prevented from reusing their previous password.
- Idle Timeout: Automatically log users out of the SoftMax Pro GxP application after a defined period of inactivity.
- Guest Access: The administrator can specify whether or not guest users are allowed to launch the SoftMax Pro GxP application. If enabled, guests can only open, view, and print existing SoftMax Pro data files.
- Forced Password Change: When set globally, all users are required to change their passwords when they first log on to SoftMax Pro GxP software.
- Lock Software after three log on attempts: inactivate specific user accounts within a User Accounts file after three failed login attempts. The administrator can reactivate users remotely by using GxP Admin software.
- Offline Use: When specified by the administrator, SoftMax Pro GxP software automatically copies an encrypted read-only copy of the User Accounts file for each user, after each user has successfully logged on at least once. This local copy of the User Accounts file (under administrative control) then allows users to work without a persistent network connection, thereby allowing users to continue using the SoftMax Pro GxP application when no network connection is available.

For information on the specific use and configuration of GxP Admin features, please consult the GxP Admin User Manual.

## Using SoftMax Pro GxP Software

### User Log On

#### Authorized Users

Once the SoftMax Pro GxP application has been installed, all authorized users can log on to the SoftMax Pro GxP application and begin working.

Simply launch SoftMax Pro GxP software and all users will be prompted to enter their UserID and password.

### **Guests**

New users who have not been added to the User Accounts file may use the SoftMax Pro GxP application as a Guest, if the administrator has enabled Guest access to the User Accounts file using GxP Admin. Guests can open, view, and print any existing Experiment data files or protocols. They are not able to modify or save changes to any existing Experiment data file or protocol.

### **Files**

Files created in the SoftMax Pro GxP application use the file name extensions .epr for protocol files and .eda for data files, and can be opened only by the SoftMax Pro GxP application.

Files saved from other editions of the SoftMax Pro application (.ppr, .pda, .dat, .pro) can be opened by the SoftMax Pro GxP application but have no associated audit trail.

### **GxP Menu**

The GxP menu in the SoftMax Pro GxP application provides access to most of the security and GxP management features in the SoftMax Pro GxP application. The available commands depend on the user's permissions.

### **Reset User Accounts**

Displays a dialog box in which the user (or administrator) links to an existing User Accounts file. The link can be made using a direct file path or a TCP/IP connection; see the *GxP Admin User Guide* for more information on configuring connections to User Accounts files.

### **Account Info**

Displays the permissions available for the currently logged on user.

### **Log On**

Prompts a user to enter a user ID and password. All users must log on if they want to read new microplates or perform any data analysis.

**Log Off**

Terminates the user's connection to the SoftMax Pro GxP application.

**Change Password**

Allows the currently logged on user to change his/her password.

**Add Note to Audit Trail**

Opens a dialog box in which the user can securely append a note to the permanent audit trail for the current file.

**Show Audit Trail**

Opens a file showing a list of all user activities recorded for the current file.

**Statements**

Within the SoftMax Pro GxP application a user with sufficient permissions can add a Statements section to an Experiment data file. Electronic statements provide a means of designating that an Experiment has been approved, reviewed, or verified. Users with appropriate permissions can also apply electronic signatures to these statements. Once a statement has been signed, the file cannot be modified.

Electronic statements are saved with Experiment data files and protocol files, but signatures are not saved with protocol files. Only one electronic statement section is allowed per file and is always the first section displayed for each Experiment data file.

**Add Statements**

**Experiment > New Statements** allows a user to enter a title for each statement or comment, and to enter a brief note for all other users who may be using or accessing a particular Experiment data file.

After the first statement has been created, it appears as the Statements section at the top of the Experiment data file.

**Statements Menu**

Once a statement has been created, the Statements menu is displayed when the statement is active (it is similar to the other section menus such as Plate, and Notes). The Statements menu

allows the logged in user to create or modify statements in the current Experiment data file.

### **Add Statement**

Allows the user to create a new statement (if that permission is available to the user).

### **Edit Statement**

Opens the Electronic Statement Setup dialog box, allowing the current user to change the title and text of the active statement.

### **Remove All Signatures**

If a statement has been signed, a user with this permission can remove all signatures from the active statement. When all signatures have been removed, the file is editable again.

### **Copy Statement**

Copies the contents of the statement onto the clipboard.

### **Paste Statement**

Inserts the clipboard contents into a Statement section of the current Experiment.

### **Sign Statement**

Allows the user to add their electronic signature to a Statement section.

Once a Statement is signed, the Experiment is locked and cannot be edited unless signatures are removed.

Multiple signatures may appear on one Statement.

Signatures are saved with data files but are not saved with protocols.

Add A Note: Allows the user to add a note to the statement prior to signing it.

Agree: Prompts the user for a password and signs the document with the user name, date, and time.

## **Locking and Unlocking Sections**

Sections (Notes, Plate, Group, Graph, CuvetteSet) can be individually locked if the user has the necessary permissions enabled for their account. Newly created sections are unlocked by default, and can be

locked and unlocked by clicking the lock/unlock button on the right side of the tool bar for each section (it is a toggle, so the icon changes from locked to unlocked).

Changing the lock setting for the Experiment changes the lock state of all sections within that Experiment.

Changing the lock setting for one or more sections within an Experiment causes the lock setting button for the Experiment to change to mixed, reflecting the fact that the Experiment contains both locked and unlocked sections. Toggling the Experiment's lock setting button overrides any changes made to individual sections.





## Introduction

In order to integrate MDS Analytical Technologies readers and the SoftMax® Pro software with robotic systems from other manufacturers, or to automate the export of data from the SoftMax Pro application to a desired LIMS package, the SoftMax Pro application supports a remote control interface. MDS Analytical Technologies has tested the interface but does not provide technical support for specific integration needs. MDS Analytical Technologies strongly recommends that customers consult a third-party automation company if internal software resources or expertise are not available.

All leading automation vendors have integrated the SoftMax Pro software and MDS Analytical Technologies' instruments with their robotics systems. A latest list of approved robotics partners can be found on the MDS Analytical Technologies website.

## Sample Source Code and Applications

Automation sample source code and applications are included as a part of the SoftMax Pro software installation package. The sample source code is available as Visual Basic, Visual C++, or Excel Macro projects, and can be found in the Automation Examples folder in the SoftMax Pro application directory.

## Remote Control

SoftMax Pro software can be controlled by other Windows applications with simple ASCII string-based messages. Using the Microsoft Windows Win32 SendMessage function, the SoftMax Pro application remote command strings such as ReturnStatus, Read or SaveAs:myFile can be sent as a Windows message to the SoftMax Pro application for remote process execution.

## Send/Receive Remote Commands

The user can choose between sending remote commands using Windows message WM\_SETTEXT or WM\_COPYDATA.

## The WM\_SETTEXT Approach

Call the Win32 RegisterWindowMessage function to register a unique message tag. The SoftMax Pro application uses this tag information when the command is received to distinguish between a Windows generated message and an external command. The tag should be a non-zero value:

```
UINT tag = RegisterWindowMessage("SOFTMaxPROMsg");
```

All messages must be sent to the SoftMax Pro main window. Note that if the SoftMax Pro application is not currently running, this returns a null value, requiring your program to discontinue until the SoftMax Pro application is restarted:

```
HWND hwnd = FindWindow("SOFTMaxPROMainWnd", "SoftMax Pro");
```

If the target SoftMax Pro application is the GxP version, the Window name should be "SoftMax Pro GxP".

```
HWND hwnd = FindWindow("SOFTMaxPRORemote", "SoftMax Pro GxP");
```

A complete list of remote commands can be found in SoftMax Pro Commands, below. Do not forget that the tag value, from RegisterWindowMessage above, is sent with the command string. (See Microsoft documentation regarding the SendMessage function for more information.)

SendMessage(hwnd, WM\_SETTEXT, tag, (LPARAM)remoteCommandString) For example, to send SoftMax Pro a Read command:

```
SendMessage(hwnd, WM_SETTEXT, tag, (LPARAM) "Read")
```

If return data is expected, such as with remote commands ReturnStatus, ReturnTiming, and ReturnData, the data is returned to the clipboard. The advantage of this approach is that it may be universally applied to Visual C++ and Visual Basic modules, and Excel macros. An example of retrieving clipboard data is shown below, where the string data retrieved is finally stored into a Microsoft CString object.

```
if ( !::OpenClipboard(NULL) )  
    return FALSE;  
  
// Error opening clipboard  
HANDLE handle = ::GetClipboardData(CF_TEXT);
```

```

if (handle == NULL) {
    ::CloseClipboard();
    return FALSE; // Error, no data in clipboard
}
// get clipboard data, store in MFC CString object
CString returnStr = CString ((char *) handle );
// close the clipboard
::CloseClipboard();

```

## The WM\_COPYDATA Approach

Call the Win32 RegisterWindowMessage function to register a unique message tag. SoftMax Pro uses this tag information when the command is received to distinguish between a Windows generated message and an external command. The value returned should be a non-zero value:

```
UINT tag = RegisterWindowMessage("SOFTMaxPROMsg");
```

Call the Win32 RegisterWindowMessage function to register another unique message tag. Your software can use this tag information when the asynchronous windows message is returned, and can distinguish between a Windows generated message and an external command. The value returned should be a non-zero value:

```
UINT replyTag =
RegisterWindowMessage("SOFTMaxPROReplyMsg");
```

All messages must be sent to the SoftMax Pro main window. If SoftMax Pro software is not currently running this returns a null value, thus, requiring your program to discontinue until SoftMax Pro software is restarted:

```
HWND hwnd = FindWindow("SOFTMaxPROMainWnd", "SoftMax
Pro");
```

If the target SoftMax Pro software is the GxP version, the Window name should be "SoftMax Pro GxP".

HWND hwnd = FindWindow("SOFTMaxPRORemote", "SoftMax Pro GxP"); Get your window handle. This is the window to receive the return message:

```
HWND MyWnd = GetSafeHwnd()
```

Assign the command to send. Place the command in heap space, which appears to be necessary with older Windows operating systems:

```
char *cmdStr = "Read";// Example remote command
// to send
char *msgStr = _strdup(cmdStr);// String duplication, copy
// command to heap space.
```

Create the Win32 COPYDATASTRUCT message that is sent. See Microsoft documentation regarding this structure for details:

```
COPYDATASTRUCT cd; // This struct is defined in
// Microsoft include file
cd.dwData = (DWORD) MyWnd; // Handle to a window
// receiving return
// information
cd.cbData = strlen(msgStr) + 1;// Length of message string,
// adding 1 for null
// terminator
cd.lpData = msgStr;// String pointer to remote
// command
```

Finally, send the message to SoftMax Pro software. A complete list of remote commands can be found in SoftMax Pro Commands, below. Don't forget that the tag value, from RegisterWindowMessage above, is sent with the command string. (See Microsoft documentation regarding the SendMessage function for more information.)

```
SendMessage(hwnd, WM_COPYDATA, tag, (LPARAM) &cd)
```

Either a completion message or return data is returned. Remote commands ReturnStatus, ReturnTiming, and ReturnData return data. In either case, data is received through an asynchronous windows message inside Win32 COPYDATASTRUCT type data packet.

For example, a typical OnCopyData window callback is shown below, where the string data retrieved is finally stored into a Microsoft CString object. Note the use of variable replyTag, discussed above, which is used to isolate the correct windows message returned.

```
BOOL CUserDlg::OnCopyData(CWnd* pWnd,
COPYDATASTRUCT* cd) {
....
if (cd->dwData == replyTag)
{
/* String pointing to status */
CString retStatus = (char*) cd->lpData;
.... }
```

## Summarizing Remote Commanding

SoftMax Pro software has two groups of remote command instructions: those that do not return data, and those that do.

### Remote commands that DO NOT return data

Commands that do not receive data are further broken into:

- The user wants to receive a completion message after the command has been sent

After receiving a remote command that used the WM\_COPYDATA message, SoftMax Pro software sends a completion message once the command has been executed. An asynchronous window message is returned as a window callback, typically an OnCopyData handler, inside the Win32 COPYDATASTRUCT types data packet. For example, if the Read command is sent, an "OKRead" text string is returned after the read is completed.

- The user does not want to receive a completion message

Use SendMessage combined with WM\_SETTEXT for messages that do not require a completion message. The advantage of this approach is that it may be universally applied to Visual C++ and Visual Basic modules, and Excel macros.

### Remote commands that DO return data

There are three remote commands that return data: ReturnStatus, ReturnTiming, and ReturnData.

If the caller has enabled queue operation, the command ReturnStatus is handled immediately, and not queued, as are other remote commands. Data is returned in one of two ways:

- After sending the remote command using the WM\_COPYDATA message, the caller can look for return data through an asynchronous Windows message inside a Win32 COPYDATASTRUCT type data packet. An asynchronous window message is returned as a window callback, typically an OnCopyData handler.
- After sending the remote command using the WM\_SETTEXT message, the return data is available immediately on the clipboard. The advantage of this approach is that it may be universally applied to Visual C++ and Visual Basic modules, and Excel macros.

### Programmer Notes

- COM ports 1–9 can be used for robotics purposes. MDS Analytical Technologies recommends using COM1 or COM2. These are typically the most stable.
- An Autosaved file from the robotics run can be saved as either a text (.txt) or SoftMax Pro experiment data file (.pda). This selection is made in the SoftMax Pro Edit > Preferences dialog box.
- More random access memory (RAM) is required to run SoftMax Pro software using remote commands. Depending on the robotics used and the type of file created with SoftMax Pro software, a computer configured with the minimum amount of RAM may restrict the size of the file that can be generated.
- It is important to put a short pause between remote commands.

An Experiment with more than one plate of data could adversely affect SoftMax Pro software efficiency. If increasing the computer's memory is not an option, MDS Analytical Technologies recommends using only one plate per file and also ensuring that the computer is not being used for other purposes while SoftMax Pro software is running.

## Remote Control in SoftMax Pro GxP Software

If the target SoftMax Pro application is the GxP version, /R is required in the program command line (for example, "SoftMaxProGxP.exe /R") when launching the SoftMax Pro GxP application in order to avoid the Log On dialog box which requires manual entry of user ID and password. Instead, the Logon remote command can be sent to log on the user through the robotic interface.

## SoftMax Pro Commands

SoftMax Pro commands are single words with no spaces. If parameters are required with a command, a colon is used between the command and the parameters.

Commands may be sent using either the WM\_SETTEXT or WM\_COPYDATA approach.

Completion messages or data are returned directly when commands are sent using the WM\_COPYDATA method.

Completion messages or data are placed in the system clipboard when commands are sent using the WM\_SETTEXT method.

### **AppendCurSectionTitle**

Append text to the title of the current Plate section.

### **CDrawer**

(FlexStation® Instrument Only)

This command returns the status [Open or Closed] of the compound or reagent plate drawer.

### **Close**

Closes the current document.

If the data has not been saved, the document is closed anyway with no user warning.

### **CloseCDrawer**

(FlexStation Instrument Only)

Tell SoftMax Pro software to close the compound drawer.

### **CloseDrawer**

Tell SoftMax Pro software to close the instrument drawer.

### **CloseTDrawer**

(FlexStation Instrument Only)

Tell SoftMax Pro software to close the tips drawer.

### **Copy**

This is equivalent to pressing CTRL+C or selecting **Edit > Copy**.

If a Plate section has just been read or is the only selected section, the plate data is copied to the clipboard.

The data copied follows the display settings for the current plate (normally Raw Data).

Other sections such as Groups can also be copied. The SelectAll command should be used first.

### **Drawer**

This command returns the status [Open or Closed] of the assay plate drawer.

### **ExportAs: -a filepath**

Tell the SoftMax Pro application to export data in the standard SoftMax Pro format.

The ExportAs command without the -a attribute overwrites any existing file without warning. However, you can choose to append export data to the end of an existing file by specifying the -a attribute. In addition, this exported file can then be imported and all data can be viewed as a single document file.

Please note that, when using the attribute -a, a single space is required between the attribute and the following directory and file string.

#### **Example**

```
ExportAs:c:\SoftMax Pro\exportfile.txt
```

```
ExportAs: -a c:\SoftMax Pro\exportfile.txt
```

### **GetAutosaveState**

Returns the AutoSave state. Return value is one of following:

1 = AutoSave is turned on

2 = AutoSave is turned off

3 = AutoSave is enabled, but the specified format is not SoftMax Pro file (.pda) format.

### **GetNumPlateSections**

Returns the number of Plate sections in active document.

### **ImportTemplate:filepath**

Have the SoftMax Pro application import a template file.



- The file should comply with SoftMax Pro's template file format.
- A Plate section should be selected prior to importing a template.

### **Logoff**

(SoftMax Pro GxP Software only)

This logs off the currently logged on user in the SoftMax Pro GxP application.

### **Logon:userid password**

(SoftMax Pro GxP Software only)

This enables the remote logon of a user with remote logon permission to the SoftMax Pro GxP application. The logon ID and password are separate by single space.

#### **Example**

Logon:mylogonID myPassword

### **NameCurSection**

This renames the current selected section. If no section is selected, nothing happens.

- This command is used typically to identify different plates in an automated environment.
- The data copied follows the display settings for the current plate (normally Raw Data).

#### **Example**

NameCurSection:Plate#002345

### **New**

Have the SoftMax Pro application send a **File > New** command. This creates a new document and reads in the values contained in the "Default Protocol.ppr" protocol file.

### **NewExperiment**

Create a new Experiment section.

### **NewNotes**

Tell the SoftMax Pro application to send an **Experiment > New Notes** command. This creates a new Notes section within the current Experiment.

### **NewPlate**

Tell the SoftMax Pro application to send an **Experiment > New Plate** command. This creates a new Plate section within the current Experiment.

### **OpenAssay:filepath**

Tell the SoftMax Pro application to open a protocol file. The file should be in SoftMax Pro's Protocols folder if the full file path is not specified. The Protocol folder can be set using the Protocols menu. Protocols can also be stored in subdirectories of the Protocols folder.

- Only protocol files can be opened.
- SoftMax Pro software appends .ppr to any file name passed in.
- If the file is not found, the SoftMax Pro application does nothing.

#### **Example**

OpenAssay:myFile

OpenAssay:\mySubDirectory\myFile

OpenAssay:c:\temp\testprotocol

### **OpenCDrawer [FlexStation only]**

Tell SoftMax Pro to open the compound drawer.

### **OpenDocument**

Tell the SoftMax Pro application to open a data file. The file should be in SoftMax Pro's application folder unless the full file path is specified.

If the file is not found, the SoftMax Pro application does nothing.

#### **Example**

OpenDocument:myDataFile

### **OpenDocument:c:\temp\testdataOpenDrawer**

Tell SoftMax Pro software to open the instrument drawer.

**OpenTDrawer**

(FlexStation only)

Tell the SoftMax Pro application to open the tips drawer.

**Quit**

Tell the SoftMax Pro application to exit the application.

**Read**

Tell the SoftMax Pro application to read the current Plate or CuvetteSet section. If the current section is neither a Plate nor CuvetteSet section, next Plate or CuvetteSet section will read.

**ReturnCalc:formulaName@sectionName@experimentName**

This command will return the result from specified formula. This command requires the name of the formula, name of either Notes or Group section, and the name of the Experiment. If the name of the Experiment is not specified, it will search from the first Experiment of the document.

**Example**

ReturnCalc:MyFormula@MyGroupSection@MyExperimentSection

**ReturnData**

This command asks the SoftMax Pro application to send data back from the currently selected plate. Generally this is used to send the data directly to the caller instead of using the clipboard. If data is to be returned to the clipboard, this command functions like the SoftMax Pro **Edit > Copy** command when a plate is the currently selected section.

**ReturnStatus**

This command returns the current instrument status. This command can, for example, be sent to check whether the instrument is busy or reading or staying idle.

The status is an ASCII string that has:

- Information about the instrument type and version.
- Whether the instrument is busy or idle.
- The instrument's current temperature.
- The state of its drawer.

The exact format of this response depends on the type of instrument in use.

If the SoftMax Pro application failed to place data in the clipboard, then no return status is available.

**Example:**

```
ReturnStatus:<tab><cr>
```

```
ReaderType<tab>Version#<cr> State:<tab>[Offline, Busy, or Idle]
```

```
Temperature:<tab>XXX
```

```
Drawer:<tab>[Open or Closed]
```

**ReturnTiming**

This command sends back the number of seconds that it would take to read the current plate. For multiplate documents, the plate to be used is chosen in the same way as described in the Read command.

**SaveAs :filepath**

Tell the SoftMax Pro application to save the current document file. The file is saved in the current SoftMax Pro AutoSave directory, which is set in the Edit > Preferences dialog box. Data files can also be saved to subdirectories of the Autosave directory.

The SoftMax Pro application appends .pda to any file name passed in to it.

If a file with the same name already exists, the SoftMax Pro application automatically overwrites the file with no warning.

If the subdirectory does not exist, the file is not saved.

**Example**

```
SaveAs:myFile
```

```
SaveAs:\mySubDirectory\myFile
```

**SelectAll**

This is equivalent to pressing CTRL+A or choosing **Edit > Select All**. SelectAll can be used to select data in group tables for copying (not needed for copying plate data).

**SelectNextPlateSection**

Selects next Plate section from the current Plate section.

**SelectSect**

Select a section by name. SelectSect can be useful for multi-plate protocols or for selecting Group tables for copying. In version 5 and above, this command can also be used to specify sections in different Experiments. If no Experiment is specified, SelectSect defaults to the active Experiment.

**Example**

```
SelectSect:Plate#0023499
```

```
SelectSect:Group#1 @Experiment#2
```

**SelectSectNum :X**

Select a section by its order in the document. SelectSectNum can be useful for multi-plate protocols or for selecting Group tables for copying.

**Example**

```
SelectSectNum:3
```

**SetFolderAs :folderpath**

Set the current protocol directory. You must specify the root directory.

**Example**

```
SetFolderAs:c:\SoftMaxPRO\Basic Protocols
```

**SetPort :X**

Tell the SoftMax Pro application to set a serial COM port.

**SetTemp :XX.X**

Send a command to set the current instrument incubator temperature.

Setting temperature to zero turns off the incubator.

**Shake:On/Off**

This command, if set to 'On', starts shaking the microplate tray until the Shake:Off command is sent. By default, the shake will stop after 30 seconds in most instruments.

**Example**

```
Shake:On
```

## Stop

Send the SoftMax Pro application a command to stop reading. If queue operation is enabled, this command is not queued, the queue is cleared, and the command is processed immediately. This is similar to clicking the Stop button on the SoftMax Pro tool bar except the Stop button does not affect remote command processing or its queue.

This command should be sent before any Close or Quit command if there is a possibility that the instrument is still reading.

If the data has not been saved, the application quits with no user warning.

## TDrawer

(FlexStation Instrument Only)

This command returns the status [Open or Closed] of the tip rack drawer.

## TrayLock: On/Off

When sent before a read, this command keeps the microplate tray locked inside the instrument after a read if the parameter is set to 'On'. If set to 'Off', the tray comes out.

### Example

```
TrayLock:On
```

## Version

Returns the version of the (active) SoftMax Pro application.

## Visual Basic (Excel Macro) Example

```
' Declarations
Private Declare Function FindWindow Lib "user32" Alias
    "FindWindowA" _ (ByVal IpClassName As String, ByVal
    IpWindowName As String) As Long
Private Declare Function SendMessage Lib "user32" Alias
    "SendMessageA" _ (ByVal hWnd As Long, ByVal wParam As Long,
    ByVal lParam As String) As Long
Private Declare Function RegisterWindowMessage Lib "user32" Alias _
    "RegisterWindowMessageA" (ByVal msgName As String) As Long _
'-----

Const WM_SETTEXT = 12
```

```
' No. of seconds that GetMsgFromPro() will wait before timing out
Const mcsngMsgTimeout As Single = 600

Public Sub Start()
    ' Start the main form
    MainForm.Show
End Sub

Public Function SendMsgToPRO(ByVal msgStr As String) As Long
    ' Function to send windows command to SoftMaxPro application
    ' This is the primary method of communication between the two
    processes, on the same PC
    Dim lVal As Long
    hWnd = FindWindow("SOFTMaxPROMainWnd", "SOFTmax Pro")
    If hWnd = 0 Then
        Beep
    Else
        softmaxMsg = RegisterWindowMessage("SOFTMaxPROMsg")
        If softmaxMsg = 0 Then
            Beep
            Beep
        Else
            lVal = SendMessage(hWnd, WM_SETTEXT, softmaxMsg, msgStr)
        End If
    End If
End Function

Public Function GetMsgFromPro(ByVal strMessage As String) As String
    ' First send a request message to the instrument and then get
    ' back a text message from the instrument via the clipboard.
    ' Wait for a message for up to (mcsngMsgTimeout) seconds
    before giving up.

    Dim sngStart As Single

    Dim Clipboard As DataObject
    Set Clipboard = New DataObject
```

```
Clipboard.Clear ' clear the clipboard for the data that we are
about to get.

SendMsgToPRO (strMessage)

sngStart = Timer ' get the time that we started looking at
clipboard

' wait for an answer for up to (mcsngMsgTimeout) seconds
Do While Timer < sngStart + mcsngMsgTimeout
  'GetMsgFromPro = Clipboard.GetText() ' get text from the
  clipboard.

  Clipboard.GetFromClipboard ' Retrieve text from clipboard
  GetMsgFromPro = Clipboard.GetText(1)

  If Len(GetMsgFromPro) > 0 Then
    ' if we got message back from the clipboard then exit
    with it in hand
    Clipboard.Clear
    Exit Do
  End If
DoEvents ' Yield to other processes while we wait.
' make sure that we didn't just span midnight with our timer
check
If Timer < sngStart Then
  ' Current timer is less than start time so reset the
  starting time.
  ' In the worst case, this routine might take twice as long
  to timeout
  ' if the message fetch spanned midnight. This is unlikely
  and even if ' it does occur, it doesn't hurt anything.
  sngStart = Timer
End If
Loop

End Function
```



## MFC C++ Interface to SoftMax Pro

Some SoftMax Pro software IPC commands return values. By default they put the data in the clipboard, but optionally they use the WM\_COPYDATA mechanism to return data directly to the caller.

In order to have the SoftMax Pro application response returned in this fashion, the commands must be sent using WM\_COPYDATA so that SoftMax Pro knows what to send the response to.

SoftMax Pro then replies with another WM\_COPYDATA message that contains the response.

### MFC Example Code

The sample MFC code is taken from the \Visual C++\RemoteCmdTest v2 project in the Automation Examples folder, which is installed in the SoftMax Pro application directory by default.

#### To send a command to SoftMax Pro

```

/*****
    SendCmdToSMP (Send a single command to SoftMax Pro)

    1.Retrieve the handle to the SoftMax Pro window
    2.Register the window message
    3.Construct the COPYDATASTRUCT
    4.Send the command to SoftMax Pro
*****/
void CMySampleClass::SendCmdToSMP( CString cmdStr )
{

    HWND hwnd = ::FindWindow( "SOFTMaxPROMainWnd", "SoftMax Pro");
    UINT smpMsg = RegisterWindowMessage("SOFTMaxPROMsg");

    ...

    // use WM_COPYDATA
    char temp[255];
    strcpy( temp, (LPCSTR)cmdStr );

    COPYDATASTRUCT cds;
    cds.dwData = (ULONG_PTR)this->m_hWnd;

```

```
    cds.cbData = (DWORD)strlen( temp ) + 1;
    cds.lpData = &temp;

    ::SendMessage( hwnd, WM_COPYDATA, (WPARAM) smpMsg, (LPARAM)
&cds );

    ...
}
```

If the command is the Read remote command, the status can be periodically checked by sending ReturnStatus remote command.

### To parse the response from SoftMax Pro

```

/*****
    OnCopyData (Message handler for WM_COPYDATA)
    1.Receive data from SoftMax Pro.
    2.Retrieve the response.
    3.If the response is for ReturnStatus, parse it accordingly.
*****/
BOOL CMySampleClass::OnCopyData(CWnd* pWnd, COPYDATASTRUCT*
pCopyDataStruct)
{
    CString aStr;

    char* temp = (char*)pCopyDataStruct->lpData;
    aStr.SetString(temp, pCopyDataStruct->cbData - 1);
    aStr += '\n';

    if( aStr.Find("ReturnStatus:") != -1 )

        CheckInstrumentState(aStr);

    ... \\or do your own parsing.
    return TRUE;
}

```

### To add a short pause between remote commands

```

const DWORD kMaxTimeOut = 1000; // 1 second
for( int i = 0; i < noOfCmdsToSend; i++ )
{
    ...
    SendCmdToSMP( cmdStr );

    // Put 1 second buffer
    DWORD start = GetTickCount();
    while( ( GetTickCount() - start ) < kMaxTimeOut ) { }
}

```



# Glossary of Terms

A

## **%Transmittance, T**

Transmittance is the ratio of transmitted light to the incident light (for absorbance readings).

$$T = (I/I_0) \\ \%T = 100T$$

where  $I_0$  is incident light, and  $I$  is transmitted light.

## **Absorbance, A**

Absorbance is the amount of light absorbed by a solution. To measure absorbance accurately, it is necessary to eliminate light scatter. In the absence of turbidity, absorbance = optical density.

$$A = \log(I_0/I)$$

where  $I_0$  is incident light, and  $I$  is transmitted light.

In this manual, we use the terms absorbance and optical density interchangeably.

## **Area Under Curve**

Reduction formula for a Kinetic or Spectrum scan that determines the area under the Kinetic or Spectrum plot, or under a plot in a Graph section.

## **AutoCalibrate**

For Absorbance instruments, an air reference reading is taken before runs (and sometimes between Kinetic readings during a run) as determined by the type of instrument and read settings. For Fluorescence instruments, the measurement occurs before reads (and sometimes between Kinetic reads if the interval is long enough).

Calibration data is stored in memory and is used by SoftMax® Pro software until the instrument is powered down or the wavelength is changed. AutoCalibration can be turned off.

### **Automix**

The Automix function determines how often, if at all, automated shaking of the micro- plate is performed during a reading.

### **AutoRead**

This feature enables automatic reading of subsequent Plate sections in the order in which they appear within an Experiment. You can set intervals (delay times) between the plate readings, if desired.

### **Baud**

A data transmission rate (measured in bits/second) for communication devices.

### **Bottom Read**

Available in the Instrument Settings, this selection causes instruments with bottom read capability to read up through the bottom of the microplate rather than down from the top.

### **Compound Source**

Compounds are transferred from the compound source plate to the assay plate.

### **Cubic Spline (Curve Fit)**

This curve fitting method generates a fit to a cubic equation between each pair of data points. The cubic equations are computed such that these equations and their first two derivatives are continuous everywhere.

### **Cutoff**

A filter used to condition the light entering or exciting the monochromators. In automatic mode, the instrument sets the cutoffs automatically based upon the wavelengths chosen for reading. With some read modes and types, you can choose a different filter wavelength (manual setting) for the emission monochromator.

### **CuvetteSet Section**

CuvetteSet sections determine how the cuvettes will be read, display data as it is acquired from the instrument, and define how the data will be displayed and used in reduction. CuvetteSet sections can display samples in three different ways: three samples per row, one sample per row, and in a grid of cells corresponding to a microplate format.

**Data Mode**

Reduction option to view absorbance data as OD (absorbance) or %Transmittance, fluorescence data as RFU (relative fluorescence units), or luminescence data as RLU (relative luminescence units).

**Default Protocol**

A protocol file, included with the SoftMax Pro program or created by a user, that resides in the default folder. The Default Protocol file defines the initial information that appears when you open SoftMax Pro software or when you create a new file from within the program. Both Mac and Windows versions of this file are called "Default Protocol" and both versions append the 3-character ".ppr" (standard) or ".epf" (GxP) extension.

**Emission Spectrum Scan**

Measures fluorescence or luminescence across a spectrum of wavelengths for emitted light for a fixed excitation wavelength. The default value reported for each well is the wave-length of maximum emission (either fluorescence or luminescence).

**End Time**

Reduction parameter used to omit the end of a Kinetic, Spectrum, or Flex run from data reduction.

**Endpoint**

A single reading made at one or more wavelengths.

**Excitation Spectrum Scan**

Measures fluorescence at a single emission wavelength across a spectrum of excitation wavelengths. The default value reported for each well is the excitation wavelength of maximum fluorescence.

**Exponential (Curve Fit)**

The exponential function used to generate this curve fit is:

$$y = A + B * (1 - \exp^{(-x / C)})$$

**Flex**

A read type using fluorescence that is available with the FlexStation® instrument. You can specify compound additions and trituration for individual wells. Data is collected over time, with readings taken at regular

intervals (Flex runs are measured in seconds and fractions of seconds). Flex readings can be single-wavelength or multiple-wavelength (up to four).

**FLIPR®** Reader

MDS Analytical Technologies' FLIPR® (Fluorometric Imaging Plate Reader) system was developed to perform cell-based, high-throughput screening (HTS) assays.

**Fluorescence**

The light emitted by certain substances resulting from the absorption of incident radiation. To measure fluorescence accurately, it is necessary to reduce light scatter.

The governing equation for fluorescence is:

Fluorescence = extinction coefficient \* concentration \* quantum yield \* excitation intensity \* pathlength \* emission collection efficiency

**Fluorescence Polarization**

By using a fluorescent dye to label a small molecule, its binding to another molecule of equal or greater size can be monitored through its speed of rotation.

SpectraMax® M5 and M5e Fluorescence Polarization read mode returns three different raw data types: Raw (S&P), Polarization (mP), and Anisotropy (r).

**Fluorophore**

A material that absorbs light energy of a characteristic wavelength, undergoes an electronic state change, and instantaneously emits light of a longer wavelength.

**Four-Parameter Logistic (Curve Fit)**

The equation used to generate this curve fit is:

$$y = ((A - D) / (1 + (x / C)^B)) + D$$

**G-Factor**

G-factor (Grating factor) is used in fluorescence polarization to correct polarization data for optical artifacts, converting relative mP data to theoretical mP data. Optical systems, particularly with reflective components, pass light of different polarization with different efficiency.



G-factor corrects this instrumental bias.

### **Gain**

The amount of increase in signal power expressed as the ratio of output to input for a photomultiplier tube.

### **Grayscale Data Display**

Raw or reduced data is displayed proportionally in a plate format using seven shades of gray, ranging from light shading at the low end to dark shading at the high end, based on user-defined high and low limits.

### **Group**

Wells can be assigned to group types using the Template Editor. Depending on the default protocol used, certain group types (such as Standard or Unknown) may be created automatically; you may create others as required.

### **Incubator**

Choosing Incubator from the Control menu or clicking the Incubator button opens a dialog box allowing you to start or stop temperature regulation and to select an elevated temperature for the microplate chamber on instruments that have temperature control capability.

### **Instrument Icon**

Located in the Status Bar, the instrument icon shows the status of the connection between the computer and the instrument. If the icon has a "?" through it, the connection is not functioning or no instrument is connected; if the instrument icon appears without a "?," the connection is working properly.

### **Interleaved Display**

A display choice in which the wells are shown in a format that skips every other well as follows: all odd columns and rows begin in the upper left corner of the plate display and are followed by all even columns and rows. This display is most useful when the 384-well plate comprises four daughter plates of 96 wells each.

### **Kinetic**

During Kinetic readings, data is collected over time, with multiple readings made at regular intervals (measured in seconds). The values

calculated based on raw Kinetic data are VMax®, VMax per Sec, Time to VMax, and Onset Time. Kinetic readings can be single- or multiple-wavelength readings (up to six, if using a SpectraMax® instrument other than the Gemini or Gemini XS which have a maximum of four, or the FlexStation which also has four).

### **Lag Time**

Lag time in a Kinetic protocol is the period of very slow growth (microorganisms) or the rate of reaction that may precede the rapid/linear phase of reaction. When a lag time is specified in SoftMax Pro software, data collected prior to the lag time is not included in data reduction.

### **Large Display**

A display choice available with 384-well plates. This setting is useful when looking at large numbers if you are using a monitor that is 17-inches or larger.

### **Linear (Curve Fit)**

The linear function fits the best straight line to the data. The equation for this fit has the form of:

$$y = A + Bx$$

where A is the y-intercept of the line and B is the slope. A linear fit should be used whenever the values appear to lie on or scattered around a straight line.

### **Log-Logit (Curve Fit)**

The Log-Logit is also called a two-parameter curve fit. The equation used to generate this curve fit is:

$$y = ((A - D) / (1 + (X / C)B)) + D$$

### **Luminescence**

The emission of light by processes that derive energy from essentially non-thermal changes, the motion of subatomic particles, or the excitation of an atomic system by radiation.

### **Mask**

Lets you hide selected data so that they are not used for calculations and are not reported. The Mask function is commonly used to suppress outliers.

**Maximum**

Reduction for a Spectrum scan that reports the maximum signal within the reduction parameters.

**MaxOD, MaxRFU, or MaxRLU**

The limit for the maximum value you want to use for displaying and analyzing Kinetic or Spectrum data. Any values above these limits are not shown and are excluded from data reduction.

**Minimum**

Reduction for a Spectrum scan that reports the minimum signal within the reduction parameters.

**MinOD, MinRFU, or MinRLU**

The limit for the minimum value you want to use for display and analysis of Kinetic or Spectrum scan data. Any values that are under this limit are not shown and are excluded from data reduction. The default is 0. To display negative Kinetics, the value should be set below 0 (zero).

**Normal Display**

The default display for 384-well plates in the Plate section.

**Onset OD, Onset RFU, or Onset RLU**

The change in signal required to compute the onset time (for Kinetic readings).

**Onset Time**

The time it takes for a given increase in signal, called the onset OD (or RFU or RLU) to occur.

**Optical Density, OD**

The amount of light passing through a sample to a detector relative to the total amount of light available. Optical Density includes absorbance of the sample plus light scatter from turbidity.

**PathCheck® Software**

Instrument Settings option within SoftMax Pro software for the SpectraMax M2, M2e, M5, M5e, Plus, Plus384, 190, or 340PC readers that allows data collected in a Plate section to be normalized to a 1-cm pathlength.

### **Photomultiplier Tube (PMT)**

A vacuum tube that detects light especially from dim sources through the use of photo emission and successive instances of secondary emission to produce enough electrons to generate a useful current.

### **Pipette Tips (FlexStation Instrument only)**

The FlexStation instrument requires the use of pipette tips. MDS Analytical Technologies suggests that you load a complete set of pipette tips each time you make a run to ensure that sufficient tips are always available. See the instrument manual for the FlexStation instrument regarding the types of tips that can be used and how they should be loaded.

### **Plate Section**

The Plate section is used to display data, to specify how microplates are read, and to define how the data received from the instrument should be reduced. The data display in this section is a matrix that corresponds to the well format of the microplate in use.

### **Plate Type**

A choice in the Instrument Settings dialog box that determines the type of microplate being read. The specification of the plate type depends on the well format and manufacturer of the microplate.

### **Point-to-Point**

This curve-fitting option fits a linear equation to each pair of data points.

### **Point Count/Number of Reads per Well**

The number of integrations performed on a single well in a Fast Kinetic read.

### **Protocol File**

Contains the template(s), instrument settings, and reduction parameters. Protocol files are useful if you repeat a particular type of protocol frequently.

### **Quadratic (Curve Fit)**

A mathematical description of a parabola:

$$y = A + Bx + Cx^2$$

**Ranged Data Display**

Raw or reduced data is assigned proportionally to integer values from 0 through 9 based on user-defined limits. Values above the high limit are displayed as plus (+) and values below the low limit are displayed as minus (-).

**Raw Data**

Signal reported from the instrument with no alteration. This is reported as OD, RFU, or RLU, depending on the instrument type.

**Read Interval**

The interval of time between the start of one reading and the start of the next reading during a Kinetic or Flex run.

**Read Type**

The method used to read the microplate or cuvette: Endpoint, Kinetic, Spectrum, Well Scan, or Flex. Not all types are available with all instruments.

**Read Mode**

The type of reading performed in the instrument: Fluorescence, Time Resolved Fluorescence, Luminescence, Absorbance or Fluorescence Polarization.

**Readings per Well**

The number of times (programmable by the user) that readings are taken on a well in Fluorescence and Time Resolved Fluorescence modes. This setting also determines the amount of time that data is collected in Luminescence mode.

**Reduced Data**

Data reduction causes the raw signal values reported by the instrument to be calculated and displayed according to user-defined formula settings.

**Reduction Limits**

Sets limits on data to be included in reduction for Kinetic, Spectrum, Well Scan, or Flex readings.

### **Reference**

A reading of buffer or water taken in the cuvette that is used as IO to calculate Absorbance or %Transmittance of sample read in the cuvette port.

### **Semi-Log (Curve Fit)**

The semi-log function fits the best straight line to a set of data for  $\log(X)$  plotted against Y. The resulting curve displayed will be a straight line with the X-axis drawn in logarithmic scale.

### **Series**

The Series function allows you to work with groups of wells in the Template so that the standard value or dilution factor increases or decreases in specified steps. The name can be incremented as well.

### **Slope**

A reduction option that determines the slope of the combined plot (for example, the slope of the line using linear regression after the wavelength combination reduction). This reduction uses all time points visible in the reduction window. Slope is the same as VMax Rate when VMax Rate is set to the same number of points as the run (for example, the default). Slope is not the same as VMax Rate if you have modified the number of VMax Points.

### **Spectrum**

A read type that allows you collect absorbance or fluorescence data across a spectrum of wavelengths. This data is displayed as OD, RFU, or RLU versus wavelength. The default reduction displays the wavelength that corresponds to the maximum signal.

### **Standard Curve**

A mathematically idealized representation of the relationship between the concentration of a standard calibrator (X-value) and the OD/RFU/RLU (Y-value). The standard curve may be set to a linear, semi-log, log-log, quadratic, 4-parameter logistic, log-logit (2-parameter logistic), exponential, cubic spline, or point-to-point curve fits.

### **Stokes Shift**

The difference between the wavelengths of the excitation and emission peaks.

### **Template Editor**

A representation of the microplate or a set of cuvettes shown as a grid of wells that can be used to designate the location of blanks, standards, controls, unknowns, empty wells, or to assign wells to other groups you create. The template is a map of the microplate. It tells the software what is in the various wells of the microplate or cuvette.

### **Threshold Data Display**

Data is displayed as plus (+) for values above, asterisk (\*) for values within, and minus (-) for values below user-defined limits.

### **Time Resolved Fluorescence**

Most fluorescent substances are not suitable for this type of reading. However, the fluorescence emitted by lanthanide dyes is delayed long enough to measure fluorescence after the lamp is turned off. Time resolved fluorescence is used to reduce the amount of background noise interfering with fluorescence. The excitation lamp flashes and, after it is off, the delayed emission is collected before the lamp flashes again.

### **Time at 1/2 Maximum**

Reduction for a Kinetic or Flex run that reports the time at half the maximum signal within the reduction limits as follows: SoftMax Pro software first determines the point (within the reduction limits) that has the maximum signal level. It then scans the plot from left to right until it finds two points that have signals bracketing one-half of that signal value. A linear interpolation between these two points is used to estimate the Time at Half Maximum.

### **Time at Maximum**

Reduction for a Kinetic or Flex run that reports the time at the maximum signal within the reduction limits.

### **Time at Minimum**

Reduction for a Kinetic or Flex run that reports the time at the minimum signal within the reduction limits.

### **Time to VMax**

The elapsed time until the maximum reaction rate of a Kinetic or Flex reading is reached, reported in seconds.

**Triturate (FlexStation Instrument only)**

To mix a solution by means of aspirate and dispense cycles.

**User Name**

The User Name is the name of the person who reads data into the shielded file. The User name is entered after the Read button is pressed but before the reading begins.

**VMax Rate**

The rate reported as signal/min (milli-OD, RFU, or RLU units per minute) for a Kinetic or Flex reading. It is calculated using a linear curve fit,  $y = Ax + B$ . A creeping iteration is performed using Vmax Points and the slope of the steepest line segment is reported as Vmax Rate. It may also be reported as units per second (the default for Fluorescence and Luminescence modes).

**VMax Points**

The number of points in the Kinetic or Flex curve included in each of a series of regressions to calculate the maximum rate, VMax, or the time segment used to calculate the Time to VMax.

**Well Scan**

Some applications involve the detection of whole cells in large-area tissue culture plates. Well Scan reads can be used with such microplates to allow maximum surface area detection in whole-cell protocols. Since many cell lines tend to grow as clumps or in the corners of microplate wells, you can choose from several patterns and define the number of points to be scanned to work best with your particular application.



# Index

## A

Absolute Values 93  
Absorbance 9, 44, 58, 97, 173  
reads 97  
add  
summary 101  
adding  
columns 100  
analysis  
parallel line 111  
application window 25  
Area  
under curve 92, 94, 173  
Assay Plate Fluid 62  
assign  
a compound 65  
a tip to an assay plate well 65  
button 72  
Autocalibrate 58, 173  
for Luminescence 59  
Automix 28, 174  
feature 57  
AutoRead 68, 174  
Autosize 102

## B

Baseline Options  
absolute 95  
with Multiplier 96  
Zero 96  
Baud 174  
blanking 57, 74  
blanks  
group-associated 76  
Bottom Read 174

## buttons

assign 72  
Automix 28  
clear 72  
Drawer 28  
edit 72  
help 27  
Incubator 27  
Read 57, 79  
Reference 27  
Series 73  
StakMax 27  
X-Axis 39  
Y-Axis 39

## C

Calibrating  
a plate 81  
Calibration 81  
change  
column formulas 101  
tip assignment 66  
Clear  
button 72  
Close Drawer command 29  
column  
adding 100  
copy 101  
format 71  
formulas 101  
paste 101  
Wavelength Priority 67  
commands  
Close Drawer 29  
Open Drawer 29  
Ref 29

---

**Compound**  
  Source 62, 174  
  Transfer 62

**Control** 28

**copy**  
  column 101  
  data 35  
  experiment 77  
  group 38  
  section 32  
  template 36, 77

**correlation coefficient** 113

**creating**  
  default protocol 40

**Cubic Spline** 174

**cubic spline** 108

**curve**  
  fit weighting 110  
  fitting 105  
  point-to-point 108  
  sigmoidal 107  
  standard 107

**Curve Fit** 176  
  Cubic Spline 174  
  Exponential 175  
  Four-Parameter Logistic 176  
  Linear 178  
  Log-Logit 178  
  Quadratic 180  
  Semi-Log 182

**custom formulas** 97

**Cutoffs** 46, 174

**cuvette**  
  data collection 79  
  reading 78

**CuvetteSet** 26, 53  
  section 38, 76, 89, 174

---

## D

**data** 175  
  collection 78  
  collection from a Cuvette 79  
  copy 35  
  display 82, 83, 177  
  Grayscale 84  
  paste 35  
  Ranged 84  
  ranged display 181  
  raw 83, 181  
  reduced 84, 181  
  reduction 88  
  scan 179  
  show 84  
  Threshold 84  
  threshold 183  
  transmission rate 174

**Data Mode** 97

**default**  
  protocol 40, 175

**defining**  
  a series 73

**Delay** 48

**delete**  
  experiment 31  
  group 38  
  section 32

**deselect**  
  tip 66

**dialog**  
  display 83  
  graph options 103  
  group settings 37  
  plot 103  
  reduction 35

display  
  data 83  
  data during reading 82  
  interleaved 177  
  large 178  
  normal 179  
  polarization 85  
  temperature 26  
  wells 85  
Drawer 28  
Dual Read 43  
duplicating  
  experiment 30

---

## E

edit  
  button 72  
  plate type 61  
editor  
  template 34, 68, 183  
Emission Spectrum 175  
End Time 175  
Endpoint 42, 175  
  readings 46  
  reads 89  
error bars 105  
Excitation Spectrum 175  
Experiment  
  delete 31  
  duplicate 30  
  manipulating 30  
  paste 77  
  rename 30  
  reorder 31  
  section 31, 77  
  selecting 30  
  window 31  
exponential 175  
  function 109

export  
  graph 40  
  template 78

---

## F

fit  
  settings 110  
fitting  
  curve 105  
Flex 43, 175  
  reads 95  
Flex Only  
  Pipette Tips Layout 64  
Flex only  
  compound source 62  
  compound transfer 62  
Fluorescence 9, 44, 176  
Fluorescence Polarization 45, 67, 176  
Fluorometric Imaging Plate Reader 176  
Fluorophore 176  
formulas  
  show 101  
Four-Parameter Logistic 176  
function 105  
  delay 48  
  exponential 109  
  injection 48  
  integration 47  
  linear 105  
  log-log 106  
  quadratic 106  
  semi-log 106  
  Shake 47

---

## G

Gain 177  
G-Factor 176

Good Curve Fit  
  judging 113  
Graph 35  
  export 40  
  options 39  
  section 39, 103  
graphing  
  wells 86  
Grayscale Data Display 177  
Group 177  
group  
  copy 38  
  delete 38  
  lists 69  
  names 37  
  settings 37, 69  
  show formulas 101  
Group-Associated Blanks 76

---

## H

Hide 102  
HTS 176

---

## I

Icons  
  instrument 26  
import  
  template 37, 78  
Incubator 177  
Injection 48  
  Wells 48  
instrument  
  controls 26  
  icon 177  
Integration 47  
Interleaved Display 177

---

Interpolate  
  raw data 96

---

## K

Kinetic 43, 177  
  limits 92  
  readings 46  
  reduction 90  
  timing 49

---

## L

Lag Time 178  
Lambda  
  Maximum 94  
  Minimum 94  
Large Display 178  
limits  
  Kinetic 92  
linear 105, 178  
lists  
  groups 69  
log-log  
  function 106  
Log-Logit 178  
Luminescence 9, 45, 178

---

## M

manipulating  
  experiment 30  
  sections 31  
Mask 35, 178  
Maximum 179  
MaxOD 179

---

---

MaxRFU 179  
menus  
    active section 31  
    Control 28  
method  
    Pre-Read 55  
microplate  
    reading 78  
Minimum 179  
    Time 63  
MinOD 179  
MinRFU 179  
MinRLU 179  
modes  
    data 97  
    Fluorescence Polarization 67

---

## N

Normal Display 179  
notes section 33

---

## O

OD (optical density) 179  
ODs 44  
Onset  
    OD 179  
    RFU 179  
    RLU 179  
    Time 91, 179  
open  
    active section 33  
Open Drawer command 29  
Optical Density 179  
options  
    graph 39

---

## P

Parallel Line analysis 111  
parameters  
    manual setting 110  
paste  
    column 101  
    data 35  
    experiment 77  
    template 36, 77  
    template regions 77  
PathCheck 50, 89, 179  
    calculation 54  
    feature 51  
    measurements 56  
pathlength 50  
Photomultiplier Tube 180  
Pipette Height 62  
Pipette Tips 180  
Pipette Tips Layout 64  
plate  
    blank 55  
    calibration 81  
    section 34, 88, 100  
    sections 75  
    type 61  
    types 60  
Plate Background Constants 55  
Plate Blank 55, 75  
Plate Section 180  
Plate Types 60, 180  
plots  
    create multiple 104  
PMT 49  
    sensitivity 49  
PMT (photomultiplier tube) 180  
Point Count 180  
Point-to-Point 180  
point-to-point curve 108  
Polarization  
    display 85  
Pre-Read method 55

---

print  
  section 33  
protocol  
  creating 40  
  default 175  
  file 180  
Protocol File 180

---

## Q

Quadratic 180  
quadratic  
  function 106

---

## R

Ranged Data Display 181  
Raw Data 181  
Raw data 83  
  interpolate 96  
Read 57  
  button 79  
read  
  mode 181  
  type 181  
Read Interval  
  read  
    interval 181  
Read Mode 181  
read mode 44  
  Fluorescence 44  
  Fluorescence Polarization 45  
  Luminescence 45  
  Time Resolved Fluorescence 45  
Read Type 181

read type 42, 61  
  Dual Read 43  
  Endpoint 42  
  FLEX 43  
  Kinetic 43  
  Spectrum 43  
  Well Scan 43  
reading  
  cuvette 78  
  data display during 82  
  Endpoint 46  
  Kinetic 46  
  microplate 78  
  reference 80  
  sample 79  
  Spectrum 46  
Readings per Well 181  
reads 58  
  absorbance 58, 97  
  EndPoint 89  
  Flex 95  
  Spectrum 94  
  Well Scan 49  
Recalculation 99  
reduced  
  data 84, 181  
reduction  
  dialog 35  
  formulas 97  
  limits 181  
  Spectrum 94  
  Well Scan 95  
Ref command 29  
reference 182  
  reading 80  
renaming  
  experiment 30  
reordering  
  experiments 31  
RFUs 44, 45, 92  
RLUs 45, 92

---

**S****Sample**

- area 72
- Descriptor 70
- Descriptor field 72
- list 72
- reading 79

**Scale**

- to Data 87
- to Limits 87

**section 31**

- active 31
- collapse 32
- copying 32
- CuvetteSet 26, 38, 76, 89
- delete 32
- expand 32
- Experiment 77
- Graph 103
- graph 39
- Group 37, 100
- manipulating 31
- name 32
- new 31
- notes 33
- open active 33
- plate 34, 41, 100
- print 33
- reorder 33
- view 32

**selecting**

- an experiment 30
- cuvettes 69
- wells 69

**semi-log 182**

- function 106

**Sensitivity 49**

- PMT 49

**Series 182**

- button 73
- defining 73

**settings 18, 63**

- Axis 113
- Compound & Tip Columns 63
- fit 110
- group 37, 69
- instrument 41
- manual 110
- X-Axis 87
- Y-Axis 87

**Settling Time 67****Shake 47****show**

- command 84
- group 101

**Slope 91, 182****Smoothing 95****spectrophotometers 9****Spectrum 43, 182**

- limits 94
- readings 46
- reads 94

**Speed Read 67****StakMax 27****Standard Curve 182****standards**

- minimum 109

**Status Bar 26****Stokes Shift 182****Summary**

- adding 101
- edit 34
- new 34

---

**T****Target Calibration 49****temperature 26**

---

## template

- copy 36
- copying 77
- editor 34, 68, 183
- export 37
- exporting 78
- import 37
- importing 78
- paste 36, 77
- regions 77

## Threshold Data Display 183

### Time

- at 1/2 Maximum 183
- at Maximum 183
- at Minimum 183
- end 93
- lag 93
- Maximum 91
- onset 91
- Point 63
- Resolved Fluorescence 45, 183
- to VMax 91, 183

### Timing 49

### tips

- transfers 65

### transfers 62

- compounds 65
- tips 65

### Transmittance 97, 173

### TRF 183

### Triturate 64, 184

---

## U

### User Name 184

---

## V

### View menu

- menus 32

### VMax 90, 184

- points 90, 184
- time to 91

### Volume 63

---

## W

### Water Constant 52

### Wavelength 46

- combination 89, 90, 95
- priority 67
- target calibration 49

### Well Display 85

### Well Scan 43, 95, 184

- Editor 59
- reads 49
- reduction 95

### wells 61

- graphing 86

### WYSIWYG 25

---

## X

### X-Axis 87, 104

- button 39

---

## Y

### Y-Axis 87, 105

### Y-Axis button 39



