

and even some tap water, even with the high grade of stainless steel used in the manufacture of this unit.

When the bath is used under certain conditions, algal growth can occur which can be mistaken for rust. This can easily be cleaned off using warm soapy water.

Boekel recommends the use of a bath algaecide, item # B1906001 to help maintain the bath water. Apply treatment at a dose of 20 drops per 1 gallon of distilled water, free of visible algae.

## 5. Specifications

Temperature Range	Ambient +5 to 100°C		
Temperature Setting Range	10 to 100°C		
Stability	±0.2°		
Temp Resolution	0.1°C		
Heat up to 65°C -	6 liter unit	40 minutes	
	14 liter unit	40 minutes	
	28 liter unit	45 minutes	
Timer	Settable from 1 min. to 99 hrs. 59 min.		
Power Rating:	115VAC 60Hz.	6L - 300w	14L - 575w 28L - 1100w
	230VAC 50/60Hz.	6L - 300w	14L - 575w 28L - 1100w
Amperage Rating :	115VAC	6L - 2.6a	14L - 5.0a 28L - 9.6a
	230VAC	6L - 1.3a	14L - 2.5a 28L - 4.8a