

Multi-Purpose Tube Rotator

Catalog Number:
88861049
88861050





Important before using this product, read this entire operation manual carefully. Users should follow all of the operational guidelines contained in this manual and take all necessary safety precautions while using this product. Failure to follow these guidelines could result in potentially irreparable bodily harm and/or property damage.

Caution all internal adjustments and maintenance must be performed by qualified service personnel.

Material in this manual is for information purposes only. Fisher Scientific committed to a continuing program of product development and improvement, and reserves the right to change information, such as specifications, appearance, and dimensions, described in this document without notice. Fisher Scientific makes no representations or warranties with respect to this manual. In no event shall Fisher Scientific be held liable for any damages, direct or incidental, arising out of or related to the use of this manual.

No part of this manual may be reproduced or transmitted in any form or by any means, including photocopying, recording, or using information storage and retrieval systems, for any purpose other than the purchaser's own use, without the express written permission of the manufacture.

Any other product names and services identified in this manual are trademarks or registered trademarks of their respective owners. No such use, or the use of any trade name, is intended to convey endorsement or other affiliation with Fisher Scientific.

@2015 Fisher Scientific. All rights reserved.

Table of Contents



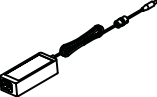

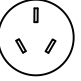
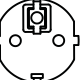
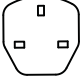

Section 1	Inspection and Installation	1-1
	Packing List	1-2
	Connections	1-3
	Structure Diagram	1-3
Section 2	Overview	2-1
	Specifications	2-1
	Environmental Conditions	2-2
	Safety Instructions	2-3
Section 3	Operation	3-1
	Operations of the Control Panel	3-1
	Settings	3-2
	Installation of Accessories	3-4
Section 4	Safety Tips and Maintenance	4-1
Section 5	TroubleShooting	5-1
Section 6	Optional Accessories/Spare parts	6-1
Section 7	Warranty	7-1

Section 1 Inspection and Installation


1. Inspect package and contents upon receipt of the instruments. If the package is severely damaged or if there are any missing pieces, please contact the manufacturer immediately.
2. Unpack the instrument, ensure all parts of the instrument and accessories are not missing or damaged. Make sure to take out all the components before discarding the packing. If there are any missing or damaged pieces, please contact the manufacturer immediately.
3. Place the instrument on a level and firm surface to avoid vibration and noise.

1.1 Packing List


Table -1 Packing List

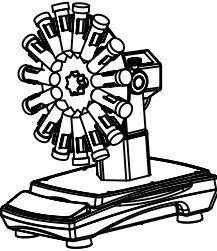Cat. No	88861049	88861050	
Model	Multi-Purpose Tube Rotator US	Multi-Purpose Tube Rotator International	
50x1.5/2.0ml Rotator Drum	1	1	
General Power Adaptor	1	1	
US Plug	1	N/A	
CN Plug	N/A	1	
EU Plug	N/A	1	
UK Plug	N/A	1	
Screw Driver	1	1	

1.2 Connections

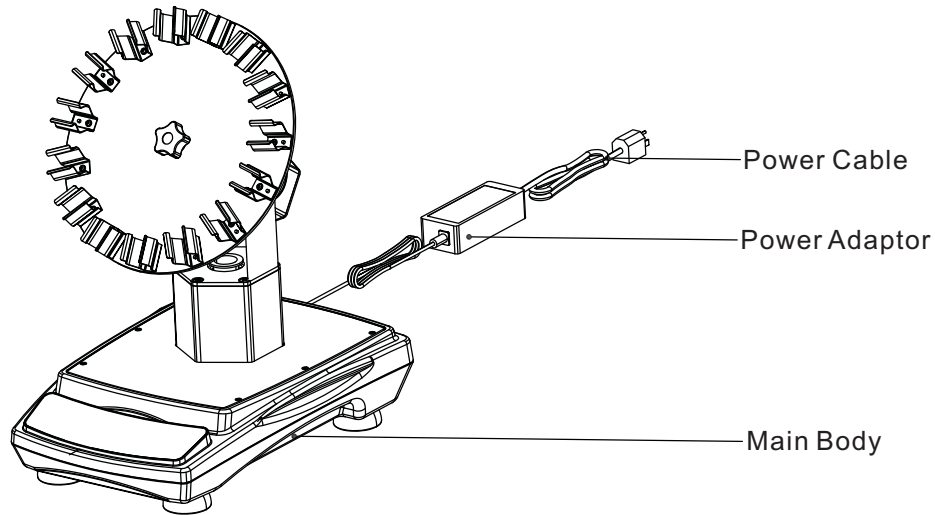

1.3 Structure Diagram

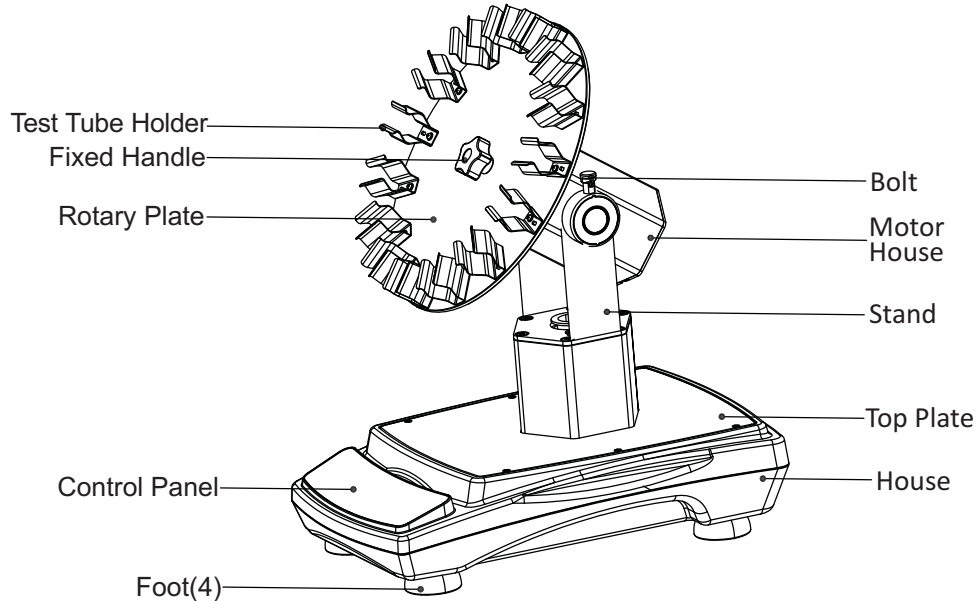

Section 2 Overview

2.1 Specifications

Speed

Speed Range	5~80rpm
Display Accuracy	1rpm
Speed Accuracy	±1rpm
Display	LED

Note: Maximum speed may vary with heavy or unbalanced loads.

Angle

Angle Adjustment Range	95° ~180°
Step Angle	17°
Adjustment Method	Manual Adjustment

Load

Maximum Capacity	64×Φ18.5mm culture tubes (drum rotor)
Maximum Load (Centered on tray)	5 kg (Clamps Included)

Time

Range	0, 1 min. ~ 99h59min.
Display	LED

Size

Overall Dimensions	415×293×464mm(16.3×11.5×18.3in)
Standard Match Dimensions	15×50ml rotary plate (Diameter Φ290mm/11.4 in)
Package Dimensions	536×502×505mm(21.1×19.8×19.9in)

Weight

Net Weight	9.4Kg(20.7lb)
Gross Weight	13.2Kg(29.0lbs)

Power Supply

Power Supply	AC100~240V, 50Hz/60Hz, 0.2A
--------------	-----------------------------

Others

Certificate	ROHS, WEEE, cCSAus, CE Mark
-------------	-----------------------------

2.2 Environmental Conditions

Application Environmental Conditions: indoor use


Temperature	5 to 40 °C (41 to 104 °F)
Humidity	20% to 85%,
Altitude	≤2,000 m
Voltage Fluctuation	±10% of the nominal voltage

Storage Environmental Conditions

Temperature	0 to 60 °C (36 to 124 °F)
Humidity	20% to 90%, non-condensing

2.3 Safety Instructions

Please read the entire instruction manual before operating the Multi-Purpose Tube Rotator.

 **WARNING DO NOT** use the Multi-Purpose Tube Rotator in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if accessories used are not provided or recommended by the manufacturer, or are used in a manner not specified by the manufacturer.

CAUTION ! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet.

Disconnect unit from the power supply prior to maintenance and service. Any spills should be removed promptly. Biohazard spills should be cleaned using approved liquid promptly. Solvent spills are a fire hazard. Stop the unit immediately, and **DO NOT** operate until clean up is complete and vapors have dissipated.

DO NOT immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.


Note: The rotary plate maybe have a little vibration below 11 rpm, but the vibration will relieve or disappear when add some loads.

Various independent clamps are available with this Multi-Purpose Tube Rotator for convenient combination use. It is applicable for 1.5 to 50 ml microtubes to mix liquid efficiently and gently. It has an angle range of 95 to 180 degrees and can maintain suspension state of biological samples such as blood mixing. Applications include blood coagulation prevention, emulsion diagnosis, immunoprecipitation, and other related areas.

Section 3 Operation

This Chapter covers the control panel and its operation.

3.1 Operations of the Control Panel

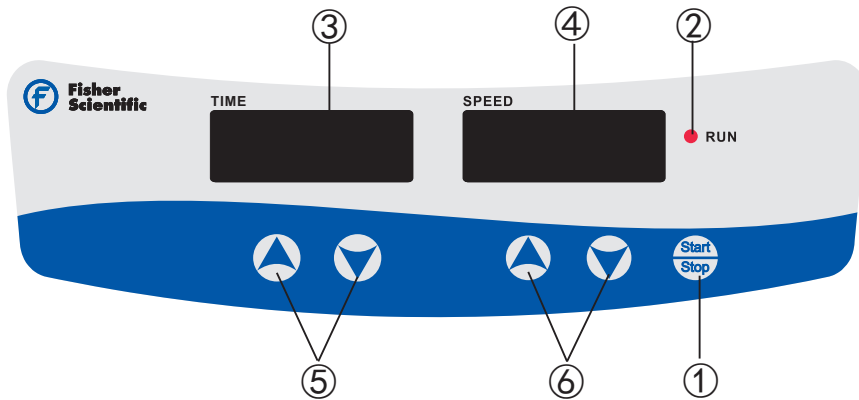

The front panel of the Multi-Purpose Tube Rotator contains all the controls needed to operate the unit.

1. Start/Stop button: Start or stop the instrument.
2. RUN indicator light: The light is on when the instrument is running and off when the instrument is in standby.
3. TIME display window: The window shows cumulative time (in continuous mode) or remaining time (in timer mode). The range of time displayed is 0 to 99 hours and 59 minutes. The accuracy is 1 minute.
4. SPEED display window: The window shows set speed (when the instrument is in standby) or current speed (when the instrument is running).
5. Set Time Buttons: UP/DOWN Arrow buttons are used to increase/decrease the set time of the instrument.
6. Set Speed Buttons: UP/DOWN Arrow buttons are used to increase/decrease the set speed of the instrument.

Preparation

1. Connect all the components according to the figures shown on page 1-3 of this manual. Use grounded power outlet.
2. Press power button on the back right side of the instrument and put it into the "I" state and then the instrument is in standby.

3.2 Settings

Time Settings

1. Continuous mode

Press the "▲" or "▼" arrow button below the TIME display window. When the number shown on the display window starts flashing, press "▼" arrow button to decrease the time to 00:00 and then release the button. The time setting is finished after the number shown on the display window has flashed twice.

2. Timer mode

Press the "▲" or "▼" arrow button below the TIME display window. When the number shown on the display window starts flashing, press "▲" or "▼" arrow button to increase or decrease the time value. Release the button when the time shown on the display window reaches the set value. The time setting is finished after the number shown on the display window has flashed twice.

Speed Settings

Press the "▲" or "▼" arrow button below the SPEED display window. When the number shown on the display window starts flashing, press "▲" or "▼" arrow button to increase or decrease the speed value. Release the button when the speed shown on the display window reaches the set value. The speed setting is finished after the number shown on the display window has flashed twice.



Note: press the "▲" or "▼" arrow button for a longer time to accelerate the setting.

Run and Stop

1. Continuous Mode

Press "Start/Stop" button and the instrument will start running with the specified settings and the RUN indicator light will be on. The TIME display window will show the cumulative time and the SPEED display window will show the current speed. Press "Start/Stop" button again and the instrument will slow down until it stops. The instrument will then be in standby and the two display windows will show the set values.

2. Timer Mode

Press “” button and the instrument will start running with the specified settings and the RUN indicator light will be on. The TIME display window will show the remaining time and the SPEED display window will show the current speed. Press “” button again and the instrument will slow down until it stops. The instrument will then be in standby and the two display windows will show the set values.

Finish Operation

After the operation is finished, please press the power switch at the back right side the instrument and put it into the “O” state. Unplug the instrument and store the instrument according to the storage guide.

Note:

To ensure shaking operation smooth and steady, it may take 1 minute for the microprocessor control system to accelerate the tray to the set speed.

Alarm System

In Timer mode, the instrument will stop operation automatically and the alarm will sound when the timer goes off.


Power Recovery

If the power supply is cut off suddenly while the instrument is in operation, the unit will automatically run at the previously set parameter upon power restoration. The display window will flash. Press any button to stop flashing.

3.3 Tray Installation and Tilt Angle Adjustment

Angle Adjustment Method


1. Pull out the black bolt

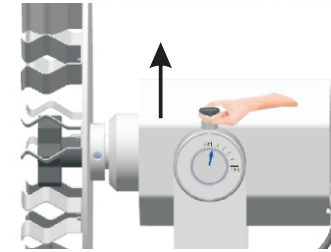

Pull bolt upward

Note: If bolt does not pull out, please operate according to the following instructions

- 1) Turn motor house back and forth gently by hand and pull bolt upward at the same time.
 - 2) If still cannot pull out bolt, please use vise grip plier to pull bolt upward and turn motor house at the same time.
 - 3) Please contact us immediately if the problem still exists.
2. Rotate the motor house gently by hand to adjust the angle.

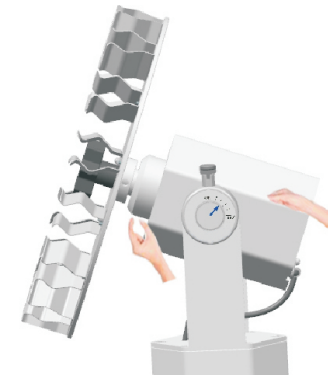

3. Pull out bolt and turn 90 degrees and insert the top of the bolt into the long slot of the hexagon block. Push the motor house gently until the bolt fits well.

Section 3
Operation



4. Angle adjustment range is shown in the figures below.



Mixing angle 95°



Mixing angle 135°



Mixing angle 180°

Single Tray Installation

1. Insert the stud on the rotation shaft into the center hole of the tray and make sure the tray is firmly touching the fixing block of the rotary plate, as shown in figure 1 (indicate on next page 3-6).
2. Screw in the fixed handle until the tray fastens, as shown in figure 2 (indicate on next page 3-6).
3. Align the tip of the screw driver with the hole at the side of the fixing block of the rotary plate. Make sure the fixing block of the rotary plate is not moving and continue screwing the fixing handle until the tray is fastened tightly (shown in figure 3 which indicate on next page 3-6).

Section 3
Operation

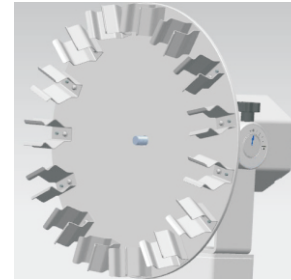


figure 1.

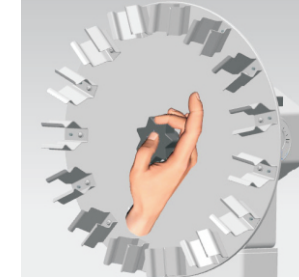


figure 2.



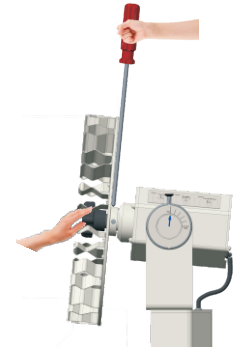


figure 3.

4. Insert the test tube from the outside of the tube holder until it reaches the appropriate position.

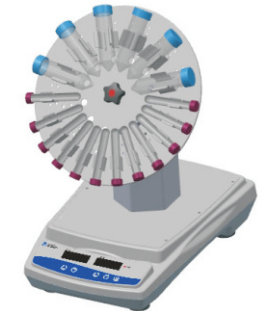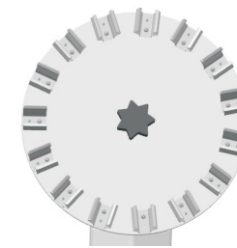

Double Tray Installation

1. Insert the stud on the rotation shaft into the center hole of the tray and make sure the tray is firmly touching the fixing block of the rotary plate, as shown in figure 4.
2. Insert the special-purpose handle into the center hole and screw in the stud until the tray is fastened (shown in figure 5).

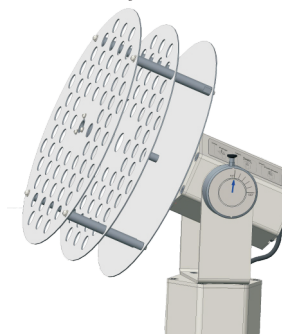


figure 4

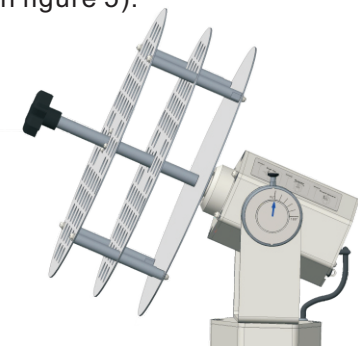


figure 5

3. Fasten the screw on the tray with the screw driver to keep the tray is fixed well, as shown in figure 6.


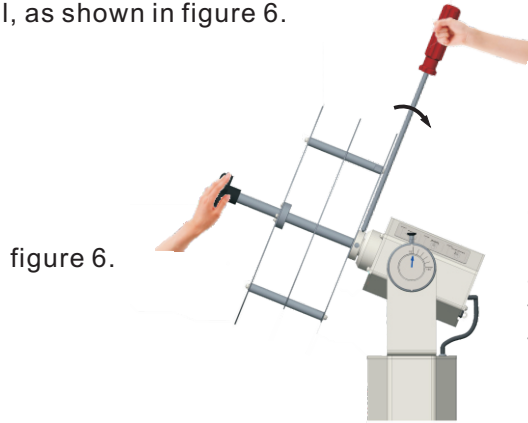


figure 6.

Caution

Failure to load or unload the tray as listed above may result in uneven tray rotation or possible stalling of the tray

4. Insert the test tube into the hole. Make sure to place the test tubes close to the center or symmetric around the axis.

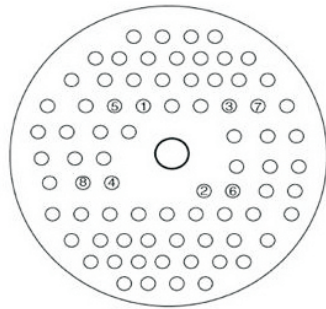

Tray, Front Facing

Note

After continued and extended operation, it is possible for the tray to exhibit uneven or erratic rotation. This does not necessarily indicate equipment malfunction and a cause for returning the unit to the manufacturer.

**Loading and Unloading the Tray:
Maintaining Tray Balance**

Because the rotator is a finely balance unit, loading and unloading the tray with culture tubes or other containers needs to maintain the overall balance of the tray, especially when it is moving.

Starting from the center and working to the outside edge of the tray, insert or remove tubes that are diametrically opposed—that is, from the top, bottom, left, right. This way the container weight is distributed evenly, allowing optimum tray rotation.

Section 4 Safety Tips and Maintenance

Safety Tips

1. Use independent power supply.
2. Check if the local power supply voltage is suitable for use.
3. Do not drag the power supply cable when unplugging.
4. Do not use non-specified power cable nor damage cable.
5. Service should only be performed by a qualified professional.
6. The power supply must be unplugged under the following situations:
 - a. when the unit is moved
 - b. when the electrical cabinet or the moving component is opened
 - c. when the equipment is malfunctioning
 - d. when the equipment is not in use
7. Wear appropriate protective equipment during operation to prevent spillage of liquid or leakage of flammable or poisonous gas.

- Reagent must be sealed inside the test tube during operation to avoid spillage while rotating.
- Avoid vibration ensure precise operation of the instrument.
- Balance loads if possible to improve service life of the instrument. (It is recommended to reduce load or rotation speed if the loads are unbalanced.)
- Make sure to pull out the test tube horizontally from the holder to avoid side flip of the instrument.
- Fixed handle must be fastened to prevent rotary plate from falling during operation.

Maintenance

- a. This instrument uses a high-accuracy motor. It is maintenance free and has a long service time, high quality, and low noise level.
- b. Surfaces can be cleaned with a mild detergent and water.

Warning: Avoid dripping detergent or water into the inside of the instrument during cleaning.

Clean Spill

If accidental spillage of liquids caused by mishandling or contained break-age occurs on the surface of the instrument, please shut down the instrument and clean up the liquid immediately.

If the liquid has already spilled into the unit, please cut off the power supply first and immediately clean up the liquid at the surface of the instrument. Then place the instrument in a ventilated and dry environment for 24 hours before reuse. If the instrument is still not functioning after drying for 24 hours, please contact the manufacturer.

Warning: Disassembling/Assembling without a qualified professional's guidance may cause malfunctioning of the instrument.




Section 5 Troubleshooting

Please refer to the following table to troubleshoot if any malfunction occurs. If the problem still exists, contact your local sales representative.

Error	Cause	Solution
Cannot start machine, LED display window off	Power disconnected	Connect the power
	Power adaptor failure	Replace power adaptor
	Switch off	Switch on
No shaking of the tray	Electrical malfunction	Contact Fisher Scientific
	Mechanical malfunction	Contact Fisher Scientific
	Over-weighted or unbalanced load	Adjust the weight and position of load, decrease rotation speed
Loud noise	Component loose	Fasten screws
	Rotary plate loose	Fasten fixed handle
	Uneven or resonant surface	Place instrument at a level and firm surface
Other	Keep record for maintenance	

Section 6 Optional Accessories/Spare Parts

Description	Cat. No.	Dimensions	Figure
50x1.5/2.0ml rotator drum	88861121	50×1.5/2ml	
30x15ml rotator drum	88861122	30×15ml	
15x50ml rotator drum	88861123	15×50ml	
10x10/15ml + 20x5/7ml drum	88861124	10×15ml 20×5ml	
64x14ml rotator drum	88861125	64×Φ14mm	
64x18.5ml rotator drum	88861126	64×Φ18.5mm	
General Power Adaptor	88861128	AC 100~240V, 50/60HZ	
US Plug	88861129	250VAC, 10A, 1.8m	

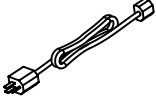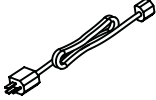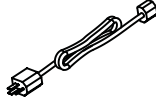Description	Cat. No.	Dimensions	Figure
CN Plug	88861130	250VAC, 10A, 1.8m	
EU Plug	88861131	250VAC, 10A, 1.8m	
UK Plug	88861132	250VAC, 10A, 1.8m	

Section 7 Warranty

When used in laboratory conditions and according to these operation instructions and maintenance, this product is warranted for 24 months against defective materials or workmanship. The 24 month warranty period begins from the delivery date of this product.

For product quality or performance issues, contact Fisher Scientific Customer Service.