

Model 310 Series

Forma Direct Heat CO₂ Incubator

Operating and Maintenance Manual 7010310 Rev. 9



Model	CO2 Sensor*	Voltage**
310	T/C	115
311	T/C	230
320	IR	115
321	IR	230
350†	T/C	115
351†	T/C	230
360†	IR	115
361†	IR	230

* T/C is a thermal conductivity sensor.

IR is an infrared sensor.

** All units are 50/60 Hz.

† Copper interior

MANUAL NUMBER 7010310

--	--	7/13/10	Re-installed "Specs are based on..." on Specification page. Unintentional removal	ccs
9	25584	8/6/09	Updated RH low alarm ringback time from 30 to 15 minutes	ccs
8	25248/IN-3945	6/9/09	Updated drawer specs - sliding and stationary	ccs
7	23793/IN-3916	5/29/09	Updated 310-201-0-D - o-ring 101004	ccs
--	24891/SI-9962	5/29/09	Replaced 190793 display board with 190609 (color)	ccs
--	25499/IN-3936	5/29/09	Updated 310-203-0-D - bezel part number 190786	ccs
6	25303/IN-3902	2/16/09	UL decom warning pg 1-6	ccs
5	25210/IN-3885	1/6/09	Updated stacking instructions (screws), pg 1-5	ccs



Important Read this instruction manual. Failure to read, understand and follow the instructions in this manual may result in damage to the unit, injury to operating personnel, and poor equipment performance. ▲

Warning If this unit is not used in the manner specified in this operating manual, the protection provided by the equipment design may be impaired. ▲

Caution All internal adjustments and maintenance must be performed by qualified service personnel. ▲

Material in this manual is for information purposes only. The contents and the product it describes are subject to change without notice. Thermo Fisher Scientific makes no representations or warranties with respect to this manual. In no event shall Thermo be held liable for any damages, direct or incidental, arising out of or related to the use of this manual.

©2003 Thermo Fisher Scientific. All rights reserved.



Important operating and/or maintenance instructions. Read the accompanying text carefully.



Potential electrical hazards. Only qualified persons should perform procedures associated with this symbol.



Equipment being maintained or serviced must be turned off and locked off to prevent possible injury.



Hot surface(s) present which may cause burns to unprotected skin, or to materials which may be damaged by elevated temperatures.



Marking of electrical and electronic equipment, which applies to electrical and electronic equipment falling under the Directive 2002/96/EC (WEEE) and the equipment that has been put on the market after 13 August 2005.



This product is required to comply with the European Union's Waste Electrical & Electronic Equipment (WEEE) Directive 2002/96/EC. It is marked with the WEEE symbol. Thermo Fisher Scientific has contracted with one or more recycling/disposal companies in each EU Member State European Country, and this product should be disposed of or recycled through them. Further information on Thermo's compliance with this directive, the recyclers in your country and information on Thermo products will be available at www.thermo.com.

- ✓ Always use the proper protective equipment (clothing, gloves, goggles, etc.)
- ✓ Always dissipate extreme cold or heat and wear protective clothing.
- ✓ Always follow good hygiene practices.
- ✓ Each individual is responsible for his or her own safety.

Do You Need Information or Assistance on Thermo Scientific Products?

If you do, please contact us 8:00 a.m. to 6:00 p.m. (Eastern Time) at:

1-740-373-4763

Direct

1-800-438-4851

Toll Free, U.S. and Canada

1-877-213-8051

FAX

<http://www.thermoscientific.com>

Internet Worldwide Web Home Page

service.led.marietta@thermofisher.com

Service E-Mail Address

Our **Sales Support** staff can provide information on pricing and give you quotations. We can take your order and provide delivery information on major equipment items or make arrangements to have your local sales representative contact you. Our products are listed on the Internet and we can be contacted through our Internet home page.

Our **Service Support** staff can supply technical information about proper setup, operation or troubleshooting of your equipment. We can fill your needs for spare or replacement parts or provide you with on-site service. We can also provide you with a quotation on our Extended Warranty for your Thermo Scientific products.

Whatever Thermo Scientific products you need or use, we will be happy to discuss your applications. If you are experiencing technical problems, working together, we will help you locate the problem and, chances are, correct it yourself...over the telephone without a service call.

When more extensive service is necessary, we will assist you with direct factory trained technicians or a qualified service organization for on-the-spot repair. If your service need is covered by the warranty, we will arrange for the unit to be repaired at our expense and to your satisfaction.

Regardless of your needs, our professional telephone technicians are available to assist you Monday through Friday from 8:00 a.m. to 6:00 p.m. Eastern Time. Please contact us by telephone or fax. If you wish to write, our mailing address is:

Thermo Fisher Scientific
401 Millcreek Road, Box 649
Marietta, OH 45750

International customers, please contact your local Thermo Scientific distributor.

Warranty Notes

Information You Should Know Before Requesting Warranty Service

- **Locate the model and serial numbers.** A serial tag is located on the unit itself.
- For equipment service or maintenance, or with technical or special application inquiries, contact Technical Services at 1-800-438-4851 or 1-740-373-4763 (USA and Canada). Outside the USA, contact your local distributor.

Repairs NOT Covered Under Warranty

- **Calibration of control parameters.** Nominal calibrations are performed at the factory; typically $\pm 1^{\circ}\text{C}$ for temperature, $\pm 1\%$ for gases, and $\pm 5\%$ for humidity. Our service personnel can provide precise calibrations as a billable service at your location. Calibration after a warranty repair is covered under the warranty.
- **Damage resulting from use of improper quality water, chemicals or cleaning agents detrimental to equipment materials.**
- **Service calls for improper installation or operating instructions.** Corrections to any of the following are billable services:
 - 1) electrical service connection
 - 2) tubing connections
 - 3) gas regulators
 - 4) gas tanks
 - 5) unit leveling
 - 6) room ventilation
 - 7) adverse ambient temperature fluctuations
 - 8) any repair external to the unit
- **Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, or improper installation.**
- **Repairs to parts or systems resulting from unauthorized unit modifications.**
- **Any labor costs other than that specified during the parts and labor warranty period, which may include additional warranty on CO₂ sensors, blower motors, water jackets, etc.**

Table of Contents

Section 1	Installation and Start-Up	1-1
	Control Panel Keys, Displays & Indicators	1-2
	Keypad Operation	1-3
	Displays	1-4
	Installing the Incubator	1-4
	Incubator Start-Up	1-11
Section 2	Calibration	2-1
	Calibrating the Temperature	2-1
	Calibrating the T/C CO ₂ System	2-2
	Calibrating the Infrared CO ₂ System	2-3
	Calibrating Relative Humidity	2-4
Section 3	Configuration	3-1
	Turn the Audible Alarm ON/OFF	3-1
	Set an Access Code	3-1
	Set a Low Temp Alarm Limit	3-2
	Enable the Low Temp Alarm to Trip Contacts	3-2
	Set a Low CO ₂ Alarm Limit	3-3
	Setting a High CO ₂ Alarm Limit	3-3
	Enable CO ₂ Alarms to Trip Contacts	3-4
	Set New Zero Number for T/C CO ₂ Sensors	3-4
	Set New Span Number for T/C CO ₂ Sensors	3-5
	Set a Low RH Alarm Limit	3-5
	Enable RH Alarms to Trip Contacts	3-6
	Enable Temp/RH to be Displayed	3-6
	Select Primary Tank w/ Gas Guard Option	3-7
	Disable the Gas Guard System	3-7
	Communications Address for RS485	3-8
Section 4	Alarms	4-1
	Temp Control Failure Alarm TMP CNTR ERR	4-2
	Sensor Fault Alarms	4-2
	CO ₂ SNSR ERR	4-2
	IR AUTOZ ERR	4-2

Section 5	Routine Maintenance	5-1
	Disinfecting the Incubator Interior	5-1
	Cleaning the Cabinet Exterior	5-2
	Cleaning the Glass Doors	5-2
	Cleaning the Humidity Pan	5-3
	Reversing the Door Swing	5-3
	HEPA Filter (Factory Installed Option)	5-8
	Replacing the Power Fuses	5-8
	Major Components	5-9
	Electronics Section	5-9
	Replacing the Sample Air Filter	5-10
Section 6	Factory Options	6-1
	Connect Remote Alarm Contacts	6-1
	Connect RS485 Interface	6-2
	Connecting the Analog Output Boards	6-2
	CO ₂ Gas Guard	6-5
	Humidity Readout	6-6
	Uninterruptible Power Supply Connections	6-7
Section 7	Specifications	7-1
Section 8	Spare Parts	8-1
Section 9	Electrical Schematics	9-1
Section 10	Warranty Information	10-1

Section 1 Installation and Start-Up

Incubator components are described below.

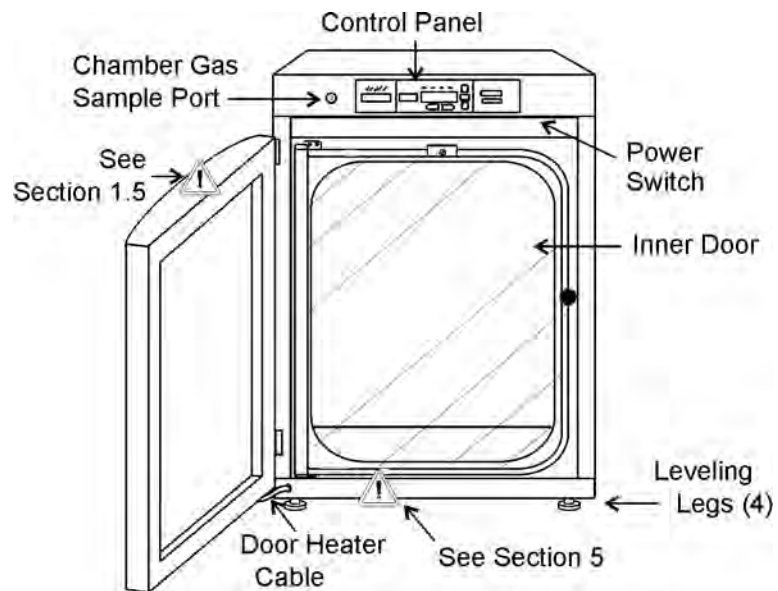


Figure 1-1. Model 310 Series Direct Heat Incubator

- Outer Door - Reversible to opposite swing, see Section 5
- Inner Door - Reversible to opposite swing, see Section 5
- Chamber Gas Sample Port - Used for sampling chamber CO₂ content using a FYRITE or similar instrument.
- Main Power Switch
- Control Panel - Keypad, Displays and indicators (See Figure 1-2)
- Leveling Legs - Used to level the unit

Note The incubators are stackable. Information follows. ▲

Control Panel Keys, Displays & Indicators

Silence - Silences the audible alarm.

Alarm Indicator - Light pulses on/off during an alarm condition in the cabinet.

Mode Select Switch - Used to select Run, Setpoints, Calibration and System Configuration Modes.

Message Center - Displays system status.

Mode Select Indicators -

Run: Run Menu

Set: Set Points Menu

Cal: Calibrate Menu

Config: Configuration Menu

Up and Down Arrows - Increases or decreases the number values, toggles between choices.

Enter - Stores the value into computer memory.

Heat Indicator - Lights when power is applied to the heaters.

Temp Display - Program to display temperature continuously, RH continuously (with RH option), or toggle between temperature and humidity (with RH option). See Section 3, Configuration.

Scroll for Parameters Arrows - Moves the operator through the choices of the selected mode.

CO₂ Inject Indicator - Lights when CO₂ is being injected into the incubator.

% CO₂ Display - Displays CO₂ percentage continuously.

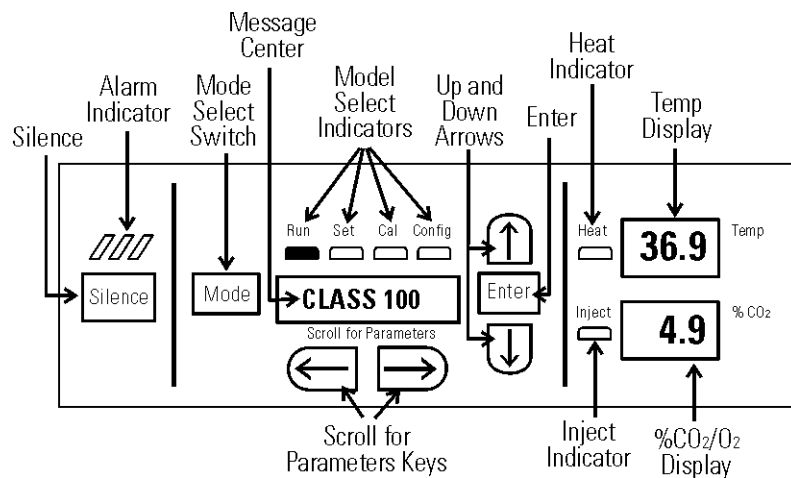


Figure 1-2. Control Panel

Keypad Operation

The Model 310 Series direct heat incubator has four basic modes which allow incubator setup: Run, Setpoints, Calibration and System Configuration.

Run is the default mode which the incubator will normally be in during operation.

Set is used to enter system setpoints for incubator operation.

Calibration is used to calibrate various system parameters.

Configuration allows for custom setup of various options.

The chart below shows the selections under each of the modes.

RUN	SETPOINT	CALIBRATION	CONFIGURATION
Default Mode	Temperature	Temp Offset	Audible
	Overtemp	CO₂ Cal¹	Access Code
	CO₂	IR Cal²	Temp Lo Alarm
		<i>RH Cal</i>	Temp Relay
			CO₂ Lo Alarm
			CO₂ Hi Alarm
			CO₂ Relay
			CO₂ Z & S #'s *
			<i>RH Lo Alarm</i>
			<i>RH Relay</i>
			<i>Display Temp</i>
			<i>Display Rh</i>
			<i>Tank Select</i>
			<i>Gas Guard</i>
			RS485 Address

¹ T/C units only

² IR units only

*T/C units only

Base Unit Displays

Option Displays

Scroll for Parameters Arrows: Steps the operator through the parameters of SET, CAL and CONFIG Modes. The right arrow goes to the next parameter, the left arrow returns to the previous parameter.

Up Arrow: Increases or toggles the parameter value that has been selected in the SET, CAL, and CONFIG Modes.

Enter: Must press Enter key to save to memory all changed values.

Down Arrow: Decreases or toggles the parameter values that have been selected in the SET, CAL and CONFIG Modes.

Silence Key: Press to silence the audible alarm. See Section 4 for alarm ringback times.

Displays

Message Center: Displays the system status (Mode) at all times. Displays SYSTEM OK during normal operation, or alarm messages if the system detects an alarm condition. See Section 4, Alarms.

Upper and Lower Displays: These 7 segment displays vary depending upon the options present and the configuration chosen. The upper display can display temp or RH, or toggle between them. The bottom display shows CO₂ continuously.

Installing the Incubator

Warning Single and stacked units must be installed against a wall or similar structure. Maintain a three-inch clearance behind the incubator for electrical and gas hook-ups. ▲

1. Locate the unit on a firm level surface capable of supporting the unit's weight of 205 lbs.
2. Locate the unit away from doors and windows and heating and air conditioning ducts.
3. Lift the unit only by the sides of the cabinet base . Do not attempt to lift it by the front and back . This places stress on the outer door hinges.

Stacking the Incubators

Caution When stacking incubators, the direct heat incubator must be the top unit. Never stack a water-jacketed incubator on top of a Model 310 Series unit. ▲

Warning With incubators in a stacked configuration, do not leave both exterior doors open at the same time. ▲

Warning If the units have been in operation, shut them both off and disconnect from the power source before beginning any service work. ▲

Note Two stacking brackets (shown at right) are included in the parts bag shipped with each incubator. ▲

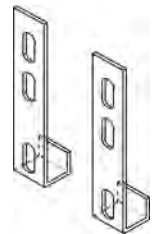


Figure 1-3.
Stacking Brackets

Stacking (continued)

1. Remove and discard the slotted head screws on the top of the bottom incubator (Figure 1-4). Replace with 5/16 HH bolts and flatwashers included in the stacking kit. Do not tighten.

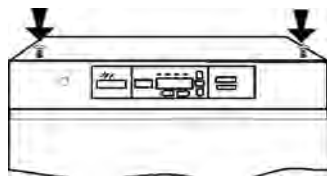


Figure 1-4. Slotted Head Screws

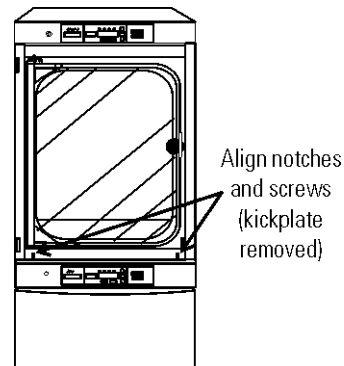


Figure 1-5. Align

2. Remove the kickplate from below the door on the top incubator (Figure 1-5). Two Phillips screws secure the plate. Note the notches at the base of the incubator frame.
3. Disconnect the door cable from the connector behind the plate. Lift the door off the hinge pins and set it aside. The hinge pins are not attached to the brackets and may fall out when the door is lifted off.
4. Unscrew and remove the leveling feet from the top unit and lift it onto the bottom unit, offsetting the base of the top unit approximately 2-3 inches behind the screws identified by the arrows in Figure 1-4.

Warning This incubator weighs 205 lbs. Have sufficient personnel available when lifting. Lift the unit only by the sides of the cabinet base. Do not attempt to lift it by the front and back as this places stress on the outer door hinge. ▲

5. Align the sides of the top unit with the bottom unit and slide the top unit forward until the notches in the base of the top unit align with the bolts in the top of the bottom unit. The flatwashers should slide over the base frame. Tighten bolts.
6. Remove the four nylon plugs on the lower back of the upper incubator.
7. Insert the stacking brackets into the slots on the back of the control panel of the bottom unit as shown in Figure 1-6.

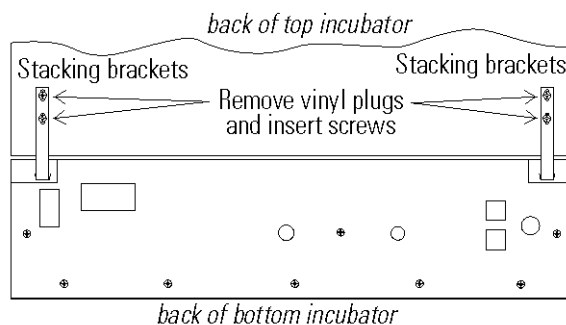


Figure 1-6. Brackets Installed

Stacking (continued)

- Align the slots in the brackets with the mounting holes on the back of the top incubator. Secure the brackets with the screws provided in the stacking kit.
- Make sure the hinge pins are in the hinge brackets. Slide the door onto the hinge pins.
- Secure the door heater cable to the connector below the door and replace the kick plate on the top unit, taking care not to crimp cable.
- The stacked incubators are ready for service.

Preliminary Cleaning and Disinfecting

Remove vinyl from shelf channels, duct sheets, and air duct, if present. Using a suitable laboratory disinfectant, disinfect all interior surfaces including shelves and shelf supports, door gaskets, blower wheel and CO₂ sensor. Refer to Section 5.

Caution Before using any cleaning or decontamination method except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment. Accidental spills of hazardous materials on or inside this unit are the responsibility of the user. ▲

Installing the Shelves

- Install the large sheet metal shelf supports with the tabs facing into the center of the chamber with their slots up. There are no right side or left side supports. Simply turn one of them to fit the opposite side. Tilt the shelf supports as they are placed in the chamber so the tops fit into the top air duct, then guide them into the vertical position. Figure 1-7 shows the support as it would be oriented for right side of chamber.

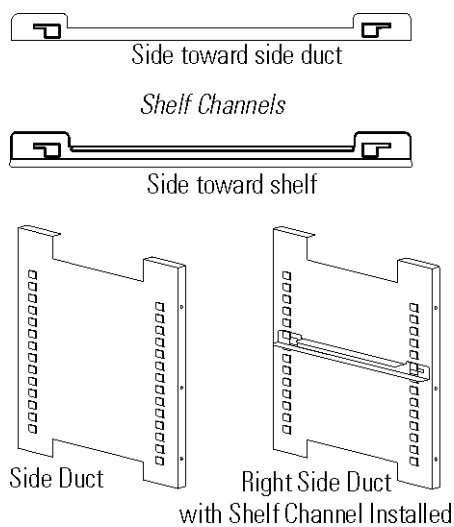


Figure 1-7. Shelf Channels and Ductwork

Installing the Shelves (continued)

2. Note there is no difference in the left side and right side shelf channels.
3. Install the shelf channels by placing the channel's rear slot over the appropriate rear tab on the shelf support. Pull the shelf channel forward and engage the channel's front slot into the shelf support's appropriate forward tab. Refer to Figure 1-8.

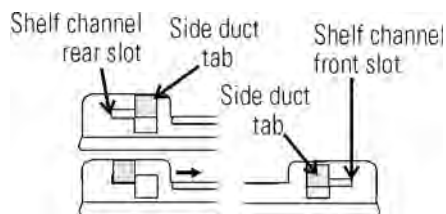


Figure 1-8. Channel Ends

Installing Access Port Filter

Locate the opening in the top left corner on the inside of the chamber. Remove the tape covering the opening on the outside of the unit. Locate the stopper with filter in the hardware bag. Install the assembly in the opening inside the chamber. See Figure 1-9.

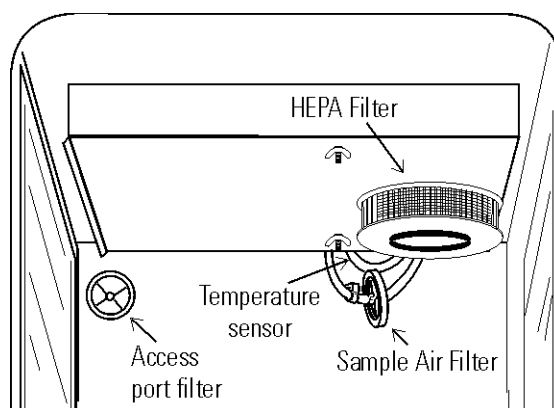


Figure 1-9. Filter and Sensor Locations

Installing (optional) HEPA Filter

Caution Be careful when handling the filter. The media can be damaged if it is mishandled.

To avoid damage to the incubator, do not operate the unit without the HEPA filter in place. ▲

1. Remove the filter from the shipping box.
2. Remove the plastic coating from the filter, using caution not to touch the filter media.
3. Install the filter as shown in Figure 1-9. Refer to Section 5 for HEPA filter maintenance.

Installing the Air Sample Filter

1. Remove the filter from the shipping bag.
2. Install the air sample filter assembly to the black hose barb behind the top duct.
3. Insert the other end of the filter assembly into the larger hole in the air intake to the blower. Refer to Figure 1-10.
4. Verify that the temperature sensor is installed into the smaller hole in the intake, as in Figure 1-10.

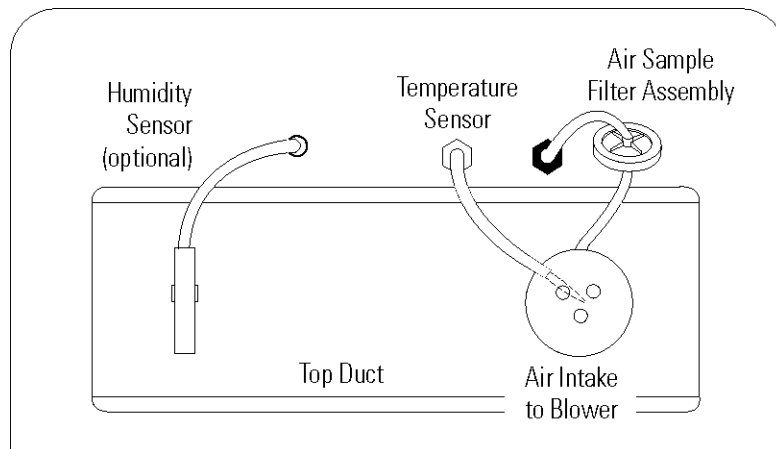


Figure 1-10. View of Chamber Ceiling

Leveling the Unit

Check for level by placing a bubble style level on one of the shelves. Turn the hex nut on the leveler counterclockwise to lengthen the leg or clockwise to shorten it. Level the unit front-to-back and left-to-right.

Connecting the Unit to Electrical Power

See the serial tag on the side of the unit for electrical specifications or refer to the electrical schematics at the end of this manual.

Caution Connect the incubator to a grounded, dedicated circuit. The power cord connector is the mains disconnect device for the incubator. Position the incubator so the unit can be easily disconnected. ▲

Plug the provided power cord into the power inlet connector on the back of the cabinet (See Figure 1-11), then into the grounded, dedicated, electrical circuit.

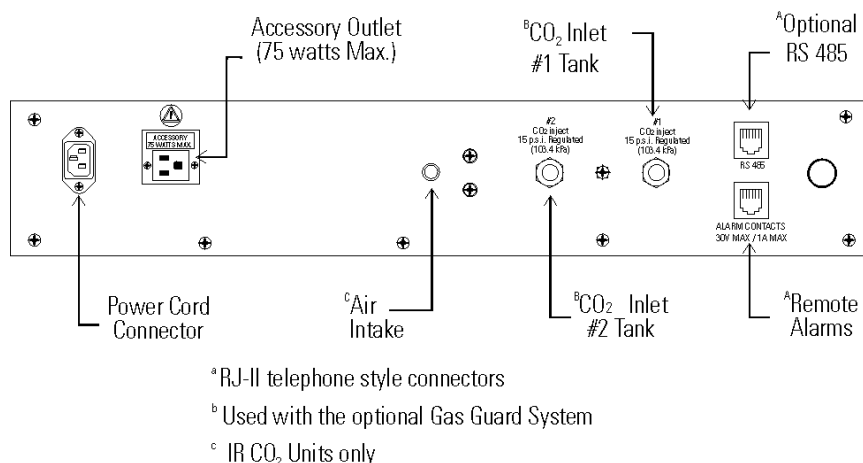


Figure 1-11. Connections on Unit Back

Filling the Humidity Pan

Fill the humidity pan with sterile pure water to within 1/2" of the top. Place the pan directly in the center of the incubator floor to ensure best humidity and temperature response.

For applications requiring abnormally high humidity, the pan should be placed against the left side wall of the incubator. The ductwork has been modified to allow the pan to be placed against the wall of the incubator. In some ambient conditions, this may cause condensation to form in the chamber.

For best operation of the incubator, sterilized distilled, demineralized or de-ionized water should be used in the humidity pan. Water purity should be in the resistance range of 50K to 1M Ohm/cm, or a conductivity range of 20.0 to 1.0 uS/cm. Refer to ASTM Standard D5391-93 or D4195-88 for measuring water purity.

Distillation systems, as well as some types of reverse osmosis water purity systems, can produce water in the quality range specified. Tap water is not recommended as it may contain chlorine, which can deteriorate the stainless steel. Tap water may also have a high mineral content, which would produce a build-up of scale in the pan. High purity or ultra pure water is not recommended as it is an extremely aggressive solvent and will deteriorate the stainless steel. High purity water has a resistance of above 1M to 18M Ohm. Even high purity water can contain bacteria and organic contaminants. Water should always be sterilized or treated with a decontaminant, safe for use with stainless steel as well as safe for the product, prior to being introduced into the humidity pan.

Caution Distilled or de-ionized water used in the humidity pan must be within a water quality resistance range of 50K to 1M Ohm/cm to protect and prolong the life of the stainless steel. Use of chlorinated water, or decontamination products containing chlorine, will deteriorate the stainless steel and cause rust, voiding the warranty. ▲

Check the level and change the water frequently to avoid contamination. Do not allow the water level to fluctuate significantly. “Dry-outs” will have an adverse effect on the humidity level and CO₂ calibration of the T/C units.

Connecting the CO₂ Gas Supply

Warning High concentrations of CO₂ gas can cause asphyxiation! OSHA Standards specify that employee exposure to carbon dioxide in any eight-hour shift of a 40-hour work week shall not exceed the eight-hour time weighted average of 5000 PPM (0.5% CO₂). The short term exposure limit for 15 minutes or less is 30,000 ppm (3% CO₂). Carbon dioxide monitors are recommended for confined areas where concentrations of carbon dioxide gas can accumulate. ▲

Warning This incubator is designed to be operated with CO₂ gas only. Connecting a flammable or toxic gas can result in a hazardous condition. Gases other than CO₂ should not be connected to this equipment. CO₂ gas cylinders have a UN1013 label on the cylinder and are equipped with a CGA 320 outlet valve. Check the gas cylinder for the proper identification labels. The CO₂ gas supply being connected to the incubator should be industrial grade, 99.5% pure. Do not use CO₂ gas cylinders equipped with siphon tubes. A siphon tube is used to extract liquid CO₂ from the cylinder which can damage the pressure regulator. Consult with your gas supplier to ensure that the CO₂ cylinder does not contain a siphon tube. Gas cylinders should also be secured to a wall or other stationary object to prevent them from tipping.

A two-stage CO₂ pressure regulator is required to be installed on the outlet valve of the gas cylinder. Input pressure to the incubator must be maintained at 15 psig (103.4 kPa), ±5 psig, for proper performance of the CO₂ control system. A single stage CO₂ pressure regulator will not maintain 15 psig (103.4 kPa) to the incubator as the pressure in the CO₂ cylinder decreases; therefore, a two stage regulator is recommended.

If higher purity CO₂ is desired inside the incubator (greater than 99.5% pure), the pressure regulator should be constructed with a stainless steel diaphragm, along with specifying the purity of the CO₂ from the gas supplier. Follow the manufacturer's instructions to ensure proper and safe installation of the pressure regulator on the gas cylinder.

Consult your facility safety officer to ensure that the equipment is installed in accordance with the codes and regulations applicable in your area. ▲

Connecting the CO₂ gas Supply (continued)

The CO₂ gas supply being connected should be industrial grade 99.5% pure and should not contain siphon tubes. Install a two-stage pressure regulator at the cylinder outlet. The high pressure gauge at the tank should have 0-2000 psig range. The low pressure gauge, at the incubator inlet, should have a 0-30 psig range. Input pressure to the incubator must be maintained at 15 psig (103.4 kPa).

The incubator has serrated fittings on the back of the cabinet to connect the gas supply. Refer to Figure 1-11. The fitting is labeled CO₂ Inlet #1 Tank. Make sure that the connections are secured with clamps. Check all fittings for leaks.

For units having the CO₂ Gas Guard option, refer to Section 6.

Incubator Start-Up

With the incubator properly installed and connected to power, the humidity pan filled, and the unit connected to gas supplies, system setpoints can be entered. The following setpoints can be entered in Set mode: Temperature, Overtemperature and CO₂. To enter Set mode, press the Mode key until the Set indicator lights. Press the right and/or left arrow keys until the proper parameter appears in the message center. See Chart 1-1 for more detail.

Setting the Operating Temperature

Model 310 Series incubators have an operating temperature range of 10°C to 50°C, depending on ambient temperature. The incubator is shipped from the factory with a temperature setpoint of 10°C. At this setting, all heaters are turned off. To change the operating temperature setpoint:

1. Press the Mode key until the Set indicator lights.
2. Press the right arrow until “Temp XX.X” is displayed in the message center.
3. Press the up/down arrow key until the desired temperature setpoint is displayed.
4. Press Enter to save the setpoint.
5. Press the Mode key until the Run indicator lights for Run mode or press the right/left arrow keys to go to next/previous parameter.

Setting the Overtemp Setpoint

Caution The independent overtemp system is designed as a safety to protect the incubator only. It is not intended to protect or limit the maximum temperature of the cell cultures or customer's equipment inside the incubator if an overtemp condition occurs. ▲

Model 310 Series incubators are equipped with a secondary temperature monitoring system to monitor the air temperature inside the cabinet. This system is designed as a safety device to turn off all heaters in the event of a temperature control failure. Temperature control in the incubator will be $\pm 1^\circ$ of the overtemp setpoint.

The factory overtemperature setpoint (default) is 40°C . However, the overtemp can be set up to 55°C in 0.5° increments.

If the incubator's operating temperature setpoint is set above the overtemp setpoint, the overtemp setpoint will automatically update to 1°C above the temperature setpoint. It is recommended that the overtemp setpoint be maintained at 1°C over the operating temperature setpoint.

To set the Overtemp setpoint:

1. Press the Mode key until the Set indicator lights
2. Press the right arrow until Otemp XX.X is displayed in the message center
3. Press the up or down arrow key until the desired Overtemp setpoint is displayed
4. Press Enter to save the setting
5. Press the Mode key until the Run indicator lights or press the right or left arrow to go to the next or previous parameter.


Setting the CO₂ Setpoint

All T/C CO₂ cells are calibrated at the factory at 37°C, high humidity, and 10% CO₂. Therefore, if a temperature setpoint of 37°C has been entered, the humidity pan has been filled and the CO₂ control is to run between 0-10% with a T/C CO₂ sensor, the CO₂ setpoint may be entered immediately. Otherwise, it is important to allow the unit 12 hours to stabilize at the temperature setpoint before entering the CO₂ setpoint.

All models of the incubator have a CO₂ setpoint range of 0.0% to 20.0%. The incubator is shipped from the factory with a CO₂ setpoint of 0.0%. At this setting, all CO₂ control and alarms are turned off. To change the CO₂ setpoint:

1. Press the Mode key until the Set indicator lights.
2. Press the right arrow until “CO₂ XX.X” is displayed in the message center.
3. Press the up/down arrows until the desired CO₂ setpoint is displayed.
4. Press Enter to save the setpoint.
5. Press the Mode key until the Run indicator lights to go to Run mode or press the right/left arrow keys to go to next/previous parameter.

Chart 1-1

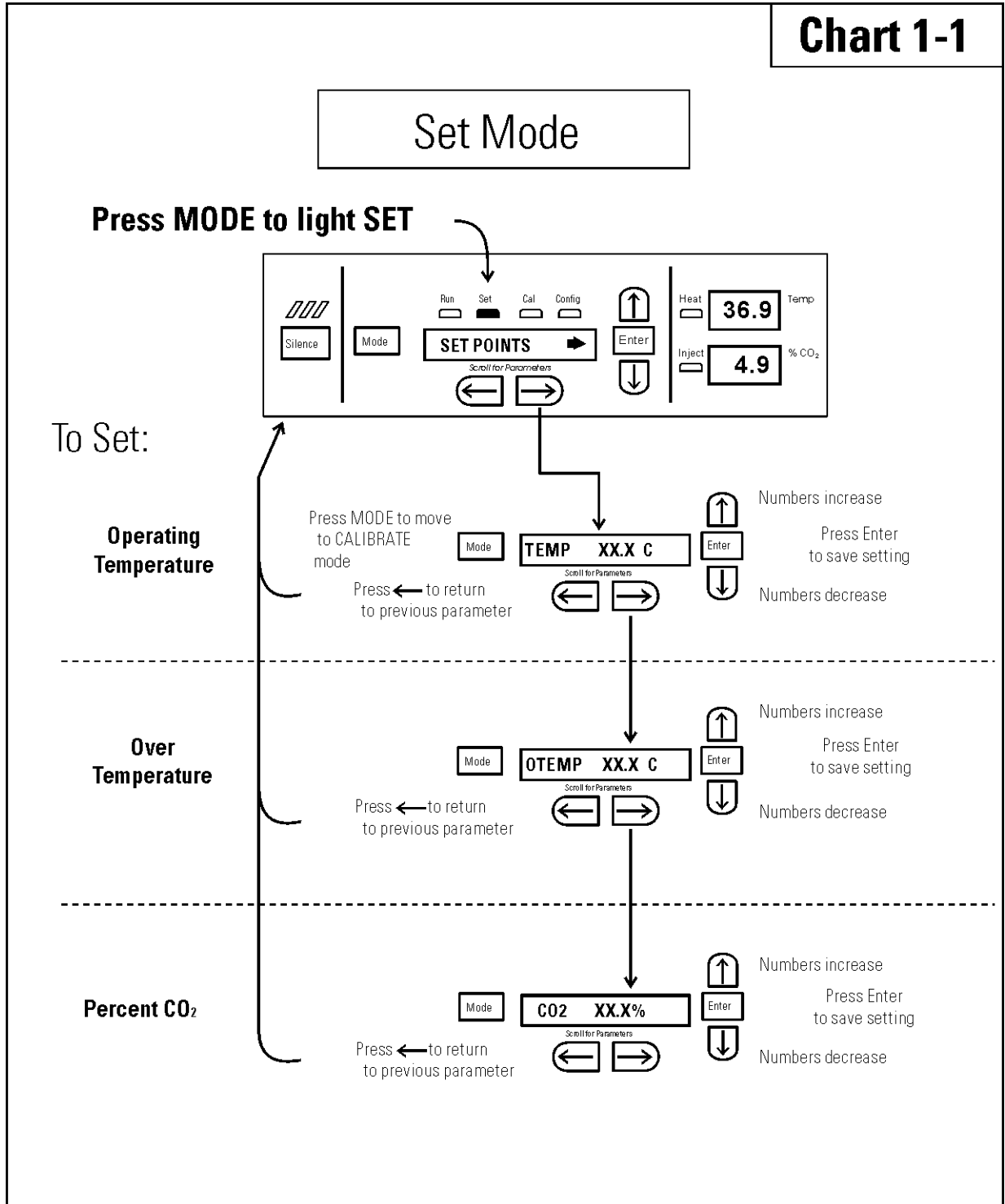

Section 2 Calibration

After the unit has stabilized, several different systems can be calibrated. In the calibration mode, the air temperature, CO₂ and RH levels can be calibrated to reference instruments. To access the calibration mode, press the Mode key until the Cal indicator lights. Press the right and/or left arrow until the proper parameter appears in the message center. See Chart 2-1 at the end of this section for more detail.

Calibration frequency is dependent on use, ambient conditions and accuracy required. A good laboratory practice would require at least an annual calibration check. On new installations, all parameters should be checked after the stabilization period.

Prior to calibration, the user should be aware of the following system functions. While the unit is in the calibration mode, all system control functions are stopped so the unit remains stable. Readout of the system being calibrated will appear on the message center. If no keys are pressed for approximately five minutes while in calibration mode, the system will reset to Run mode so control functions are reactivated.

Caution Before making any calibration or adjustments to the unit, it is imperative that all reference instruments be properly calibrated. ▲

Calibrating the Temperature

Before calibration, allow the cabinet temperature to stabilize. Place the calibrated instrument in the center of the chamber. The instrument should be in the airflow, not against the shelf.

Temperature Stabilization Periods

Startup - Allow 12 hours for the temperature in the cabinet to stabilize before proceeding.

Already Operating - Allow at least 2 hours after the display reaches setpoint for temperature to stabilize before proceeding.

Calibrating Temp (continued)

1. Press the Mode key until Cal indicator lights.
2. Press the right arrow until “TEMPCAL XX.X” appears in the message center.
3. Press up/down arrow to match display to calibrated instrument.
4. Press Enter to store calibration.
5. Press the Mode key to return to Run or the right/left arrow to go to next/previous parameter.

Calibrating the T/C CO₂ System

Models 310, 311, 350 and 351 have a thermal conductivity (T/C) CO₂ sensor. Thermal conductivity of the incubator atmosphere is not only effected by the quantity of CO₂ present, but also by the air temperature and water vapor present in the incubator atmosphere. In monitoring the effects of CO₂, air temperature and absolute humidity must be held constant so any change in thermal conductivity is caused by a change in CO₂ concentration.

Changing temperature or changing from elevated humidity to room ambient humidity levels will necessitate a re-calibration of the CO₂ control.

T/C CO₂ Sensor Stabilization Periods

Some T/C CO₂ sensors go through an aging period, especially on new installations. Calibration should be checked on a weekly basis, and adjusted as necessary. When stabilization occurs, checks may become less frequent.

Start -Up - The CO₂ sensor has been calibrated at the factory for 37°C and elevated humidity. Allow the temperature, humidity, and CO₂ levels in the chamber to stabilize at least 12 hours before checking the CO₂ concentration with an independent instrument.

Presently Operating - Make sure the chamber doors are closed. Allow at least 2 hours after the temperature and CO₂ displays reach their setpoints for chamber atmosphere stabilization.

Calibrating T/C CO₂ System (continued)

1. Make sure the stabilization periods outlined above are followed.
2. Sample the chamber atmosphere through the sample port with an independent instrument. Sample the atmosphere at least 3 times to ensure accuracy of the instrument.
3. Press the Mode key until the Cal indicator lights.
4. Press the right arrow until “CO₂ CAL XX.X” is displayed in the message center.
5. Press the up /down arrows to change the display to match the independent instrument.
6. Press Enter to store the calibration.
7. Press the Mode key to return to Run or the right or left arrows to go to the next/ previous parameter.

Calibrating the Infrared CO₂ System

Models 320, 321, 360 and 361 have an infrared (IR) CO₂ sensor. Infrared CO₂ sensors are not effected by chamber atmosphere temperature or humidity. However, the light detector in the sensor is effected by wide temperature changes. Therefore, changing temperature setpoints could necessitate a recalibration of the CO₂. Chamber temperature should be allowed to stabilize before checking CO₂ concentrations with an independent instrument, especially on start-up.

All models equipped with an IR/CO₂ sensor have an automatic calibration which occurs every 24 hours and lasts for 5 to 6 minutes. During automatic calibration, the CO₂ display is blanked out and HEPA filtered room air is pumped through the CO₂ sensor. A new CO₂ calibration value is stored in memory for use as the 0.0% CO₂ reference point. The keypad/control panel is locked during calibration, with the message center reading “CO₂ AUTO CAL”.

IR CO₂ Sensor Stabilization Times

Start-Up- Allow the temperature and the CO₂ of the cabinet to stabilize at least 12 hours before proceeding.

Presently Operating - Allow CO₂ to stabilize at least 2 hours at setpoint before proceeding.

Calibrating the IR CO₂ System (continued)

1. Measure the CO₂ concentration in the chamber through the gas sample port with a FYRITE or other independent instrument. Several readings should be taken to ensure accuracy.
2. Press the Mode key until the Cal indicator lights.
3. Press the right arrow until “IR CAL XX.X” appears in the message center.
4. Press the up/down arrow to adjust the display to match the independent instrument reading.
5. Press Enter to store the calibration.
6. After Enter is pressed, the unit will go into a calibration cycle that will last 5 to 6 minutes. The control panel is locked during this cycle.
7. Press the Mode key to return to Run mode.

Calibrating Relative Humidity

All 310 Series incubators can be equipped with an optional direct readout relative humidity (RH) sensor. This is a readout only of the chamber relative humidity level. It does not provide any control of the relative humidity in the cabinet.

Relative Humidity Stabilization Times

Start-Up - Allow 12 hours for the relative humidity and temperature in the chamber to stabilize before proceeding.

Already Operating - Allow at least 2 hours after temperature display reaches setpoint for relative humidity to stabilize before proceeding.

Calibrating Relative Humidity (continued)

1. Place an accurate independent instrument in the center of the chamber. Allow at least 30 minutes for RH to stabilize.
2. Press the Mode key until the Cal indicator lights.
3. Press the right arrow key until “RH CAL XX” appears in the message center.
4. Press the up/down arrow to match the display to the independent instrument.
5. Press Enter to store the calibration.
6. Press the Mode key to return to Run mode.

If a reliable RH measuring device is not available, the display may be calibrated to a typical level.


1. Follow the RH stabilization periods outlined previously.
2. With a full humidity pan and stable temperature, the relative humidity in the chamber will be 95%.
3. Using Step 3-5 of the relative humidity sensor adjustment, adjust the display to 95%.

This calibration method should be accurate to within 5%.

Chart 2-1

Calibrate Mode

Press Mode to light CAL

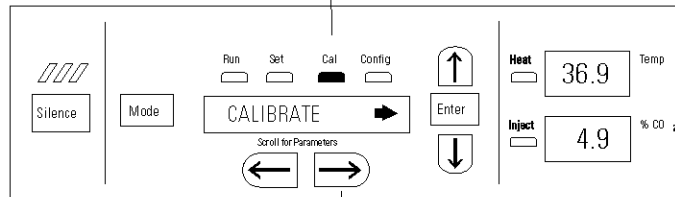

To Calibrate:

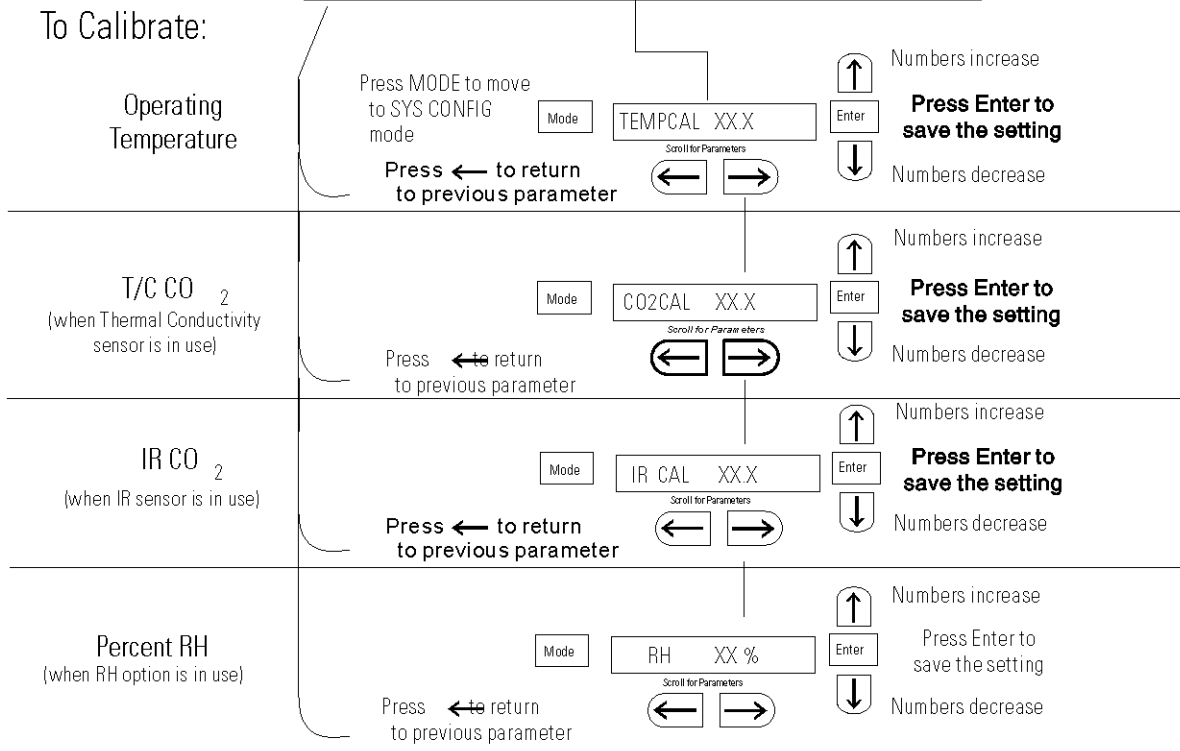


Chart2-1.edr

Section 3 Configuration

Several features available in Configuration Mode allow custom setup of the incubator. These features are listed and described below. All features may not be necessary in all applications, but are available if needed. To enter Configuration mode, press the Mode key until the Config indicator lights. Press the right and/or left arrow until the appropriate parameter appears in the message center. See Chart 3-1 for more detail.

Turn the Audible Alarm ON/OFF

The audible alarm can be turned on or off. The factory setting is ON.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Audible XXX is displayed in the message center.
3. Press up/down arrow to toggle Audible ON/OFF.
4. Press Enter to save the setting.
5. Press the Mode key to return to run mode or right/left to go to next/previous parameter.

Set an Access Code

A 3-digit Access Code can be entered to avoid unauthorized personnel from changing the setpoints, calibration, or configuration. A setting of 000 will bypass the access code. The factory setting is 000.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Acc Code XXX is displayed in the message center.
3. Press up/down arrow to change the access code.
4. Press Enter to save the access code.
5. Press the Mode key to return to the Run mode or right/left to go to next/previous parameter.

Set a Low Temp Alarm Limit

The low temp alarm limit (tracking alarm) is the deviation from the temperature setpoint which will cause a low temp alarm. The low temp alarm is variable from 0.5° below setpoint to 5° below setpoint. The factory setting is 1° below setpoint. A minus sign (-) in the display indicates that the alarm setting is below the setpoint.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Temp Lo -X.X is displayed in the message center.
3. Press up/down arrow to change the low temp alarm limit.
4. Press Enter to save the low temp alarm limit.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Enable the Low Temp Alarm to Trip Contacts

The low temperature alarm can be programmed to trip the remote alarm contacts. A setting of ON allows it, a setting of OFF will not allow the temperature alarm to trip the contacts. The factory setting is ON.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until TMP RLY XXX is displayed.
3. Press the up/down key to toggle the setting ON/OFF.
4. Press Enter to save the setting
5. Press the Mode key to return to Run or the right/left arrow key to go to next/previous parameter.

Set a Low CO₂ Alarm Limit

The low CO₂ alarm limit (tracking alarm) is the deviation from the CO₂ setpoint which will cause a low CO₂ alarm. The setpoint is variable from 0.5% CO₂ below setpoint to 5.0% CO₂ below setpoint. The factory setting is 1.0% CO₂ below setpoint. A minus sign (-) in the display indicates that the alarm setting is below the setpoint.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until CO₂ Lo -X.X is displayed in the message center.
3. Press up/down arrow to change the low CO₂ alarm limit.
4. Press Enter to save the low CO₂ alarm limit.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Setting a High CO₂ Alarm Limit

The high CO₂ alarm limit (tracking alarm) is the deviation from the CO₂ setpoint which will cause a high CO₂ alarm. The setpoint is variable from 0.5% CO₂ above setpoint to 5.0% CO₂ above setpoint. The factory setting is 1.0% CO₂ above setpoint.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until CO₂ Hi X.X is displayed in the message center.
3. Press up/down arrow to change the high CO₂ alarm limit.
4. Press Enter to save the high CO₂ alarm limit.
5. Press the Mode key to return to run mode or right/left to go to next/previous parameter.

Enable CO₂ Alarms to Trip Contacts

High and Low CO₂ alarms can be programmed to trip the remote alarm contacts. A setting of ON allows it, a setting of OFF will not allow the CO₂ alarms to trip the contacts. The factory setting is ON.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until CO₂ Rly XXX is displayed in the message center.
3. Press up/down arrow to toggle the setting ON/OFF.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run Mode or right/left to go to next/previous parameter.

Set New Zero Number for T/C CO₂ Sensors

If a new T/C CO₂ sensor is being installed, the two numbers on the factory installed sticker on the T/C cell must be entered to calibrate the CO₂ in the unit.

Note For the technician's convenience, a label containing the two numbers on the T/C cell is affixed inside the electronics drawer. ▲

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until T/CZ# XXXX is displayed in the message center.
3. Press up/down arrow to change the zero number to match the sticker.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Set New Span Number for T/C CO₂ Sensors

If a new T/C CO₂ sensor is being installed, the two numbers on the factory installed sticker on the T/C cell must be entered to calibrate the CO₂ in the unit.

Note For the technician's convenience, a label containing the two numbers on the T/C cell is affixed inside the electronics drawer. ▲

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until T/CS# XXXX is displayed in the message center.
3. Press up/down arrow to change the span number to match the sticker.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Set a Low RH Alarm Limit

On units that have the RH option installed, a low RH alarm limit may be entered. The low RH alarm limit is the %RH in the cabinet which will cause a low RH alarm. The setpoint is variable from setpoint 0 to 90% RH. The factory setting is 0% RH which disables the alarm.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until RH Lo XX is displayed in the message center.
3. Press up/down arrow to change the RH low alarm limit.
4. Press Enter to save the RH low alarm limit.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Enable RH Alarms to Trip Contacts

The low RH alarm can be programmed to trip the remote alarm contacts. A setting of ON will cause this, a setting of OFF will not allow the RH alarm to trip the contacts. The factory setting is ON.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until RH Rly XXX is displayed in the message center.
3. Press up/down arrow to toggle the setting ON/OFF.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Enable Temp/RH to be Displayed

On units that are equipped with the RH option, the upper seven segment display on the control panel can be configured to display Temp continuously, RH continuously, or toggle between Temp and RH. If the units does not have RH, the upper display will always display temperature. If temperature is set to ON and the RH is set OFF, temperature will be displayed continuously. If temperature is set to OFF and RH is set to ON, RH will be displayed continuously. If both are turned ON, the display will toggle between the two. The factory setting will default to toggle mode if the RH option is present.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Disp Tmp XXX or Disp RH XXX is displayed in the message center.
3. Press up/down arrow to toggle the setting ON/OFF.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Select Primary Tank w/ Gas Guard Option

On units equipped with the Gas Guard option, a primary tank can be selected. The primary tank will be either Tank 1 or 2. The factory setting is Tank1.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Tnk Sel X is displayed in the message center.
3. Press up/down arrow to toggle the setting between 1 and 2.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Disable the Gas Guard System

On units equipped with the Gas Guard option, the Gas Guard system may be turned ON, or OFF if not in use. The factory setting is ON.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until Gas Grd XX is displayed in the message center.
3. Press up/down arrow to toggle the setting ON/OFF.
4. Press Enter to save the setting.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.

Communications Address for RS485


On units that have the RS485 option, direct communication with the Model 1535 alarm system can be established. Each piece of equipment connected to the 1535 must have a unique address. An address of 0-24 can be entered for the incubator. A setting of 0 is an invalid address that the 1535 will ignore. The factory setting for the RS485 address is 0.

1. Press the Mode key until the Config indicator lights.
2. Press the right arrow until RS485 XX is displayed in the message center.
3. Press up/down arrow to move the RS485 address.
4. Press Enter to save the RS485 address.
5. Press the Mode key to return to Run mode or right/left to go to next/previous parameter.


Chart 3-1

Configure Mode

Press Mode to light CONFIG

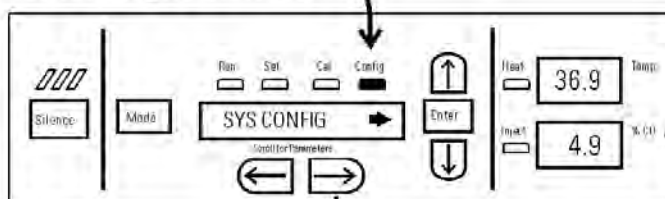

To Configure:

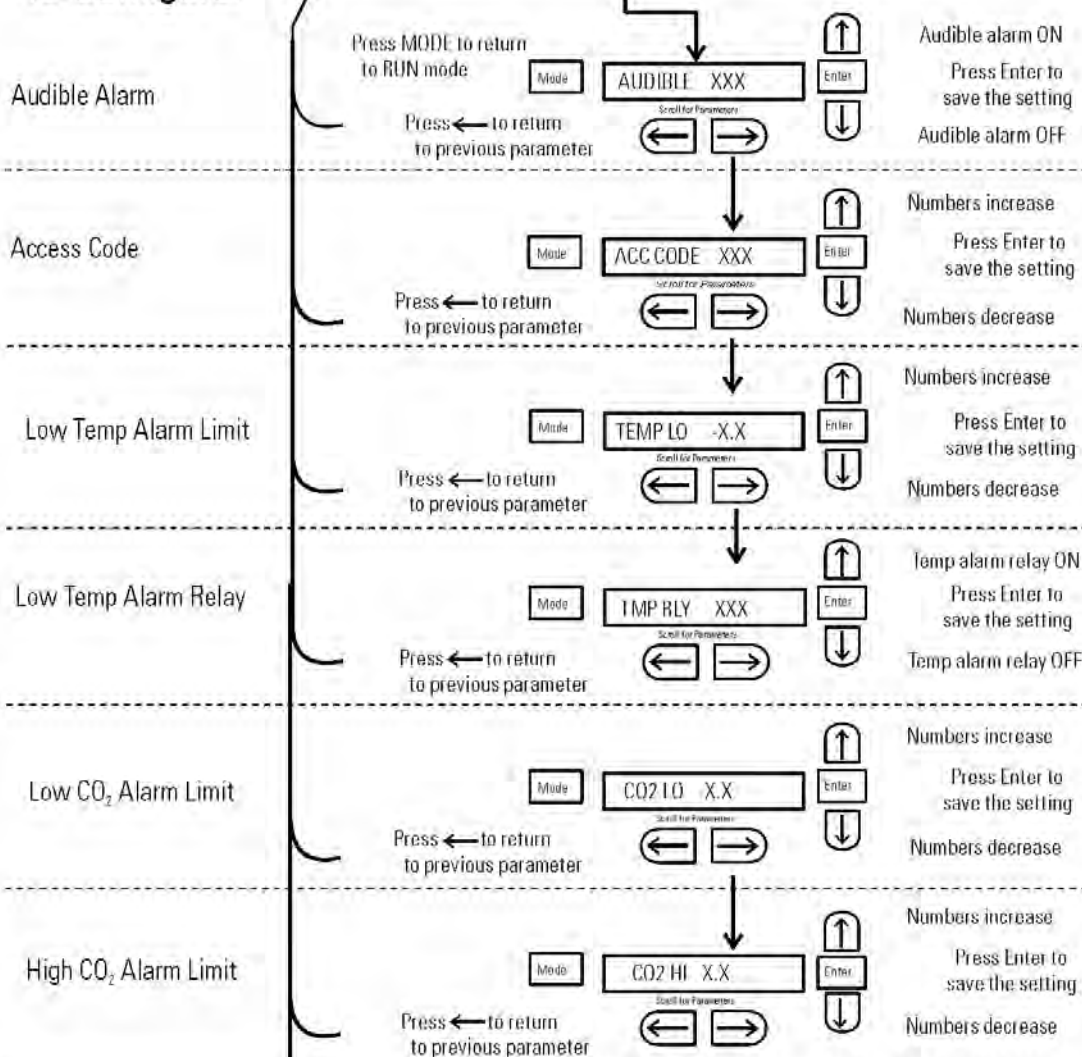

continue on next page

Chart3-1-1.cdr

Configure Mode

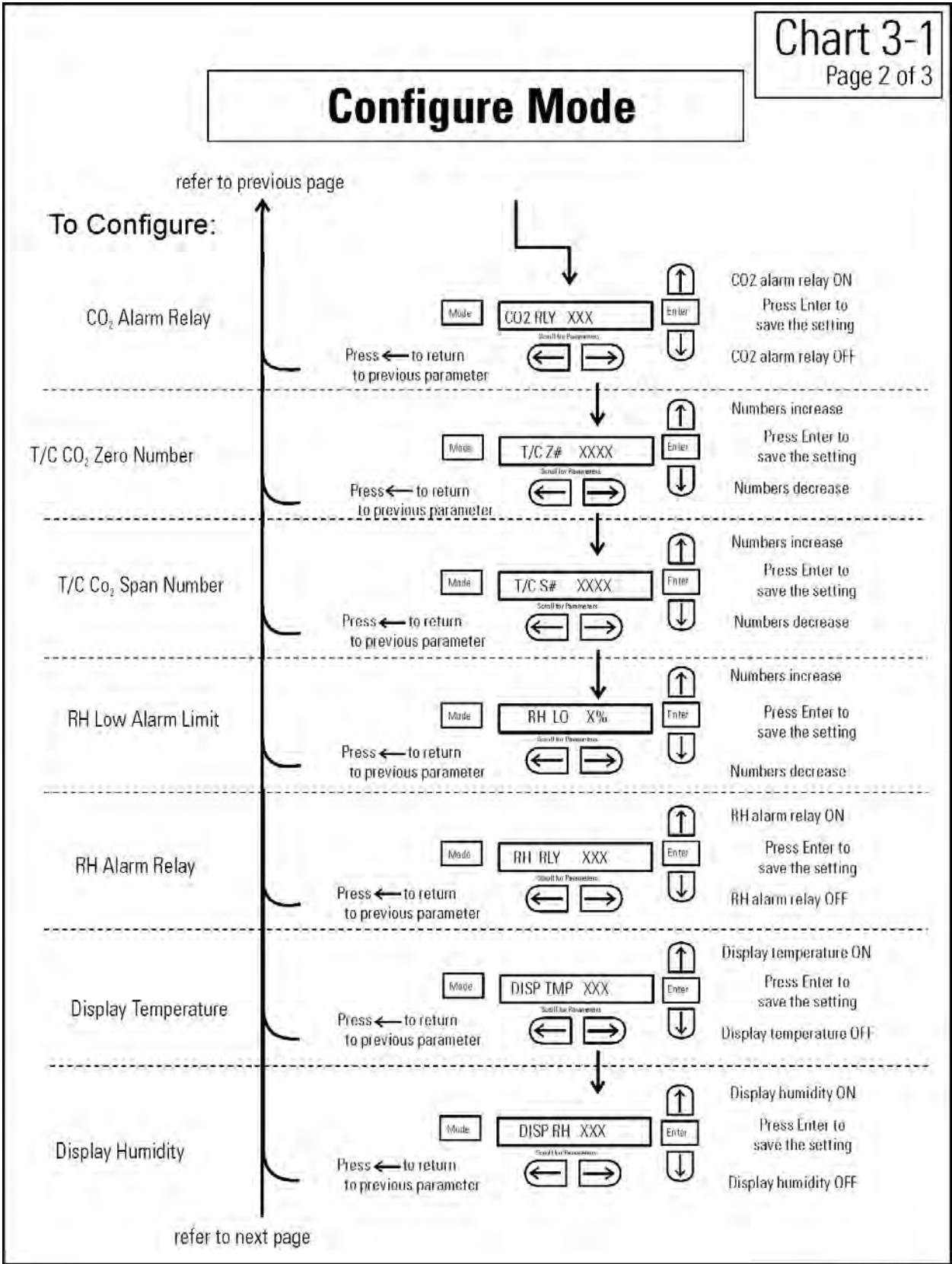



Chart 3-1
Page 3 of 3

Configure Mode

refer to previous page

To Configure:


Tank Select

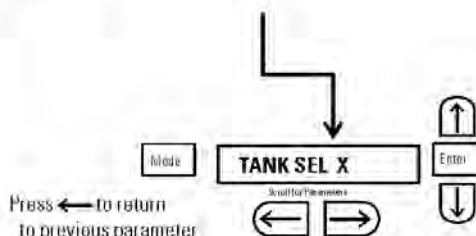

Tank 1 prime
Press Enter to save the setting
Tank 2 prime

Press ← to return to previous parameter

Gas Guard

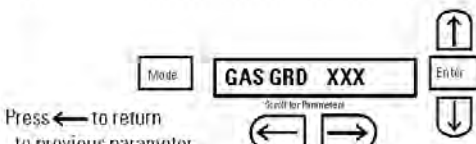

Gas Guard ON
Press Enter to save the setting
Gas Guard OFF

Press ← to return to previous parameter

RS485



Numbers increase
Press Enter to save the RS485 address
Numbers decrease

Press ← to return to previous parameter

Chan3-1 3.cdr

Section 4 Alarms

The Model 310 Series incubator alarm system is shown in the table below. When an alarm is active, the message appears in the LED message center. Pressing Silence disables the audible alarm for the ringback period. However, the visual alarm continues until the incubator returns to a normal condition. The alarms are momentary alarms only. When an alarm condition occurs and then returns to normal, the incubator automatically clears the alarm condition and the message center.

Table 4-1. Alarms and Descriptions

Description	Message	Delay	Ringback	Relay
No alarm condition exists	SYSTEM OK	----	----	----
CO2 System Auto Calibrating	CO2 AUTO CAL	----	----	----
Temp > Otemp Set point	SYS IN OTEMP	0 min.	15 min.	Yes
Air Temp Sensor Fault (See Section 4)	TMP SNSR ERR	0 min.	15 min.	No
Temperature Controller Failure (See Section 4)	TMP CTRL ERR	0 min.	15 min.	YES
CO2 Sensor Fault (See Section 4)	CO2 SNSR ERR	0 min.	15 min.	No
CO2 Auto Zero Fault (IR option only, See Section 4)	IR AUTOZ ERR	0 min.	15 min.	No
CO2 Sensor Cannot be Calibrated (IR option only, See Section 4)	REPL IR SNSR	0 min.	15 min.	No
Inner Door is Open	DOOR OPEN	15 min.	15 min.	No
CO2 is higher than CO2 High Tracking Alarm	CO2 IS HIGH	15 min.	15 min.	Programmable
CO2 is lower than CO2 Low Tracking Alarm	CO2 IS LOW	15 min.	15 min.	Programmable
TEMP is lower TEMP Low Tracking Alarm	TEMP IS LOW	15 min.	15 min.	Programmable
RH is lower than RH Low Limit Alarm (RH option)	RH IS LOW	30 min.	15 min.	Programmable
Tank 1 is low, switch to Tank 2 (Gas Guard only)	TANK1 LOW	0 min.	----	No
Tank 2 is low, switch to Tank 1 (Gas Guard only)	TANK2 LOW	0 min.	----	No
Both tanks are low (Gas Guard only)	TANK 1 and 2 LOW	0 min.	15 min.	No

- All alarm delays and ringback times are ± 30 seconds -

When multiple alarm conditions occur, active messages are displayed in the message center one at a time, updating at 5 second intervals. Pressing Silence during multiple alarms causes all active alarms to be silenced and to ring back in 15 minutes.

The temperature alarms are disabled when the Temp set point is 10°C. The CO₂ alarms are disabled when the CO₂ set point is 0.0%.

Temp Control Failure Alarm TMP CNTR ERR

In addition to other safety features designed into Model 310 Series incubators, a thermostat is also provided to monitor the cabinet's temperature. In the unlikely event of a temperature control failure, the thermostat will turn off all heaters at a cabinet temperature of 65°C, $\pm 10\%$. This is intended to be a safety feature to protect the incubator, and is not intended to protect the cell cultures or the equipment inside the chamber should a temperature control failure occur.

Sensor Fault Alarms

The microprocessor in Model 310 Series incubators continually scans all available sensors to ensure that they are operating properly. Should an error be detected, the incubator sounds an alarm and displays the appropriate message. If such an alarm occurs, contact your local distributor or the Technical Services department at 740-373-4763 or 1-800-438-4851 (USA and Canada) or fax 740-373-4189.

CO2 SNSR ERR

If the cables or connectors between the main microprocessor board and the CO₂ sensor or between the CO₂ board and the sensor head on the I/R CO₂ units become loose or disconnected, the CO₂ SENS ERR alarm occurs.

I/R Units

On I/R incubators, the red light on the I/R module (see Figure 5-8 for module circuit board location) will be lit continuously if communication is lost between the CO₂ board and the sensor head. The CO₂ display will also be locked at 00.0 without injection. Turning the incubator Off and On will not clear the alarm. Only proper connection of all components will correct the alarm.

IR AUTOZ ERR

On incubators equipped with I/R CO₂ control, calibration is done automatically using an auto zero system. Auto zero occurs once every 24 hours. If, during the auto zero cycle, a CO₂ correction of more than 0.45% is detected, the IR AUTOZ ERR alarm occurs.

IR AUTOZ ERR (continued)

Possible problems which cause this alarm are:

- Auto Zero pump, orifice, filter or tubing will not let air to the sensor.

Possibilities are:

Defective or electrically disconnected air pump

Kinked auto zero vinyl tubing

Disconnected tubing between the air pump and the sensor

Plugged filter or orifice on the auto zero assembly

Defective auto zero circuit

- Cabinet temperature has been increased significantly from a previous setpoint. For example, the unit was calibrated and operating at 30°C and the setpoint is increased to 50°C. In this example, calibrating the CO₂ will correct the alarm.
- High background CO₂ in the laboratory. This could be from leaks in the tank, regulator or vinyl CO₂ tubing.
- High CO₂ sensor calibration drift. This will require replacement of the sensor.
- Calibration of the CO₂ system to less than 3% actual CO₂. If this has occurred, contact the Technical Services Department.

PREVENTIVE MAINTENANCE

Incubators

Your equipment has been thoroughly tested and calibrated before shipment. Regular preventive maintenance is important to keep your unit functioning properly. The operator should perform routine cleaning and maintenance on a regular basis. For maximum performance and efficiency, it is recommended the unit be checked and calibrated periodically by a qualified service technician.

The following is a condensed list of preventive maintenance requirements. See the specified section of the operating manual for further details.

Thermo Electron Corporation has qualified service technicians, using NIST traceable instruments, available in many areas. For more information on Preventive Maintenance or Extended Warranties, please contact us at the number listed below.

Cleaning and calibration adjustment intervals are dependent upon use, environmental conditions and accuracy required.

Tips for all incubators:

- Do NOT use bleach or any disinfectant that has high chloros
- Use sterile, distilled or demineralized water.
- Avoid spraying cleaner on the CO₂ sensor.
- Do not use powdered gloves for tissue cultures.

Preventive Maintenance for Direct Heat Incubators

Action	Daily	Weekly	Monthly	6 Months	Yearly	2 Years
Inspect door latch, hinges and door gasket seal.					✓	
Check water level in the humidity pan, 1/2" from top. See "Filling the Humidity Pan".		✓				
Verify and document CO ₂ , O ₂ , humidity and temperature calibration, as applicable. See Calibration section					✓	
Perform a complete decontamination procedure. Wipe down interior, shelves and side ducts with disinfectant. Change or clean blower wheel and scroll. Clean top duct. Clean humidity pan. Rinse everything well with sterile distilled water. See "Cleaning the Interior, "Cleaning the Humidity Pan".	Between experiments <i>More frequent decontamination may be required, depending on use and environmental conditions</i>					
Change HEPA and gas filters, if applicable (<i>or as needed</i>). See "Installing Air Sample Filter", "Installing HEPA Filter".				✓		
Replace AquaTec water cleaning cell, if applicable.				✓		

Section 5 Routine Maintenance

Warning If the unit has been in service, turn it off and disconnect the power cord connector before proceeding with any maintenance. ▲

Disinfecting the Incubator Interior

Caution Before using any cleaning or decontamination method except those recommended by the manufacturer, users must check with the manufacturer that the proposed method will not damage the equipment. ▲

Use an appropriate disinfectant. All articles and surfaces to be disinfected must be thoroughly cleaned, rinsed with sterile water, and rough dried.

Warning Alcohol, even a 70% solution, is volatile and flammable. Use it only in a well ventilated area that is free from open flame. If any component is cleaned with alcohol, do not expose the component to open flame or other possible hazard.

Do not spray T/C sensor with flammable solutions. The internal temperature of the CO₂ sensor is approximately 150°C when the unit is in operation. Allow sufficient time for the sensor to cool before cleaning. ▲

Caution Do not use strong alkaline or caustic agents. Stainless steel is corrosion resistant, not corrosion-proof. Do not use solutions of sodium hypochlorite (bleach) as they may also cause pitting and rusting. ▲

1. Remove the shelves, HEPA filter (a factory installed option), access port filter, sample air filter, left and right duct sheets and the top duct. The top duct requires the two wing nuts to be removed.
2. Set the HEPA filter aside. Wash the shelves, ducts, wing nuts and stopper with disinfectant. Optional: The shelves, ducts and wing nuts may be autoclaved.
3. Remove the blower scroll held on by four screws. Remove the blower wheel and the access port filter. If a new blower wheel and scroll are going to be used, discard the old ones. If the old ones are being reused, wash and rinse all parts with disinfectant.

4. The CO₂ sensor may be lowered by unscrewing the retaining ring around the sensor. This will allow the sensor gasket to be removed, cleaned and replaced. WIPE the CO₂ sensor, do not spray or soak.
5. Wash the inner door gasket with disinfectant. The inner door gasket may be removed and/or replaced.
6. Wash the cabinet interior with disinfectant, starting at the top and working down. Refer to the disinfectant directions for length of time needed before rinsing. Wash the inner door both inside and out. The cabinet and door must be rinsed with sterile water until the disinfectant has been removed. After the cabinet has been rinsed, spray with 70% alcohol.
7. If the CO₂ sensor gasket was removed, spray it with 70% alcohol and install it. Install the sample filter.
8. Install the top duct, left and right ducts, inner door gasket, and access port stopper spraying each with 70% alcohol.
9. Install new HEPA Filter.
10. Install the shelves and spray with 70% alcohol.

Cleaning the Cabinet Exterior

Clean the incubator exterior with a damp sponge or soft, well-wrung cloth and mild detergent dissolved in water. Dry with a soft cloth.

Cleaning the Glass Doors

The chamber glass door and the optional independent inner glass doors may be cleaned using the same disinfectant used on the incubator interior. It is imperative that they be rinsed with sterile distilled water to remove the disinfectant residue. The doors should then be dried with a soft cloth.

Precautions in the cleaning and care of the incubator glass doors:

Moisture leaches alkaline materials (sodium, Na) from the surface of the glass. Evaporation of the moisture concentrates the alkaline and may produce a white staining or clouding of the glass surface. Cleaning chemicals with a pH above 9 and heat (autoclaving) accelerate the corrosion process. Therefore, it is very important to rinse and dry the glass doors after cleaning. Autoclaving the glass doors should be avoided.

There is no simple method for repairing corroded glass. In most cases, the glass must be replaced.


Cleaning the Humidity Pan

Clean the humidity pan with soap and water and a general use laboratory disinfectant. Rinse with sterile water and spray with 70% alcohol. The humidity pan may be autoclaved.

Reversing the Door Swing

For side-by-side operation or changing lab layouts, the inner and outer doors are field reversible. The procedure is written from the perspective of changing the door swing from a left-hand to a right-hand swing. All screw holes are pre-drilled for reversing the door. The tools required are a Phillips screwdriver, a 1/8" Allen wrench and a flatblade screwdriver.

The door reversing procedure takes about 30 minutes.

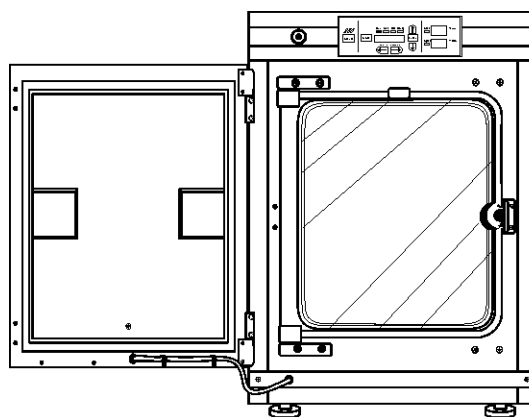

Left hand door swing

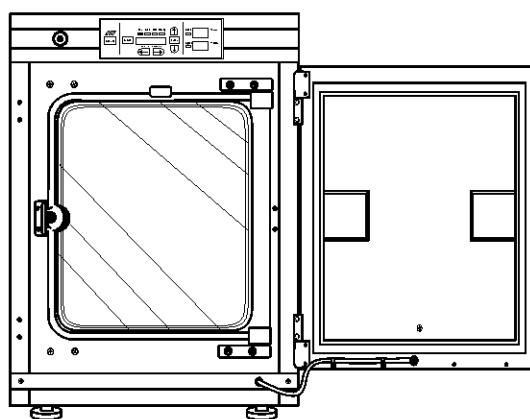


Figure 5-1. Door Swings

Reversing the Door Swing (continued)

1. Remove the kick plate at the bottom of the cabinet by removing the two Phillips screws indicated by the black arrows in the illustration. Disconnect the cable inside the cabinet.

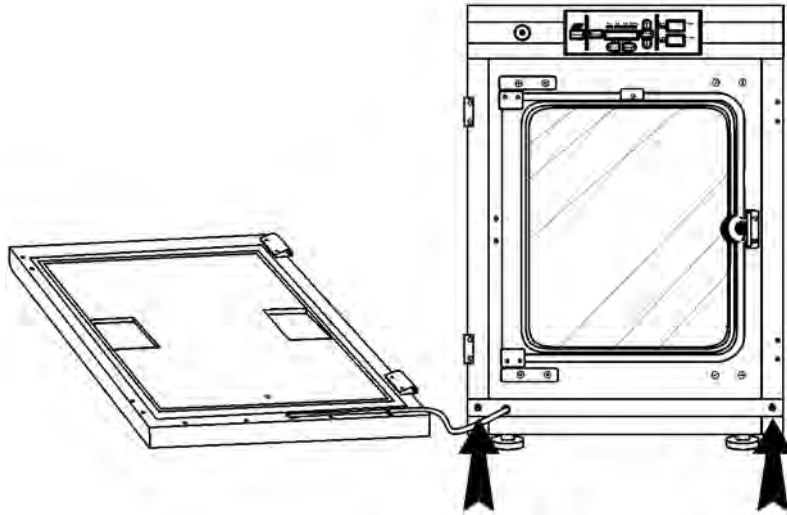


Figure 5-2. Remove Two Screws

2. Open the outer door and remove it by lifting it off its hinge pins. Place the door on a padded surface to prevent scratches (Figure 5-2).

Note The hinge pins are not connected to the brackets and may fall out when the door is lifted off. ▲

3. Refer to the illustration of the outer door in Figure 5-2. With a Phillips screwdriver, remove and save the four nylon screws from the door hinge mounting holes. The black arrows identify the screw locations.
4. Remove the door hinges, rotate them 180° and secure them on the left side of the door. Install the nylon screws into the newly exposed holes.
5. Remove the two plastic clips securing the door heater cable. These clips will be rotated 180° when installed on the left lower edge of the door frame.

6. Rotate the kick plate and attach the door heater cable with the clips as shown in Figure 5-3. The insert in the illustration shows the proper positioning of the clips.

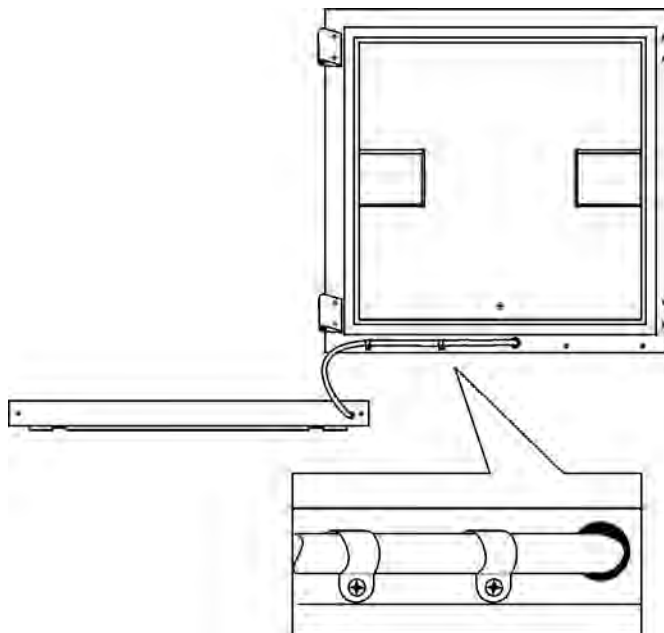


Figure 5-3. Rotate Kickplate

7. The incubator and door should look similar to Figure 5-4.

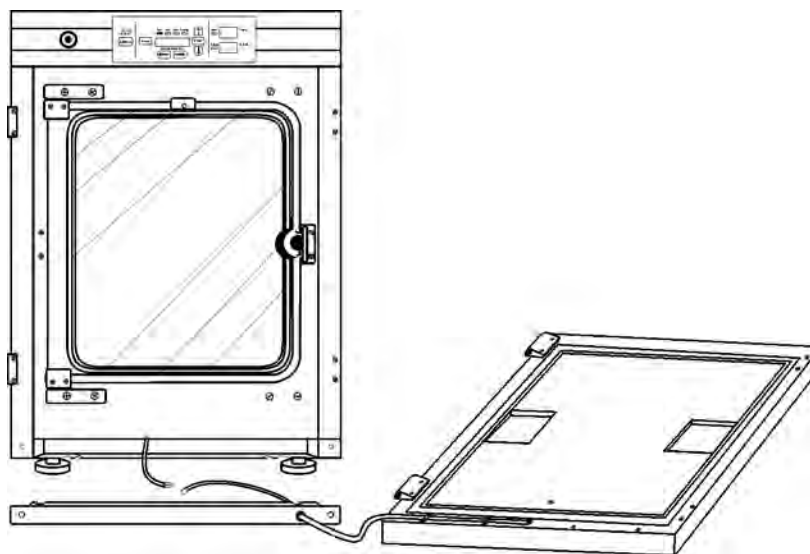


Figure 5-4. Unit with Detached Door

8. Refer to Figure 5-5 for the next steps. Remove the nylon screws along the right side of the cabinet. Remove the outer door hinges. Rotate them 180° and install them on the right side of the cabinet. Install the nylon screws in the holes vacated by the hinges.

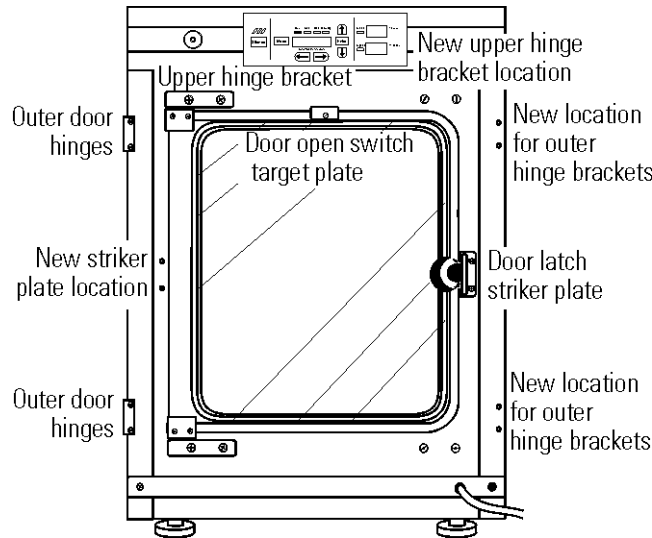


Figure 5-5. Component Locations

9. With a 1/8" Allen wrench, remove the door open switch target plate on the top of the inner door.
10. Remove the door latch striker plate.
11. Hold the door securely and remove the upper hinge bracket on the cabinet. This hinge bracket will be rotated 180° and be placed in the lower right corner of the cabinet where indicated.
12. Lift the glass door from the lower hinge and set it aside. Then remove the lower hinge from the cabinet. This hinge bracket will be rotated 180° and be placed in the upper right corner of the cabinet at where indicated.
13. Remove the screws and washers from the new striker plate location. These screws and washers can be threaded into the holes vacated by the striker plate where indicated.
14. Rotate the striker plate 180° from its original position and attach it to the cabinet where indicated.
15. Using a flatblade screwdriver, remove the threaded nylon plugs from the new hinge bracket locations. Install the lower hinge bracket where indicated.

16. Rotate the door 180° and place it in the lower hinge bracket. Install the upper hinge bracket where indicated.
17. Screw the threaded plugs into the holes vacated by the door hinge brackets where indicated in Figure 5-5.

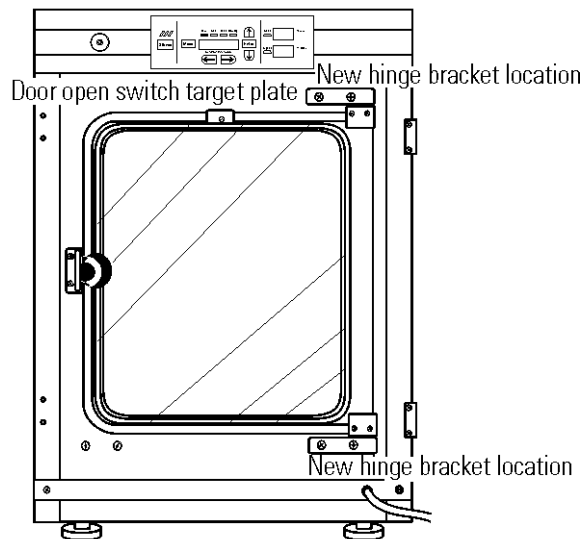


Figure 5-6. Reversed Components

18. Install the door open switch target plate. Refer to Figure 5-6 and 5-7. The plate must be positioned exactly 12.5 inches from the outside edge of the cabinet to the leading edge of the plate. Do not overtighten.

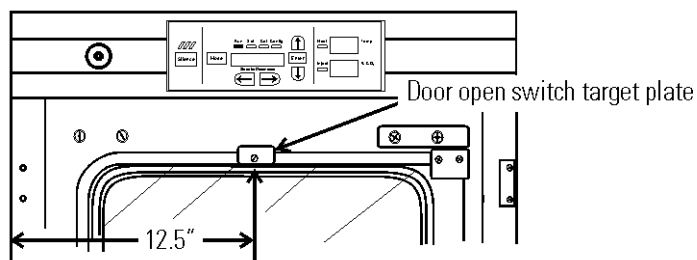


Figure 5-7. Plate Position

19. Connect the door heater cable and install the kick plate with the door heater cable on the right side.
20. Installing the outer door onto the hinge pins completes the door reversal procedure.

HEPA Filter (Factory Installed Option)

Replace the HEPA filter routinely on an annual basis, or more often if it changes color. The filter should also be changed when the chamber is cleaned.

Warning The electronics section contains hazardous voltages. Only qualified service personnel should access this area. ▲

Replacing the Power Fuses

There are only two replaceable fuses in the incubator. To access the fuses:

1. Turn off the incubator's power switch and unplug the power cord.
2. Open the outer door to the chamber.
3. Remove the two screws located below the control panel.
4. Lift up on the control panel, then pull it out.
5. Refer to Figure 5-8 for the location of the two fuses.

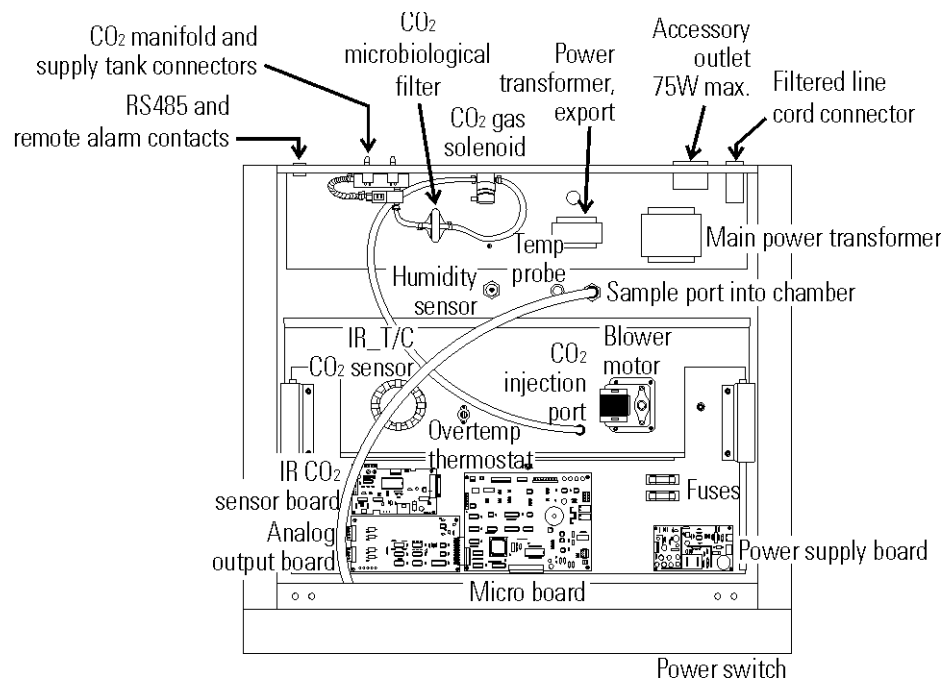


Figure 5-8. Major Components in Electronics Drawer

6. Slide the control panel back in, replace the screws and return the unit to service. If the fuse(s) blow after restoring power to the incubator, contact Technical Services Department (see Page iv near the front of this manual).

Table 5-1. Fuses

Fuse Voltage	Manufacturers Part #	Amperage Rating	Rupture Speed	IEC Letter Code
115 VAC ACC	BUSS GMC - 1.0A	1.0 Amp	Time-Lag	T
230 VAC ACC	BUSS BK-GMC-500ma	0.5 Amp	Time-Lag	T
24 VAC Door Heat	Schurter 034.3124	5.0 Amp	Time-Lag	T

Major Components

Expanded descriptions for some of the components indicated in Figure 5-8 follow.

Power Switch - Located beneath the control panel. The outer door must be open to access the switch.

RS 485 and Remote Alarm Contacts - Connectors and auxiliary wire through port

CO₂ manifold and supply tank connectors - Part of the Gas Guard system, a factory installed option.

Humidity sensor (optional)

Analog output board - 4-20 mA, 0-1 Volt, 0-5 Volt

Electronics Section

1. Turn off the incubator's power switch and unplug the power cord.
2. Remove the two screws from the front of the cabinet top and the two screws in the uppermost corners of the back panel.
3. Slide the top about an inch to the rear and lift it off.

Note There are small locking tabs in the rear which will be bent if the top is lifted off without sliding it first to the rear. ▲

Replacing the Sample Air Filter

1. Connect one end of the sample port filter to the ceiling port.
2. Connect the other end of the filter to the back of the blower fan inlet.

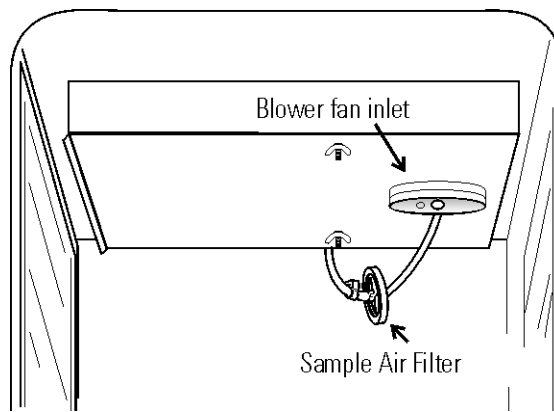


Figure 5-9. Air Sample Filter

Section 6 Factory Options

Descriptions for connections to external equipment follow.

Connect Remote Alarm Contacts

A set of relay contacts is provided to monitor alarms through an RJ11 telephone style connector on the back of the cabinet. Refer to Figure 6-3 for the location of the alarm connector. The 12-foot telephone cord (P/N 190388) and RJ11-to-screw terminal conversion box (190392) are available through our service department.

The remote alarm provides a NO (normally open) output, an NC (normally closed) output and COM (common). Refer to Figure 6-1.

The contacts will trip on a power outage or an overtemperature condition. The contacts may also be programmed to trip or not trip on temperature alarms, CO₂ alarms, and RH alarms. See Section 3, Configuration Mode.

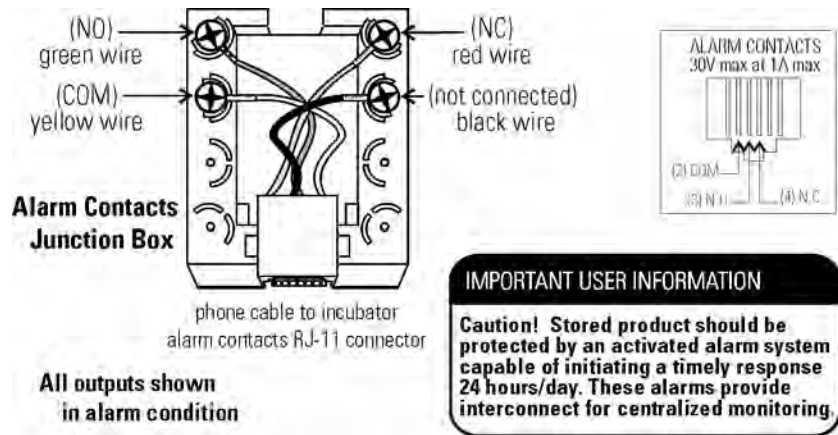


Figure 6-1. Remote Alarm Output



Connect RS485 Interface

All incubator models can be purchased with the RS485 communications option (P/N 190523). This option allows the incubator to be directly connected to a Model 1535 alarm system without the use of a communications module. A junction box is provided with each RS485 option. See Figure 6-2 for wiring details. Figure 6-3 shows the RS485 connector location on the back of the incubator cabinet.

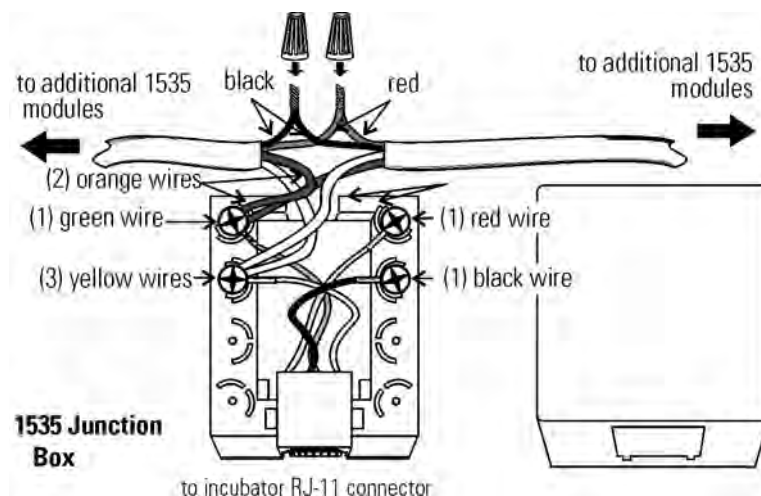


Figure 6-2. RS485 Wiring

To allow the incubator and the 1535 to communicate, an address must be allocated on the 1535. Refer to Section 5.8 of the 1535 operating manual. The same address number must be assigned to the incubator. Refer to Section 3 of this manual.

Connecting the Analog Output Boards

Warning The electronics section contains hazardous voltages. Only qualified personnel should access this area. ▲

The analog output board option (P/N 190512, 190543, 190544) allows the incubator to output analog signals representing the air temperature of the cabinet, CO₂ content, and relative humidity, depending upon which systems are in the incubator. There are three different analog output board options available : 0-1V, 0-5V, or 4-20mA signals. Negative display readings will output 0V. The outputs do not have isolated grounds. Refer to Table 6-1 for output specifications of the three boards.

For the 0-1V and 0-5V boards, the recording device must supply a load ≥ 1000 ohm. For the 4-20mA board, the recording device must supply a load of ≤ 100 ohm.

Table 6-1. Analog Output Boards

	190512 4-20 mA Output Scaling 4-20 mA Equals	190544 0-1V Output Scaling 0-1 V Equals	190543 0-5V Output Scaling 0-5V Equals
Temperature	0.0-100.0°C	0.0-100.0°C	0.0-100.0°C
RH	0-100 %RH	0-100 %RH	0-100 %RH
CO2	0.0-100.0 %CO2	0-100.0 %CO2	0-100.0 %CO2

To wire in the analog output board, a shielded 22 gauge, 3-conductor wire (P/N 73041) is recommended, with the maximum length being 50 feet (15.2m). This wire is readily available from other vendors including Alpha P/N 2403 and Deerborn P/N 972203.

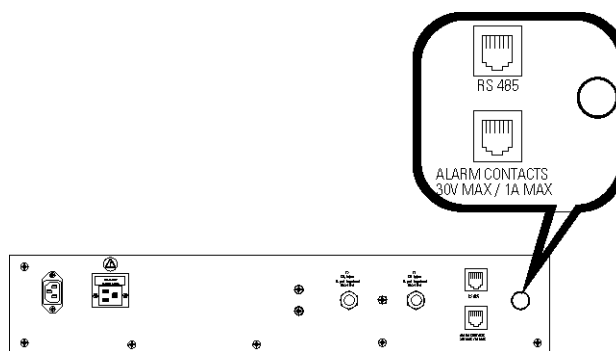


Figure 6-3. Auxiliary Through-Port on Back Panel

Caution Accuracy of the output at the board terminal strip to the incubator display is ± 1 unit. There is no calibration from the incubator. Calibration to the incubator display must be at the instrument connected to the output board. ▲

To install the analog board:

1. Turn off the incubator and unplug it from the wall outlet.
2. Remove the top of the electronics section. Refer to Section 5 of this manual.*
3. Locate the Analog Output board, Figure 6-4 and Figure 5-8.

* Alternate access to the electronics is to open the outer door and remove the two screws in the upper corners of the cabinet just under the display top. Lifting up the top section, the electronics drawer can be pulled out to the limit of the internal wiring.

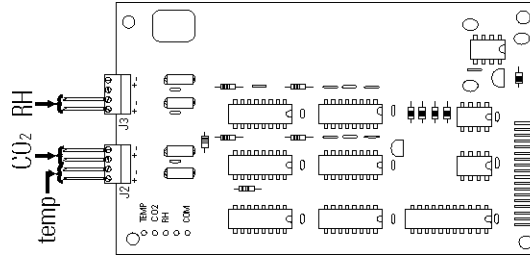


Figure 6-4. Connector Terminals

4. Each system monitored (Temp, CO₂, RH) requires two conductors (refer to Figure 6-4). Feed the wires through the auxiliary wire through-port on the back panel. This port is a circular fitting adjacent to the remote alarm and RS485 connectors as illustrated in Figure 6-3. The wire routing is shown by the black arrows in Figure 6-5.

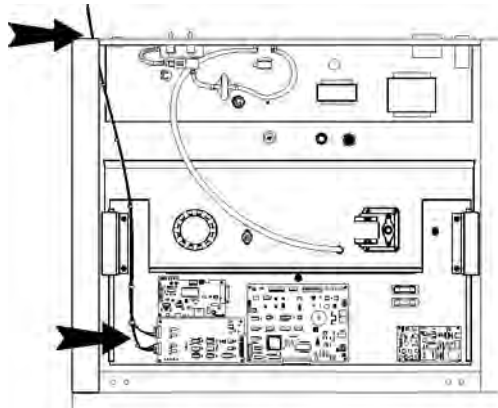


Figure 6-5. Wire Routing

5. Strip the ends of each conductor and wire it to the appropriate terminals of connectors J2 and/or J3 on the analog board. Refer to Figure 6-4.
6. When wiring is completed, slide in the electronics drawer or the replace the sheet metal cabinet top.
7. Replace the screws removed earlier and return the unit to service.

CO₂ Gas Guard

Warning CO₂ can be lethal in high concentrations. Refer to warnings in Section 1 of this manual. ▲

The Model 310 Series incubators can be equipped with a built-in Gas Guard system (P/N 190640) that will operate with a CO₂ gas supply. The Gas Guard uses two pressure switches to continuously monitor the pressures of two independent CO₂ supplies and automatically switches from one supply to the other when the supply pressure drops below 10 psig (0.690 bar). The Gas Guard's design does not facilitate use by multiple incubators.

The CO₂ gas supplies must be equipped with two-stage pressure regulators with gauges. The high pressure gauge at the tank should have a 0-2000 psig range and the low pressure gauge should have a 0-30 psig range. The gas supply to the incubator must be maintained at 15 psig (1.034 bar), ±5 psig. Gas pressures below 15 psig will cause nuisance alarms to occur on incubators equipped with the built-in Gas Guard.

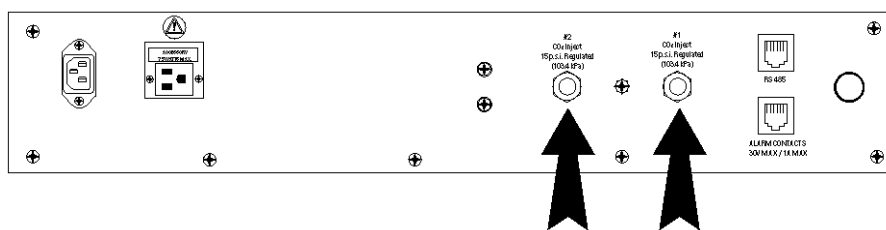


Figure 6-6. Connect Gas Supplies

Connecting the CO₂ Gas Supplies

The CO₂ inlets for the Gas Guard are located on the rear of the cabinet. Using 1/4" ID tubing, connect one of the CO₂ supply tanks to the fitting labeled CO₂ Inlet #1 Tank. Connect the second CO₂ supply tank to the fitting labeled CO₂ Inlet #2 Tank. Install 3/8" hose clamps to secure the 1/4" ID tubing to the fittings on the rear of the drawer. (Refer to Figure 6-6.)

De-activating the Gas Guard

The built-in Gas Guard is turned ON when shipped from the factory. In addition, the Tank Sel for the Gas Guard is specified as Tank 1 when shipped. Refer to Section 3, Configuration, to de-activate the Gas Guard or change the Tank Sel from #1 to #2. If the Gas Guard system is not used, the incubator will function normally by supplying CO₂ from the supply tank connected to Inlet #1.

Operation of the CO₂ Gas Guard

With the Gas Guard in operation, the incubator will use the gas supplied by the tank connected to Inlet #1 until the pressure drops below 10 psig (0.690 bars). At this time, the Gas Guard automatically switches to gas supplied through CO₂ Inlet #2.

In addition, the incubator automatically changes the Tank Sel in Configuration Mode from 1 to 2 to indicate that the incubator is now using gas supplied through Inlet # 2. If the gas supply to Inlet #1 is replenished, the incubator will continue to operate using the gas supplied through Inlet #2 unless the operator changes the Tank Sel from #2 to #1 through Configuration Mode. Refer to Section 3, Configuration.

Audible and visual alarms occur on the control panel when the gas guard switches from one supply to the other. The audible alarm sounds until the operator presses the Silence key on the control panel. The visual alarm in the Message Center reads Tank 1 Low while the audible alarm is sounding, but the message is removed when the operator presses the Silence key. However, the 'Tank Low' indicator on the control panel stays lighted until the condition is resolved. The unit operates normally.

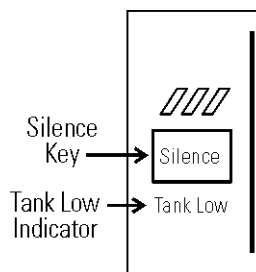


Figure 6-7. Silence Key

Caution Both the audible and visual alarms described above do not ring back once the Silence key is pressed. ▲

If the Gas Guard system does not detect an adequate gas supply at CO₂ Inlet #1 or Inlet #2, a visual and audible alarm will again occur on the control panel. The visual alarm in the Message Center will read Tank 1&2 Low. The audible alarm will continue to ring until the Silence key is pressed. The audible alarm will ring back every 15 minutes after the alarm is silenced if the Gas Guard continues to detect that both gas supply pressures are below 10 psig (0.690 bars).

Humidity Readout

The 310 Series incubators can be equipped with a humidity sensor (P/N 190643) to monitor the relative humidity inside the chamber. The sensor is mounted to the top air duct and provides a signal that is displayed in 1% increments on the control panel. The humidity readout can be displayed continuously or toggles with the temperature readout. In addition, a low alarm limit can be set on the humidity readout, which will detect when the humidity pan runs dry. Refer to Section 3, Configuration.

Humidity Level Factors

- Water level in the humidity pan
- Frequency of door openings
- The humidity pan located on bottom of chamber versus a shelf.

Caution Incubators equipped with a Thermal Conductivity CO₂ sensor rely on a constant level of relative humidity in order to accurately measure and control the CO₂ concentration in the incubator. ▲

Humidity Readout Accuracy

The sensor is capable of measuring relative humidity from 10% to 100% with an accuracy of $\pm 5\%$ above 90%. See Section 2, Calibration, for details on calibrating the humidity readout.

Uninterruptible Power Supply Connections

Warning The Uninterruptible Power Supply is not rated for life-support applications. ▲

The optional Uninterruptible Power Supply ([UPS] P/N 270078 and 270082) provides a back-up power source for the incubator. When the batteries are fully charged and the incubator operating at 37°C in an 18°C ambient temperature, the UPS provides approximately one hour of uninterrupted power during a power failure. A higher operating temperature decreases this back-up time, a higher ambient temperature increases it. This estimated time assumes that the incubator will be already at its operating temperature and that the incubator door will not be opened during the back-up period.

The remote alarm contacts will not trip during a power failure while connected to the UPS until battery power is exhausted. There is however, an audible and visual alarm when the unit switches to battery operation.

The UPS batteries are fully charged when the unit has been connected to utility power for a minimum of four hours. The UPS power switch does not have to be on for the batteries to charge. When battery replacement becomes necessary, refer to the UPS user's manual.

Caution Use only factory recommended UPS units which produce a pure sine wave output. Other UPS units could damage the incubator and void the warranty. Typical facility back-up systems will not effect the efficiency of the incubator. ▲

Connecting UPS
P/N 270078, 115V, 50/60Hz

1. Plug the incubator line cord into one of the receptacles on the back of the UPS. Then plug the UPS line cord into a suitable wall outlet.
2. Turn the UPS power switch on.

Connecting UPS
P/N 270082, 230V, 50/60Hz

Use the line cord supplied with the incubator to connect the UPS to a suitable wall outlet, and the UPS coupler to connect the incubator to the UPS.

Warning The Uninterruptible Power Supply contains potentially hazardous voltages even when not connected to line voltage.
Do not attempt to disassemble UPS as it contains no user-serviceable parts. Repairs must be performed by factory trained service personnel. ▲

Section 7 Specifications

** Specifications are based on nominal voltages of 115V or 230V in ambient of 22°C to 25°C.*

Temperature

Control ±0.1°C Microprocessor PID Control
Setpoint Digital - Touch pad, 0.1°C
Range +5°C above ambient to 50°C
Uniformity ±0.2°C @ +37°C
Tracking alarm . . User programmable (low) indicator
Overtemp . . Tracking, user programmable, action, and indicator
Display Digital, LED, 0.1°C increments

Temperature Safety

Type . . Extreme temperature safety, action, and indicator
Sensor . . Thermostat, independent of temperature control system
Indicator . . Message center, audible and visual alarms

Relative Humidity

Control Humidity pan - natural vaporization
Humidity with pan 95% RH at 37°C
Display Optional in 1% increments
Alarm Low RH with optional RH monitor

CO₂

Control ±0.1% microprocessor PID control
Sensor T/C or IR
Readability 0.1%
Range 0 to 20%
Inlet pressure 15 psig, ±5 psig
Display Digital LED, 0.1% increments

Shelves

Dimensions18.5" x 18.5" (47cm x 47cm)
Construction . . .Stainless steel (belt sanded, both sides)
Surface area2.4 sq. ft. (0.22 sq. m) per shelf
Max. per chamber38.4 sq ft (3.6 sq. m)
Loading35 lbs (16kg) slide in and out, 50 lbs (23kg) stationary
Standard4
Maximum16

Construction

Interior volume6.5 cu. ft.
InteriorType 304 stainless steel shiny finish
Exterior18 gauge cold roll steel
Outer door gasket . . .Four-sided molded, magnetic vinyl
Inner door gasketFeather-edge, removable
InsulationFiberglass

Fittings

Access port1-1/4 inch removable silicone plug
CO₂ inlet1/4" barbed
Sample portFront mounted barbed

Electrical

115 Volt models -
90-125VAC, 50/60 Hz, 1PH, 2.4 FLA
230 Volt models -
180-250, 50/60 Hz, 1 PH, 2.0 FLA
Power switch/circuit breaker2 Pole, 6.0 Amp
Accessory outlet . . . Voltage equal to the cabinet input,
75 Watts maximum, 0.5ma leakage current
Alarm contacts . . . Deviation of temperature, CO₂, power, NO and NC

Unit BTU Output

115/230V344 BTUH (100 watt)

Dimensions

Exterior26.0" W x 38.5" H x 25.0" F-B
Interior21.4" W x 26.8" H x 20.0" F-B

Weight

Net Weight Stainless Steel205 lbs

Net Weight Copper230 lbs.

Options/Accessories

HEPA filter . . Factory installed: Stock no. 190858

Battery back-up . . Free Standing: Stock no. 270078, 120VAC, 60Hz

Stock no. 270082, 220VAC, 50Hz

Humidity . . Factory installed: Stock no. 190463 - provides humidity readout and low alarm

Inner door kit . . Customer installed: Stock no. 190650 - independent eight separate glass doors

Built-in Gas Guard . . Factory installed: Stock no. 190640 - CO₂ only, automatically switches supply tanks

Recorders . . Customer installed: Standard recorders used on water-jacket incubators

Decontamination kits

HEPA filter, blower assembly, inner door gasket, etc. -

Stock no. 190868 (with HEPA filter),

Stock no. 190869 (without HEPA filter)

Copper Options

Interior ductwork .Solid copper: Stock no. 190725

Perforated shelfCustomer Installed: Stock no. 224166 - solid copper

Humidity PanCustomer Installed: Stock no. 237015 - solid copper

Copper interior kitFactory installed: Replace stainless steel duct work, shelves, and

humidity pan with solid copper Stock no. 190870

Optional Data Outputs

Analog Factory installed: Stock no. 190544 - 0-1 volt

Analog Factory installed: Stock no. 190543 - 0-5 volt

Analog Factory installed: Stock no. 190512 - 4-20mA

Digital Factory installed: Stock no. 190523 - RS-485

(Compatible with Model 1535 Alarm Monitor only)

Certifications

Refer to the Declarations of Conformity at the back of this manual

Safety Specifications

Indoor Use Only

Altitude - 2,000 meters

Temperature - 5°C to 40°C

Humidity - 80% RH at or below 31°C, decreasing linearly to 50% RH at 40°C

Mains Supply Fluctuations - Mains supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage

Installation Category II ¹

Pollution Degree 2 ²

Class of Equipment I

¹ Installation Category (overvoltage category) defines the level of transient overvoltage which the instrument is designed to withstand safely. It depends on the nature of the electricity supply and its overvoltage protection means. For example, in CAT II which is the category used for instruments in installations supplied from a supply comparable to public mains such as hospital and research laboratories and most industrial laboratories, the expected transient overvoltage is 2500V for a 230V supply and 1500V for a 120V supply.




² Pollution Degree describes the amount of conductive pollution present in the operating environment. Pollution Degree 2 assumes that normally only non-conductive pollution such as dust occurs with the exception of occasional conductivity caused by condensation.

Section 8 Spare Parts

Part Number	Description
190699	Removable feather gasket
190670	Duct sheet and shelf channel kit
190772	Micro board, (main control)
230135	Fuse, 1A, TD, 5 x 20mm
230120	0.5A fuse, TD, 5 x 20mm (accessory outlet)
230161	Fuse, 5A, 5 x 20mm
156126	Motor, 2-pole, 115VAC, 50/60 Hz
190793	Display board
290184	Probe, 2252 Ohm/25°C, 1/8 x2
770001	Disposable filter, 99.97 (CO ₂ line, air sample)
760175	HEPA filter, 5.5" diameter
760199	HEPA2 VOC Filter Replacement Kit (includes 760200 HEPA2, 117036 silicone plug, 101018 o-ring)
1900067	Filter Replacement Kit (includes [2] 770001 filter and 760175 HEPA)
1900094	HEPA2 VOC Filter Replacement Kit (760200 HEPA2, 190985 access port filter, 770001 filter)
224175	Shelf
190868	Decontamination kit with HEPA filter
190869	Decontamination kit without HEPA filter
130077	Stopper with 3/8" hole
290090	Thermal conductivity type CO ₂ sensor
190885	Infrared type CO ₂ sensor w/ auto zero
250087	CO ₂ valve with fittings
290154	RH sensor assembly (RH display option)
190883	Sample port filter assembly
190661	IR sensor filter
190985	Filtered air exchange assembly
230153	6A, DPDT circuit breaker switch (power)
400199	Switcher, 40W, 12, 5, -12V
420096	130VA transformer, int'l, 12/24V S
194021	34 position control to display ribbon cable
420097	43VA transformer (230VAC units only)
360213	Pressure switch (Gas Guard option)
250121	Gas valve (Gas Guard option)
190512	4-20mA output board
190543	0-5V output board
190544	0-1V output board

Section 8
Spare Parts


ID#	STK #	DESCRIPTION
1	310-201-0-0	IR/TC SENSOR ASSEMBLY
2	310-202-0-0	MOTOR/BLOWER ASSEMBLY
3	116022	HINGE, 1/4" GLASS
4	105060	DOOR, TEMPERED GLASS
5	180174	FACE FRAME, PVC
6	190611	HINGE PIN, STN STEEL
7	190612	HINGE BLOCK, ALUMINUM 2.0"
8	190643	TEMPERATURE PROBE ASSEMBLY
9	191534	REMOVEABLE FEATHER GASKET
10	113002	ADJUSTABLE FOOT
11	190788	OUTER DOOR W/ HEATER
12	190795	MAGNET MOUNT, DOOR
13	190833	LATCH MOUNT, INNER DOOR
14	180210	KNOB, BLACK
15	XXXXX	GAS INJECT PORT
16	360230	SWITCH, MAGNETIC ACTUATED
17	360231	MAGNET, DOOR
18	400133	THERMOSTAT, SNAP ACTION
19	990033	MAGNETIC GASKET, MOLDED
20	290184	"OPTIONAL" RH SENSOR
21	840020	SAMPLE PORT BULKHEAD FITTING
22	132054	LEFT HEATER, 13 WATTS
23	132050	BACK HEATER, 30 WATTS
24	132051	BOTTOM HEATER, 24 WATTS
25	132055	RIGHT HEATER, 22 WATTS
26	132053	COLLAR HEATER, 9 WATTS
27	NA	DOOR HEATER, 69 WATTS
28	132056	FACE HEATER, 27 WATTS
29	1900252	MOTOR GUARD

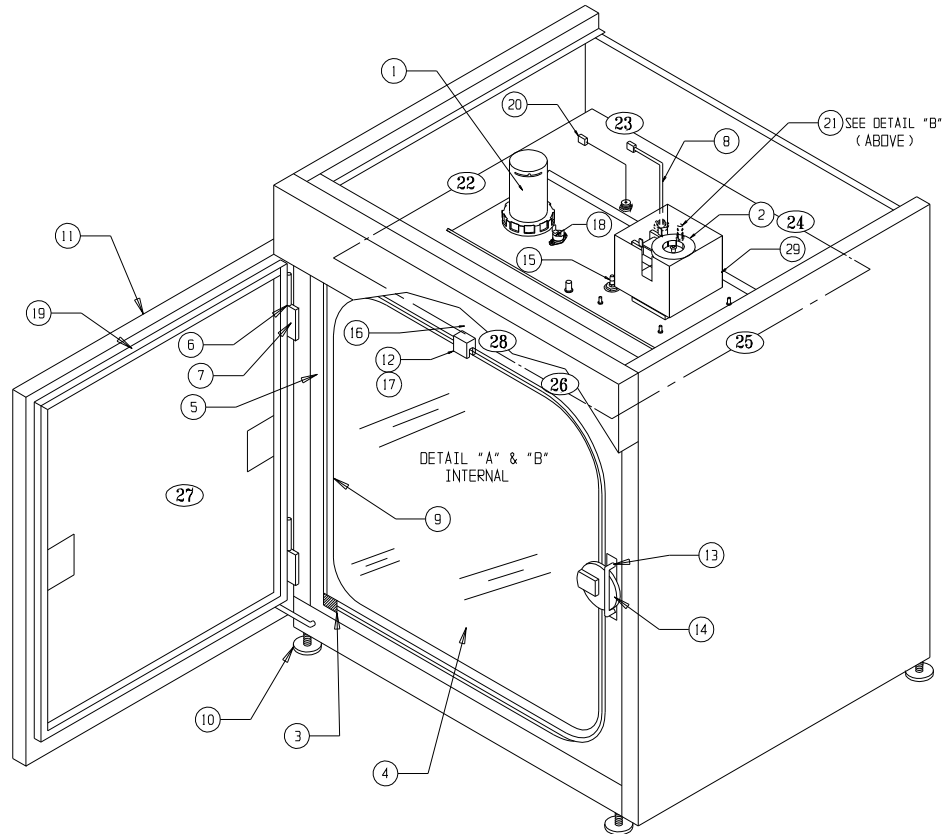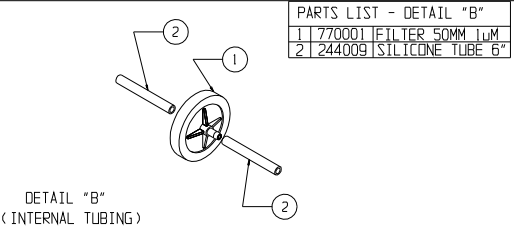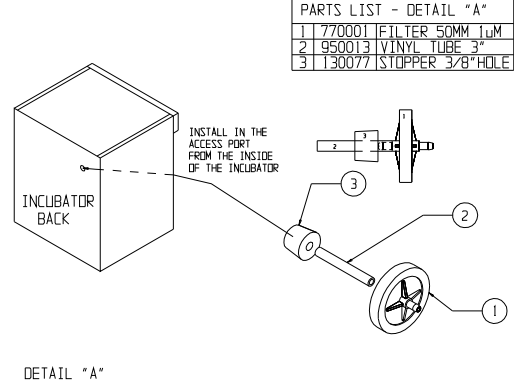

310-200-0-0



CUSTOMER APPROVAL/REFERENCE	5	SI-9227/09-27-04	REVISED	ENG 10568 GLASS DOOR TO 10569
APPROVED BY	4	DN-3180/02-24-04	GLG/SLJ/LON	REVISED MOTOR & GUARD
DATE OF APPROVAL	3	DN-3124/09-04-03	JAL/ROJ/LON	ENG: 290137 PROBE TO 290184
THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM THERMO SCIENTIFIC.				
REVISION NO.	DATE	BY	EAQ/APPD	DESCRIPTION OF REVISION
1	11-02-06	GLG/SLJ/LON	ENG	10569 GASKET TO 191534
2	07-24-00	GLG/SLJ/LON	REVISED PART NUMBERS	
3	09-04-03	JAL/ROJ/LON	REVISED MOTOR & GUARD	
4	02-24-04	GLG/SLJ/LON	REVISED MOTOR & GUARD	
5	09-27-04	SI-9227	REVISED	ENG 10568 GLASS DOOR TO 10569
DATE 11-9-98				
CUSTOMER				
JOB TITLE MAIN CABINET				
ENG TITLE ILLUSTRATED PARTS DRAWING				
LOCATION	JOB NUMBER	DRAWING NUMBER		
INCUBATR		310-200-0-0		

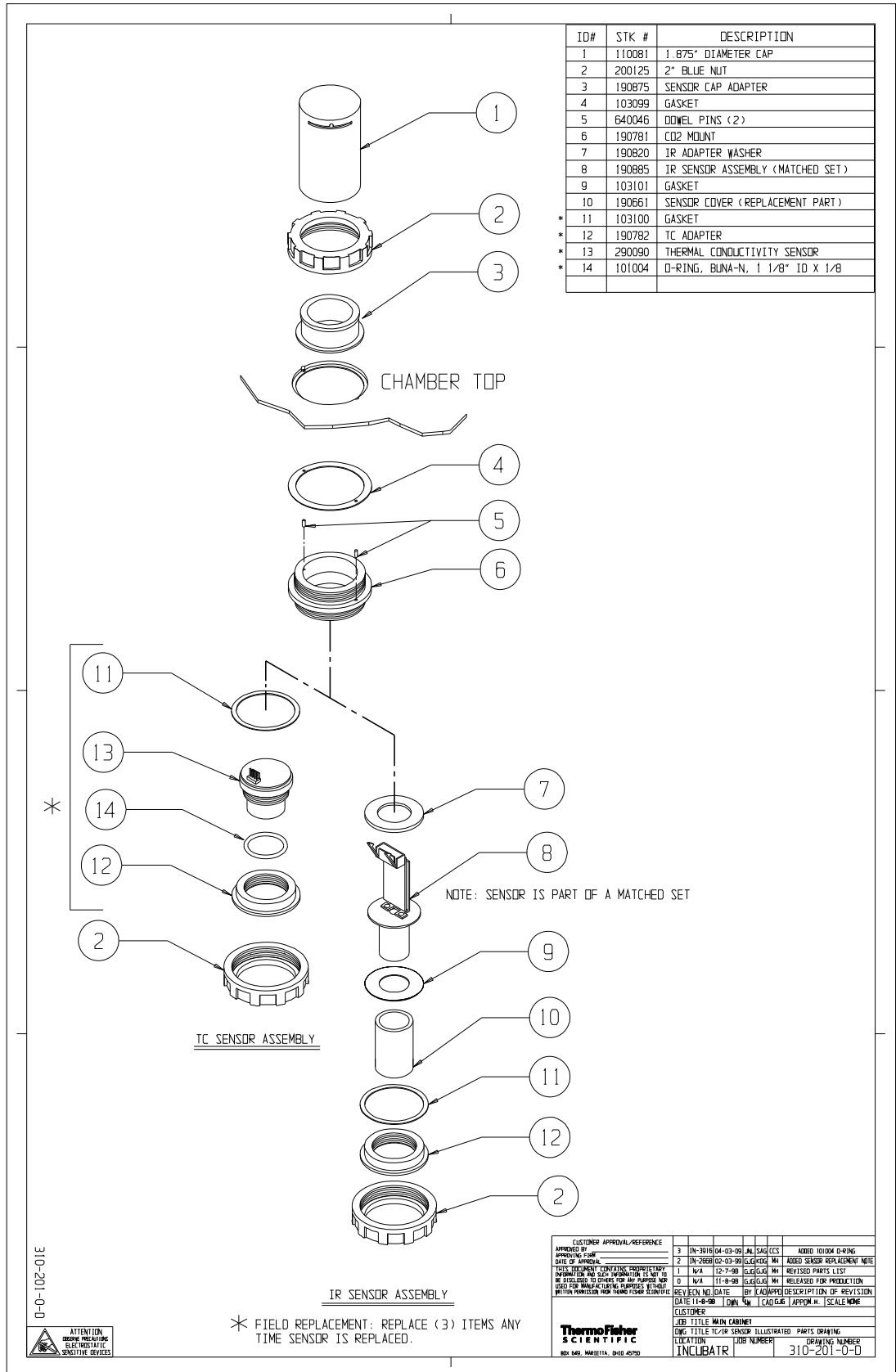


Section 8
Spare Parts

BLOWER ASSEMBLY

BILL OF MATERIALS

ITEM NO.	PART NO.	PART DESCRIPTION	QTY
1	23052	#10-32 S.S. WING NUT	1
2	22002	#4-40 X 1/2 SS PHP SCREW	4
3	640047	#8-32 x .750" LONG STN STL SELF-CLINCHING STUD	5
4	22142	#8-32 x 3/4 FHP SCREW	4
5	23002	#8-32 ZP LKWASH HEX NUT	4
6	730068	.310 ID TEFLON FLAT WASHER	1
7	100113	3.500 X 1.500 BLOWER WHEEL POLY.	1
8	190790-31-	310 INCUBATOR BLOWER SUPPORT	1
9	1900252	BLOWER MOTOR SHIELD & GUARD	1
10	190846	BLOWER SCROLL -RESTRICTED, 310/3110	1
11	190831	BLOWER SCROLL- NO HEPA, 310	1
12	156126	MOTOR	1
13	190832	MOTOR SPACER	1
14	730034	NYLON WASHER 1.500"OD X .500"ID	1
15	730069	SILICONE WASHER	1

310-202-1-D Rev. 5

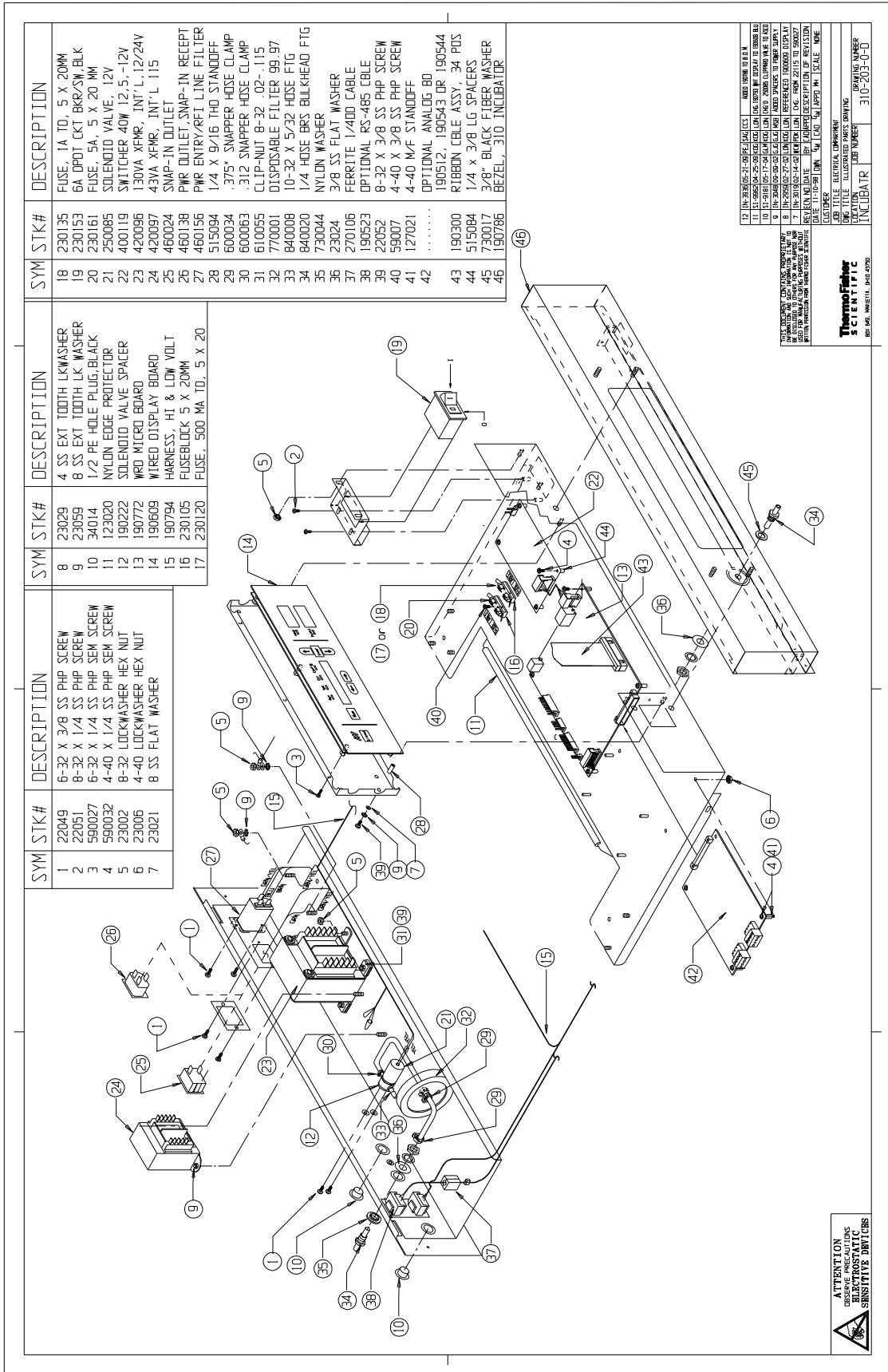



Section 8
Spare Parts

SYM	STK#	DESCRIPTION
1	22052	8-32 X 3/8" SS PHP SCREW
2	22115	6-32 X 1/4" SS PHP SCREW
3	110081	CAP, MODIFIED
4	190856	IR SENSOR HARNESS (2PC)
5	190885	IR CO2 SYSTEM (SENSOR & BOARD)
6	246008	VINYL TUBING 1/16" ID X 1/8" OD X 30"
7	246019	WIRE RNF TUBING 1/4" OD X 3-3/4"
8	250121	SOLENOID VALVE, 3 WAY 12VDC
9	270140	FERRITE PLATE
10	270141	FERRITE END CLIP
11	350006	BRASS ADAPTER, 1/8 HOSE X 10-32
12	360213	PRESSURE SWITCH
13	640057	1/4" HOSE TO 1/8" NPT BRASS 90°
14	440030	FLAT CABLE MOUNT, ADHESIVE
15	600034	HOSE CLAMP .375"
16	600044	SWAPPER HOSE CLAMP 1/2"
17	640027	NYLON FITTING 1/16 TUBE X 10-32
18	190522	HARNESS, GAS GUARD

ATTENTION
 OBSERVE PRECAUTIONS
 FOR SENSITIVE DEVICES

THIS EQUIPMENT CONTAINS PROPRIETARY INFORMATION. IT IS THE PROPERTY OF THERMO SCIENTIFIC AND IS NOT TO BE REPRODUCED, COPIED, OR TRANSMITTED IN ANY FORM OR BY ANY MEANS, ELECTRONIC OR MECHANICAL, INCLUDING PHOTOCOPYING, RECORDING, OR BY ANY INFORMATION STORAGE AND RETRIEVAL SYSTEM.	THE TITLE ILLUSTRATED PARTS DRAWING IS THE PROPERTY OF THERMO SCIENTIFIC.
THERMO SCIENTIFIC 1000 UNIVERSITY AVENUE BOSTON, MA 02116-5050	THERMO SCIENTIFIC 310-204-0000

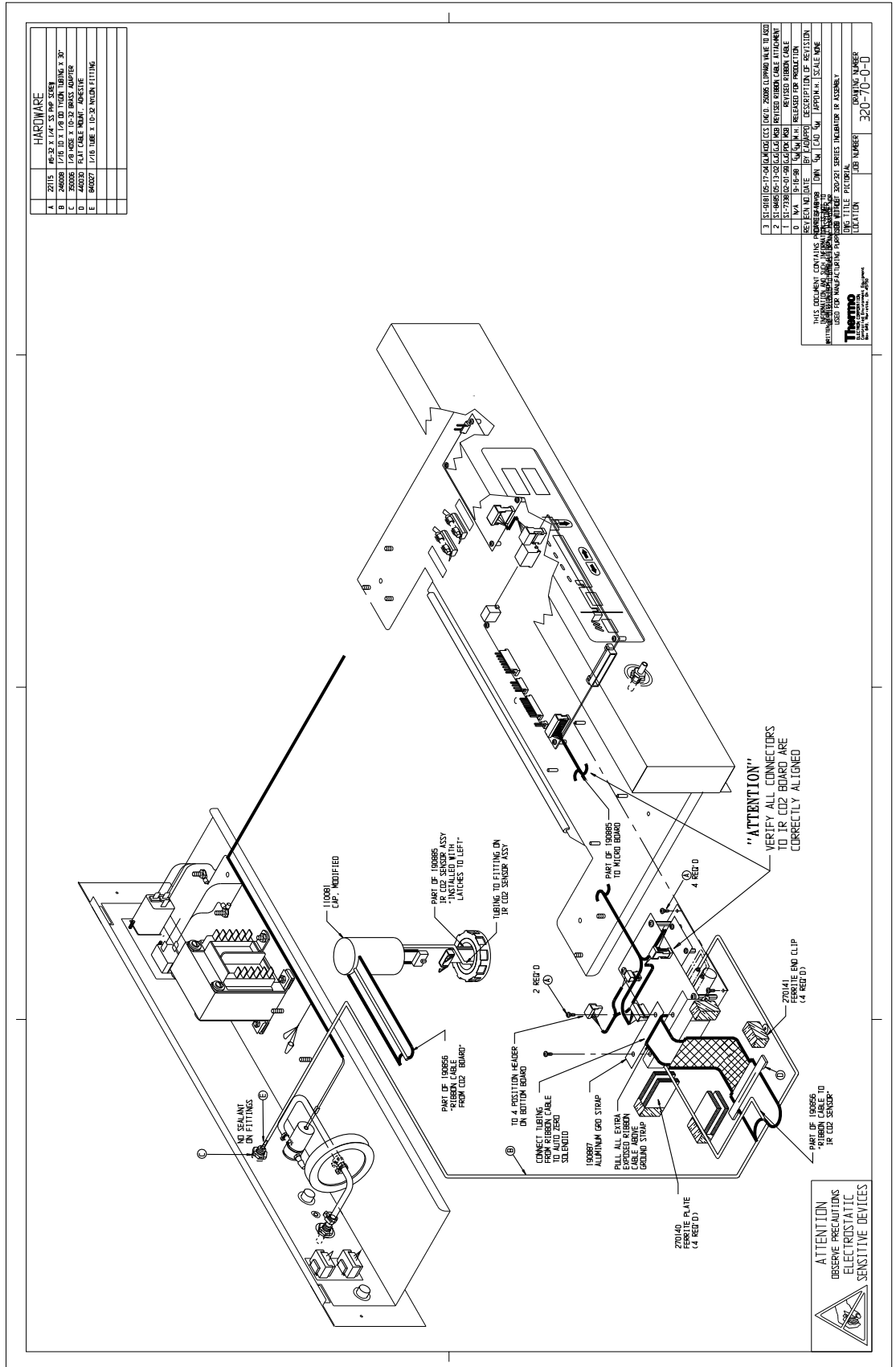

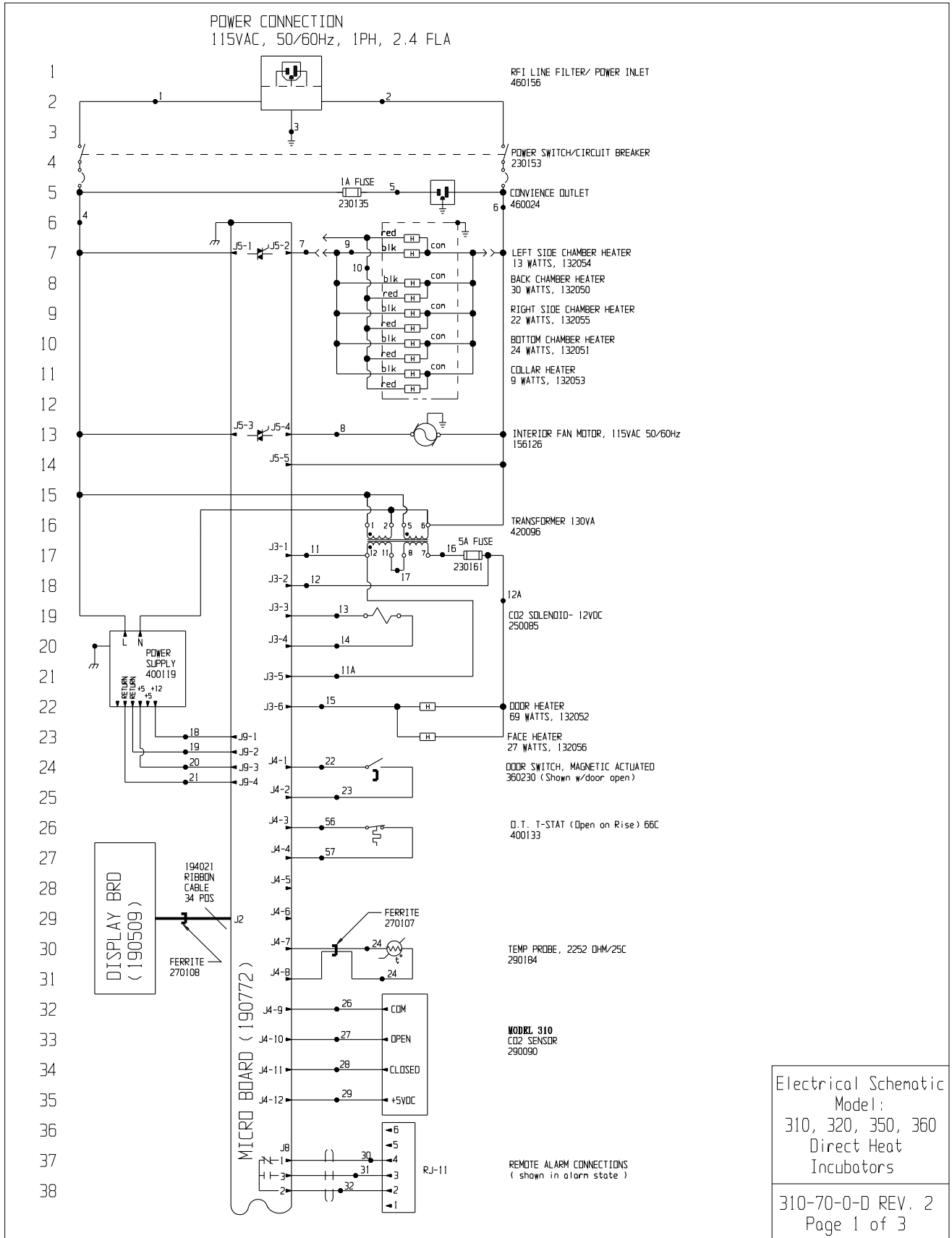


Electrical Schematic
Model:
310, 320, 350, 360
Direct Heat
Incubators

310-70-0-D REV. 2
Page 1 of 3

Section 9
Electrical Schematics

39
40
41
42
43
44
45
46
47
48
49
50
51
52
53
54
55
56
57
58
59
60
61
62
63
64
65
66
67
68
69
70
71
72
73
74
75
76

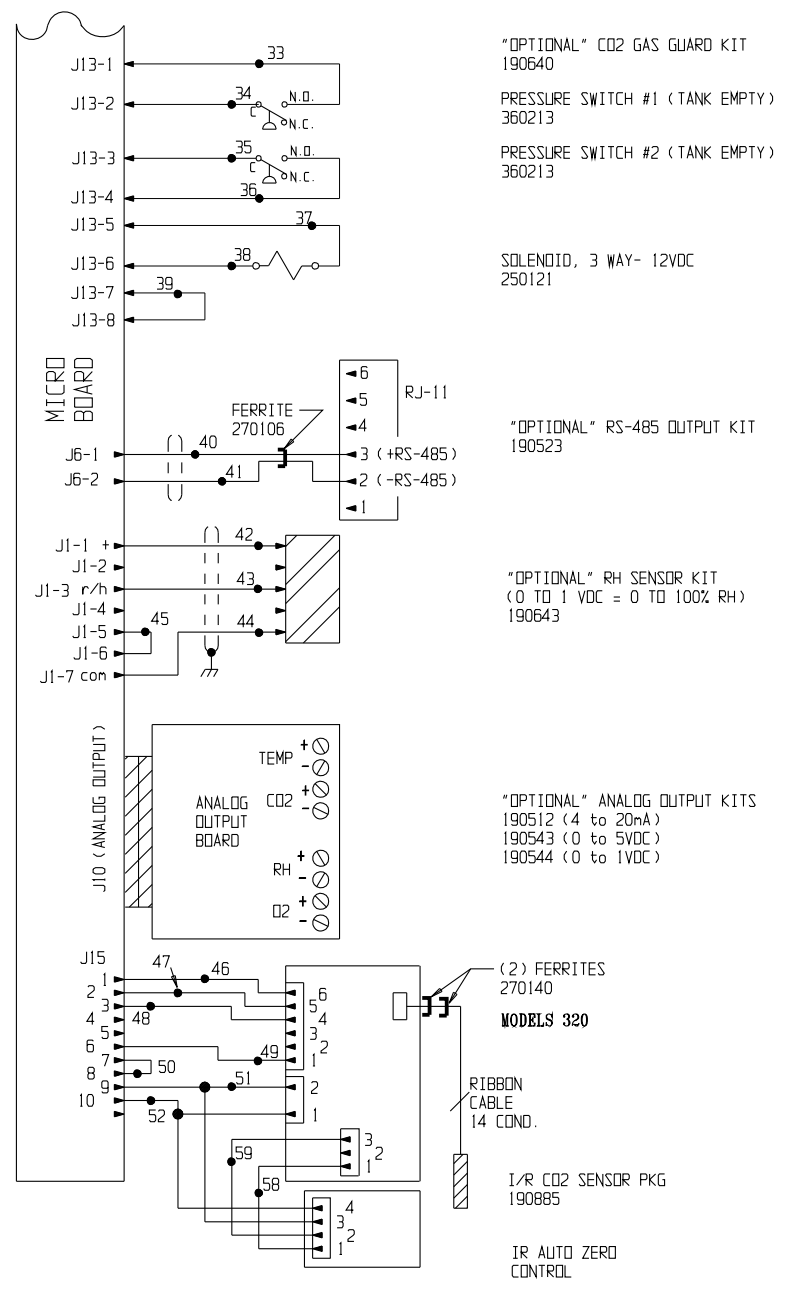

Electrical Schematic
Model:
310, 320, 350, 360
Direct Heat
Incubators


310-70-0-D REV. 2
Page 2 of 3

77
78
79
80
81
82
83
84
85
86
87
88
89
90
91
92
93
94
95
96
97
98
99
100
101
102
103
104
105
106
107


WIRE REFERENCE CHART								
NO.	GA.	COLOR	NO.	GA.	COLOR	NO.	GA.	COLOR
1	16	BROWN	22	22	BLUE	47	22	ORANGE
2	16	BLUE	23	22	BLUE	48	22	YELLOW
3	16	GRN/YEL	24	22	RED	49	22	BROWN
3B	16	GREEN	25	22	RED	50	22	RED
4	18	BLACK	26	22	GREEN	51	22	GREEN
4A	18	BLACK	27	22	GRAY	52	22	BLACK
5	18	BLACK	28	22	PURPLE	53	18	RED
6	18	WHITE	29	22	ORANGE			
7	18	ORANGE	30	22/3	BLACK			
8	18	BLUE	31	22/3	RED	56	22	BROWN
9	18	BLACK	32	22/3	WHITE	57	22	BROWN
10	18	RED	33	20	RED	58	22	BROWN
11	18	BROWN	34	20	RED	59	22	WHITE
11A	18	BROWN	35	20	BLUE			
12	18	RED	36	20	BLUE			
12A	18	RED	37	20	ORANGE			
13	18	YELLOW	38	20	ORANGE			
14	18	YELLOW	39	20	BLACK			
15	18	BLUE	40	22/2	RED			
16	18	RED	41	22/2	BLACK			
17	18	ORANGE	42	22/3	BLACK			
18	18	RED	43	22/3	RED			
19	18	GREEN	44	22/3	GREEN			
20	18	ORANGE	45	22	BLACK			
21	22	GREEN	46	22	BLUE			

SENSOR REFERENCE VALUES

CO2 (290090) DIFFERENCE VOLTAGE OF 3-6MV X CO2 J4-9 & J4-10 TO J4-9 & J4-11
RH (190643) J1-7 & J1-1 = 12VDC J1-7 & J1-3 = 10MV X RH


NOTES:		CUSTOMER APPROVAL/REFERENCE						Electrical Schematic Model: 310, 320, 350, 360 Direct Heat Incubators		
<input checked="" type="checkbox"/> Denotes Terminal Strip Connection	Parts List Reference Number	APPROVED BY _____								
Last Relay Number	<input type="checkbox"/> Assembly	APPROVING FIRM _____	2	IN-3180	03-01-04	GJG	GJG		CCS	REVISED MOTOR PART NUMBER
Last Terminal Number	<input type="checkbox"/> Panel	THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM THERMO ELECTRON.	1	IN-3124	08-04-03	JNL	KDG		MAH	CHG. 290137 TEMP. PROBE TO 290184
Last Wire Number	<input type="checkbox"/> Refrigeration		0	N/A	6/03/98	GJG	GJG		MAH	RELEASED FOR PRODUCTION
	<input type="checkbox"/> Wiring		REV	ECN NO.	DATE	BY	CAD		APPD	DESCRIPTION OF REVISION
			DATE	3-18-98	DWN	GJG	CAD	GJG	APPM.H. SCALE	
			CUSTOMER MODELS 310, 320							
			JOB TITLE AIR JACKET INCUBATOR			115 VOLT (ODM.)				
			DWG TITLE ELECTRICAL SCHEMATIC							
			LOCATION INCUBATR	JOB NUMBER	DRAWING NUMBER					
				310-70-0-D				310-70-0-D REV. 2 Page 3 of 3		

Section 9
Electrical Schematics

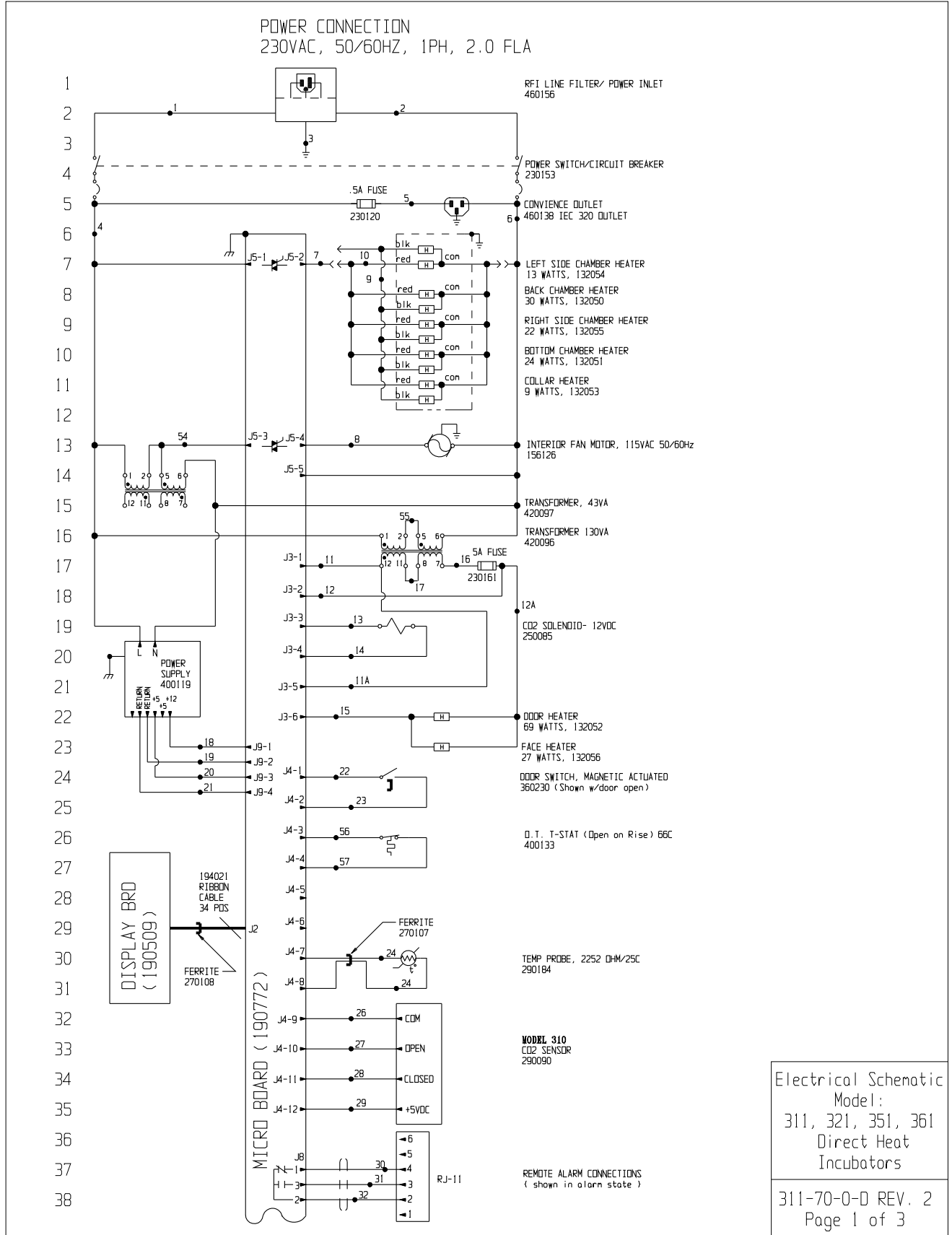

Electrical Schematic
Model:
311, 321, 351, 361
Direct Heat
Incubators

311-70-0-D REV. 2
Page 1 of 3

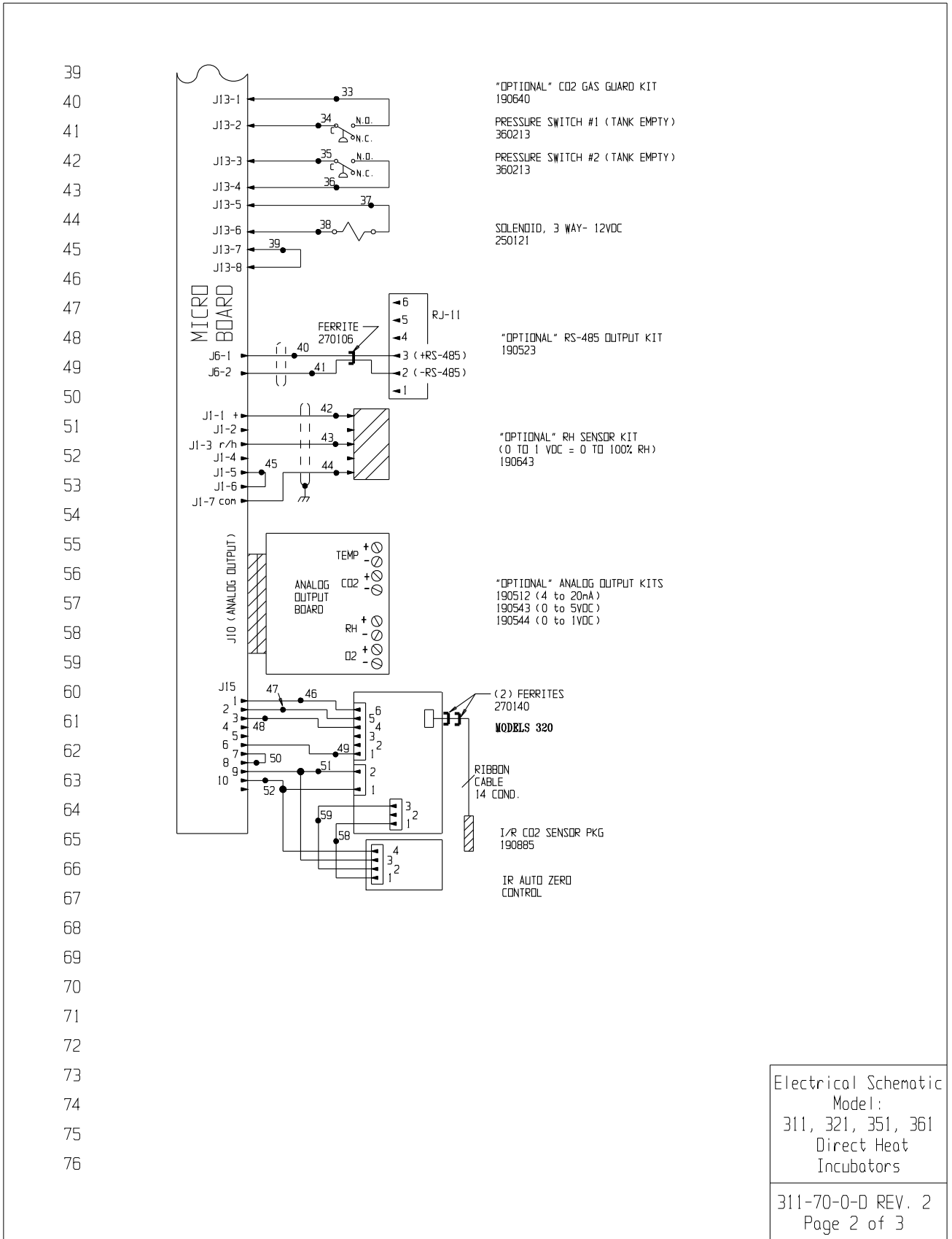

Electrical Schematic
Model:
311, 321, 351, 361
Direct Heat
Incubators

Section 9
Electrical Schematics

77
78
79
80
81
82
83
84
85
86
87
88
89
90
91
92
93
94
95
96
97
98
99
100
101
102
103
104
105
106
107

WIRE REFERENCE CHART								
NO.	GA.	COLOR	NO.	GA.	COLOR	NO.	GA.	COLOR
1	16	BROWN	22	22	BLUE	47	22	ORANGE
2	16	BLUE	23	22	BLUE	48	22	YELLOW
3	16	GRN/YEL	24	22	RED	49	22	BROWN
3B	16	GREEN	25	22	RED	50	22	RED
4	18	BLACK	26	22	GREEN	51	22	GREEN
4A	18	BLACK	27	22	GRAY	52	22	BLACK
5	18	BLACK	28	22	PURPLE	53	18	RED
6	18	WHITE	29	22	ORANGE	54	18	YELLOW
7	18	ORANGE	30	22/3	BLACK	55	18	YELLOW
8	18	BLUE	31	22/3	RED	56	22	BROWN
9	18	BLACK	32	22/3	WHITE	57	22	BROWN
10	18	RED	33	20	RED	58	22	BROWN
11	18	BROWN	34	20	RED	59	22	WHITE
11A	18	BROWN	35	20	BLUE			
12	18	RED	36	20	BLUE			
12A	18	RED	37	20	ORANGE			
13	18	YELLOW	38	20	ORANGE			
14	18	YELLOW	39	20	BLACK			
15	18	BLUE	40	22/2	RED			
16	18	RED	41	22/2	BLACK			
17	18	ORANGE	42	22/3	BLACK			
18	18	RED	43	22/3	RED			
19	18	GREEN	44	22/3	GREEN			
20	18	ORANGE	45	22	BLACK			
21	22	GREEN	46	22	BLUE			

SENSOR REFERENCE VALUES

CO2 (290090) DIFFERENCE VOLTAGE OF 3-6MV/%CO2
J4-9 & J4-10 TO J4-9 & J4-11

RH (190643) J1-7 & J1-1 = 12VDC
J1-7 & J1-3 = 10MV/%RH

NOTES:		CUSTOMER APPROVAL/REFERENCE								Electrical Schematic Model: 311, 321, 351, 361 Direct Heat Incubators	
<input checked="" type="checkbox"/> Denotes Terminal Strip Connection	Parts List Reference Number	APPROVED BY	APPROVING FIRM	2	IN-3180	03-01-04	GJG	GJG	CCS		
Last Relay Number	<input type="checkbox"/> Assembly	DATE OF APPROVAL		1	IN-3124	08-04-03	JNL	KDG	LON	CHG. 290137 TEMP. PROBE TO 290184	
Last Terminal Number	<input checked="" type="checkbox"/> Panel	THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM THERMO ELECTRON		0	N/A	6/03/98	GJG	GJG	MAH	RELEASED FOR PRODUCTION	
Last Wire Number	<input type="checkbox"/> Refrigeration	REV	ECN NO.	DATE	BY	CAD	APPD	DESCRIPTION OF REVISION			
	<input type="checkbox"/> Wiring	DATE	3-18-98	DWN	GJG	GJG	APPDM.H.	SCALE			
		CUSTOMER	MODELS 311, 321								
		JOB TITLE	AIR JACKET INCUBATOR		230 VOLT (EXP.)						
		DWG TITLE	ELECTRICAL SCHEMATIC								
		LOCATION	JOB NUMBER	DRAWING NUMBER							
		INCUBATR		311-70-0-D						311-70-0-D REV. 2 Page 3 of 3	

THERMO FISHER SCIENTIFIC DIRECT HEAT INCUBATOR WARRANTY USA*

The Warranty Period starts two weeks from the date your equipment is shipped from our facility. This allows shipping time so the warranty will go into effect at approximately the same time your equipment is delivered. The warranty protection extends to any subsequent owner during the first year warranty period.

During the first year, component parts proven to be non-conforming in materials or workmanship will be repaired or replaced at Thermo's expense, including labor. In addition, after the initial warranty period of one year the CO₂ sensor (IR) will be warranted for a second year, parts only. The CO₂ sensor (T/C) will be warranted for four additional years for a total of five years, parts only. Installation and calibration is not covered by this warranty agreement. The Technical Services Department must be contacted for warranty determination and direction prior to performance of any repairs. Expendable items, glass, filters and gaskets are excluded from this warranty.

Replacement or repair of components parts or equipment under this warranty shall not extend the warranty to either the equipment or to the component part beyond the original warranty period. The Technical Services Department must give prior approval for return of any components or equipment. At Thermo's option, all non-conforming parts must be returned to Thermo postage paid and replacement parts are shipped FOB destination.

THIS WARRANTY IS EXCLUSIVE AND IN LIEU OF ALL OTHER WARRANTIES, WHETHER WRITTEN, ORAL OR IMPLIED. NO WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE SHALL APPLY. Thermo shall not be liable for any indirect or consequential damages including, without limitation, damages relating to lost profits or loss of products.

Your local Thermo Sales Office is ready to help with comprehensive site preparation information before your equipment arrives. Printed instruction manuals carefully detail equipment installation, operation and preventative maintenance.

If equipment service is required, please call your Technical Service Department at 1-800-438-4851 (USA and Canada) or 1-740-373-4763. We're ready to answer your questions on equipment warranty, operation, maintenance, service and special applications. Outside the USA, contact your local distributor for warranty information.



Rev. 4 4/09

THERMO FISHER SCIENTIFIC INTERNATIONAL DIRECT HEAT INCUBATOR WARRANTY*

The Warranty Period starts two months from the date your equipment is shipped from our facility. This allows shipping time so the warranty will go into effect at approximately the same time your equipment is delivered. The warranty protection extends to any subsequent owner during the first year warranty period.

During the first year component parts proven to be non-conforming in materials or workmanship will be repaired or replaced at Thermo's expense, excepting labor. In addition, after the initial warranty period of one year the CO₂ sensor (IR) will be warranted for a second year, parts only. The CO₂ sensor (T/C) will be warranted for four additional years for a total of five years, parts only. Installation and calibration is not covered by this warranty agreement. The Technical Services Department must be contacted for warranty determination and direction prior to performance of any repairs. Expendable items, glass, filters and gaskets are excluded from this warranty.

Replacement or repair of components parts or equipment under this warranty shall not extend the warranty to either the equipment or to the component part beyond the original warranty period. The Technical Services Department must give prior approval for return of any components or equipment. At Thermo's option, all non-conforming parts must be returned to Thermo postage paid and replacement parts are shipped FOB destination.

THIS WARRANTY IS EXCLUSIVE AND IN LIEU OF ALL OTHER WARRANTIES, WHETHER WRITTEN, ORAL OR IMPLIED. NO WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE SHALL APPLY. Thermo shall not be liable for any indirect or consequential damages including, without limitation, damages relating to lost profits or loss of products.

Your local Thermo Sales Office is ready to help with comprehensive site preparation information before your equipment arrives. Printed instruction manuals carefully detail equipment installation, operation and preventive maintenance.

If equipment service is required, please call your Technical Services Department at 1-800-438-4851 (USA and Canada) or 1-740-373-4763. We're ready to answer your questions on equipment warranty, operation, maintenance, service and special application. Outside the USA, contact your local distributor for warranty information.



Rev. 4 4/09

Declaration of Conformity

Manufacturer's Name: Thermo Electron Corp.

Manufacturer's Address: 401 Millcreek Road
Marietta, Ohio 45750
U.S.A.

Product Description: Forma[®] Laboratory & IVF Incubators

Product Designations: 310, 311, 320 & 321

Year of Initial Marking (CE): 1998

Affected Serial Numbers: Release 2
(Release Level [REL#] shown on Serial Tag)

These products conform to the following European Union Directive(s):

EMC: 89/336/EEC

MDD: 93/42/EEC

These products conform to the following Harmonized, International and National Standards:

EMC:

EN 50081-1 & EN 50082-1

EN 50022-1, Class B

EN 61326-1

EN 61000-3-2 and -3

EN 61000-4-2, -3, -4, -6, -8 and -11


MDD:

EN 61010-1:1993

IEC 1010-1 Amendment 2

CSA C22.2 No. 1010.1-92

UL 61010A-1

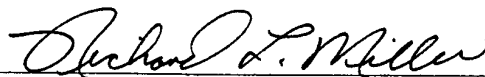

Richard L. Miller, CQE
Regulatory Compliance Manager

Thermo
ELECTRON CORPORATION

03 February 2004

Declaration of Conformity

Manufacturer's Name: Thermo Electron Corp

Manufacturer's Address: 401 Millcreek Road
Marietta, Ohio 45750
U.S.A.

Product Description: Forma[®] Laboratory Incubators

Product Designations: 350, 351, 360 & 361

Year of Initial Marking (CE): 1999


Affected Serial Numbers: Release 2
(Release Level [REL#] shown on Serial Tag)

These products conform to the following European Union Directive(s):

EMC: 89/336/EEC
LVD: 73/23/EEC

These products conform to the following Harmonized, International and National Standards:

EMC:	LVD:
EN 50081-1 & EN 50082-1	EN 61010-1:1993
EN 50022-1, Class B	IEC 1010-1 Amendment 2
EN 61326-1	CSA C22.2 No. 1010.1-92
EN 61000-3-2 and -3	UL 61010A-1
EN 61000-4-2, -3, -4, -6, -8 and -11	

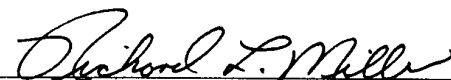

Richard L. Miller, CQE
Regulatory Compliance Manager

Thermo
ELECTRON CORPORATION

03 February 2004

Thermo Fisher Scientific
Controlled Environment Equipment
401 Millcreek Road
Marietta, Ohio 45750
United States

www.thermo.com