

Specifications

Shaking speed	4 to 160 tilts per minute
Timer	1 minute to 2 hours or "hold"
Operating Environment	4° to 65°C
Platform Dimensions	360 x 280mm
Rocking Motion	Adjustable, $\pm 20^\circ$ tilt
Maximum Load	4.6kg
Clearance between platforms (Double platform unit only)	82mm
Voltage Requirements	120V/60Hz
Power/Current	60W/0.5amps

Unpacking and Installation

Upon receipt of the Model 35/35D Benchtop Rocker, examine the exterior carton for damages. In the event of shipping damage, a claim must be filed with the transportation carrier within 10 days.

Carefully unpack the unit and check for completeness. Included in the carton should be the following:

- Rocker with platform(s)
- Securing hardware for second platform (35D only)
- Instruction Manual
- Warranty Card

If any of these items are missing, contact Labnet International's Customer Service Department immediately.

Save the original packaging until the unit has been shown to be operating properly.

Verify that the unit is rated for the proper voltage by checking the panel on the back. Check that the proper power cord has been received. The serial number is also found on the back panel. Make note of the serial number in the space provided in the Warranty section of this manual.

The Model 35/35D Benchtop Rocker should be located on a level, stable surface near a grounded electrical outlet. Allow sufficient clearance on all sides of the unit, at least 3 inches, for proper ventilation. Ensure that the platform will not touch other objects while in operation.

If the unit is supplied with an optional upper platform, it should be