
Technical Specifications

Overall Dimensions: 12 5/8" W x 7 1/4" H x 12" D

Platform Working Dimensions: 10 5/8" x 12 5/8"

Platform Angle Variation (Pitch): 0 to ± 24 angle

rpm (No Load Condition): 5-30 rpm \pm 2 rpm

Maximum Load:

Centered: 15 lbs.

On Double-Tier Platform: 15 lbs.

Ambient Operating Range: 0°C to 65°C

Relative Humidity Operating Range:

20-80% non-condensing

Timer Range:

0-2 hrs. (60 Hz.)

0-3 hrs. (50 Hz.)



Warning

Use a properly grounded electrical outlet of correct voltage and current handling capacity.

Installation

Remove your Platform Vari-Mix Rocker from the carton. Inspect to ensure that the unit has not been damaged during shipment. Set the rocker on a flat, stable surface and plug into an appropriate outlet. Remove the white silicone rubber top from the protective plastic packaging and place the top on the metal platform.

Set the power toggle switch into the "ON" position. The green power switch will illuminate to indicate that the unit is "ON." Set the pitch and the speed and the timer control knobs.