

## TROUBLESHOOTING

Observation	Suggestion(s)
The display does not light	There is no power to the unit. Secure the plug into the back of the unit and into the wall socket, and then depress the circuit breaker on the back panel
The display does not count down after <i>Start</i> is pushed	The <i>Energy</i> or <i>Time</i> modes have not been selected. Select the desired mode ( <i>Autocross-link</i> selects <i>Energy</i> automatically)
	The door switch is not engaged. Close the door completely
	The UV sensor is not completely pushed into the connector on the interior sidewall. Ensure connections are secure
Countdown is slower than normal (e.g., <i>Autocrosslink</i> takes more than one minute)	UV bulbs are burned out. Check for burned-out bulbs and replace as necessary
	The starter is out. Replace the starter. Rotate the defective starter one-quarter turn counterclockwise and pull out
	The UV sensor is blocked and may need to be replaced. Contact the Stratagene Technical Services Department

## SPECIFICATIONS

### Weight

Model 1800: 9.7 kg  
 Model 2400: 15.5 kg  
 Model 2400 (220V): 18.9 kg

### Dimensions (in cm)

Model 1800, internal: 33.7 W × 18.1 D × 16.8 H  
 Model 1800, external: 46.2 W × 22.4 D × 26.7 H  
 Model 2400, internal: 47.7 W × 34.3 D × 16.8 H  
 Model 2400, external: 58.9 W × 38.9 D × 26.7 H

### Bulbs

Model 1800: 8 watts each  
 Model 2400: 15 watts each

### Power Delivered

Model 1800: ~3000 μwatts/cm<sup>2</sup>  
 Model 2400: ~4000 μwatts/cm<sup>2</sup>

### Starters

Sylvania starters (FS-5 for model 1800 and FS-2 for model 2400) may be used as replacement starters.